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50 bp-DNA-Leiter (ready-to-use)

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50 bp DNA-ladder

(ready-to-use)

This marker was generated by PCR amplification based on Lambda phage as template DNA. Following PCR, the DNA was purified, spectroscopically measured and used for marker mixture in the amounts given below.

Fragment sizes ¹ (in bp)	DNA-mass (ng/ μ l marker solution)
500	10
450	10
400	10
350	10
300	10
250	15
200	10
150	10
100	10
50	10

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Application:

Our 50 bp DNA Ladder is solubilised in 1X concentrated sample buffer (Tris -Cl, Ph 7.5, 10 mM, EDTA 10 mM, glycerol 10 %, orange G 0.4 mg/ml) and may be applied directly to the gel. The concentration is 1 μ g DNA / 10 μ l marker solution.

This gel loading buffer resembles a 6x diluted solution of gel loading buffer Roti®-Load DNA orange 1 (Art.-No.: HP04.1) with EDTA concentration enhanced to 10 mM, in order to provide optimal stabilisation of the PCR fragments.

Sample application:

The standard loading of MINI to MIDI gels per lane is:

- with bands visible in UV-light after Et.-Br.-staining: 0.5 – 1.0 μ g
- with detection after Et.-Br.-staining with signal-integrated camera systems: 0.1– 0.5 μ g

Storage:

Optimal storage temperature is -20 °C.

Repeated (>10 times) thawing and freezing will damage the marker and should be avoided. If necessary aliquote the marker in portions. Short term storage (few weeks) may be done at 4 °C.

Please note:

The 50 bp DNA-ladder is compatible with the 100 bp DNA-ladder, *equimolar* (T834, see figure, right lane). Mixed in equal shares (f. e. 5 μ l 20 bp DNA-ladder and 5 μ l 100 bp DNA-ladder, *equimolar*) they result in a versatile range of DNA-marker bands for analysis of DNA-fragments from 50 to 1000 bp.

¹ Taking the sequencing accuracy as a basis, we can guarantee an error rate of <1 % when indicating fragment sizes.

50 bp DNA-ladder

2810.1

500 μ l

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Carl Roth GmbH + Co. KG