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Roti®-Load 1

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(4x-concentrated)

Introduction:

Roti-Load 1 is a well tested and proven sample buffer for protein gel electrophoresis that has been modified and further developed based on the original formulation acc. to Laemmli (1970) (1).

Roti®-Load 1 prevents degradation of proteins while heating during sample preparation (2,3).

The phosphate-buffer system keeps the sample pH value constant over a wide temperature range, additionally stabilizing the sensitive aspartyl-prolyl peptide bonds (4). The ready-to-use solution Roti®-Load 1 already contains denaturing and reducing reagents, a dye and glycerol as density increasing reagent.

For a sample preparation under non reducing conditions we recommend our Roti®-Load 2 (Art. No. K930.1).

Use:

1. Prepare your polyacrylamide gel.
2. Determine sample volume.
3. Add Roti®-Load 1 (one third of sample volume).

Example:

60 µl sample + 20 µl Roti®-Load 1 (4x-concentrated.)

4. Mix by pipetting.

5. Heat samples for 3-5 min to 80-90 °C.
6. Directly load samples to the gel.

Tip:

While using Roti®-Load 1 and Roti®-Load 2 for comparative studies please leave at least one free gel lane between the differently treated samples.

Store at +20°C

Carefully close the flask after use. We recommend additional sealing using parafilm, in order to prevent evaporation of β-mercapto ethanol.

Please note: Roti®-Load 1 contains SDS and may form precipitates when stored below 25 °C. Dissolve precipitates completely before use (approx. 50 °C).

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Carl Roth GmbH+Co. KG