Produkt-Datenblatt

D/E-NEUTRALISIERUNGSPUFFER

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Product Data Sheet

D/E NEUTRALIZING BUFFER

For isolation and enumeration of microorganisms in presence of disinfectants or preservatives.

1L8A

Approximate formulation in g/l:

Dextrose	10
Lecithin	7
Sodium thiosulfate	6
Casein peptone.....	5
Polysorbate 80.....	5
Yeast extract.....	2.5
Sodium bisulfite	2.5
Sodium thioglycolate.....	1
Bromocresol purple.....	0.02
pH value.....	7.6 ±0.2

USES

The Neutralizing Agar is recommended for growing a large span of microorganisms.

The medium is especially suitable for hygiene monitoring, containing 5 disinhibitors: Sodium bisulfite, sodium thioglycolate, sodium thiosulfate, lecithin and polysorbate 80.

These neutralizers inactivate bacteriostatic disinfectants and preservatives like aldehydes, mercurial, iodine and chlorine, quaternary ammonium compounds and substituted phenolics. As the growth inhibition is inactivated, it is possible to detect present living microorganisms after disinfection. Performing a before and after examination, the efficacy of a disinfectant can be checked.

ISO standards mention this medium as an alternative to culture enrichment in the detection of aerobic mesophilic bacteria (ISO 21149), in the detection of *Escherichia coli* (ISO 21150) and for checking the presence of *Pseudomonas aeruginosa* (ISO 22717) and *Staphylococcus aureus* (ISO 22718).

MICROBIOLOGICAL TEST

The following results were obtained in the performance of the given medium from type cultures after incubation at a temperature of 30-35 °C and observed after 18 - 72 hours.

Mikroorganisms	Growth
<i>Escherichia coli</i> ATCC 8739	Good
<i>Bacillus subtilis</i> ATCC 6633	Good
<i>Pseudomonas aeruginosa</i> ATCC 9027	Good
<i>Staphylococcus aureus</i> ATCC 6538	Good
<i>Candida albicans</i> ATCC 10231	Good

D/E NEUTRALIZING AGAR	10 x 100 ml	1L8A.1
------------------------------	--------------------	---------------

Carl Roth GmbH + Co. KG

Schoemperlenstraße 3-5 • 76185 Karlsruhe • P.O. Box 100121 • 76231 Karlsruhe
Phone: +49 (0) 721/ 5606-0 • Fax: +49 (0) 721/ 5606-149 • info@carlroth.com • www.carlroth.com

The company is a limited partnership with headquarters in Karlsruhe, reg. court Mannheim HRA 100055. Roth Chemie GmbH, with headquarters in Karlsruhe, reg. court Mannheim HRB 100428, is the personally liable partner. Managing Director: André Houdelet. Sales tax identification number: DE 143621073.

jh 08/2021