



# Instructions for use

## ROTIPHORESE® Ready-to-Use Gel Solutions

3029, 3030, A515, T802, A516, A121, A124

Hazard and precautionary statements see page 3 of this instructions for use.

### A) SDS-PAGE

**Resolving gel** (data apply to 20 ml gel solution)

|                        |                           |            |            |             |             |             |
|------------------------|---------------------------|------------|------------|-------------|-------------|-------------|
| 30 %                   | <b>Gel concentration</b>  | <b>6 %</b> | <b>8 %</b> | <b>10 %</b> | <b>12 %</b> | <b>15 %</b> |
| acryl-<br>amide<br>mix | Aqua dest. (ml)           | 10.6       | 9.3        | 7.9         | 6.6         | 4.6         |
|                        | 30 % acrylamide mix (ml)  | 4          | 5.3        | 6.7         | 8           | 10          |
|                        | Tris (1.5 M, pH 8.8) (ml) | 5          | 5          | 5           | 5           | 5           |
| 40 %                   | <b>Gel concentration</b>  | <b>6 %</b> | <b>8 %</b> | <b>10 %</b> | <b>12 %</b> | <b>15 %</b> |
| acryl-<br>amide<br>mix | Aqua dest. (ml)           | 11.6       | 10.6       | 9.6         | 8.6         | 7.1         |
|                        | 40 % acrylamide mix (ml)  | 3          | 4          | 5           | 6           | 7.5         |
|                        | Tris (1.5 M, pH 8.8) (ml) | 5          | 5          | 5           | 5           | 5           |

Add in this order:

200 µl 10 % SDS solution (mix carefully, avoid bubbles)

200 µl 10 % ammonia persulphate solution (prepare freshly)

20 µl TEMED (mix carefully, avoid bubbles)

Pour gel immediately and overlay with isopropanol

**Stacking gel** (data apply to 5 % gels)

|                        |                           |             |             |             |             |              |
|------------------------|---------------------------|-------------|-------------|-------------|-------------|--------------|
| 30 %                   | <b>Gel volume</b>         | <b>1 ml</b> | <b>3 ml</b> | <b>5 ml</b> | <b>8 ml</b> | <b>10 ml</b> |
| acryl-<br>amide<br>mix | Aqua dest. (ml)           | 0.68        | 2.1         | 3.4         | 5.5         | 6.8          |
|                        | 30 % acrylamide mix (ml)  | 0.17        | 0.5         | 0.83        | 1.3         | 1.7          |
|                        | Tris (1.0 M, pH 6.8) (ml) | 0.13        | 0.38        | 0.63        | 1           | 1.25         |
|                        | SDS (10 % solution) (µl)  | 10          | 30          | 50          | 80          | 100          |
|                        | APS (10 % solution*) (µl) | 10          | 30          | 50          | 80          | 100          |
|                        | TEMED (µl)                | 1           | 3           | 5           | 8           | 10           |

|                        |                           |             |             |             |             |              |
|------------------------|---------------------------|-------------|-------------|-------------|-------------|--------------|
| 40 %                   | <b>Gel volume</b>         | <b>1 ml</b> | <b>3 ml</b> | <b>5 ml</b> | <b>8 ml</b> | <b>10 ml</b> |
| acryl-<br>amide<br>mix | Aqua dest. (ml)           | 0.725       | 2.185       | 3.645       | 5.84        | 6.3          |
|                        | 40 % acrylamide mix (ml)  | 0.125       | 0.375       | 0.625       | 1           | 1.25         |
|                        | Tris (1.0 M, pH 6.8) (ml) | 0.13        | 0.38        | 0.63        | 1           | 1.25         |
|                        | SDS (10 % solution) (µl)  | 10          | 30          | 50          | 80          | 100          |
|                        | APS (10 % solution*) (µl) | 10          | 30          | 50          | 80          | 100          |
|                        | TEMED (µl)                | 1           | 3           | 5           | 8           | 10           |

\* prepare freshly!

Be careful to mix the solution thoroughly before and after addition of SDS and TEMED.

Avoid bubbles. Pour the stacking gel immediately and insert the comb carefully.

## **B) Separation of nucleic acids**

### ***Denaturing TBE gels for separation of single stranded nucleic acids (e.g. sequencing gels)***

(data apply to 100 ml gel solution)

|   |                                      |            |            |            |
|---|--------------------------------------|------------|------------|------------|
| 25 % sequencing gel concentrate with urea | <b>Gel concentration</b>             | <b>4 %</b> | <b>6 %</b> | <b>8 %</b> |
|   | Sequencing gel diluent (ml)*         | 74         | 66         | 58         |
|   | 25 % sequencing gel concentrate (ml) | 16         | 24         | 32         |
| 30 % acrylamide mix (29:1)                | <b>Gel concentration</b>             | <b>4 %</b> | <b>6 %</b> | <b>8 %</b> |
|   | Aqua dest. (ad 90 ml) (ml)*          | app. 52    | app. 45    | app. 39    |
|   | 30 % acrylamide mix (ml)             | 13.3       | 20         | 26.5       |
|   | Urea (g)**                           | 42         | 42         | 42         |
| 40 % acrylamide mix (19:1 or 29:1)        | <b>Gel concentration</b>             | <b>4 %</b> | <b>6 %</b> | <b>8 %</b> |
|   | Aqua dest. (ad 90 ml) (ml)*          | app. 55    | app. 50    | app. 45    |
|   | 40 % acrylamide mix (ml)             | 10         | 15         | 20         |
|   | Urea (g)**                           | 42         | 42         | 42         |

\*If required for resolution of secondary structures the gel may be supplemented with formaldehyde by replacing 25 ml aqua dest. with 25 ml formaldehyde.

\*\*Results in gels with 42 % urea (7 M)

#### Add in this order:

10 ml 10 x TBE buffer\*\*\* (mix carefully, avoid bubbles, degas if required)

400 µl 10 % ammonia persulphate solution (prepare freshly)

50 µl TEMED (mix carefully, avoid bubbles)

Pour gel immediately and insert the comb carefully

\*\*\*Results in 45 % urea if sequencing gel concentrate and sequencing gel diluent are used. If 50% urea are required replace 10 x TBE by the ready-to-use sequencing gel buffer concentrate with 50 % urea (Ord. No. 3050.1).

### ***TBE gels for electrophoresis of ds nucleic acid***

(data apply to 100 ml gel solution)

|                                    |                          |            |             |             |
|------------------------------------|--------------------------|------------|-------------|-------------|
| 30 % acrylamide mix (29:1)         | <b>Gel concentration</b> | <b>6 %</b> | <b>10 %</b> | <b>15 %</b> |
|                                    | Aqua dest. (ml)          | 69         | 56          | 39          |
|                                    | 30 % acrylamide mix (ml) | 20         | 33          | 50          |
| 40 % acrylamide mix (19:1 or 29:1) | <b>Gel concentration</b> | <b>6 %</b> | <b>10 %</b> | <b>15 %</b> |
|                                    | Aqua dest. (ml)          | 74         | 64          | 51.5        |
|                                    | 40 % acrylamide mix (ml) | 15         | 25          | 37.5        |

#### Add in this order:

10 ml 10 x TBE buffer (mix carefully, avoid bubbles, degas if required)

1 ml 10 % ammonia persulphate solution (prepare freshly)

60 µl TEMED (mix carefully, avoid bubbles)

Pour gel immediately and insert the comb carefully

**Detailed instructions for use and further information can be found in the info brochure **PAGE Instructions** (see [www.carlroth.de](http://www.carlroth.de) next to the product description).**


## Hazard and precautionary statements

Please wear protective gloves and eye protection when handling the solutions.


Waste (solution residues, poured gels) must be disposed of as hazardous waste in accordance with the relevant national or regional provisions.

Please note safety data given on label and MSDS.

- **ROTIPHORESE® Ready-to-Use Gel Solutions: 3029, A124**

 **Danger** H302-H315-H319-H317-H340-H350-H361f-H372  
P201-P280-P301+P312-P302+P352-P305+P351+P338-P308+P313

- **ROTIPHORESE® Ready-to-Use Gel Solutions: 3030, A516, A121, A515, T802**

 **Danger** H301-H315-H319-H317-H332-H340-H350-H361f-H372  
P201-P280-P301+P310-P302+P352-P305+P351+P338-P308+P313

| Art. No.  | Product                                 | Conc. | Mixing ratio |
|---|---|-------|--------------|
| <b>Acrylamide/bisacrylamide mixtures, ready-to-use</b>                                |   |       |              |
| 3029  | ROTIPHORESE® Gel 30 (37.5:1):           | 30 %  | 37.5:1       |
| 3030  | ROTIPHORESE® Gel 40 (19:1):             | 40 %  | 19:1         |
| A515  | ROTIPHORESE® Gel 40 (29:1):             | 40 %  | 29:1         |
| T802  | ROTIPHORESE® Gel 40 (37.5:1):           | 40 %  | 37.5:1       |
| <b>Acrylamide/bisacrylamide mixtures for automated sequencing (fluorescence free)</b> |   |       |              |
| A516  | ROTIPHORESE® NF acrylamide/bis-solution | 40 %  | 19:1         |
| A121  | ROTIPHORESE® NF acrylamide/bis-solution | 40 %  | 29:1         |
| A124  | ROTIPHORESE® NF acrylamide/bis-solution | 30 %  | 29:1         |

## Further ROTIPHORESE® Gel Solutions

| Art. No.  | Product   | Conc.        | Mixing ratio     |
|---|---|--------------|------------------|
| <b>Acrylamide- and bisacrylamide solution, ready-to-mix</b> |   |              |                  |
| 3037  | ROTIPHORESE® Gel A acrylamide solution  | 30 %         | x                |
| 7748  | ROTIPHORESE® Gel A-40 acrylamide solution                                     | 40 %         | x                |
| 3039  | ROTIPHORESE® Gel B bisacrylamide solution                                     | 2 %          | x                |
| <b>Sequencing gel solutions, ready-to-use</b>               |   |              |                  |
| A431  | ROTIPHORESE® DNA sequencing system  | Kit          | 3043. 3047. 3050 |
| 3043  | ROTIPHORESE® sequencing gel concentrate:<br>urea and acrylamide/bisacrylamide | 50 %<br>25 % | 19:1             |
| 3047  | ROTIPHORESE® sequencing gel diluent: urea                                     | 50 %         |                  |
| 3050  | ROTIPHORESE® sequencing gel buffer concentrate<br>urea in 10x TBE             | 50 %.        |                  |

|  |       |        |      |
|--|-------|--------|------|
| <b>ROTIPHORESE®Gel 30 (37,5:1)</b>                       | glass | 1 L    | 3029 |
| <b>ROTIPHORESE®Gel 40 (19:1)</b>                         | glass | 1 L    | 3030 |
| <b>ROTIPHORESE®Gel 40 (29:1)</b>                         | glass | 1 L    | A515 |
| <b>ROTIPHORESE®Gel 40 (37,5:1)</b>                       | glass | 1 L    | T802 |
| <b>ROTIPHORESE®NF-Acrylamid/Bis-solution 40 %(19:1)</b>  | glass | 250 ml | A516 |
| <b>ROTIPHORESE®NF-Acrylamid/Bis-solution 40 % (29:1)</b> | glass | 250 ml | A121 |
| <b>ROTIPHORESE®NF-Acrylamid/Bis-solution 30 % (29:1)</b> | glass | 1 L    | A124 |

### Carl Roth GmbH + Co. KG

Schoemperlenstraße 3-5 • 76185 Karlsruhe

P.O. Box 100121 • 76231 Karlsruhe

Phone: +49 (0) 721/ 5606-0

Fax: +49 (0) 721/ 5606-149

info@carlroth.com • www.carlroth.com

sse 06/2021

The company is a limited partnership with headquarters in Karlsruhe, reg. court Mannheim HRA 100055. Roth Chemie GmbH, with headquarters in Karlsruhe, reg. court Mannheim HRB 100428, is the personally liable partner. Managing Director: André Houdelet. Sales tax identification number: DE 143621073.