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Neisser-Lösungen I-III

Best.-Nr.: 3463; 3464; 3466

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Neisser's Solutions I-III

Art. No.: 3463; 3464; 3466

Microbiological staining method for visualising diphtheria bacteria

The Neisser staining is used for detection of special ingredients of cytoplasm – so-called polar nuclei (metachromatic granules) - typically for some Gram positive bacteria, especially diphtheria bacteria (*Corynebacterium diphtheriae*). Therefore, Neisser staining is very important for diagnosis of diphtheria.

For use with air-dried, heat-fixed smears of various cytological materials such as sputum, pus etc.

The staining is carried out in two steps:

First the polar nuclei are stained with a mixture of Neisser's solutions I and II.

Then the bacterium body is counterstained with Neisser's solution III.

Reagents:

- *Neisser's solution I* 3463.1 100 ml
 (Methylen blue solution) 3463.2 500 ml
- *Neisser's solution II* 3464.1 100 ml
 (Crystal violet solution) 3464.2 500 ml

Working solution

Methylen blue Crystal violet solution - mix prior to use:

- 2 parts Neisser's solution I
- 1 part Neisser's solution II

- *Neisser's solution III* 3466.1 100 ml
 (Chrysoidine solution) 3466.2 500 ml

Instructions:

For air-dried, heat-fixed smears.

1. Stain with Methylen blue Crystal violet solution. 30 sec	2. Wash briefly with aqua dest.
3. Counterstain with Neisser's solution III. 1 min	4. Wash briefly with aqua dest. Dry between sheets of filter paper.

Result: Polar nuclei: dark brown to black; bacterium body: brownish

Note: The colour intensity depends on the pre-treatment and the composition of the samples to be stained. It may initially be necessary to adapt the method to the respective conditions.

 **Warning** H371-H412

Neisser's solution I	100 ml	3463.1
	500 ml	3463.2

Neisser's solution II	100 ml	3464.1
	500 ml	3464.2

Neisser's solution III	100 ml	3466.1
	500 ml	3466.2

Carl Roth GmbH + Co. KG

Schoemperlenstraße 3-5 • 76185 Karlsruhe
P.O. Box 100121 • 76231 Karlsruhe
Phone: +49 (0) 721/ 5606-0
Fax: +49 (0) 721/ 5606-149
info@carlroth.com • www.carlroth.com

sse 06/2021

The company is a limited partnership with headquarters in Karlsruhe, reg. court Mannheim HRA 100055. Roth Chemie GmbH, with headquarters in Karlsruhe, reg. court Mannheim HRB 100428, is the personally liable partner. Managing Director: André Houdelet. Sales tax identification number: DE 143621073.