



Guter Rat ist Roth.

Information zu Ameisensäure

Empfohlene Lagerungsbedingungen:

Zersetzung unter Bildung gasförmiger Produkte möglich, besonders bei längerer Lagerung.

Daher Gefäße so verschließen, dass Innendruck entweichen kann (z.B. Überdrucksicherung).

Trocken und lichtgeschützt, an einem gut belüfteten Ort aufbewahren. Von Hitze- und Zündquellen fernhalten. Unter Verschluss oder nur für Sachkundige oder deren Beauftragte zugänglich aufbewahren.

Bitte ersetzen Sie beim Einlagern den Transportverschluss durch den beiliegenden Entgasungsverschluss. Gebinde mit Entgasungsverschluss müssen unter einem Abzug oder in einem gut belüfteten Lager gelagert werden. Der Transport von Ameisensäuregebinden mit Entgasungsverschluss ist nicht erlaubt.

Information on formic acid

Recommended storage conditions:

May decompose forming gaseous products, especially when stored over long periods.

Close vessels such that the internal pressure can escape (e.g. overpressure protection).

Store in a dry place that is protected from light and well-ventilated. Keep away from heat and ignition sources. Keep under lock and key or only accessible to competent personnel or their agents.

Please replace the transport cap with the degassing cap included when stored.

Packages with degassing cap must be stored beneath a vent or in a well-ventilated storage location.

The transport of formic acid packages with degassing cap is not permitted.