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## H&E-Schnellfärbekit

Best.-Nr. 9194

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## H&E Fast Staining Kit

Art. No. 9194

### Modified Hematoxylin-Eosin Staining (H&E Staining)

The H&E staining is a standard staining method in histology. It gives an overview of the different tissue structures.

The kit contains a modified hematoxylin solution for staining of cell nuclei and a modified eosin y solution 1 % for counterstaining of cell structures.

*The staining process is completed after ca. 14 minutes.*

#### Kit includes:

- H&E staining solution 1 (modified hematoxylin solution, Art. No. 8947) 250 ml / 500 ml
- H&E staining solution 2 (modified eosin y solution, Art. No. 8950) 250 ml / 500 ml

*The staining solutions should be filtrated before use! Solutions cannot be bought separately.*

#### Further Reagent:

Hydrochloric acid 0.1 %

#### Material:

Fixed, dewaxed tissue sections, last rinsing step: Aqua dest.  
All common fixing media can be used (no restrictions).

#### Instructions:

1. Incubate in H&E solution 1	6 min	4. Rinse in flowing tap water (blueing)	6 min
2. Rinse in tap water	10 sec	5. Incubate in H&E solution 2	30 sec
3. Differentiate with hydrochloric acid 0.1 %	10 sec	6. Rinse in flowing tap water	30 sec

#### Result:

- Cell nuclei: blue
- All other structures: reddish with different intensity  
e.g. cytoplasm: pink / erythrocytes: orange

#### Mounting:

All common mounting media are suitable. Before mounting with water free media please dehydrate by ascending alcohol series.

**Please note:** The colour intensity depends on the pre-treatment and the composition of the samples to be stained. It may initially be necessary to adapt the method to the respective conditions.

 **Danger** H225-H301+H311+H331-H318-H317-H370-H373

#### Carl Roth GmbH + Co. KG

Schoemperlenstraße 3-5 • 76185 Karlsruhe  
P.O. Box 100121 • 76231 Karlsruhe  
Phone: +49 (0) 721/ 5606-0  
Fax: +49 (0) 721/ 5606-149  
info@carlroth.com • www.carlroth.com

sse 06/2021

The company is a limited partnership with headquarters in Karlsruhe, reg. court Mannheim HRA 100055. Roth Chemie GmbH, with headquarters in Karlsruhe, reg. court Mannheim HRB 100428, is the personally liable partner. Managing Director: André Houdelet. Sales tax identification number: DE 143621073.

#### H&E Fast Staining Kit

<b>1 Kit</b>	<b>2 x 250 ml</b>	<b>9194.1</b>
<b>1 Kit</b>	<b>2 x 500 ml</b>	<b>9194.2</b>