## ROTICLEAR®

Best.-Nr. A538

### I. Eigenschaften

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Art. No. A538

### I. Characteristics

ROTICLEAR® is an organic solvent consisting of different hydrocarbons - similar to petroleum benzene. It is **nearly free from aromatic compounds (≤0.01 %)**. For it is flammable (flp 25 °C), it should not be used near open flame.

ROTICLEAR® has only **a slight inherent smell** and can be mixed with IPA, butanol, ethanol and acetone. It evaporates easily. **The vapours are - in contrary to xylene- not harmful.** Nevertheless, we recommend working in a fume cupboard, because high concentrations of vapour can cause drowsiness. Like most solvents ROTICLEAR® has a strong degreasing effect, therefore, contact with skin and eyes should be avoided.

It is totally non-polar and does not absorb water.

### II. Application

ROTICLEAR® can be used like xylene

- **for dewaxing histological sections**
- **as a clearing agent**
  - **during paraffin-embedding**
  - **prior to coverslipping with hydrophobic mounting media**

ROTICLEAR® produces non-sticky cuts and, due to a short evaporation period, enables fast and effective working.

ROTICLEAR® **can be used in automates without any difficulties.** The absorption capacity for paraffin is even better than that known from xylene. Therefore, water baths for removing the paraffin need not to be changed so frequently.

**Roti®- Mount (Art. No. HP68) is an appropriate mounting medium.** You can start coverslipping directly after clearing with ROTICLEAR®.

ROTICLEAR® is suitable for immunohistochemical assays (e.g. for working with antisera and monoclonal antibodies).

### III. Storage

Store ROTICLEAR® **in tightly closed glass bottles** to avoid evaporation. It does not cause harm to plastics.

**nlg Danger** H304-H226-H336-H412-EUH066  
P210-P261-P280-P301+P310-P331

<b>ROTICLEAR®</b>	<b>A538.1</b>	<b>1 l</b>
	<b>A538.5</b>	<b>2,5 l</b>
	<b>A538.6</b>	<b>5 l</b>
	<b>A538.2</b>	<b>10 l</b>
	<b>A538.3</b>	<b>25 l</b>
	<b>A538.4</b>	<b>200 l</b>

#### Carl Roth GmbH + Co. KG

Schoemperlenstraße 3-5 • 76185 Karlsruhe  
P.O. Box 100121 • 76231 Karlsruhe  
Phone: +49 (0) 721/ 5606-0  
Fax: +49 (0) 721/ 5606-149  
info@carlroth.com • www.carlroth.com ip 07/2021

The company is a limited partnership with headquarters in Karlsruhe, reg. court Mannheim HRA 100055. Roth Chemie GmbH, with headquarters in Karlsruhe, reg. court Mannheim HRB 100428, is the personally liable partner. Managing Director: André Houdelet. Sales tax identification number: DE 143621073.