



## Instruction Components

Push-button, suction tip with sucker.

## Application

aspirette is an accu-operated vacuum tweezer for picking-up and placing flat objects up to a maximum weight of 70 grams.

## Preparation

Before use charge the batteries with the enclosed charger. Plug the recharger into the aspirette and the recharger into the socket (7 hours are sufficient though overloading is harmless).

## Operation

Select the suction tip most suited to the object's size. Using only the suction jet - with the suction tip removed - very small and light objects may be picked up. The vacuum pump is operated by depressing the button and simultaneously the airtight channel is closed by the finger. After picking up and placing the object, the vacuum pump is switched off when the finger is lifted from the push button, destroying the vacuum.

## Air Filter

Between suction jet insert and the apparatus there is a particle filter, which guarantees the removal of dust particles from the pump system.

### Attention:

Working without the air filter will lead to the pump breaking down after only a very short time!

## Maintenance

With ordinary use the aspirette is maintenance free!

## Ni / Cd batteries

The built-in batteries are rechargeable up to 1000 times for long years use. Due to the integral Ni / Cd batteries after 1st lifetime the unit should be sent for special waste management or be returned to us for disposal.

## Spare Parts

- 990 16 04 air filter set (5 pcs. filter paper)
- 990 16 05 suction tip set  
(3 suction tips with suckers of 3, 2, 6 and 10 mm Ø).

## Exchange of filter paper

