# HIRSCHMANN®

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## Instruction Manual

Please read before use!

# pipetus® junior

Manuelle Pipettierhilfe  
Manual pipet filler

Hirschmann Laborgeräte GmbH & Co. KG  
Hauptstraße 7-15 · 74246 Eberstadt Germany  
Fon +49 7134 511-0 · Fax +49 7134 511-990  
[www.hirschmannlab.com](http://www.hirschmannlab.com) · [info@hirschmannlab.com](mailto:info@hirschmannlab.com)

X2120472001

## Inhalt

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We are delighted that you have chosen our pipet filler. The unit is a manual pipet filler which will facilitate your work considerably.

The pipet filler is designed to the latest technical standards and is operationally safe. However, residual dangers can be associated with use of the unit, so it is imperative that the instruction manual be read carefully and all instructions therein heeded.

All the same, the instruction manual cannot illustrate every safety problem which may arise. Each user bears individual responsibility for adhering to safety and health regulations and determining and heeding limitations prior to using the unit.

Please observe the safety instructions prior to start up in all cases!

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## 1. Safety Instructions

- Observe all general safety practices and guidelines for laboratories, such as the use of protective clothing, safety goggles and gloves when handling hazardous chemicals.
- Never point the pipette in anyone's direction when in use!
- Observe the guidelines and procedures of the chemical manufacturers (see also point 5. and 6.).

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---

### **Attention!**

If the unit is opened, the warranty is invalidated and any resultant claims void.

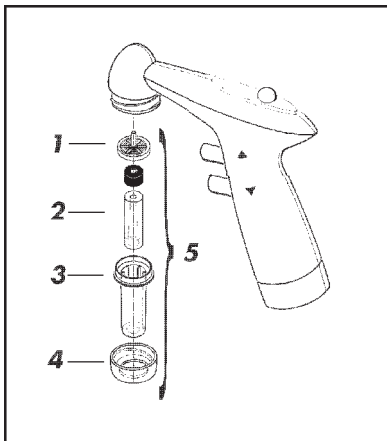
---

- The filter supplied contains a hydrophobic 0.2 µm membrane for air filtration in either direction.
  - Only have the unit professionally serviced.
  - Never use undue force!
- 

### **Attention!**

Only ever use the original accessories and replacement parts to avoid damaging the unit.

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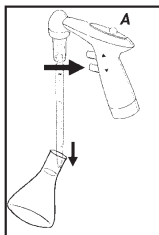
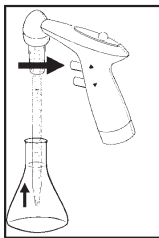
### 3. Functional principle

- The unit is a manual pipet filler for use with all standard pipettes from 0.5 ml to 100 ml.

The medium is suctioned into the pipette by pressing the PUMP button. Complete depression conveys approximately 2 ml of water, whereas partial or repetitive pressing conveys a respective volume.

The pipette is emptied through drainage (EX) when the OUT button is pressed. Drainage speed can be regulated by reducing pressure on the OUT button.

Residual liquid in the tip of BLOW-OUT pipettes is expelled by pressing the rubber ball (A) once.

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#### 4. Scope of delivery

The following components should be contained in the original carton:

- Pipet filler
- 1 hydrophobic membrane filter 0.2  $\mu\text{m}$  (yellow)
- 1 instruction manual

Please check that all parts have been delivered and are undamaged!

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---

### **The unit can not be used with:**

liquids of which vapours attack the materials silicone, PTFE, Novodur, PVDF, NBR, PCPA, PP and PA6.

---

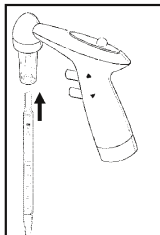
## 6. Restricted operations

Liquids which emit copious volumes of gas, as vapours are deposited inside the pipetting aid and can corrode the unit.

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## 7. Pipetting

### 7.1 Place the pipette

Place the pipette into the holder to ensure a  
secure and airtight fit!

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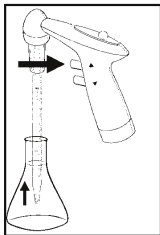
### **Attention!**

Danger of glass breakage, especially with thin pipettes.

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## 7.2 Fill the pipette

Press the PUMP button to fill the pipette, filling the pipette to a level just above the desired volume. The volume can be set precisely by pressing the OUT button lightly.

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---

### Attention!

No liquid should get into the unit as this can adversely affect the operating functions.

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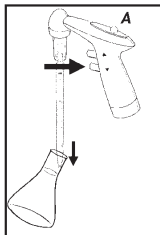
### 7.3 Empty the pipette

The pipette is emptied through drainage (EX) when the OUT button is pressed.

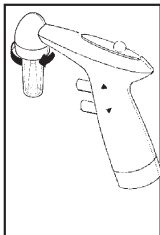
Drainage speed can be regulated by reducing pressure on the OUT button.

Residual liquid in the tip of BLOW-OUT pipettes is expelled by pressing the rubber ball (A) once.

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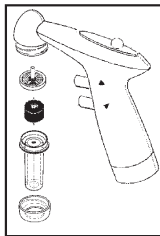
## 8. Spare parts replacement

### 8.1 Replacement of pipetholder complete

- Screw off the nut.
- Replace the pipetholder by a new one.

### 8.2 Replacement of filter

Screw off the nut.  
Remove the old filter and place the new one in the pipetholder case so that the pivot shows upwards. Screw the pipetholder onto the unit, by turning the nut clockwise.

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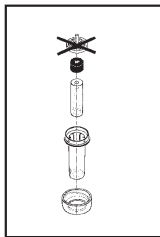
## 9. Maintenance, Sterilisation and Cleaning

### 9.1 Maintenance

Under normal conditions of use, the unit is maintenance free.

In the event, nevertheless, that they do become dirty or the suction becomes impaired, the membrane filter should be replaced.

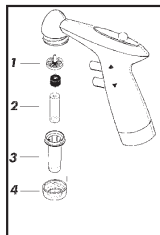
### 9.2. Sterilisation

The pipetholder, except the membrane filter, can be sterilised at 121°C.

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### 9.3 Cleaning

The pipetting aid handles can be cleaned with a damp cloth.

Do not spray the unit, as liquid can penetrate to the interior.

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---

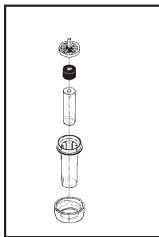
#### **Attention!**

When cleaning and exchanging the safety valve and membrane filter, protective gloves and goggles must be worn (danger of infection)!

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## 10. Test of tightness

Once the required volume has been aspirated into the pipette, the liquid has to be kept in it unless the buttons are pressed. If the pipette however is dripping, disassemble the pipetholder and reassemble it again carefully.

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## 11. Repair at the manufacturer

Please observe the following points should it prove necessary to return the unit for repair (despite having handled it with all due care):

- Before sending the unit back, the user must ensure that it is carefully cleaned and decontaminated. The healthwise harmlessness must be confirmed, send the confirmation together with the unit to the manufacturer or the dealer. (A copy of the clearance certificate can be found on page 22).

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#### 11. Repair at the manufacturer

- Units will only be repaired, when they have been cleaned and decontaminated by the user.
- For safety reasons dirty units will be returned to the sender without being repaired.
- Please describe the malfunction exactly.
- Please quote the pipetted liquid.
- In the event of the unit being tampered with or being used in any improper way, (e.g. the application of undue force) any claims under the warranty are void.

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## 12. Technical data

CE:

This mark is the confirmation that the unit conforms to the established requirements of the EU guidelines and has been tested according to these criteria.

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### 13. Accessories and spare parts

The spare parts for your pipet filler listed below can be obtained from your specialised laboratory retailer:

|   |            |
|---|------------|
| Silicone adapter  | 990 32 01  |
| Pipetholder complete                                      | 990 24 00  |
| Pipetholder case  | 990 38 03  |
| Nut magenta   | 990 24 02  |
| Round filter set<br>(consisting of 5 replacement filters) | 990 37 04  |
| O-ring for cone   | P990 32 05 |
| Wall bracket (incl. dowels and screws)                    | 990 56 00  |
| Stand   | 990 57 00  |

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for unit .....(Serial number)

Medium used: .....

|   | YES                      | NO                       | Please define in case of YES |
|---|--------------------------|--------------------------|------------------------------|
| Blood, bodily fluids, patholog. samples                 | <input type="checkbox"/> | <input type="checkbox"/> | _____                        |
| Other hazardous biologic substances                     | <input type="checkbox"/> | <input type="checkbox"/> | _____                        |
| Biologically-degradable materials which can pose a risk | <input type="checkbox"/> | <input type="checkbox"/> | _____                        |
| Chemicals containing substances harmful to health       | <input type="checkbox"/> | <input type="checkbox"/> | _____                        |
| Radioactive materials                                   | <input type="checkbox"/> | <input type="checkbox"/> | _____                        |
| Other hazardous and toxic substances                    | <input type="checkbox"/> | <input type="checkbox"/> | _____                        |

### We hereby declare that

- the unit has been carefully cleaned and decontaminated prior to shipment,
- no risks resulting from contamination exist,
- we are aware that liability claims can be made against us in the event of damage being caused by contaminated units.

### Address of sender:

Name: \_\_\_\_\_

Date: \_\_\_\_\_

Signature: \_\_\_\_\_