





Blender

Dear Customer

Before using the appliance, the user should read the following instructions carefully.

Connection to the mains supply

Make sure that the supply voltage corresponds with the voltage marked on the rating label of the appliance. This product complies with all binding CE labelling directives.

Familiarisation

1. Feeder opening stopper
2. Blender jug
3. Motor housing
4. Push switch
5. Control switch
6. Safety switch
7. Feeder opening
8. Lid
9. Contact pin
10. Power cord with plug
11. Rating label
12. Blade
13. Drive shaft coupling

Important safety instructions

- Before the appliance is used, the main body as well as any attachments should be checked thoroughly for any defects. If the appliance, for instance, has been dropped onto a hard surface, it must no longer be used: even invisible damage may have adverse effects on the operational safety of the appliance.
- Do not leave the appliance unattended while in use and keep children away.
- This appliance is not intended for use by any person (including children) with reduced physical, sensory or mental capabilities, or lacking experience and knowledge, unless they have been given supervision or instruction concerning the use of the appliance by a person

responsible for their safety.

- Children should be supervised to ensure that they do not play with the appliance.
- **Caution:** the blade is extremely sharp. To prevent injuries, take special care when cleaning the appliance.
- Do not touch any moving parts during operation - this can be extremely dangerous.
- Do not allow the appliance or its power cord to touch hot surfaces or to come into contact with any heat sources.
- **Always switch the appliance off and remove the plug from the wall socket,**
 - before opening the lid,
 - before removing or fitting the blender jug,
 - before cleaning,
 - in case of any malfunction during use,
 - after use.
- When removing the plug from the wall socket, never pull on the power cord; always grip the plug itself.
- No responsibility will be accepted if damage results from improper use, or if these instructions are not complied with.
- This appliance is only intended for domestic use and not for commercial applications.
- In order to comply with safety regulations and to avoid hazards, repairs of electrical appliances must be carried out by qualified personnel. If repairs are needed, please send the appliance to one of our customer service departments. The address can be found in the appendix to this manual.

Before using for the first time

Before the appliance is used for the first time, the blender jug and the lid should be thoroughly cleaned as described in the section *General Care and Cleaning*.

Switch settings

The following settings can be selected:

- P** The appliance remains switched on as

- long as the switch is depressed
- 0 The appliance is switched off
 - 1 Low power
 - 2 Medium power
 - 3 Maximum power

Short-term operation

The appliance is designed to process average domestic quantities of food; it can be operated without interruption for a maximum of 45 seconds, after which it should be allowed to cool down for about a minute.

Safety switch

A special safety switch in the motor housing ensures that the appliance can only be operated once the blender jug and lid are properly locked in position.

Before removing the blender jug, make sure that the control switch is in position **0**. This also applies when fitting the blender jug to the motor housing. If the switch is not in its off-position (**0**), the motor will start immediately.

Operation

- Fit the blender jug to the motor housing with its handle on the right and lock it by turning anti-clockwise into the operation position.
- Put the liquid ingredients into the blender jug. The jug has a maximum capacity of 1500 ml.
- Put the lid on and lock it by turning anti-clockwise. To add ingredients during operation, the stopper in the centre of the lid can be removed by turning it in an anti-clockwise direction and then lifting it off. Do not put any objects into the blender jug while the appliance is operating.
- Insert the plug into the wall outlet and use the control switch to select the required power setting. Depending on the food to be processed, the appliance may also be operated for short periods of time with the pulse setting "**P**". Do not

operate when the blender jug is empty.

- After using the blender, switch off and remove the plug from the wall socket. Wait until the motor has come to a complete standstill before removing the blender jug. To remove the jug and the lid, turn them in the opposite direction from that described in the assembly instructions.

General Care and Cleaning

- Always remove the plug from the wall socket before cleaning the appliance.
- To avoid the risk of electric shock, do not clean the motor housing or the power cord with water and do not immerse them in water. The motor housing may be wiped with a slightly damp, lint-free cloth.
- Do not use abrasives, harsh cleaning solutions or hard brushes for cleaning, and do not clean in a dish-washer.
- For cleaning after use, fill the blender jug with approximately 1250 ml warm water and a mild cleaning detergent. Let the appliance run for a short period as described in the section *Operation*. Then rinse the blender jug, lid and stopper with fresh water. All parts must be thoroughly dried afterwards.

Disposal



Do not dispose of old or defective appliances in domestic garbage; this should only be done through public collection points.

Guarantee

This product is guaranteed against defects in materials and workmanship for a period of two years from the date of purchase. Under this guarantee the manufacturer undertakes to repair or replace any parts found to be defective, providing the product is returned to one of our authorised service centres. This guarantee is only valid if the appliance has been used in accordance with the instructions, and provided that it has not

been modified, repaired or interfered with by any unauthorised person, or damaged through misuse.

This guarantee naturally does not cover wear and tear, nor breakables such as glass and ceramic items, bulbs etc. If the product fails to operate and needs to be returned, pack it carefully, enclosing your name and address and the reason for return. If within the guarantee period, please also provide the guarantee card and proof of purchase.