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**Scaled magnifier with illumination**

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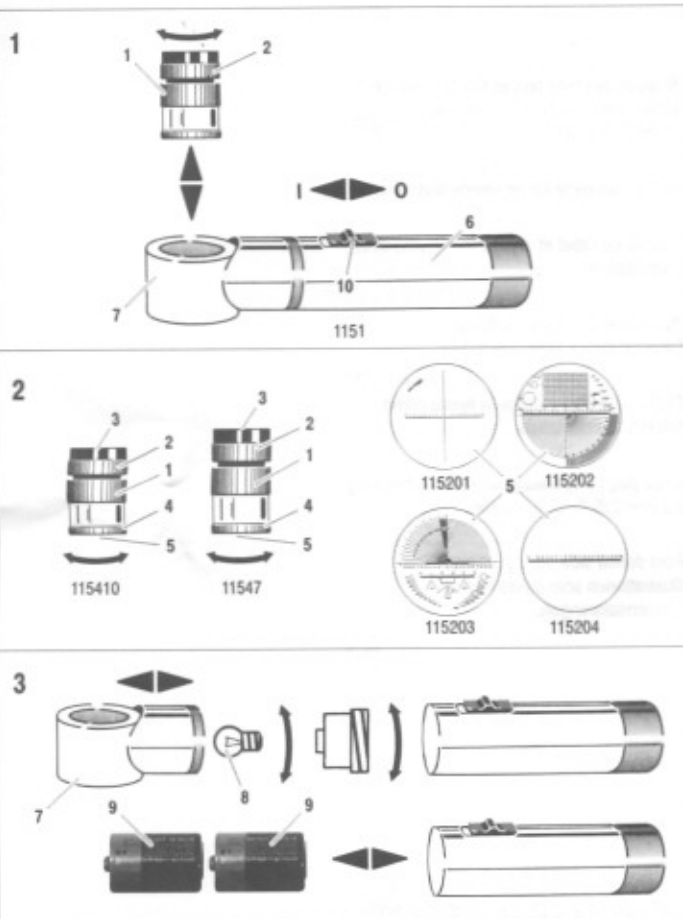
**Batterie- / Lampenwechsel**

(→ Abb. 3)

- [1] Precision scaled magnifier (lenses ø 23.5 mm)
  - No. 11547 (7x magnification)
  - No. 115410 (10x magnification)
- [2] Focusing
- [3] Achromatic lens system
- [4] Locking ring for measuring scale
- [5] Measuring scales
  - No. 115201 (standard scale)
  - No. 115202, 115203, 115204
- [6] Illumination unit
- [7] Retainer for scaled magnifier
- [8] Light bulb, round 2.5 V/200 mA (Order No. 1545)
- [9] Batteries (2x), Baby type, each 1.5 V (Order No. 1547)
- [10] Switch: I = ON; O = OFF

## Lenses

The precision scaled magnifier has a coated achromatic (colour defect correcting) lens system which provides a sharp brilliant image.

## Procedure (→ Fig. 1)

- Place the illumination unit [6] on a level support and push the scaled magnifier [1] into the retainer [7] until it touches the support.
- Spectacle wearers do not have to remove their spectacles to use the scaled magnifier.

## Changing the measuring scale (→ Fig. 2)

1. To change the measuring scale, unscrew it [5] from the locking ring [4].
2. Insert the required measuring scale into the scaled magnifier so that the numbers and measuring sections on the measuring scale can be seen on the side through the lens [3].
  - To avoid damaging the measuring scale, only hold it by the sides and store it in a safe place!

## Battery / bulb replacement (→ Fig. 3)

Do not touch the lens with your fingers. The lens is coated with a special anti-reflective coating. Do not use alcohol or other solvents to clean the lens. Use only a soft, lint-free cloth to clean the lens.

## ⚠ Safety instructions

- **Danger of fire!**  
Lenses used in optical equipment may cause considerable damage through their „burning glass effect“ if used or stored incorrectly. Ensure that optical lenses are never left uncovered in the sunlight.
- **Danger of blinding or injury!**  
Never look at the sun through any optical equipment.

## Instructions for care

- Do not clean the scaled magnifier with alcoholic solvents or with cleaners that contain softeners!

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