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## Description:

QUANTOFIX® Chlorine is a test kit for semi-quantitative determination of free chlorine in solutions.

## Pack content:

1 aluminum container with 100 test strips  
 1 bottle Cl<sub>2</sub>-1 1 measuring vessel with 5 mL mark  
 1 bottle Cl<sub>2</sub>-2 1 small measuring spoon

## Measuring range:

1–100 mg/L

## Color gradation:

0 · 1 · 3 · 10 · 30 · 100 mg/L

## Safety precautions:

Reagent Cl<sub>2</sub>-2 contains pyridine 32–44% CAS 110-86-1 and potassium cyanide 0.1–1% CAS 151-50-8. **WARNING** Flammable liquid and vapor. Harmful if swallowed. Harmful in contact with skin. Harmful if inhaled. Contact with acids liberates toxic gas. Keep away from heat/sparks/open flames/hot surfaces. – No smoking. Keep container tightly closed. Do not breathe vapors. **IF SWALLOWED:** Call a POISON CENTER or doctor/physician if you feel unwell. **IF ON SKIN:** Wash with plenty of soap and water. **IF INHALED:** Remove to fresh air and keep at rest in a position comfortable for breathing. Rinse mouth. Store in a well-ventilated place. Keep cool. With reduced labeling acc. CLP Annex I chapter 1.5.2.1. For further information, please ask for safety data sheet (see [www.mn-net.com/MSDS](http://www.mn-net.com/MSDS)).

## General indications:

Remove only as many test strips as are required. Close the container immediately after removing a strip. Do not touch the test field.

## Specific indications:

To obtain correct results with this test, the pH of the sample must be < 10. Therefore, the pH value of the sample must first be measured (for example with pH-Fix 0–14, REF 92110). If the solution has a pH > 10, add some drops of diluted sulfuric acid to lower it to 6–7.

## Instructions for use:

1. Rinse the measuring vessel with the test solution and fill it to the 5 mL mark.
2. Add 1 spoon Cl<sub>2</sub>-1 (buffer solution) to the sample.
3. Add 5 drops Cl<sub>2</sub>-2.
4. Shake carefully.
5. Immediately dip the test strip into the sample.
6. Wait 45 s.
7. Shake off excess of liquid.
8. Compare within 10 s with the color scale.

## Interferences:

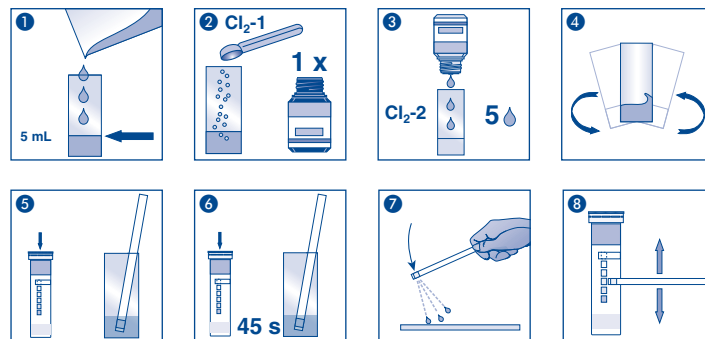
Bromide and iodide in concentrations of 5 mg/L and above lead to a fading of the color reaction. Major concentrations may inhibit the color reaction. Nitrite can also cause a decrease in color intensity with corresponding low test results. If the pH value of the sample is > 10, then the capacity of the buffer solution will not be sufficient to lower the pH to the optimal value of 6–7. In this case, prior to the test, the pH value of the sample should be adjusted to 6–7 by dropwise addition of diluted sulfuric acid.

## Disposal:

The used analysis specimens can be flushed down the drain with tap water and channeled off to the local sewage treatment works. Used test strips can be disposed in household waste.

## Storage:

Avoid exposing the strips to sunlight and moisture. Keep container cool and dry (storage temperature 4–30 °C). If correctly stored, the test strips may be used until the use-by-date printed on the packaging.

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