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Operating instructions electronic probe thermometer **with max.-min.-function art.no. 30.1018**

With this probe thermometer you have bought an electronic precision device for measuring the current temperature as well as the maximum and minimum temperature.

measuring range: -40 °C to +200 °C (-40 °F to 392 °F)

operating temperature: -20 °C to +50 °C (-4 °F to 122 °F)

precision: +/- 1 °C (+/- 1.8 °F) from -20°C to + 100°C (-4°F to 212°F), +/- 2°C (+/- 3.6°F) otherwise

resolution: 0.1°C (0.1 °F)

battery 1.5 V button cell

auto switching-off

function: after 10 minutes

Use:

1. Open the battery cover with a coin in an anti-clockwise direction, remove the battery stripe and replace the cover. Remove the protection foil of the display. The thermometer is now ready for use.
2. The display shows the currently measured temperature. The sensor is located in the point of the probe. The temperature unit can be changed from °C to °F by the °C/°F-key.
3. Press the MODE-key once in normal mode and the hold-function is activated (indicated by HOLD).
4. Press the MODE-key again and the maximum temperature is shown since the last reset (indicated by MAX).
5. Press the MODE-key again and the minimum temperature is shown since the last reset (indicated by MIN).
6. Press the MODE-key again to return to normal mode.
7. To clear the maximum and minimum memory, press and hold the MODE-key for 3 seconds in the corresponding mode while MAX or MIN is displayed (indicated by ---). Press the MODE-key again to return to normal mode.

Attention: Used batteries have to be put in the specially reserved collecting receptacles.

Source of error: Please change the battery before complaining.

No guarantee if the instrument is handled or opened improperly.