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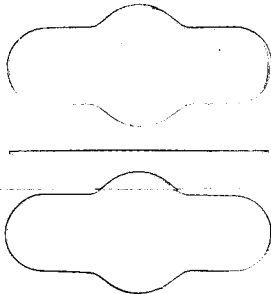
### OPERATING INSTRUCTIONS:

Open the battery cover in the direction of the arrow and insert the AAA-battery, +/- polarity as indicated. Close the battery cover and remove the protection foil from the display.

### OPERATION:

#### 1. Timer:

- Set the desired time with the buttons 0-9.
- Push the "START/STOP" button to begin the countdown.
- To stop the countdown process, press the "START/STOP" button. Push the "START/STOP" button again to continue the countdown process.
- When the timer counts down to "0.00.00", a beeping alarm will sound. The timer is now counting the elapsed time until the alarm is turned off by the START/STOP button. The timer will recall the last selected time automatically.
- Press the CLEAR button to reset the display.

#### 2. Stopwatch:

- Push the "START/STOP" button when 0.00.00 is indicated to begin the counting up.
- To stop the count-up process, press the "START/STOP" button. Push the "START/STOP" button again to continue the count-up process.
- Press the CLEAR button to reset the display.

**Source of error:** Please change the battery before complaining.

**Caution:** Used batteries have to be put in the specially reserved collecting receptacles!

No guarantee if the instrument is handled or opened improperly.