**MIN/SEC 3 KEYS  
COUNTDOWN TIMER**

**MODEL: TR112**

*User Manual*

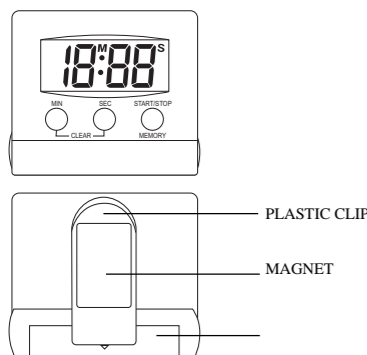
FIG. 1

**ENGLISH**

**TIME FEATURE**

- LCD count-down timer.
- 4 digits display showing minute time ("M" mark) and second time ("S" mark).
- Timer can be programmed up to a maximum of 99 minutes, 59 seconds and counts down at 1 second intervals.
- Buzzer alarm output when timer counts down at 1 second intervals.
- Timer memory recall function.
- Individual buttons for minute and second setting.

**TIMER TIME SETTING [FIG. 1]**

- Press **MIN** and **SEC** buttons at the same time to reset timer to zero.
- Press **MIN** button to advance minute digits (sound can be heard). Press and hold **MIN** button to advance minute digits more quickly.
- Press **SEC** button to advance second digits (beep sound can be heard). Press and hold **SEC** button to advance second digits more quickly.

**TIME START/STOP**

- After time setting is ready, press **START/STOP** button once and timer will start to count down, "M" and "S" marks will flash when timer is running.
- When timer is counting down, press **START/STOP** button once and timer will stop. "M" and "S" marks will stop flashing and remain on the display.
- Press **START/STOP** button once and timer will resume counting again.

**TIME BUZZER ALARM**

- When timer counts down to 00M and 00S, timer buzzer alarm will sound for 30 seconds.
- Timer buzzer alarm can be stopped by either pressing **MIN**, **SEC** or **START/STOP** buttons.

**TIMER MEMORY RECALL**

- After timer buzzer alarm stops, press **START/STOP** button once to recall pre-set timer and timer will count down for another lap.
- Press **START/STOP** button second time can start timer and timer will count down for another lap.

**BATTERY REPLACEMENT**

Follow the arrow direction to open the battery cover at back side of timer.

Remove old battery, insert a new 1.5V "AAA" battery as indicated by the polarity symbols (+&-) marked and close battery cover.

**CLIP ATTACHMENT**

The timer can clip onto a thin board or shirt pocket, using the plastic clip at rear of timer.

**MAGNETIC ATTACHMENT**

The timer can attach to iron or steel surfaces, using the magnet at rear of timer.

**TABLE TOP USE**

The timer can stand on table surface, flip out the metal stand bar at rear of timer.

**PRECAUTION**

When handled properly, this unit is engineered to give you years of satisfactory service. Here are a few product care instructions:

- Do not immerse the unit in water. If the unit comes in contact with water, dry it immediately with a soft lint-free cloth.
- Do not clean the unit with alcohol containing detergent, abrasive or corrosive materials. Abrasive cleaning agents may scratch the plastic parts and corrode the electronic circuit.
- Do not subject the unit to excessive: force, shock, dust, temperature, or humidity. Such treatment may result in malfunction, a shorter electronic life span, damaged batteries, or distorted parts.
- Do not tamper with the unit's internal components. Doing so will terminate the unit's warranty and may cause damage. The unit contains no user-serviceable parts.
- Only use new batteries as specified in this instruction manual. Do not mix new and old batteries as the old batteries may leak.
- Read this instruction manual thoroughly before operating the unit.

**CAUTION**

- The content of this manual and the product specifications are subject to change without further notice.
- The technical specifications of this product are subject to change without notice.
- The contents of this manual may not be reproduced without the permission of the manufacturer.

**ABOUT OREGON SCIENTIFIC**

Visit our website ([www.oregonscientific.com](http://www.oregonscientific.com)) to learn more about Oregon Scientific products such as digital cameras; MP3 players; children's electronic learning products and games; projection clocks; health and fitness gear; weather stations; and digital and conference phones. The website also includes contact information for our customer care department in case you need to reach us, as well as frequently asked questions and customer downloads.

We hope you will find all the information you need on our website, however if you'd like to contact the Oregon Scientific Customer Care department directly, please visit:

[www2.oregonscientific.com/service/support](http://www2.oregonscientific.com/service/support) OR Call 949-608-2848 in the US.

For international enquiries, please visit: <http://www2.oregonscientific.com/about/international/default.asp>

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