MODE D'EMPLOI

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A - IMPORTANT : ALWAYS BE CAREFUL WHEN USING GAS !

The purpose of these instructions is to enable you to use your Campinggaz® LABOGAZ 206 correctly and in complete safety.
 Please read them carefully to familiarise yourself with the equipment prior to assembling the gas container.
 Please observe these instructions and the safety measures printed on the Campinggaz® C 206 cartridges. Failure to observe these instructions may result in the user and people in the immediate vicinity being put at risk.
 Keep these instructions in a safe place so that you can refer to them if needed.
 This equipment must be used exclusively with Campinggaz® C 206 cartridges. Using other gas containers may be dangerous. The Société Application Des Gaz accepts no responsibility when cartridges of any other make are used.
 This equipment must only be used in adequately ventilated areas (minimum 2 m³/h/kW) and at a good distance from any flammable material.
 Do not use equipment which leaks, functions incorrectly or is damaged. Return it to your retailer who will inform you of your nearest after-sales service.
 Never modify this equipment or use it for purposes for which it is not intended.

B - PUTTING INTO OPERATION

INSTALLING THE CAMPINGGAZ® C 206 CARTRIDGE (fig.2)

(When the existing cartridge is empty, read paragraph D : "Removing the cartridge"). To install or remove a cartridge, always work in a well-ventilated area, preferably outside and never in the presence of a flame, heat source or spark (lit cigarette, electrical equipment, etc.) at a good distance from other people and flammable material.

- a) Remove the equipment from its protective box.
- b) Check the seal (4) is present and in good condition (fig.1). Do not use the appliance if the seal is missing or damaged, return it to your retailer.
- c) Check that the end of the piercing pin does not protrude beyond the seal. It must be just level (see detail fig. 1).
- d) Open the cartridge holder (6) following arrows A then B (fig.2). Check that the area where the C 206 cartridge is pierced is totally clean (no label, or adhesive). Place the C 206 cartridge inside the cartridge holder (6) following arrow C and close following arrows D. Check that both plastic parts are locked together properly following arrows E and F.
- e) Check that the valve is properly closed by turning the knob (3) in a clockwise direction (following arrow A) (fig. 3).
- f) Stand the cartridge holder upright and tightly screw the valve/burner unit following arrow B (fig. 3). The bottom of the valve then rests on the threaded bush of the cartridge holder.
- g) The equipment is ready for use.

In case of leakage (gas smell before opening the valve), put immediately the equipment outside, in a well-ventilated area, without inflammation source, where it will be possible to locate and stop the leak. If you wish to check the tightness of the equipment, do it outside. Do not try to locate a leak with a flame, use gas leak detection liquid.

C - USING THE APPLIANCE

1. Additional instructions:

- Never use your equipment at a distance of less than 25 cm from a wall or combustible object. Position the equipment on a stable, horizontal surface and do not move when in use so as to avoid the appearance of large flames resulting from the combustion of liquid butane rather than vaporised butane. Should this occur, position the equipment and turn the tap off.
- 2. Lighting the burner (fig.4):**
 Bring a flame close to the burner (1) then gradually turn on the gas by turning the knob (3) anti-clockwise (see detail in fig. 4).
- 3. Adjusting the flame (fig. 4):**
 1/ Power: To increase or decrease the power of the flame, turn the knob (3) towards the "•" or "••" symbols on the arrow as applicable.
 2/ Temperature (fig. 4): To change the flame temperature, slide the washer (2) along the burner tube:
 - Primary air inlet open = very hot blue flame
 - Primary air inlet closed = oxidising yellow flame

4. Extinguishing:

Turn off the gas supply by turning the wheel (3) clockwise to its full extent ("•" direction of the arrow head).

D - CHANGING THE CARTRIDGE

- Never change a cartridge which is not empty (shake the appliance to check there are no liquid movements).
 a) Check that the gas connection is tight by turning the knob (3) fully in a clockwise direction.
 b) Unscrew the valve/burner unit and check that the seal (4) is present and in good condition (fig. 1).
 c) Open the cartridge holder (6) following arrows A then B (fig.2) and remove the empty cartridge.
 d) Insert a new cartridge, following the instructions given in paragraph B "Installing a CAMPINGGAZ® C 206 cartridge".

E - STORAGE AND TROUBLESHOOTING

In case of burn back (ignition under the burner cap) switch the product off, allow it to cool down (approximately five minutes), then relight the appliance. If the problem of burn back persists, then please contact your local Campinggaz® representative.

- Once your equipment has fully cooled**
 a) Store the equipment and the cartridge in a cool, dry and well-ventilated area, out of reach of children. Never store in a basement or cellar.
 b) The burner does not light when the valve is opened
 1/ If no sound of liquid can be heard when the unit is shaken, the cartridge is empty. Replace it with a new cartridge, following the instructions in paragraph D.
 2/ If the sound of liquid can be heard when the unit is shaken, there is some gas left but the injector is blocked. Check that the gas inlet is properly closed (turn the knob (3) as far as it will go in a clockwise direction). While holding the valve firmly, unscrew the burner (1) as indicated by arrow E (fig. 3) then the injector, as shown by arrow B (fig. 5). Replace it with a new or clean injector. Then screw the burner back in place (1). The defective injector can normally be easily unlocked simply by ejecting the impurity with a jet of compressed air directed against the external face of the injector after removing it from the unit. (Never use a needle or any other metal implement).
 c) Whenever the valve burner unit except to change the cartridge and then only if it is empty. Never work with the flame lit.
 d) If you use this equipment indoors, the area must observe elementary ventilation conditions ensuring the supply of air required for combustion and avoiding any build-up of a dangerous mixture of non-burnt gas (2 m³/h/kW minimum).

F - WARRANTY APPLICATION TERMS AND CONDITIONS

The product is covered by a total parts and labor warranty for a period of 2 (two) years from its purchase date. The warranty applies to products that do not conform to the order that are defective. Within the guarantee period, the product shall be either repaired, replaced or reimbursed - in full or in part. Liability of Buyer under a claim shall in no event exceed the price of the product. The warranty is null and void and does not apply should the damage arise (i) abuse of the product, (ii) failure to operate and maintain the products in accordance with the instructions of use (iii) repair, service, alteration or modification of the product by unauthorized third parties (iv) original parts are not used. The warranty is excluded in case of professional use. Proof of the purchase date (ie invoice, bill receipt) and claim report are required to obtain a free warranty service. Do not return the appliance with the gas container (cylinder or cartidge) fitted; ensure it is dismantled before sending back. Service under the guarantee does not affect the expiry date of the warranty. All other claims including for damages resulting from the use of the product, will be handled on a case-by-case basis. (The conditions of service clients 1604 and suivants et 1386.1 et suivants du Code Civil relatifs à la garantie légale. Consulter notre service consommateurs pour toute réclamation.

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