


Instruction Guide for the **MINI-CENTRIFUGE**


Congratulations on your purchase of a Carl Roth Personal Mini-Centrifuge!
Following are some recommended operating procedures:

Operation


- After unpacking, place the Mini-Centrifuge on a flat, clear surface.
- To open, press firmly down on release tab while lifting dome.
- Insert tubes into rotor completely. Be sure that each tube is securely closed and of equal volume. If necessary, add dummy-tubes to balance.
- Load rotor as indicated below. Properly balanced loads are essential for safe and quiet operation.

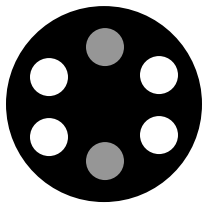


2 Tubes

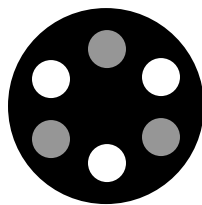

3 Tubes

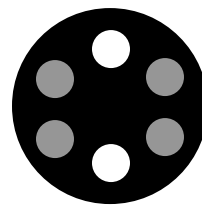

4 Tubes

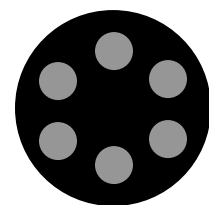

6 Tubes

- Close rotor securely and activate side switch. Device will not operate with dome open.
- To stop device, deactivate side switch. Wait until device has completely stopped to release dome. Remove samples.

Safeguard

- Do not immerse device in water. Injury due to shock or fire may result.
- Do not operate device with dome open, or open dome before rotor has ceased spinning.
- Do not insert non-standard tubes or other foreign articles into rotor.
- Do not operate under uneven loads. Destabilized operation may result in damage to the device or injury to the user.
- Store item at room temperature in a dry area. Do not expose to sun, moisture or extreme temperatures for prolonged periods of time.



Our one-year warranty assures you reliable quality at an unbeatable value!

**Carl Roth GmbH + Co. • Address: Schoemperlenstr. 3-5
Phone: (07 21) 56 06 - 0 • Fax: (07 21) 56 06 - 253 • www.carlroth.de**