## Kühl-Gefrierschrank-Thermometer

(D)

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## Digital freezer-fridge-thermometer

(GB)

- 1. Before you start using it**
  - Please make sure to read the instruction manual carefully.
  - Following and respecting the instructions in your manual will prevent damage to your instrument and loss of your statutory rights arising from defects due to incorrect use.
  - We shall not be liable for any damage occurring as a result of non-following of these instructions. Likewise, we take no responsibility for any incorrect readings and for any consequences which may result from them.
  - Please take particular note of the safety advice!
  - Please keep this instruction manual for future reference.
- 2. Field of operation and all the benefits of your new instrument at a glance**
  - Temperature measurement in rooms, refrigerator, etc.
  - Maximum-and minimum values
  - With stand and frame for mounting on surfaces or behind glass (display fridge, vegetable drawer)
  - According to EN 13485 Norm
- 3. For your safety**
  - This product is exclusively intended for the field of application described above. It should only be used as described within these instructions.
  - Unauthorized repairs, modifications or changes to the product are prohibited.
  - This product is not to be used for medical purpose or for public information, it is only intended for home use. The product is not recommended for taking body temperature.

### Caution! Risk of injury:

- Keep this instrument and the battery out of reach of children.
- Batteries must not be thrown into a fire, short-circuited, taken apart or recharged. Risk of explosion!
- Batteries can be fatal if swallowed. If a battery has been swallowed, get medical assistance immediately.
- Batteries contain harmful acids. Low batteries should be changed as soon as possible to prevent damage caused by leaking.
- Remove the battery if you do not use the product for a long period of time.
- Wear chemical-resistant protective gloves and glasses when handling leaking batteries.

### Important information on product safety!

- Do not expose the instrument to extreme temperatures, vibrations or shocks.
- Clean it with a soft damp cloth. Do not use solvents or scouring agents.
- Protect it from moisture.

### 4. Elements:

#### LCD (Fig. 1):

- |                       |                                   |
|-----------------------|-----------------------------------|
| ① Temperature (°C/°F) | ② Maximum and minimum temperature |
| ③ Battery symbol      | ④ Freezer symbols                 |
| ⑤ Fridge symbol       |                                   |

#### Housing (Fig. 2):

- |                    |                                      |
|--------------------|--------------------------------------|
| ⑥ RESET button     | ⑦ Battery compartment                |
| ⑧ Stand (fold out) | ⑨ Frame (removable, adhesives incl.) |

### 5. Getting started

- Pull the protective foil off the display.
- Open the battery compartment using a coin and remove the insulation strip.
- The display shows the current temperature, the maximum and minimum temperature, and should the occasion arise, the corresponding freezer-fridge symbol (Fig. 3).
- The unit is now ready to use.

### 6. Operation

#### 6.1 Maximum/Minimum function

- Press RESET button to clear the recorded MAX "Λ" / MIN "V" reading.
- "-.-" appears on the display.
- The maxima and minimum values are automatically reset after 3 seconds.

#### 6.2 Temperature display

- Press and hold RESET button to change between °C (Celsius) or °F (Fahrenheit) as temperature unit.

#### 6.3 Positioning of the thermometer

- With the stand (fold out) at the back of the unit, the thermometer (without frame) can be placed onto any flat surface.

#### 6.4 Fixation of the thermometer

- You can fasten the frame on a surface or behind a glass surface.
- Please clean the surface before.
- Pull off the protection foil of both adhesive tapes.
  - a) Installation behind glass: The holder for the thermometer is facing backwards.
  - b) Installation on a surface: The holder for the thermometer is facing forward.
- Fix the frame.
- Now you can easily pull in and out the thermometer.
- **Note:** The thermometer is locked in the frame when you hear a "click" during the inserting.

### 7. Battery replacement

- Change the battery CR 2032 button cell when the battery symbol appears on the display
- Open the battery compartment using a coin, remove the battery using a pin and insert the new battery, observe correct polarity (+ pole above).

### 8. Troubleshooting

Problems	Solution
No display / Incorrect display	→ Ensure battery polarity is correct (+ pole above) → Change the battery

### 9. Waste disposal

- Never dispose of empty batteries and rechargeable batteries in household waste. As a consumer, you are legally required to take them to your retail store or to appropriate collection sites depending on national or local regulations in order to protect the environment. The symbols for the heavy metals contained are: Cd=cadmium, Hg=mercury, Pb=lead
- This instrument is labeled in accordance with the EU Waste Electrical and Electronic Equipment Directive (WEEE). Please do not dispose of this instrument in household waste. The user is obligated to take end-of-life devices to a designated collection point for the disposal of electrical and electronic equipment, in order to ensure environmentally-compatible disposal.

### 10. Technical data

Measuring range temperature:	-30 °C... +50 °C (-22 °F...+122 °F)
Accuracy:	±0,8 °C (±1,4 °F) at -20...+10 °C (-4...+50 °F); otherwise: ±1 °C (±1,6 °F)
Resolution:	±0,1 °C (±0,1 °F)
Power consumption:	1 x CR2032 button cell battery (included)
Housing dimension:	91 x 15 (35) x 72 mm
Weight:	36 g (instrument only)

### This product fulfills the guidelines according to EN 13485.

- Suitability: S, T (Stockage, Transport)  
Location: C  
Accuracy class: 1  
Measuring range: -30°C...+50°C  
In accordance with EN 13485, this instrument is subject to regular inspections as per EN 13486 (recommendation: yearly).  
TFA Dostmann GmbH & Co.KG, Zum Ottersberg 12, D-97877 Wertheim.  
No part of this manual may be reproduced without written consent of TFA Dostmann. The technical data are correct at the time of going to print and may change without prior notice.  
www.tfa-dostmann.de 08/2013

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