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Operating instructions

electronic probe thermometer with max-min-function art. no. 30.1021

With this probe thermometer you have bought an electronic precision device for measuring the current temperature as well as the maximum and minimum temperature.

measuring range: -40 °C to +200 °C (-40 °F to 392 °F)
operating temperature: -20 °C to +50 °C (-4 °F to 122 °F)
precision: ± 1°C (± 1.8°F) from -20°C to +100°C (-4°F to 212°F), ± 2°C (± 3.6°F) otherwise
resolution: 0.1 °C (0.1 °F)
battery: 1.5V button cell
Pivoting head (180°)
auto switching-off
function: after 5 minutes

Use:

- Open the battery cover with a coin in an anti-clockwise direction, remove the battery stripe and replace the cover. Remove the protection foil of the display. The thermometer is now ready for use.
- The display shows the currently measured temperature. The sensor is located in the point of the probe. The temperature unit can be changed from °C to °F by the °C/F-key.
- Press the MODE-key once in normal mode and hold-function is activated (indicated by HOLD).
- Press the MODE-key again and the maximum temperature is shown since the last reset (indicated by MAX).
- Press the MODE-key again and the minimum temperature is shown since the last reset (indicated by MIN).
- Press the MODE-key again to return to normal mode.
- To clear the maximum and minimum memory, press and hold the MODE-key for 3 seconds in the corresponding mode while MAX or MIN is displayed (indicated by ---). Press the MODE-key again to return to normal mode.

Do not dispose of empty batteries in household waste. Take them to special local collection sites.

Maintenance:

- Do not expose the instrument to extreme temperatures, vibration or shock
- Clean it with a soft damp cloth. Do not use solvents or scouring agents.
- Please do not try to repair the unit. Contact the original point of purchase. Please change the battery before complaining. No guarantee if the instrument is handled or opened improperly.

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