Digital-Refraktometer

DR301-95

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Digital Refractometer

DR301-95

Operating Manual

Inhalt / Index

DEUTSCH

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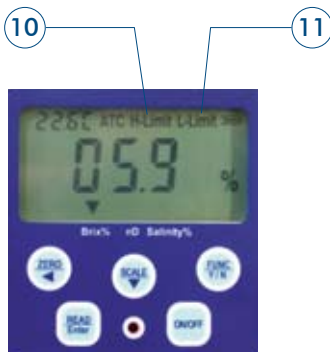
Email: service@kruess.com

1. Explanation DR301-95

1. Alarm lamp
2. Read | Enter Key
3. Zero | Tab Key
4. Scale | Setting Key
5. Function and Yes/No Key
6. ON/OFF Key
7. Prism and prism plate under lid

8. Temperature
9. Temperature compensation
10. Alarm High Limit
11. Alarm Low Limit
12. Battery
13. Measured value
14. Shows the active unit
15. Units

2. Before starting

- Inspect the prism window and prism to ensure it is completely clean and undamaged. Before every measurement clean the prism with a soft cloth and if necessary a glass cleaner.
- During the measurement keep the instrument in horizontal position.
- Calibration (see point 4)
 - At least once a day before starting measuring.
 - After moving the instrument between places with large ambient temperature difference.
 - After changing battery.

3. ON/OFF

Press ON/OFF to switch on or off DR301-95.

4. Calibration

- Put some drops of distilled water (approx. 1 ml) onto the prism and close the lid.
- The distilled water and the DR301-95 should be at the same temperature of 18–20 °C.
- Press **ZERO** for 3 seconds until **CAL** flashes on the display. The calibration is completed when **END** is displayed.

- When this symbol ⊕ is flashing, the calibration has to be restarted.

5. Measuring

- Put the sample onto the prism (approx. 1 ml).
- Close the lid.
- Press **READ | ENTER**.
- The measuring result is shown.

6. Measurement units

- Press **SCALE** to change the measurement unit (% Brix, nD or % Salinity).
- The arrow shows the current unit.

7. Functions

- By pressing **FUNC | Y/N** you can change the following functions:
 1. Alarm setting (High Limit H.L. and Low Limit L.L.)
 2. Switching on and off the temperature compensation
 3. Changing temperature unit °C or °F

To change the setting of the functions menu 1–3, always start with function menu 1 and continue with function menus 2 and 3.

To enter the functions menus start with the start display.

7.1 Function menu 1: Activation of alarm

- If you do not want to activate the alarm, press 3x **FUNC | Y/N**, to go to functions menu 2.

- To enter the alarm, press 1x **FUNC | Y/N**

- Press **READ | ENTER**
- Enter the high limit with the **SCALE** and **ZERO** button. Confirm with **READ | ENTER**

- Press **READ | ENTER**
- Enter the low limit with **SCALE** and **ZERO** button.
- Confirming with **READ | ENTER** will take you to the function menu 2

7.2 Function menu 2: Activating temperature compensation

- Now you are in the menu **Automatic Temperature Compensation ATC**.

- Press **SCALE** to switch temperature compensation on (y) or off (n).
- Confirming with **READ | ENTER** will take you to the function menu 3

7.3 Functions menu 3: Temperature units

- Now you are in the menu **temperature unit t.P**.

- By pressing **SCALE** you can change the temperature unit to °C or °F.
- Confirm with **READ | ENTER**.
- Now you are back to the start display and can continue measuring.

8. Deactivation of alarm

- The entered high and low limits can be switched off. The settings will be stored

- Start with the start display and press **FUNC | Y/N**

- Now you are in the high limit menu, which is activated. It shows **y**.
- Press **FUNC | Y/N** to deactivate the high limit. It will show **n**.
- Confirm with **READ | ENTER**.
- Now you are in the low limit menu. It shows **y**.
- press **FUNC | Y/N** to deactivate the low limit. It will show **n**.
- Confirm with 3 x **READ | ENTER** to return to the start display.

9. Replacing battery

- Replace the battery as soon as the battery voltage display (12) disappears.
- The battery case is at the bottom of the instrument.
- Open the battery case gently and replace the battery.
- Close the battery case

10. Interface

- Connect interface cable to serial port of the instrument and PC.
- Data transfer happens through the program. The instrument can only be switched on and off.

11. Specifications

	nD	% Brix	% Salinity
Measuring range	1.3330–1.5318	0–95	0–30
Accuracy	± 0.00015	± 0.1	± 0.1
Resolution	0.0001	0.1	0.1

	°C	°F
Temperature compensation	5–40	41–104
Temperature range	0–40	32–104
Temperature accuracy	± 1	± 2
Temperature resolution	0.1	0.1

Alarm Limit	A range of alarm is set. If the measured results are higher or lower than the set points, the DR301-95 indicates alarm and the red alarm light (1) flashes.
Automatic shut-off	Automatic shut off after 3 minutes
Working voltage	9 V block battery Battery-voltage display
Weight	900 g
Dimensions	180 X 100 X 60 mm

Warranty Conditions

A. KRUESS Optronic warrants that the instruments of the digital handrefractometer family will be free from defects in materials and workmanship after the date of delivery for a time period of 24 months.

During this warranty period A. KRUESS Optronic will repair or replace products which fall under the warranty conditions. For warranty repairs or service the instrument must be returned to A. KRUESS Optronic. In the case of warranty repairs A. KRUESS Optronic will bear the delivery costs. For all other deliveries the customer will bear the delivery costs. A. KRUESS Optronic warrants that the hardware which A. KRUESS Optronic has determined for use with this instrument will operate free from defects when applied according our manufacturer's instructions.

A. KRUESS Optronic does not warrant error-free and uninterrupted operation of the instrument or that this user manual is free of error. A. KRUESS Optronic will not be liable for consequential damages.

Limited Warranty:

The preceding warranty does not apply to errors and defects caused by improper handling, software not supplied by A. KRUESS Optronic, modification, abuse or by operation outside of the stated environment or by unauthorized maintenance.

Further claims are excluded and will not be accepted. A. KRUESS Optronic explicitly does not warrant the applicability or the economic use for certain applications.

A. KRUESS Optronic reserves the right to change this user manual and the technical specifications of the described instrument at any time.

This digital refractometer is only then ready for shipment when it is appropriately packed in the complete original packaging. If necessary, please order a replacement packaging from your supplier.

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A.KRÜSS Optronic GmbH
A.KRÜSS Optronik GmbH