North America

Whatman Inc.
200 Park Avenue, Suite 210
Florham Park, NJ 07932 USA
Tel.: 1-800-WHATMAN (US and Canada)
Fax: 1-973-245-8329
Email: info@whatman.com

Europe

Whatman International Ltd.
Springfield Mill, James Whatman Way
Maidstone
Kent ME 14 2LE UK
Tel.: +44 (0) 1622 676670
Fax: +44 (0) 1622 677011
Email: information@whatman.com

Whatman GmbH
Hahnstraße 3
D-37586 Dassel
Germany
Tel.: +49 (0) 5564 204 100
Fax: +49 (0) 5564 204 533
Email: information@whatman.com

Japan

Whatman Japan KK
JPR Ichigaya Building 6F
4-7-15 Kudan-Minami
Chiyoda-ku. Tokyo 102-0074 Japan
Tel.: +81 (0) 3 5215 1240
Fax. +81 (0) 3 5215 1245
Email: japaninfo@whatman.com

Asia Pacific

Whatman Asia Pacific Pte Ltd.
1 Kim Seng Promenade # 15-12
Great World City, West Tower
Singapore 237994
Tel.: +65 (0) 6534 0138
Fax: +65 (0) 6534 2166
Email: wap@whatman.com

Whatman[®]
Part of GE Healthcare

Whatman[®]
Part of GE Healthcare

Bedienungsanleitung Operating Instruction Instrucciones de uso

MV 050/0

Vakuum-Filtrationsgerät aus Edelstahl
Vacuum Filtration Apparatus made of Stainless Steel
Aparato de filtración por vacío de acero especial

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1 General

This unit can be used in combination with dangerous materials, operating procedures and apparatuses. However the Operating Instruction cannot indicate all safety problems, which may arise. It is the responsibility of the user to ensure that all safety and health related regulations are obeyed and to establish the necessary limitations prior to use.

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2 Preparing the Filtration Apparatus

Prior to the first operation the filtration unit must be cleaned (see 5).

- a. Push a fitting rubber stopper (see 8.2) onto the lower section (7) outlet pipe and insert lower section (7) into SF 100 suction flask (if the filter residue is to be examined) or a WT 100 Witt's bottle (if it is the filtrate which is required).
- b. Place PTFE sealing (6) with support sieve (5) (alternatively steel frit) onto the lower section (7).
- c. Using tweezers PZ 001, place an appropriate membrane filter on the filter support. For very cloudy fluids place a pre-filter exactly in the centre on top of the membrane filter.
- d. Fix spring-action clamp (4) to funnel (3) using the attached screws (4a). Put funnel (3) onto the lower section (7) and close filtration apparatus with spring-action clamp (4).
- e. Plug opening in lid (1) with cotton wool to prevent airborne micro organisms from entering during filtration.

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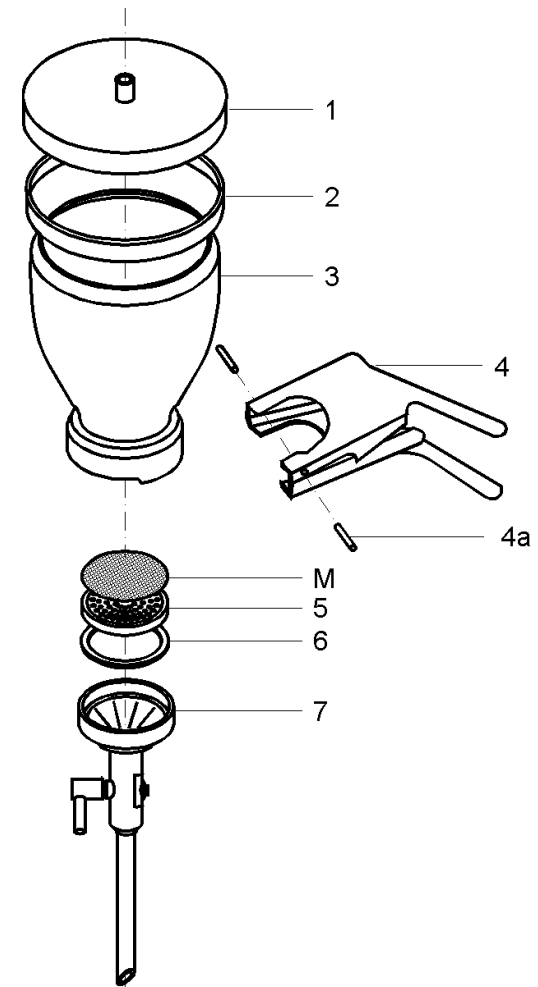
Pre-filter for MV 050/0

Material	Format mm	Type	Thick- ness µm	Ref. No.
Glass fibre	37	GF/A	260	1820-037

3 Filtration

For sterile filtration please refer to point 4.

- Using a vacuum tubing (SV 006), connect the assembled apparatus with filter in place - the whole sterilized beforehand where necessary - via a Wouffe bottle to a water jet pump or an electric pump (VP 003).
- Close tap and switch on pump.
- Lift the lid (1) slightly on one side, pour in the liquid to be filtered, shut the lid (1) and open the tap. If it is necessary to add more liquid this must be done before the funnel (3) is empty.
- When carrying out bacteriological control tests, thoroughly rinse the funnel (3) from top to bottom with sterile water in order to bring all micro-organisms to the surface of the membrane filter.

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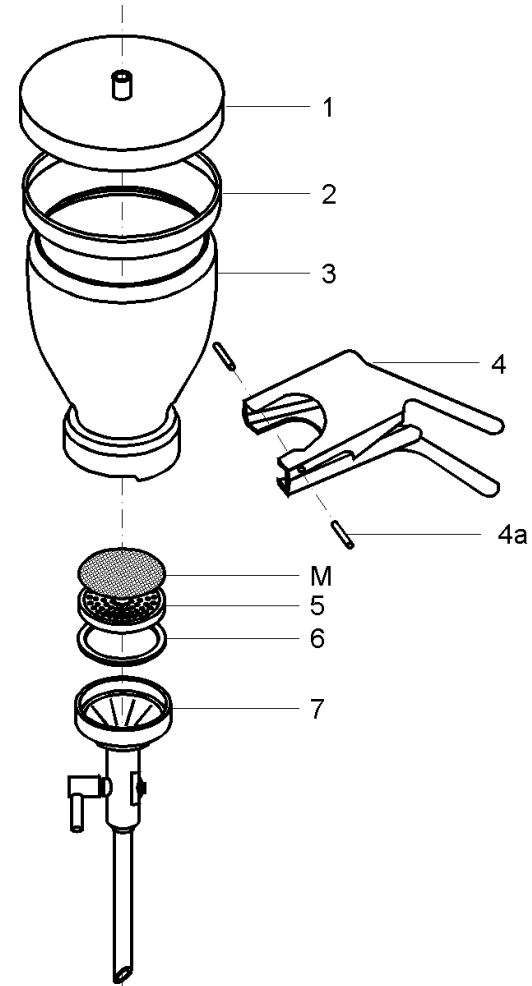
8 Spare Parts List/Accessories

8.1 Spare Parts

Pos.	Designation	Ref. No.
1	Cover with silicone sealing	10440005
2	Silicone sealing	10440006
3	Funnel (500 ml)	10440001
4	Clamp	10440004
4a	Screw	10440007
5	Support sieve	10440003
6	PTFE sealing	10440102
7	Lower part (for rubber stopper)	-

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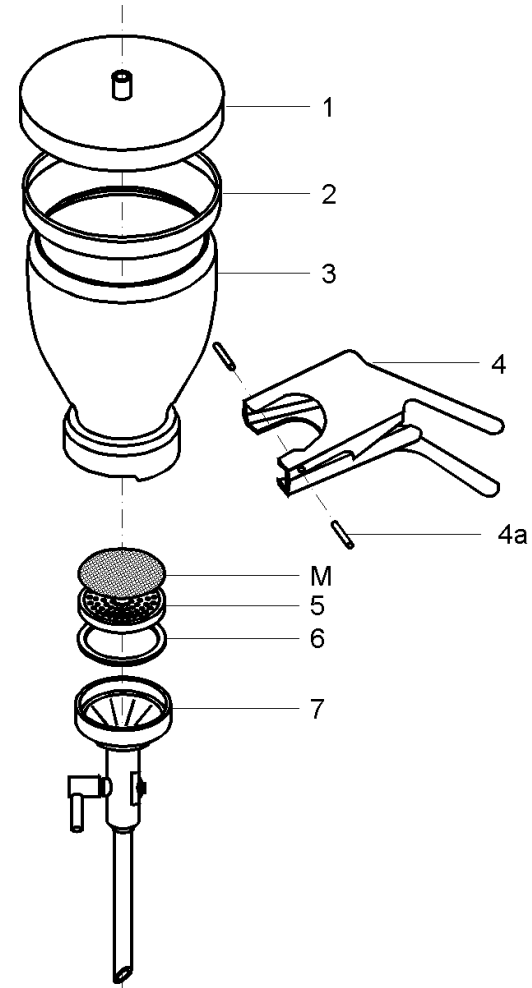
4 Sterilisation

When using the apparatus for sterile filtration, we recommend sterilization in an autoclave.

- Prepare filtration apparatus as described under point 2.
- Wrap equipment in sterilization paper and hold in place with adhesive fabric tape or rubber bands.
- After heating, hold the autoclave for 15 - 20 min. with an atmospheric pressure of 1.1 bar at a temperature of 121°C.

Temperature above 121°C can cause damage to membrane filters produced from cellulose nitrate or cellulose mixed ester.

The water or steam must not contain any additives, because their action can make the membrane filters susceptible to tearing.

- Afterwards, gradually vent the autoclave or, even better, allow the pressure to equalize by cooling down. Too rapid release of steam causes pressure shocks which may damage the filter.
- Allow filtration apparatus to cool down to "lukewarm", if possible in the autoclave.
- After sterilisation in the autoclave check for damages and replace damaged parts if necessary.

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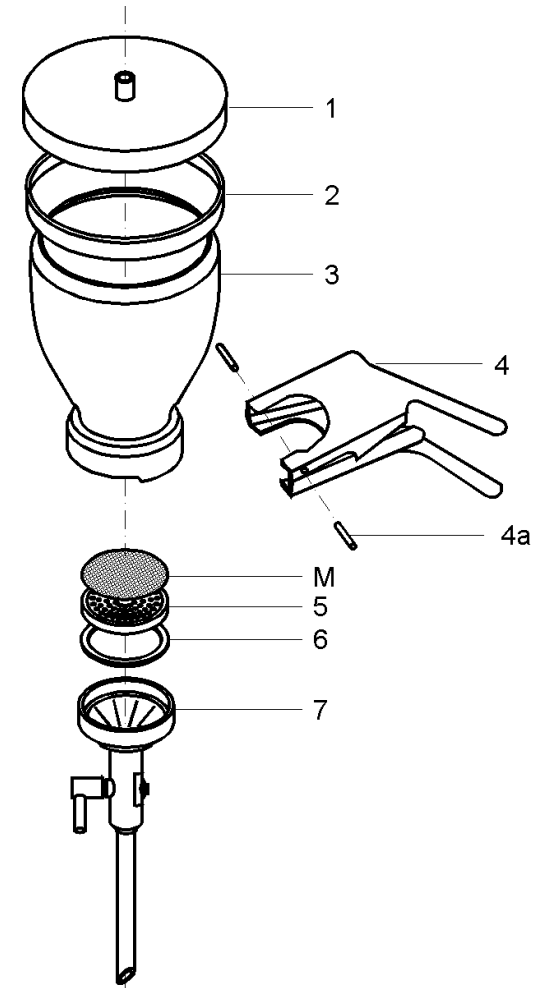
6 Technical Data

Materials		
Upper/Lower part	Stainless Steel	
Support sieve	Stainless Steel	
Sealing	PTFE and silicone	
Clamp	Aluminium	
Funnel volume	500 ml	
Filter size	50 mm	
Filter area	12,5 cm ²	
Pre-filter size	max. 40 mm	
Dimensions		
Height	320 mm	
Diameter	110 mm *	
	* without clamp	

7 Resistance

As the surface finish is important for the resistance of stainless steels to corrosion, filtration apparatus have electrolytically polished surfaces. Therefore, careful handling is necessary to avoid damaging the surface.

There is, however, no absolute protection against corrosion. Even a very low HCl content in the room air can soon lead to rusting. Solutions containing e.g. chloride or hypochlorite also damage the steel surface.

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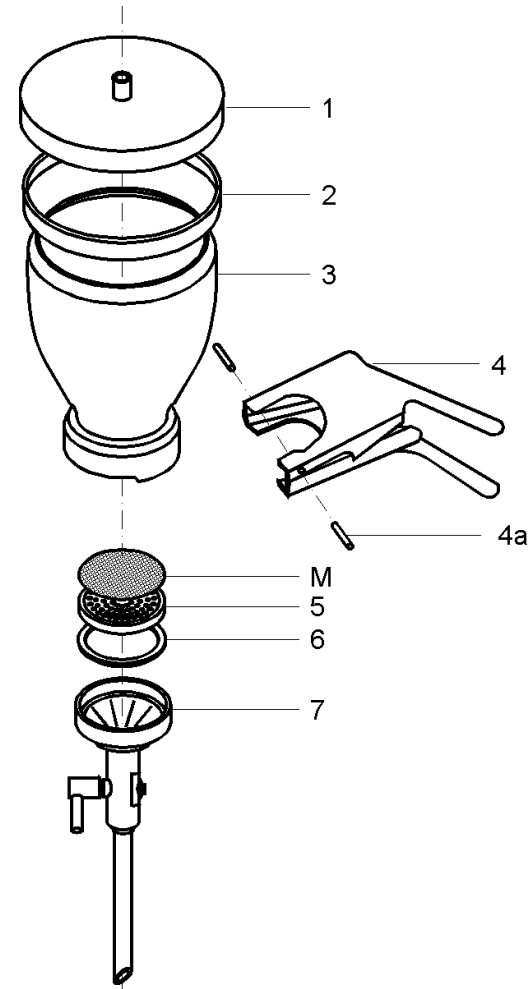
5 Cleaning

 **The equipment must be carefully cleaned before being used for the first time and always after every filtration.**

- Dismantle the filtration apparatus into its separate parts.
- Check surfaces, PTFE sealing (6) and support sieve (5) (alternatively steel frit) regularly for damage. Replace damaged parts.
- Rinse each part in hot water (if necessary, with a soft brush and a small amount of detergent).
- Any residue left in the pores of the steel frit (accessory) after filtration will require removal by chemical means. The solvent chosen will depend on the type of deposit to be removed.
- Also rinse the other parts several times with clean water. The final rinse must be carried out with distilled water, particularly before carrying out critical filtrations.
- Air-dry the parts or put them in a drying oven at 105 °C. To avoid any fibres or particles contaminating the filtrate, cloths may not be used for drying any parts.

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