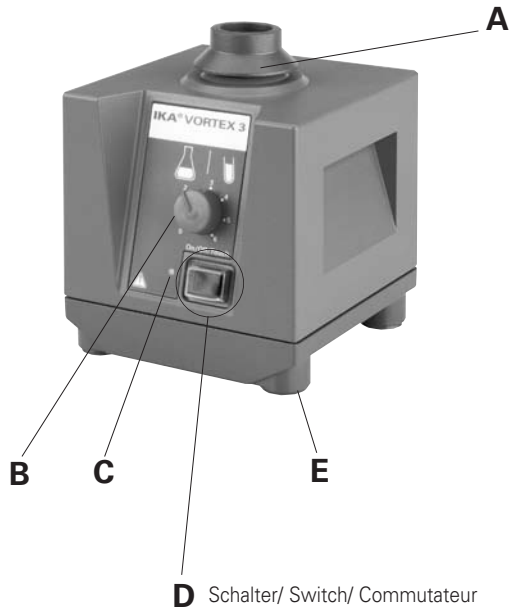
# IKA® Vortex 3

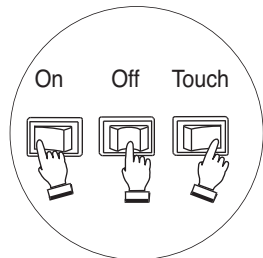
Betriebsanleitung	DE	3
Operating instructions	EN	8
Mode d'emploi	FR	13
Indicaciones de seguridad	ES	18
Veiligheidsinstructies	NL	18
Norme di sicurezza	IT	19
Säkerhetsanvisningar	SV	20
Sikkerhedshenvisninger	DA	21
Sikkerhedshenvisninger	NO	22
Turvallisuusohjeet	FI	22
Instruções de serviço	PT	23
Wskazówki bezpieczeństwa	PL	24
Bezpečnostní upozornění	CS	25
Biztonsági utasítások	HU	25
Varnostna navodila	SL	26
Bezpečnostné pokyny	SK	27
Ohutusjuhised	ET	28
Drošības norādes	LV	28
Saugos reikalavimai	LT	29

**D** Schalter/ Switch/ Commutateur

**Fig. 1**

Pos. Bezeichnung	Item Designation	Pos. Désignation
A Standardaufsatz VG 3.1	A Standard attachment VG 3.1	A Support standard VG 3.1
B Drehknopf	B Rotating knob	B Bouton rotatif
C Kontrollleuchte	C Control light	C Lumière cōntrol
D Geräteschalter	D Switch	D Commutateur
E Gerätefuß	E Base	E Pied

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**DE**

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	Page
CE-Declaration of conformity	2
Safety instructions	8
Unpack	9
Correct use	9
Commissioning	9
Accessories	10
Maintenance	11
Warranty	11
Technical data	12

## Safety instructions

### For your protection

• **Read the operating instructions in full before starting up and follow the safety instructions.**

- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Socket must be earthed (protective ground contact).
- Wear your personal protective equipment in accordance with the hazard category of the medium to be processed. Otherwise there is a risk of:
  - splashing liquids
  - projectile parts
  - body parts, hair, clothing and jewellery getting caught.
- Keep hand/arm vibration to a minimum (Directive 2002/44/EC) when operated manually:
  - Keep the sample container as vertical as possible.
  - Only exert the minimum pressure necessary for the mixing process.
  - Use attachments for multiple samples when there is a large

number of samples.

- Set up the appliance in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- The feet of the appliance must be clean and undamaged.
- Position the knob (B) at the left stop before starting up. Gradually increase the speed.
- If the shaking movement of the appliance increases (resonance), reduce the speed or pass through the critical phase as quickly as possible.
- Reduce the speed if
  - the medium splashes out of the vessel because the speed is too high
  - the appliance is not running smoothly
  - the appliance begins to move around because of dynamic forces.
- Firmly secure the accessories and vessels in place, otherwise shaking vessels could be damaged or projected out.
- Check the appliance and accessories beforehand for damage each time you use them. Do not use damaged components.
- Position one single shaking vessel in the centre and several shaking vessels so that they are evenly spread out.
- The appliance is not suitable for manual operation.
- Sharp edged vessels cause wear debris on the attachments.
- Beware of the risk of
  - flammable materials,
  - glass breakage as a result of mechanical shaking power.
- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
- **Do not** operate the appliance in explosive atmospheres, with hazardous substances or under water.
- Safe operation is only guaranteed with the accessories described in the "Accessories" chapter.
- Always disconnect the plug before fitting accessories.
- The appliance starts up again automatically following a cut in the power supply.
- The appliance may heat up when in use.

### For protection of the equipment

- The voltage stated on the nameplate must correspond to the

- mains voltage.
- Protect the appliance and accessories from bumps and impacts.
- The appliance may only be opened by experts.

## Unpack

- **Unpack**
  - Please unpack the device carefully
  - In the case of any damage a fact report must be set immediately (post, rail or forwarder)
- **Delivery scope**
  - Shaker **Vortex Genius 3**
  - Standard attachment VG 3.1
  - Operating instructions

## Correct use

- **Use**
  - ▶ For mixing liquids
  - Touch mode for single test tubes
  - Continuous mode for single or several vessels
- **Range of use**
  - Laboratories
  - Pharmacies
  - Schools

This device is suitable for use in all areas except:

- Residential areas
- Areas that are connected directly to a low-voltage supply network that also supplies residential areas.

The safety of the user cannot be guaranteed if the appliance is operated with accessories that are not supplied or recommended by the manufacturer or if the appliance is operated improperly contrary to the manufacturer's specifications.

## Commissioning

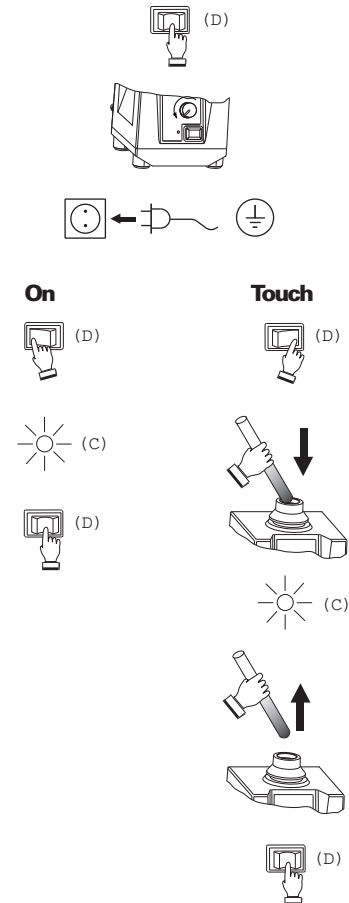
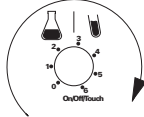
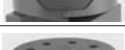
Fig. 2

## Accessories

Possible uses and permitted speed ranges of attachments:

Designation	Description	Id.- No.	Touch Mode	Continuous Mode	Range of speed
 VG 3.1 Standard attachment	For test tubes and small vessels upto diameter 30 mm	33 412 00	X	X	
 VG 3.2 One-hand attachment	88 mm-plate attachment with rubber pad	33 423 00	X	X	
 VG 3.3 Universal attachment	150 mm-plate attachment with rubber pad	33 424 00	-	X	
 VG 3.31 Test tube insert	For 54 Eppendorf tubes for VG 3.3	33 443 00	-	X	
 VG 3.32 Test tube insert	For 18 test tubes 10 mm, for VG 3.3	33 439 00	-	X	
 VG 3.33 Test tube insert	For 12 test tubes 12 mm, for VG 3.3	33 440 00	-	X	
 VG 3.34 Test tube insert	For 8 test tubes 16 mm, for VG 3.3	33 441 00	-	X	
 VG 3.35 Test tube insert	For 8 test tubes 20 mm, for VG 3.3	33 442 00	-	X	
 VG 3.36 Erlenmeyer flask attachment	For 1 Erlenmeyer/ round-bottomed flask 100-250 ml, for VG 3.3	33 421 00	-	X	
 VG 3.37 Microtiter plate insert	For 1 microtiter plate, für VG 3.3	33 444 00	-	X	

### Changing attachments

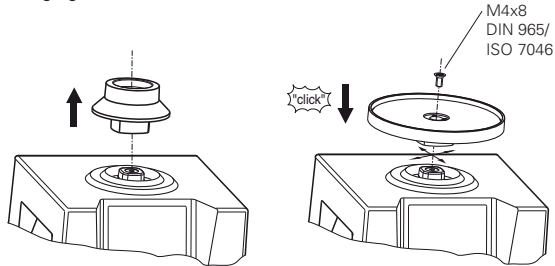
Fig. 3

### Using inserts

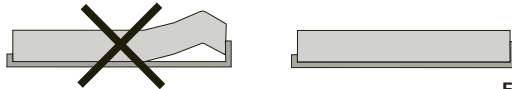
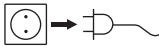
Fig. 4

## Maintenance

The appliance is maintenance-free.

### Cleaning

Only use cleansing agents which have been recommended by **IKA**

Use to remove:

Dyes	isopropyl alcohol
Construction materials	water containing tenside/ isopropyl alcohol
Cosmetics	water containing tenside/ isopropyl alcohol
Foodstuffs	water containing tenside
Fuels	water containing tenside

- Do not allow moisture to get into the appliance when cleaning
- Wear protective gloves during cleaning the devices.
- Before using another than the recommended method for cleaning or decontamination, the user must ascertain with **IKA** that this method does not destroy the instrument.
- The rubber of the attachments' pads (VG 3.2 and VG 3.3) is a natural product, so talcum can discharge. Clean the rubber pad with glycerin.

### Spare parts order

When ordering spare parts, please give:

- Machine type
- Manufacturing number, see type plate
- Item and designation of the spare part, see [www.ika.de](http://www.ika.de), spare parts diagram and spare parts list

### Repair

In case of repair the device has to be cleaned and free from any materials which may constitute a health hazard.

If you require servicing, return the appliance in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

## Warranty

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

## Technical data

Operating voltage	VAC	100 – 120 ± 10%
Design frequency	Hz	50/60
Nominal voltage	V/Hz	115/60
Input power	W	60
Output power	W	10
Speed	1/min	0/ 500 – 2500 infinitely adjustable
Speed adjustment		analog (infinitely) rotating knob in front of device
Speed display		Scale 0 – 6
Drive		Shaded pole motor
Agitation stroke	mm	4
Shaking motion		horizontal approximately- circular
Perm duration of operation	%	100
Perm ambient temperature	°C	+5 bis +40
Perm. relative humidity	%	80
Protection type acc. to DIN EN 60529		IP 21
Protection class		I (protection earth)
Overvoltage category		II
Contamination level		2
Operation at a terrestrial altitude	m	max. 2000
Dimensions total W x D x H	mm	127 x 149 x 136
Weight without supported load	kg	4.5
Max. supportes load incl. attachment	kg	0.5

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