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English

Operating Instructions Safety Showers and Eye Washes

1. Attention! Important Note

The water pipelines have to be rinsed before connecting the shower to avoid damage to the valve caused by dirt accumulation where new installations are concerned. Claims for defects that result from water pipes that had not been rinsed prior to connection cannot be honoured.

2. Introduction

Safety showers and eye washes are important first aid installations. Burns, scalds and alkali burns are among the injuries for which immediate first aid is decisive to reduce damage. The quicker one of these injuries is treated with cold water, the more likely it is that damage to the deeper tissue is avoided.

3. Standarts and Directives

All BREUELL Safety showers and eye washes comply with all the points of DIN EN 15154-1:2006, the DIN EN 15154-2:2006, the ANSI Z 358.1 - 1998, the TRGS 526 standards and the directives of the chemical employers' liability insurance BGR 120. These operating instructions are to be considered a supplement to the applicable directives and regulations. Please also heed the above named standards and directives.

4. Operating Pressure

The flow pressure must be at least 1 bar to guarantee perfect operation. All B-SAFETY safety showers and eye washes have been designed for a maximum operating pressure of 10 bar (PN10).

Hand eye washes with rose head and eye washes:

Volume flow is 7 to 9 litres/minute per rose head. This volume flow required to comply with the standard is automatically set, no matter what the operating pressure is, by a flow limiting device integrated in the shower.

Hand eye washes with two rose heads:

The flow pressure must be between 1.8 and 3.5 bar to guarantee perfect operation. At this flow pressure the water jet is between 15 and 30 cm high (measured at an inclination angle of 10° according to DIN EN 15154-2:2006). If the pressure applied is higher than 3.5 bar, we recommend the additional installation of the ball valve included in the supplies. The jet spray height should be set at 15 to 30 cm according to DIN EN 15154-2:2006.

Showers:

The volume flow is approx. 80 litres/minute at 1 bar flow pressure.

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5. Location

The location is chosen to fulfil the directives of the chemical employers' liability insurance BGR 120. Free access is to be provided at all times. Particular care must be taken that a location is chosen that can be reached quickly and accessed from the danger zones. The injured person must be able to reach the shower facilities within a maximum of 10 seconds, they must therefore be installed at a maximum of 10 to 15 metres from the workplace.

6. Water Supply

The B-SAFETY safety showers and eye washes should be connected to clear, drinking water supply lines. The drinking water lines must comply with DIN 1988. Cold water has an anaesthesiological effect and achieves the quickest results in cases of alkali burns. Besides, the problem that may occur when heating the water (too hot water, germ occurrence) is avoided.

Hand eye washes:

The B-SAFETY hand eye washes are fitted with a return flow inhibitor acc. to DIN 1988 (protection class 2). In addition, a further return flow inhibitor can be obtained from specialised trade to achieve a higher protection class.

7. Installation

To avoid damage to furniture and floors the shower should be installed in the drain area enabling the water from the shower to flow off into the canal. The attached safety signs are to be applied to the location at which the B-SAFETY safety shower and eye wash is installed.

8. Operation

The B-SAFETY safety showers and eye washes are started-up by actuating the red or green handle for the hand eye washes and compact eye washes, by actuating the large push lever for the eye washes and by pulling down the large pull bar for the showers. The trigger mechanism locks in position each time and must then be released again.

9. Maintenance

The operation of the B-SAFETY safety showers and eye washes has to be checked at least once a month as is also laid down in the directives of the chemical employers' liability insurance BGR 120. The shower has to be actuated for approx. 10 seconds for this test. The purpose is not only to check operability, but the available water in the shower is replaced by fresh water (protection against germs). In the scope of this function check, the shower is visually checked for wear and tear, possible leakage and soilage such as for instance dust or scale deposits. Defective parts must be replaced immediately.

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