

**C-MAG HS4**    **C-MAG HS7**    **C-MAG HS10**  
**C-MAG HP4**    **C-MAG HP7**    **C-MAG HP10**  
**C-MAG MS4**    **C-MAG MS4**    **C-MAG MS10**

OPERATING INSTRUCTIONS	EN	3
BETRIEBSANLEITUNG	DE	10
MODE D'EMPLOI	FR	17
INDICACIONES DE SEGURIDAD	ES	24

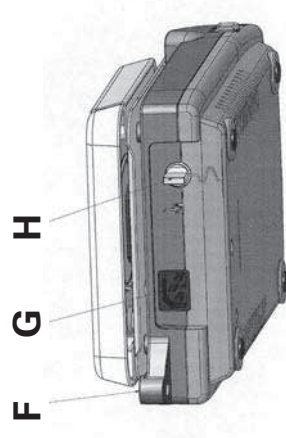Ref. No. 4380.01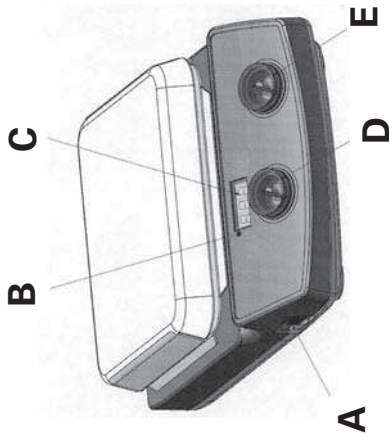

Item Designation	Pos. Bezeichnung	Pos. Désignation
A Switch	A Geräteschalter	A Commutateur
B LED	B LED	B DEL
C Control light	C Display	C Affichage
D Rotating knob	D Bedienknopf	D Bouton rotatif
E Rotating knob	E Bedienknopf	E Bouton rotatif
F Threaded support bore	F Stativgewindebohrung	F Alésage fileté du statif
G Mains socket	G Netzbuchse	G Prise secteur
H Contact-thermometer-jack	H Kontakt-thermometer-Buchse	H douille de thermomètre de contact

**CE-DECLARATION OF CONFORMITY**

**EN**

We declare under our sole responsibility that this product corresponds to the regulations 73/023/EG and 89/336/EG and conforms with the standards or standardized documents EN 61010-1 and EN 61326-1.

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Fig. 1

## Contents

	Page
CE - Declaration of conformity	2
Safety instructions	3
Unpack	4
Correct use	4
Commissioning	5
Error codes	6
Accessories	6
Maintenance	6
Warranty	7
Technical data	8
List of spare parts	9
Spare parts diagram	26/27

hazard category of the medium to be processed. Otherwise there is a risk of:

- splashing liquids
- projectile parts
- release any toxic or combustible gases.
- Set up the appliance in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- The feet of the appliance must be clean and undamaged.
- Position the knob at the left stop before starting up. Gradually increase the speed.
- Reduce the speed if
  - the medium splashes out of the vessel because the speed is too high
  - the appliance is not running smoothly
  - the container moves on the set-up surface.
- Never heat up any medium with fire point higher (min. 50°C) than the adjusted temperature
- When using PTFE-coated magnetic bars, the following have to be noted: *Chemical reactions of PTFE occur in contact with molten or dissolved alkaline and alkaline - earth metals, as well as with fine-particled powders of metals of the 2. and 3. group of the periodical system at temperatures above 300-400°C. Only elementary fluorine, chlorine trifluoride and alkaline metals do attack PTFE, halogen hydrocarbons have a reversibly swelling effect.*
- Source: *Römpfs Chemie-Lexikon and „Ullmann“ Bd.19*
- Check the appliance and accessories beforehand for damage each time you use them. Do not use damaged components.
- Only replace damaged parts with spare parts identical to the original in function and quality.
- Do not use the device if the ceramic set-up surface is damaged e.g. scratches, splinters or corrosion. A damaged set-up surface could break if used.
- Beware of the risk of
  - flammable materials
  - glass breakage as a result of mechanical shaking power
  - incorrect container size
  - too much medium

## Safety instructions


### To your protection

- **Read the operating instructions in full before starting up and follow the safety instructions.**
- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the appliance.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- Socket must be earthed (protective ground contact).
- **Attention - Magnetism!** Effects of the magnetic field have to be taken into account (e.g. data card, carriers, pacemakers...).
- **Risk of burns!** The heating plate can reach temperatures in excess of 500 °C. Pay attention the residual heat after switching off..  
Please make sure that the mains cable does not contact the heating plate
- Wear your personal protective equipment in accordance with the

- unsafe condition of container
- Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.
- **Do not** operate the appliance in explosive atmospheres, with hazardous substances or under water.
- A separation from the line is made with the equipment only by pulling net and/or device plug.
- Safe operation is only guaranteed with the accessories described in the "Accessories" chapter.
- Always disconnect the plug before fitting accessories.
- Accessories must be securely attached to the device and cannot come off by themselves. The centre of gravity of the assembly must lie within the set-up surface.
- The appliance starts up again automatically following a cut in the power supply.
- The appliance may heat up when in use.
- Abrasion of the dispersion equipment or the rotating accessories can get into the medium you are working on.

**To the protection of the equipment:**

- The voltage stated on the nameplate must correspond to the mains voltage.
- Do not cover the device, even partially e.g. with metallic plates or film. This results in overheating.
- Protect the appliance and accessories from bumps and impacts.

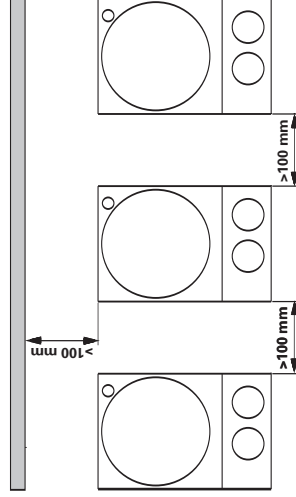- Observe the minimum distances between devices, between the device and the wall and above the assembly (min. 800mm)

**Unpack**

- **Unpack**
  - Please unpack the device carefully
  - In the case of any damage a fact report must be set immediately (post, rail or forwarder)

• **Delivery scope**

- Heating magnetic stirrer or Magnetic stirrer or Heating device
- Mains cable
- Operating instructions

**Correct use**

- **Use**
  - For mixing and/or heating liquids
- **Range of use**
  - Laboratories
  - Chemical industry
  - Schools
  - Pharmacies

## Commissioning

	MS 4	MS 7	MS 10	HS 4	HS 7	HS 10	HP 4	HP 7	HP 10
<b>Commissioning</b>	Put device switch (A) in the OFF position								
	Plug in (G) mains cable								
	Once connected to the power supply the device is in "stand-by" mode The right decimal point on the display (C) is lit								
<b>Stirring</b>	Put device switch (A) in the ON position								
	Any set values are retained when device is switched off and even after the device is disconnected from the power!								
	Set the engine speed with the operating button (E) on the right								
<b>Heating</b>	Put device switch (A) in the ON position								
	Any set values are retained when device is switched off and even after the device is disconnected from the power! Set the target temperature for the heating plate using operating button (D) [for VHP (E)]								
	The set value is indicated on the display (C). If energy is being supplied to the heating plate, the red LED (B) lights up								
	In the stir and stand-by modes, "hot" flashes on the display after the heat has been switched off for as long as the temperature of the set-up surface exceeds 50°C								
<b>Contact-thermometer connection</b>	Put device switch (A) in the OFF position, unplug contact plug								
	Safety contact thermometer acc. to DIN 12878 class 2 connected with jack (H) Put device switch (A) in the ON position								
	Beware the instruction manual of the contact thermometer Important: Display (C) also indicates the target temperature of the heating plate when the contact thermometer is connected								

## Error codes

Footstep during the enterprise a disturbance up, is indicated these with the devices to **HS** and **HP** by an error message in the display (C). Whenever an error message appears switch the device off to cool down.

Error code	Cause	Solution
E1	Inner temperature too high	<ul style="list-style-type: none"> <li>switch off the unit and allow it cool down</li> </ul>
E6	Motor seized	<ul style="list-style-type: none"> <li>Stirring bars ungeeignet</li> <li>Reduce the viscosity</li> <li>Foreign bodies in the equipment remove</li> </ul>
E9	Safety circuit ETC	<ul style="list-style-type: none"> <li>Plug the contact plug (G)</li> <li>Plug the contact thermometer</li> <li>Connecting cable, plug or contact thermometer defect, exchange</li> </ul>

Have the device repaired if the error is not corrected using the measures described or if another error code is displayed.

## Accessories

### Stirring organs

stirring bars:  
 Ø 6 mm, lengths to 15mm  
 Ø 8 mm, lengths to 50mm  
 Ø10mm, length to 80mm


### Any other accessories

RSE stirring bar remover  
 H15 bath top  
 H28 bath top  
 H36 holding rod  
 H16V support rod

H44 cross sleeve  
 ETS-D4 electrical contact thermometer

### Installing the support rod

The support rod is attached using the threaded support bore (F).

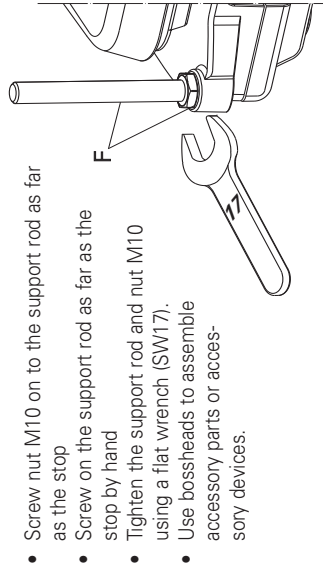- Screw nut M10 on to the support rod as far as the stop
- Screw on the support rod as far as the stop by hand
- Tighten the support rod and nut M10 using a flat wrench (SW17).
- Use bossheads to assemble accessory parts or accessory devices.

Fig. 3

## Maintenance

The appliance is maintenance-free.

### Cleaning

Only use cleansing agents which have been recommended by **IKA**.

Use to remove:

Dyes	isopropyl alcohol
Construction materials	water containing tenside/ isopropyl alcohol
Cosmetics	water containing tenside/ isopropyl alcohol
Foodstuffs	water containing tenside
Fuels	water containing tenside

- Do not allow moisture to get into the appliance when cleaning
- Wear protective gloves during cleaning the devices.
- Before using another than the recommended method for cleaning or decontamination, the user must ascertain with **IKA** that this method does not destroy the instrument.
- The rubber of the attachments' pads (VG 3.2 and VG 3.3) is a natural product, so talcum can discharge. Clean the rubber pad with glycerin.

#### Spare parts order

When ordering spare parts, please give:

- Machine type
- Manufacturing number, see type plate
- Item and designation of the spare part, see spare parts list and diagram

#### Repair

In case of repair the device has to be cleaned and free from any materials which may constitute a health hazard.

If you require servicing, return the appliance in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

### **Warranty**

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

## Technical data

	MS 4	HS 4	HP 4	MS 7	HS 7	HP 7	MS 10	HS 10	HP10
Operating voltage	VAC								
	VAC				230 ±10%				
	VAC				120 ±10%				
	VAC				100 ±10%				
Nominal voltage	VAC				230 / 50Hz				
	VAC				120/ 60 Hz				
	VAC				100/ 60 Hz				
Design frequency	Hz				50/60				
Input power max. at 230 and 120 VAC	W	30	270	255	30	1020	1005	30	1520
100 VAC		30	270	255	30	1020	1005	30	1070
Power consumption in stand-by mode	W				2.5				
Perm. duration of operation	%				100				
Perm. ambient temperature	°C				+5 bis +40				
Perm. relative humidity	%				80				
Protection type acc. to DIN EN 60529					IP 21				
Protection class					I				
Overvoltage categorie					II				
Contamination level					2				
Operation at a terrestrial altitude	m				max. 2000				
Dimensions (W x D x H)	mm	150 x 260 x 105			220 x 335 x 105			300 x 415 x 105	
Weight	kg	3			5			6	
<b>Motor</b>									
Speed range (indefinitely)	1/min	100-1500	-	100-1500	100-1500	-	100-1500	100-1500	-
Speed display	Scale	Scale	-	Scale	Scale	-	Scale	Scale	-
Power consumption	W	15	15	15	15	15	15	15	15
Power output	W	1,5	1,5	1,5	1,5	1,5	1,5	1,5	1,5
Max. stirring quantity (water)	ltr	5	5	10	10	10	15	15	15
<b>Heating plate</b>									
Heating plate dimension	mm	120x120	120x120	200x200	200x200	200x200	280x280	280x280	280x280
Heating power at 230 and 120 VAC	W	-	250	250	-	1000	1000	1500	1500
at 100 VAC		-	250	250	-	1000	1000	1050	1050
Surface temperature min.	°C	-	50	50	-	50	50	50	50
Surface temperature max.	°C	-	500	500	-	500	500	500	500
Temperature fluctuation	°C	-	±5	±5	-	±5	±5	±5	±5
Limit of safety temperature	°C	-	550	550	-	550	550	550	550

## List of spare parts

Pos.	Designation	MS-C4/C7/C10	HP-C4/C7/C10	HS-C4/C7/C10
7	Conter sunk screw	X	X	X
8	Connect distributor	X	X	X
12	Insulating		X	X
13	Danger sign		X	X
17	Distance holder LP	X	X	X
18	Clip-Triac fastening	X	X	X
19	Angel connector	X	X	X
20	Cover	X	X	X
21	Mounting plug	X	X	X
31	Placing plate	X	X	X
32	Heating foil		X	X
33	Insulating		X	X
35	Ceramic bushing		X	X
37	Holding bracket	X	X	X
38	Protection sheet	X	X	X
40	Distance bushing	X	X	X
41	Distance bushing	X	X	X
45	Mylar-foil	X	X	X
46	Color filter foil		X	X
47	Front foil	X	X	X
51	Control knob	X	X	X
53	Contact plug		X	X
54	Magnet	X		X
59	Shaded pole asynchron motor	X		X
65	Device pedestal	X	X	X
66	Impulse disk	X		X
70	Distance bushing	X		X
71	Magnett holder	X		X
75	Temperature sensor		X	X
76	Pressure sheet	X	X	X
80	Contact thermometer bushing		X	X
80	Cap (Contact-thermometer)	X		
81	Connecting cable	X	X	X
100	Label	X	X	X
2001	PCB compl.	X	X	X
2002	PCB accessories consisting of: Poti (motor); Poti (heating); Opto- coupler and LED-display	X	X	X

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**IKAWorks, Inc.**

LABORATORY TECHNOLOGY  
ANALYZING TECHNOLOGY  
PROCESSING EQUIPMENT

**IKAWorks (Asia) Sdn Bhd**

LABORATORY TECHNOLOGY  
ANALYZING TECHNOLOGY  
PROCESSING EQUIPMENT

Europe - Middle East - Africa

**IKAWERKE GMBH & CO.KG**

Janke & Kunkel-Str. 10  
D-79219 Staufen  
Germany  
TEL. +49 7633 831-0  
FAX +49 7633 831-98  
E-mail: [sales@ika.de](mailto:sales@ika.de)  
<http://www.ika.net>

North America

**IKAWorks, Inc.**

2635 North Chase Pkwy SE  
Wilmington, NC 28405-7419  
USA  
TEL. +1 800 733-3037  
TEL. +1 910 452-7059  
FAX +1 910 452-7693  
E-mail: [usa@ika.net](mailto:usa@ika.net)

Asia - Australia

**IKAWorks (Asia) Sdn Bhd**

No. 17 & 19, Jalan PJU 3/50  
Sunway Damansara Technology Park  
47810 Petaling Jaya  
Selangor, Malaysia  
TEL. +60 3 7804-3322  
FAX +60 3 7804-8940  
E-mail: [sales@ika.com.my](mailto:sales@ika.com.my)

**IKAJapan Y.K.**

LABORATORY TECHNOLOGY  
ANALYZING TECHNOLOGY  
PROCESSING EQUIPMENT

Japan

**IKAJapan Y.K.**  
293-1 Kobayashi-cho  
Yamato Koriyama Shi  
639-1026 Japan  
TEL. +81 74358-4611  
FAX +81 74358-4612  
E-mail: [japan@ika.de](mailto:japan@ika.de)

**IKAWorks Guangzhou**

LABORATORY TECHNOLOGY  
ANALYZING TECHNOLOGY  
PROCESSING EQUIPMENT

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## CE-DECLARATION OF CONFORMITY

**EN**  
We declare under our sole responsibility that this product corresponds to the regulations 89/336/EG and 73/23/EG and conforms with the standards or standardized documents DIN EN IEC 61010-1 and DIN EN IEC 61326-1.

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## Warranty

You have purchased an original IKA laboratory machine which meets the highest engineering and quality standards.

In accordance with IKA guarantee conditions, the guarantee period is 24 months. For claims under the guarantee please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

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In order to be able to use the appliance properly and safely, every user must first read the operating instructions and observe the safety instructions contained therein. Take care of these operating instructions and keep them in a place where they can be accessed by everyone.

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## Contents

	Page
Safety instructions	9
Correct use	10
Unpacking	10
Factory settings	10
Installation	10
Operating modes	11
Setting the HI-TEMP	11
Adjusting the set-point temperature	11
Replacing extension cable and sensor	11
Error messages	12
Accessories	12
Technical data	12

## Safety instructions

**Caution!** The supply voltage (Limited Power Supply) for the **ETS-D5** must be provided using an IEC 61558 compliant transformer with double or reinforced insulation. Alternatively, laboratory equipment that meets the requirements of IEC 61010 or NEC Class 2 may be used.

Do not touch the temperature sensor while measurements are being taken. This will prevent incorrect results.

Always use the extension cable H 70 when the media being processed produces vapour. This ensures that the control unit does not come into contact with the vapour.

**Note:** This equipment must only be used with media that have a flash point higher than the temperature safety limit set for the magnet agitator and/or heating system used.

The equipment must not be operated in areas where there is a risk of explosion.

Only use glass encapsulated temperature sensors for electrolysis

procedures.

The stainless steel temperature sensor must not be used with aggressive media such as acids, caustic solutions or distilled water, due to the risk of corrosion. The H66 glass sensor should be used in such cases.

Please ensure that the spiral cable does not come into contact with the heating plate.

Take care when touching the temperature sensor!

**Burn hazard!** The temperature sensor may heat up to 450°C when inserted in the media!

Please observe all accident prevention regulations applicable to the activity and the work station. Personal protective equipment should be worn when operating the equipment.

### Danger:

The user must check that all conditions have been met for the safe and correct operation of the equipment before operating the IKA-TRON® ETS-D5 with a heater that does have a DIN 12878 compliant connector, and therefore does not provide cut-out protection.

(The cut-out protection is a mechanism that ensures that the electrical heating circuit is switched off should the contact thermometer (ETS-D5) be broken or if the connector is removed.).

### H+P magnetic agitator systems (no break fuse protection included)

When the supply current for the ETS-D5 is interrupted, the electronics for the heatable magnetic agitator system will take over the temperature control. This means that the magnetic agitator system will heat up to the temperature set (e.g. 300 °C). Readjust the set-point temperature for the magnetic agitator system or switch the heating off.

**⚠ Danger when using flammable media with a low boiling temperature.**

### Corning Magnetic Agitator PC-400, PC-420, PC-600, PC-620

When using the ETS-D5 in conjunction with a Corning Magnetic Agitator it is absolutely essential that the rotary "heat" dial on the magnetic agitator is set to the "O" (off) position. The ETS-D5 will

If it  
pos  
re s  
**Wa**

only control the set temperature when it is in this position.  
**Warning:** The heating plate temperature limiter is not active in this operating mode.


If the rotary "heat" dial on the magnetic agitator is not in the "O" position then the magnetic agitator will heat up to the temperature set, regardless of the settings on the ETS-D5.

**Warning:** ETS-D5 is inactive.

The safety advice given above represents the current status of the known risks. Nonetheless, the user must check equipment for correct and safe operation before operating any system comprising the ETS-D5 and a magnetic agitator/heating system.

## Correct use

The IKATRON ETS-D5 is an instrument for the precise control of temperature. The sensor on this instrument must be inserted at least 20mm into the medium to be tempered. The instrument can be connected to any magnetic agitator or heating plate that has a contact thermometer connection, provided this conforms with the requirements listed in the technical data (see section Technical Data).

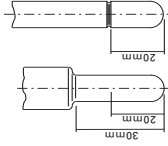## Unpacking

Please unpack the equipment carefully and inspect for damage. It is important that any transit damage should be identified at the time of unpacking. If necessary the equipment should be returned immediately for inspection (by post, rail or freight forwarder).

As supplied, the package contains: one IKATRON ETS-D5 and one set of operating instructions.

## Factory settings

The ETS-D5 is supplied with the following factory settings:

Operating mode	<b>A</b>
Set-point temperature	<b>-10 °C</b>
HI-TEMP	<b>450 °C</b>

### Restoring parameter to factory settings (RESET)

The following steps must be followed to restore factory settings:

- Switch off the equipment
- Press and hold down the membrane keys "▲" and "▼" and switch on the equipment
- The factory settings will now be restored (see above)

## Installation

Plug the instrument into the contact thermometer jack on the magnetic agitator (e.g. IKA RCT basic, IKA RET basic, RH basic, RH digital KT/C, Heidolph MR3001K, Heidolph MR3002, VWR VMS-D, VWR VMS-A, H+P Variomag, etc.).

Please observe all safety instructions in this respect.

Procedure for connecting a Corning Magnetic Agitator:

Turn the rotary "heat" dial on the magnetic agitator to the "O" (off) position. Next, connect the ETS-D5 to the magnetic agitator using the AD-C1 adapter provided.

Please observe all safety instructions in this respect.

The device will be activated when the magnetic agitator is switched on. If the supply voltage at the contact thermometer connector is not between 8V and 16V DC then it will be necessary to power the device using the H52 power supply set (H50 double connector adapter and H51 plug-in power supply unit). The power supply and/or double connector adapter will provide the necessary supply voltage for the device. An automatic self-test will be performed when the device is activated. All segments on the LCD display will be visible during this test (Figure 1).

## Operating modes

The ETS-D5 can be run in three different operating modes (**A**, **B**, **C**). The selected operating mode will be permanently displayed (Figure 4).

**Operating mode "A"** This is the factory-setting mode for this instrument. When the instrument is switched on, the set-point temperature is always set to -10°C. The set-point temperature can be adjusted using the "▲" and "▼" keys. The HI-TEMP (maximum set-point temperature) can only be set in this mode.

**Operating mode "B"** In this mode the most recent set-point temperature is used, and will become active when the instrument is switched off and on again. The settings can be adjusted using the "▲" and "▼" keys.

**Operating mode "C"** In this mode the current settings are used and will become active when the instrument is switched off and on again. The setting cannot be adjusted.

Holding down the "▲" key when switching on the instrument switches to the next operating mode in the sequence **A-B-C-A-B-C-A** etc.

## Setting the HI-TEMP

The "HI-TEMP" (maximum set-point temperature) can only be set when in operating mode "A". After the self-test, HI-TEMP will be visible on the LCD display (Fig.2). The HI-TEMP can now be set between 0 and 450°C using the "▲" and "▼" keys. The displayed HI-TEMP value will be set and stored if no key is pressed for approximately 5 seconds. The LCD display will then appear as shown Fig.3.


## Adjusting the set-point temperature

The set-point temperature can be adjusted in 0,1K steps by pressing the "▲" and "▼" keys. If the key is held down for longer than 5 seconds then the set-point temperature will change in 1K or 10K steps.

The LCD display (Figure 4) shows: the instantaneous actual temperature with physical units; the set-point temperature (underneath actual temperature);  to indicate that the heating for the connected device is switched on; the selected operating mode (top left).

## Replacing extension cable and sensor

To replace the sensor or the extension cable, pull the sensor and protective cap downwards until the plastic snap fastener is released. Then push the sensor or extension cable and protective cap over the fitting on the instrument to re-establish the connection.

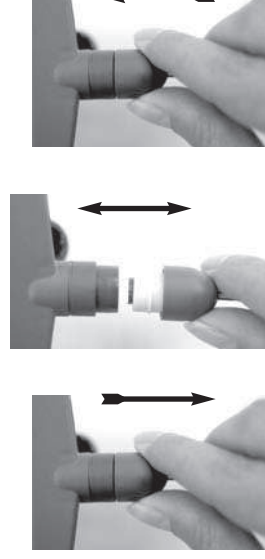## Error messages

The following error messages may be displayed (**Fig.5**) and have the causes listed below:

- Er1 Sensor not connected or defective (open circuit).
- Er5 Sensor not in media (no change in temperature of the media after heater active for 3 minutes).
- Er6 Sensor short circuit.
- Er7 Maximum temperature for the media exceeded (HI-TEMP setting).
- Er8 Calibration error (re-calibration by field service required).

## Accessories

- H 16 V** Support rod
- H 38** Retaining bar
- H 44** Cross sleeve
- H 50** Twin-jack adapter
- H 52** Power supply kit
- H 62** Standard sensor 270 mm stainless steel
- H 66** Glass sensor 270 mm glass
- H 70** Extension cable, 1 m long
- AD-C1** Adapter (for Corning magnetic stirrer)

## Technical data

Measuring range:	°C	-50 ... 450
Resolution:	K	0.1
Measuring accuracy:	K	±0.2 + sensor tolerance PT 1000 DIN IEC 751 Class A
Setting accuracy:	K	0.1
Type of control:		PID
Control deviation:	K	± 0,5
<small>(500 ml water in 600 ml glass beaker, magnetic stirring bar 40, 600 min<sup>-1</sup>, 50 °C)</small>		
Supply voltage:	V DC	8 ... 16 (Limited Power Supply)
Current consumption:	mA	8 (at 9V)
Ambient temperature:	°C	0 ... 60
Relative humidity:	%	80
Duty cycle:	%	100
Protection type:		IP-54
Cable length:	mm	350
Male connector:		6-pin DIN 45322
Socket:		DIN 12878 Class 2
Dimensions: (W x D x H)	mm	80 x 20 x 90 (without sensor)
Weight:	kg	0.2

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