# SEVERIN

---

Gebrauchsanweisung	Kühlschrank (D)
Instructions for use	Refrigerator (GB)
Mode d'emploi	Réfrigérateur (F)
Gebruiksaanwijzing	Koelkast (NL)
Instrucciones de uso	Frigorífico (E)
Manuale d'uso	Frigorifero (I)
Brugsanvisning	Køleskab (DK)
Bruksanvisning	Kylskåp (S)
Käyttöohje	Jääkaappi (FIN)
Instrukcja obsługi	Lodówka (PL)
Οδηγίες χρήσεως	Ψυγείο (GR)
Руководство по эксплуатации	Холодильник (RUS)

---

**Liebe Kundin, lieber Kunde,**

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**Dear Customer,**

Before using the appliance, the user must read the following instructions carefully.

**Connection to the mains supply**

The appliance should only be connected to an earthed socket installed in accordance with the regulations. Make sure that the supply voltage corresponds with the voltage marked on the rating label. This product complies with all binding CE labelling directives.

- Do not operate this unit by means of a transformer or, together with other electrical appliances, by means of an extension cable.
- To prevent noise vibration, ensure that the power cord does not touch any part of the rear of the unit (e.g. the heat-exchanger).

**General information**

This appliance is a compressor refrigerator, designed for short-term storage of food.

The unit is rated as Climate Class 'N' (Normal) and designed for domestic use. It complies with all relevant norms for domestic refrigeration appliances.

**Important safety instructions**

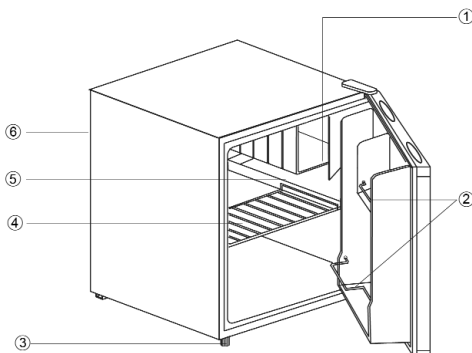
- To prevent the risk of personal injury or damage to the unit, it must only be transported while in its original packaging. For unpacking and installation, 2 persons are needed.
- The cooling circuit in this appliance contains the refrigerant isobutane (R600a), a natural gas with a high level of environmental compatibility, which is nevertheless flammable. Ensure, therefore, that none of the components of the refrigerant circuit become damaged during transport and installation of the appliance. If damage to the cooling circuit does occur, however, do not switch on or connect the appliance to the mains power. If there is an open fire or any other sources of ignition in the vicinity of the refrigerant gas, make sure it is removed immediately from the area and that the room is then thoroughly ventilated.
- **Warning:** Do not damage the cooling circuit. Any escaping refrigerant causes damage to the eyes; there is also a danger of the gas igniting.
- **Warning:** Keep ventilation openings, in the appliance enclosure or in the built-in structure, clear of obstruction. Sufficient ventilation must be ensured at all times.
- **Warning:** Do not use any external device (e.g. heaters or heating fans) to accelerate defrosting; follow only the methods recommended in this manual.
- **Warning:** Do not operate any electrical appliances (e.g. ice makers) inside the food storage compartment.
- Before it is connected to the mains power, the unit must be thoroughly checked for transport damage, including its power cord. In the event of such damage being found, the appliance must not be operated.
- To avoid the risk of fire, do not place any thermo-electric appliances on top of the unit. Do not operate electric water kettles on top of the unit. Do not place any liquid containers on top of the unit: this will prevent any leaking or escaping liquid damaging the electrical insulation.

- Only food commonly used in domestic households should be stored in the unit.
- Alcoholic substances may be stored only in properly sealed containers and in an upright position.
- To avoid the risk of food poisoning, do not consume food after its storage time has expired.
- Do not lean or put undue weight on the shelves, compartments, door etc.
- Protect the inside of the refrigerator at all times from open flames and any other sources of ignition.
- **Remove the plug from the wall socket**
  - in case of any malfunction,
  - before de-frosting,
  - during cleaning.
- When removing the plug from the wall socket, never pull on the power cord; always grip the plug itself.
- This appliance is not intended for use by any person (including children) with reduced physical, sensory or mental capabilities, or lacking experience and knowledge, unless they have been given supervision or instruction concerning the use of the appliance by a person responsible for their safety.
- Children should be supervised to ensure that they do not play with the appliance.
- In order to comply with safety regulations and to avoid hazards, any repairs or modifications to this unit must be carried out by our authorised service personnel, including the replacement of the power cord.
- We reserve the right to introduce technical modifications.
- If the unit is not used for an extended period of time, we recommend keeping the door open.
- If the unit is sold, handed over to a third party or disposed of at a suitable recycling facility, attention must be drawn to the presence of the insulation agent 'Cyclopentane' as well as to the refrigerant R 600 a.

## Familiarisation

1. Cold storage compartment
2. Bottle storage
3. Adjustable foot
4. Removable storage rack
5. Drip tray basin
6. Temperature control (on the rear)

## Before using for the first time

- Remove any exterior and interior packaging materials completely, including the adhesive strips.
- **Caution:** Keep any packaging materials well away from children, as they are a potential source of danger e.g. from suffocation.
- Before it is connected to the mains, the unit must be thoroughly checked for transport damage, including its power cord.
- Clean the unit as described in the section *De-frosting and cleaning*.
- To ensure proper functioning, wait for about 2 hours before connecting the unit and

switching it on after it has been positioned.

- When the unit is first switched on, a slight 'new' smell may be noticed. However, this will disappear once the cooling process has begun.

### Installation

The appliance should be set up in a well-ventilated, dry room. It should be operated in an ambient temperature of 16 to 32 °C and a relative humidity of max. 70 %. Do not expose the appliance to direct sunlight, and do not position it next to any heat sources (radiators, cookers etc).

- The rear and the sides should have a minimum clearance of 10 cm from the wall.  
The unit must not be installed inside a cabinet, nor directly below a hanging cupboard, shelf or similar object .
- Do not place any objects on top of the unit, and do not otherwise cover it.

If the floor is uneven, the adjustable foot can be used to compensate and ensure optimum stability. The unit should be installed at a slight angle in such a way that it leans to the rear, which will cause the door to close automatically so that it cannot be left open by mistake. Ensure, however, that both front feet are set to the same height.

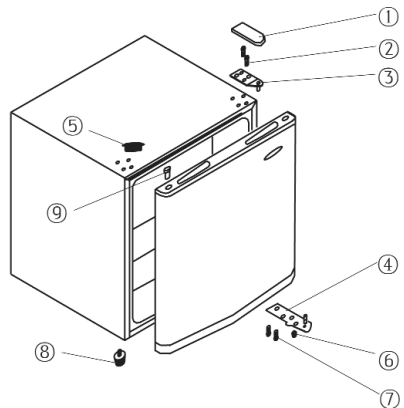
### Ventilation

The heated air around the unit must be able to circulate freely. Therefore, ensure that the circulation is not obstructed.

### Reversible door

The design provides for the door to be reversed, i.e. for right-hinged (default installation) to become left-hinged.

1. Remove the upper hinge cover (1).
2. Use a suitable pointed tool to remove the cover (5).
3. Use a screw-driver to remove the two screws (2) holding the upper hinge to the right side of the housing.
4. Carefully lift up the door and place it on a soft surface to protect it from scratches.
5. Remove the blind cover for the hinge mounting hole (9) and fit it in the corresponding place on the right-hand side of the door. Ensure that the cover is firmly pushed down into the hole.
6. Pull out the two bolts (7) which hold the lower right hinge plate (4) in position.
7. Remove the bolt on the left and insert it into the corresponding thread on the right.
8. Install the lower hinge plate on the left
9. Replace the door in its new position. Ensure that the hinge pin is inserted into the guide bushing (hole) on the lower part.
10. Install the upper hinge (3) on the left-hand side, using the screws (2). Ensure that the hinge pin is inserted into the guide bushing (hole) on the upper part.
11. Replace the cover (5) in the corresponding holes on the right.
12. Replace the hinge cover (1) on the upper hinge.

## Temperature control

The unit is activated by connecting it to the mains. It is only switched off completely when the plug is removed from the wall socket.

The temperature in the refrigerator cabinet can be adjusted by setting the temperature control.

- After switching on, the temperature control should be set to position **5**. After 20 minutes, the setting should be changed to position **3**, which is usually sufficient for home or office use.

## Storing food

- The temperatures inside the refrigerator cabinet and cold storage compartment depend mainly on the ambient temperature, the temperature control setting and the amount of food stored.

### *Cold storage compartment:*

- The cold storage compartment is designed for short-term storage of fresh meat, fish and poultry.

### *Main cabinet*

- Initially, the temperature control should be set to position **3**. If the cooling effect is too strong, lower it to position **1-3**; if it is insufficient, raise it to **3-5**.

**If you notice frost on the rear wall, the most likely cause is that the door has been left open for too long, or that warm food has been placed inside the refrigerator or that the temperature control has been set too high.**

Liquid or strong-smelling food, or food which is likely to absorb strong odours (e.g. cheese, fish, butter), should be stored separately or in hermetically sealed containers.

- **Warm food should be left to cool down to room temperature before being placed in the refrigerator.**
- Always ensure that the door is properly closed and that the seal is not obstructed by food stored inside the cabinet or door.

## Cleaning

- **Always remove the plug from the wall socket and turn the temperature control down to '0' before cleaning.**
- Do not pour water over or inside the unit.
- For frequent cleaning we recommend the use of warm water with some mild detergent. To prevent any unpleasant odour within the refrigerator, it should be cleaned about once a month.
- Any accessories should be cleaned separately with soapy water. Do not put them in a dishwasher.
- Do not use abrasives or harsh cleaning solutions or any cleaning agents that contain alcohol.
- After cleaning, thoroughly clean all surfaces with fresh water before wiping them completely dry. When re-inserting the plug into the wall socket ensure that you have **dry hands**.
- Take care not to remove or damage the rating plate inside the cabinet during cleaning.

### *De-frosting the freezer compartment:*

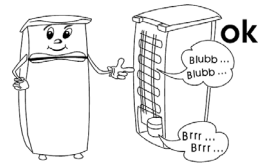
Once the build-up of ice around the door frame or on the vapouriser has reached a thickness

of 4 mm, the freezer compartment must be de-frosted. It should be cleaned at least twice a year as described below.

### Trouble-shooting

Certain typical sounds can be heard when the appliance is switched on. These sounds are:

- caused by the electrical motor within the compressor assembly; during compressor start-up the sound level will be slightly higher for a limited period of time.
- caused by the cooling agent flowing through the circuit.

The following table lists possible malfunctions and their probable causes and solutions. In the event of operational problems, check first whether a solution can be found using this table. If the problem persists, disconnect the appliance from the mains power and contact our Customer Service Department.

Problem	Possible cause and solution
The appliance is not working at all.	<ul style="list-style-type: none"> <li>- There is a power failure</li> <li>- The main fuse has blown.</li> <li>- The temperature control is set to '1'.</li> <li>- The fuse in the wall socket (if applicable) has blown. This can be checked by connecting another electrical device to the socket and checking for function.</li> </ul>
The normal operating sound changes or becomes louder.	<ul style="list-style-type: none"> <li>- Check for proper positioning (uneven floor?).</li> <li>- Are any adjacent objects affected by the running of the cooling unit (i.e. causing them to vibrate?)</li> <li>- Are there any objects on top of the unit which might be vibrating?</li> </ul>

### Transporting the appliance

The unit must only be transported in its vertical position.

Do not tilt it any more than 45°. During transport, ensure that the unit is properly secured and protected against shock, vibration and shifting.

### Disposal

 This appliance has been manufactured from recyclable materials. After removing the plug from the wall socket, the unit should be rendered unusable by cutting off the power cord before disposing of it through special waste collection points in accordance with local regulations. The refrigerant and the chemical agent in the insulation foam must be disposed of through a competent agency. Take special care not to damage the

cooling circuit before the unit is handed over for such disposal.

### Guarantee

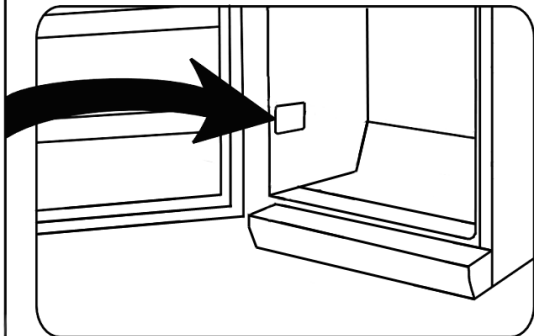
This product is guaranteed against defects in materials and workmanship for a period of two years from the date of purchase. Under this guarantee the manufacturer undertakes to repair or replace any parts found to be defective, providing the product is returned to one of our authorised service centres. This guarantee is only valid if the appliance has been used in accordance with the instructions, and provided that it has not been modified, repaired or interfered with by any unauthorised person, or damaged through misuse.

This guarantee naturally does not cover wear and tear, nor breakables such as glass and ceramic items, bulbs etc. This guarantee does not affect your statutory rights, nor any legal rights you may have as a consumer under applicable national legislation governing the purchase of goods. If the product fails to operate and needs to be returned, pack it carefully, enclosing your name and address and the reason for return. If within the guarantee period, please also provide the guarantee card and proof of purchase.

### Information needed by the Customer Service Department

Should repairs to the appliance become necessary, please contact our Customer Service Department, providing a detailed fault description and quoting the model number *KS ...* on the rating plate of the appliance (see picture). This information will help us to handle your request efficiently.

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## Technical specifications

Art.	<b>KS 9827</b>
Type of appliance	Refrigerator
Energy efficiency classification in compliance with directive 94/2 EC	<b>A+</b>
Energy consumption: <b>kWh/24 h</b> <b>kWh/ year</b>	<b>0.308</b> <b>113</b>
<b>The actual consumption depends on the pattern of use and the location of the appliance</b>	
Usable refrigerator capacity (litres)	47
Usable freezer capacity (litres)	-
Climate classification	N
Ambient temperature range, °C	16 - 32
Dimensions (mm)	525 x 440 x 490
Weight (kg)	16
Electrical specifications	See rating plate
Insulation foam	Cyclopentane

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Oy Harry Marcell Ab