Bürkle GmbH

Rheinauen 5 | D - 79415 Bad Bellingen
Telefon +49 (0) 76 35 8 27 95-0
Telefax +49 (0) 76 35 8 27 95-31
info@buerkle.de | www.buerkle.de

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood.
- ▶ When working with explosive substances, follow the technical rules for combustible liquids (in accordance with TRBF – (German) Technical Rules on Combustible Liquids) and wear appropriate protective equipment (zone 0, see annex).

⚠ WARNING!

Possible sparking due to electrostatic discharge. Danger of explosion if combustible liquids are used.

- ▶ Connect the earthing terminal before use (see accessories).
- ▶ Wear the necessary protective clothing.

- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.

Safety instructions during and after use

- ▶ The instrument generates overpressure. Possible bodily injury due to escaping liquids. Hold the shut-off valve on the discharge tube closed.
- ▶ The instrument generates overpressure. If there are differences in pressure, loud popping noises can be heard with some types of containers.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

Layout and function

- ▶ The instrument is made of conductive material (1.4301).
- ▶ The instrument is suitable for filling combustible liquids in accordance with groups IIA and IIB (see annex).
- ▶ Always use the instrument with pressure-resistant containers (up to 220 l, >1 bar) for filling into open containers (<1 l).
- ▶ When filling combustible liquids, always use an antistatic set (see accessories).
- ▶ The instrument is maintenance-free and suitable at a temperature range from 0 to 60 °C.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

⚠ WARNING!

Possible electrostatic discharge. Danger of explosion if combustible liquids are used.

- ▶ Connect the earthing terminal before use (see accessories).
- ▶ Wear the necessary protective clothing.

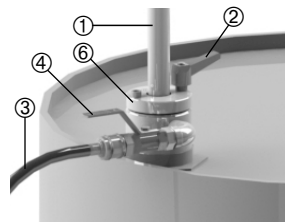
⚠ WARNING!

Some liquid will continue to run. Bodily injury due to escaping liquid.

- ▶ Close the shut-off valve on the discharge tube before mounting.

Assembly

1. Guide the barrel screw connection ⑥ and riser pipe with the discharge tube ① into the container and screw it together finger-tight.
2. If necessary open the clamping lever ② and adjust the riser pipe. Then close the clamping lever and lock the riser pipe in place.
3. Screw the discharge hose onto the riser pipe ①
4. Attach the connecting hose to the foot bellows on the instrument ③.

Safe filling

1. Open the shut-off valve ④ on the instrument.
2. Activate the foot bellows and generate overpressure.
3. Hold the discharge tube above the filling container.
4. Open the shut-off valve ⑤ on the discharge tube.
5. After filling is complete, close the shut-off valve on the discharge tube.

WARNING!

The container is under pressure. Bodily injury due to escaping liquid.

- ▶ Move the shut-off valve on the instrument to a 45° position ④.
- ▶ Allow excess air to escape completely from the container.
- ▶ After pressure decrease release any remaining liquids by the shut-off valve.

WARNING!

Some liquid will continue to run. Bodily injury due to escaping liquid.

- ▶ Close the shut-off valve on the discharge tube before dismounting.

Completing filling and dismounting

1. Close the shut-off valve ④ on the instrument.
2. Unscrew the connecting hose to the foot bellows ③.
3. Carefully unscrew the instrument from the container.
4. Rinse the instrument with a pH-neutral fluid.

Accessories

Type	Specification	Number	Item No.
Antistatic set	to prevent electrostatic charging	1	5602-1000

More accessories can be found at www.buerkle.de

Annex

Operational directive in accordance with 1999/92/EC:

Zone 0	a potentially explosive atmosphere is assumed to be continuously present
Zone 1	a potentially explosive atmosphere is assumed to be occasionally present
Zone 2	a potentially explosive atmosphere is not anticipated (and if it does occur, then only rarely and briefly)

Product directive in accordance with 94/9/EC:

Group I	for use in mining
Group II	for use in other areas subject to explosion hazard
Group IIA	Maximum permitted gap > 0.9 mm
Group IIB	0.5 mm < maximum permitted gap < 0.9 mm
Group IIC	Maximum permitted gap < 0.5 mm

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