ROTI®Store Cryoröhrchen

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Instructions for use

ROTI®Store cryo-vials

For preparing deep-freeze cultures of microorganisms

A. Product description

ROTI®Store is a product for preparing deep-freeze cultures. It enables safe storage and recultivation of microorganisms. These bind to beads of exceptional surface quality and are protected by a special cryo-medium when freezing. They may be stored at -20 °C for maximal one week. For long term-storage we recommend a minimum of -80 °C. In order to recultivate simply remove a bead and bring it into contact with desired culture medium.


Also suitable for conservation of marine microorganisms, even during field work (Siebert et al., *Appl. Environ. Microbiol.* 2004 (70) 5912-5).

B. Pack contents and storage

ROTI®Store contains 50 cryo-vials with 25±5 beads per vial in a special cryo-medium. The non-inoculated ROTI®Store cryo-vials can be stored between +4 °C to +25 °C (optimal storage: +4 to +8 °C). They should not be constantly exposed to sunlight.

C. Inoculation

- You will require a fresh overnight culture or colonies resuspended in culture medium with an OD₆₆₀ of at least 2 *
- Check the vial prior to use for possible contamination (cloudiness). Please label the cryo-vial with a waterproof marker (e.g. ROTILABO®LabMarker, Roth Art. No. 3083.2) and add approx. 0.5 ml culture solution.
- Close the vial tightly and invert it carefully several times.
- Remove the supernatant above the pellet with a sterile pipette (Roth Art.No. EA69.1) and dispose of it correctly.
- The inoculated and tightly closed ROTI®Store cryo-vial can now be frozen and stored at -20 °C for one week or at -80 °C (or lower) for years.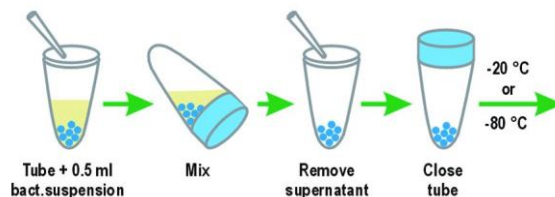

*Direct inoculation of ROTI®Store cryo-vials using picked colonies is possible, it does result, however, in much less efficient recultivation rates. In these cases, we recommend to resuspend the colony well in the cryo-medium. Slightly agitate the tube 15-20 mins. at 4 °C and remove the supernatant only then.

D. Recultivation

Remove the ROTI®Store cryo-vial from the deep-freezer and transport it in a cold cryo-block (e.g. ROTILABO®ice box, Roth Art.No. K088.1). The vial should not be allowed to thaw.

Remove a bead under sterile conditions with a sterile pincette or inoculation needle (e.g. Roth Art.No. EA88.1) and then close the cryo-vial.

The bead may now be added directly to the liquid medium or rolled out onto a solid culture medium.

Ø: 12 mm

length: 49 mm

Carl Roth GmbH + Co. KG

Schoemperlenstraße 3-5 • 76185 Karlsruhe

P.O. Box 100121 • 76231 Karlsruhe

Phone: +49 (0) 721/ 5606-0

Fax: +49 (0) 721/ 5606-149

info@carlroth.com • www.carlroth.com

jh 12/2021

The company is a limited partnership with headquarters in Karlsruhe, reg. court Mannheim HRA 100055. Roth Chemie GmbH, with headquarters in Karlsruhe, reg. court Mannheim HRB 100428, is the personally liable partner. Managing Director: André Houdelet. Sales tax identification number: DE 143621073.

ROTI®Store cryo-vials

P730.1	50 vials with 25 glass beads per vial
P730.2	250 vials with 25 glass beads per vial
P730.3	5 vials with 25 glass beads per vial