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🇬🇧 Cleaning & Maintenance

1. Clean b.safe Caps with a damp cloth and, if necessary, with detergents and warm water. **Attention:** Water must not get into the b.safe Air Valve.
2. Possible labelling or adhesive residues on b.safe Caps can be removed with isopropyl.
3. Once the b.safe Timestrip is activated, the white inspection window turns into red with increasing time of use. If the white writing "CHANGE" is completely visible on the red background, the b.safe Air Valve is dead and has to be replaced.

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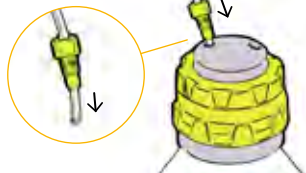
Push the capillary without frit through the b.safe Fitting and the b.safe Cap. Put the frit again on the tube.

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Put the fitted b.safe Cap on the eluent bottle and tighten it. If the frit touches the bottle base, also tighten the b.safe Fitting. If not, push the capillary further on.

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Screw in the Air Valve to the b.safe Cap.

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Push the button to activate the service life indicator.

Optional: Note the exchange date on the back side of the service life indicator.

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Fix the service life indicator on the b.safe Air Valve.

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Mark the b.safe Cap with a foil pen or an adhesive label if necessary.

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b.safe Cap with Stopcocks

Push the capillary through a suitable b.safe Fitting (length: distance from bottle neck to bottle base). Fix this capillary with the fitting as a suction pipe on the bottom side of the b.safe Cap. Push the capillary coming from your HPLC device through a b.safe Fitting and screw it into the top side. Screw the fitted cap onto the eluent bottle. If the lever of the stopcock stands horizontally, the suction way is interrupted.

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Important: Use b.safe Blind Fittings for gas tight closing of unused openings on a b.safe Cap!

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Bohlender GmbH · Waltersberg 8 · 97947 Grünsfeld