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Compatibility

All b.safe consumables are manufactured at highest quality standards. They are completely compatible with all known caps for eluent bottles and waste canisters on the market.

Please note: b.safe Air Vales require a thread UNF 1/4", b.safe Exhaust Filters are connected to thread GL14.

For optimum safety we recommend to use the complete b.safe system!

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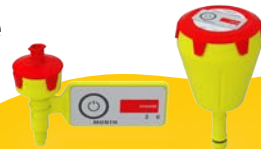
b.safe Abluftfilter / b.safe Exhaust Filter

Größe Size	Standzeit Service life	Verpackungseinheit Packing unit	Artikelnummer Cat. No.
Small	3 Monate / months	2 Stück / pieces	M 506-01
Medium	6 Monate / months	2 Stück / pieces	M 506-02
Large	6 Monate / months	2 Stück / pieces	M 506-03

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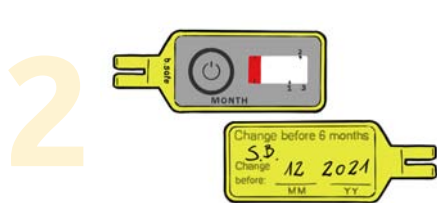
Screw in the Air Valve to the b.safe Cap.

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Push the button to activate the service life indicator.

Optional: Note the exchange date on the back side of the service life indicator.

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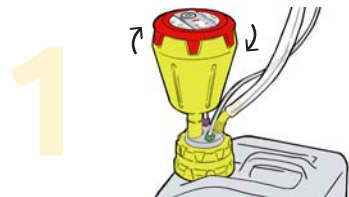
Fix the service life indicator on the b.safe Air Valve.

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Use the integrated luerlock connector for optional connection of dry tubes or insertion of protective gas.

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Screw a b.safe Exhaust Filter into the free thread connection of the b.safe Waste Cap.

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Push the button to activate the Service Life Indicator.
Optional: Note the exchange date on the sticker.

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Cleaning & Maintenance

Once the b.safe Timestrip is activated, the white inspection window turns into red with increasing time of use. If the white writing "CHANGE" is completely visible on the red background, the b.safe Air Valve is dead and has to be replaced.

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Important: Use b.safe Blind Fittings for gas tight closing of unused openings on a b.safe Waste Cap!

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Bohlender GmbH · Waltersberg 8 · 97947 Grünsfeld