

## Technical Information:

- TYPE:** Chill Checker Indikator, round, self-adhesive  
**SIZE:** 32 mm diameter at 6 mm depth  
**PRINT:** Plain-coloured. Blue on white background  
**PROCESS:** 1  
**TEMPERATURE:** in °F and °C, +/-1°C possible difference

Type	1:	-17°C
	2:	-8°C
	3:	-5°C
	4:	-2°C
	5:	+5°C
	6:	+9°C
	7:	+17°C
	8:	+20°C



Befor activation



Partial discoloration when the temperature has been reached for a short time, then drops below the measured value again.



Complete discoloration if the measured value has been permanently reached or exceeded.

### APPLICATIONS:

- To be activated by hand
- Instructions to use

1. Please make sure that the Chillcheckers are stored below their measured value for at least 1 hour.
2. Please remove the Chillchecker from the backing paper and attach it to the object to be measured.
3. To activate, please press the middle of the convex surfaces until it becomes concave. Note: Chillcheckers should ideally always be stored below their measured value. When this has been done, the Chillchecker can be removed from the cool storage area and placed and activated directly on the pre-cooled product. If the product has not been cooled, the Chillchecker can be activated.

### • Indication

After a short period of time that does not exceed five minutes, the colour changes of the white paper if it exceeds the specified dimension. If the temperature is constantly above the specified dimension, a complete discolouration will take place. If the product is frozen again, the discolouration will stop and start again when the dimension is exceeded.

- Packaging Unit: 100 pieces per box.

### • Main Application:

Difficult temperature conditions, for example during the transport of goods, medical transports, to identify and control or to better protect food during transit to warehouses or depots.

### • Durability:

In cool storage, ideally in the refrigerator, 6 months from the invoice date.