# HIRSCHMANN®

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## Instruction manual

Please read before use!

# pipetuses® standard

Elektrische Pipettierhilfe  
Electrical pipet filler

Hirschmann Laborgeräte GmbH & Co. KG  
Hauptstraße 7-15 • 74246 Eberstadt Germany  
Fon +49 7134 511-0 • Fax +49 7134 511-990  
[www.hirschmannlab.com](http://www.hirschmannlab.com) • [info@hirschmannlab.com](mailto:info@hirschmannlab.com)

X2120471000

## Inhalt

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## Contents

	Page
1. Safety instructions .....	5
2. Components .....	9
3. Functional principle .....	10
4. Start-up .....	11
4.1 Scope of delivery .....	11
4.2 Use .....	12
5. Prohibited operations .....	13
6. Restricted operations .....	14
7. Pipetting .....	15
7.1 Place the pipette .....	15
7.2 Fill the pipette .....	16
7.3 Empty the pipette .....	17
7.4 Operating speed adjustment .....	18
8. Spare parts replacement.....	19
8.1 Replacement of pipetholder complete .....	19
8.2 Replacement of filter.....	19
8.3 Replacement of safety valve.....	20
8.4 Safety valve - function check and trouble-shooting.....	21
9. Maintenance, cleaning and sterilisation.....	23
9.1 Maintenance .....	23
9.2 Sterilisation .....	23
9.2 Cleaning .....	24
10. Test of tightness .....	25
11. Repair at the manufacturer.....	25

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We are delighted that you have chosen our pipette filler. The unit is a modern, extremely accurate pipette filler which will facilitate your work considerably. The pipette filler is designed to the latest technical standards and is operationally safe. However, residual dangers can be associated with use of the unit, so it is imperative that the instruction manual be read carefully and all instructions therein heeded.

All the same, the instruction manual cannot illustrate every safety problem which may arise. Each user bears individual responsibility for adhering to safety and health regulations and determining and heeding limitations prior to using the unit.

Please observe the safety instructions prior to start-up in all cases!

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## 1. Safety instructions

- Observe all general safety practices and guidelines for laboratories, such as the use of protective clothing, safety goggles and gloves when handling hazardous chemicals.
- Never point the pipette in anyone's direction when in use!
- To avoid the risk of an electrical shock, never immerse either the unit or the recharger unit in water or any other liquid!
- Please ensure that the current of the power network matches with the voltage on the charger unit.

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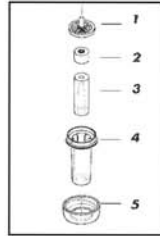
#### 1. Safety instructions

- In the event of any damage of the cable or the recharger unit, the recharger unit can not be used.
- Never touch the plug of the recharger unit with wet hands!
- Observe the guidelines and procedures of the chemical manufacturers (see also points 5 and 6).
- The unit may only be used in rooms. Protect from moisture!

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### 1. Safety instructions

- A safety valve (2) in the pipetholder case (4) is located on top of the silicone adapter (3) which prevents the liquid from getting into the handle, if the pipette is accidentally overfilled. If the safety valve has been activated, it must be removed by unscrewing the pipetholder and then cleaned and dried. Afterwards it must be inserted in the prescribed direction (see point 8). Wear safety gloves and goggles (Danger of infection)!

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---

### **Attention!**

If the unit is opened, the warranty is invalidated and any resultant claims void.

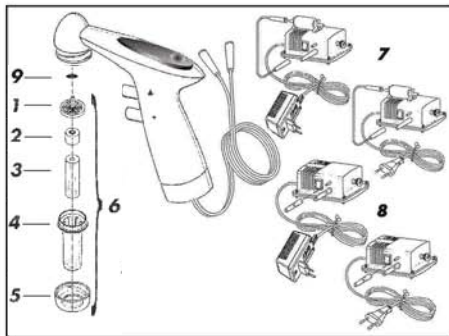
---

- The filter supplied contains a hydrophobic 0.2 µm membrane for air filtration in either direction.
  - Only have the unit professionally serviced!
  - Never use undue force!
- 

### **Attention!**

Only ever use the original recharger unit, accessories and replacement parts to avoid damaging the unit.

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## 2. Components

- 1 Membrane filter
- 2 Safety valve
- 3 Silicon adaptor
- 4 Pipetholder case
- 5 Nut
- 6 Pipetholder complete
- 7 Pump with filter
- 8 Pump without filter
- 9 O-ring

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### 3. Functional principle

- The unit is an electrical pipetting controller for use with all standard pipettes from 0.1 ml to 100 ml.
- The pipetting medium in the pipette is suctioned by pressing the In button ( $\triangle$ ) when the pump is activated, while pressing the Out button ( $\nabla$ ) empties the pipette.
- Pipetting speed is controlled by the pressure with which the button is pressed.

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## 4. Start-up

### 4.1 Scope of delivery

The following components should be contained in the original carton:

- Pipette filler
- 1 pump
- 1 hydrophobic membrane filter 0.2  $\mu\text{m}$  (yellow)
- 1 instruction manual

Please check that all parts have been delivered and are undamaged!

- The charger should not be used again if the cable or charger are damaged in any way!

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---

## 4.2 Use

- Connect the pump plug to the mains supply connection.

---

### **Attention!**

Compare the mains voltage to the voltage specification on the pump prior to connecting!

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---

### **The unit can not be used with:**

- liquids of which vapours attack the materials silicone, PTFE, Novodur, PVDF, NBR, PCPA, PP and PA6.
  - easily inflammable liquids (flame point below 0°C, for instance, ether and acetone).
  - usage in explosive areas.
-

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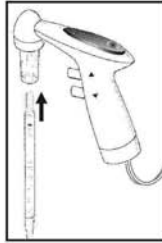
## 6. Restricted operations

- Working and charging temperatures between +10°C and +40°C.
- Usage only in areas in which a dangerous, explosive atmosphere occurs only seldom and then only short-termed.
- Do not keep unsupervised in an explosive area.
- Liquids which emit copious volumes of gas, as vapours are deposited inside the pipette filler and can corrode the unit.

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## 7. Pipetting

### 7.1 Place the pipette

- Place the pipette into the holder to ensure a secure and airtight fit!

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---

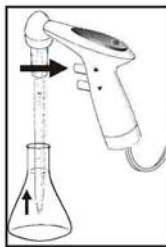
### **Attention!**

Danger of glass breakage, especially with thin pipettes.

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## 7.2 Fill the pipette

- For filling the pipette, press the In-button ( $\Delta$ ).
- Working speed is controlled by the pressure exerted on the button.

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---

### Attention!

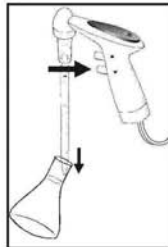
No liquid should get into the unit as this can adversely affect the operating functions.

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### 7.3 Empty the pipette

You can choose between two pipetting modes, blow-out and flow.

- When using blow-out pipettes:  
The green hose should be connected to the pump.  
The pipette is blown out when the Out button is pressed.
- When using Ex- (flow) pipettes:  
The green hose should not be connected to the pump. The pipette empties under atmospheric pressure.

⇒ To empty the pipette, press the Out-button (▽), the intensity of finger pressure on the button regulates the working speed.

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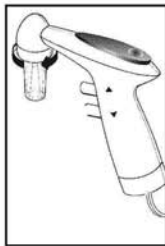
## **7.3 Operating speed adjustment**

The performance of the pump can also be influenced using the knurled nut on the pump (990 42 01).

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
## 8. Spare parts replacement

### 8.1 Replacement of pipetholder complete

- Screw off the nut.
- Replace the pipetholder by a new one.

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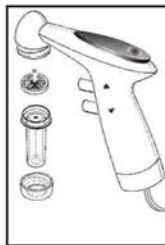### 8.2 Replacement of filter

- Screw off the nut.  
Remove the old filter and place the new one in the pipetholder case so that the pivot shows upwards. Screw the pipetholder onto the unit, by turning the nut clockwise.

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### 8.3 Replacement of safety valve!

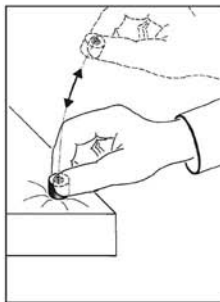
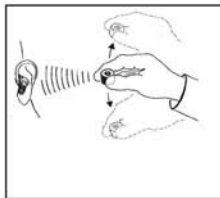
- Screw off the nut (5) counter-clockwise. Take out the filter (1) and place it aside. After that, remove the old safety valve (2) and place the new one in the pipetholder case (4), so that the black sealing ring shows upwards. Place filter (1) on the safety valve with the pivot facing up. Screw the pipetholder onto the unit by turning the nut clockwise.

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## 8.4 Safety valve - function check and trouble-shooting

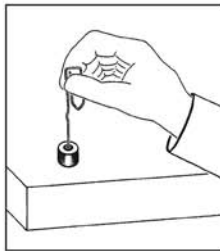
- IMPORTANT HINT FOR ASSEMBLING  
Black O-Ring of safety valve must be visible upside above silicone adapter!

Function check of the safety valve. The membrane of the valve must create a „rattling“ sound, while shaking the valve.

A sticking valve might become flexible again, by gently knocking it e. g. on the table edge. However in the event that the valve is glued, it must be soaked in a cleaning solution, rinsed with water and dried afterwards. Do not open the valve.

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
#### 8.4. Safety valve - function check and trouble-shooting

In case the function of the valve is not restored yet, try to make it work again, by gently pushing the membrane with a thin and blunt object.

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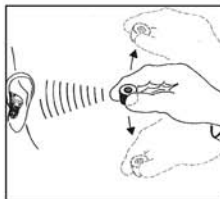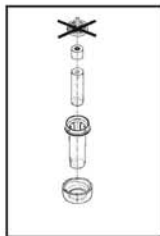
## 9. Maintenance, sterilisation and cleaning

### 9.1 Maintenance

Under normal conditions of use, the unit is maintenance-free.

However, the safety valve and membrane filter should be replaced in the case of dirt or a reduction in suction capacity (see point 8.2 and 8.3).

### 9.2. Sterilisation

The pipetholder, except the membrane filter, can be sterilised at 121°C.

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---

### 9.3 Cleaning

The safety valve (2) must be cleaned and dried before autoclaving. While reassembling the unit make sure that the side of the safety valve (2) with the black O-ring shows towards to the filter (1).

The pipette filler handles can be cleaned with a damp cloth. Do not spray the unit, as liquid can penetrate to the interior.

---

#### **Attention!**

When cleaning and exchanging the safety valve and membrane filter, protective gloves and goggles must be worn (Danger of infection)!

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## 10. Test of tightness

Once the required volume has been aspirated into the pipette, the liquid has to be kept in it unless the buttons are pressed. If the pipette however is dripping, disassemble the pipetholder and reassemble it again carefully.

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## 11. Repair at the manufacturer

Please observe the following points should it prove necessary to return the unit for repair (despite having handled it with all due care):

- Before sending the unit back, the user must ensure that it is carefully cleaned and decontaminated. The healthwise harmlessness must be confirmed, send the confirmation together with the unit to the manufacturer or the dealer (a copy of the clearance certificate can be found on page 36).

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#### 11. Repair at the manufacturer

- Units will only be repaired, when they have been cleaned and decontaminated by the user.
- For safety reasons dirty units will be returned to the sender without being repaired.
- Please describe the malfunction exactly.
- Please return the unit complete with the power supply unit in all cases, as this will facilitate checking of all components.
- Please quote the pipetted liquid.
- In the event of the unit being tampered with or being used in any improper way, (e.g. the application of undue force) any claims under the warranty are void.

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## 12. Trouble shooting

Problem	Possible cause	Solution
Liquid column in the pipette does not hold (pipette drops or runs out)	<ul style="list-style-type: none"><li>- Safety valve built-in incorrectly</li><li>- Pipetholder is untight</li></ul>	<ul style="list-style-type: none"><li>- Insert the safety valve so, that the side with the black O-ring points to the filter (see point 8).</li><li>- Test the correct assembly sequence for the pipetholder. If required, test or replace individual components.</li><li>- Silicone adapters can become worn or cracked. Change if necessary.</li><li>- Filters can be blocked. Change if necessary.</li></ul>
The pump will not blow out the pipette	<ul style="list-style-type: none"><li>- The green hose is not attached to, or not attached firmly to, the pump</li></ul>	<ul style="list-style-type: none"><li>- Push the hose on firmly.</li></ul>
The pump does not function neither filling nor blowing out	<ul style="list-style-type: none"><li>- Safety valve incorrectly positioned</li><li>- Pump not connected to net work</li><li>- Pump defective</li></ul>	<ul style="list-style-type: none"><li>- Place the safety valve so that the black O-ring shows towards the filter.</li><li>- Connect the unit plug to the mains supply.</li><li>- Return the unit to the manufacturer for repair.</li></ul>

Overfilled pipette

- Safety valve is activated immediately and prevents liquid from getting into the unit
- Remove safety valve, clean and reassemble. If required, replace the safety valve. During assembly take note of point 8.3.

---

Suction too slow or too weak

- Knurled nut has been set too loosely
- No tight, secure connection of the pipetholder
- Filter is full
- Silicone adaptor is damaged
- Knurled nut more tightly onto the handle.
- Screw the pipetholder more tightly onto the handle.
- Exchange the filter.
- Exchange the silicone adaptor.

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Input voltage: 15 V DC

Weight: handle complete 120 g  
incl. pump 420 g

#### CE:

This mark is the confirmation that the unit conforms to the established requirements of the EU guidelines and has been tested according to these criteria.

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## 14. Accessories - spare parts

The spare parts for your pipette filler listed below can be obtained from your specialised laboratory retailer:

Silicone adapter	990 32 01
Safety valve	990 32 02
Pipetholder complete	990 33 00
Pipetholder case	990 38 03
Nut green	990 39 02
Round filter set (consisting of 5 replacement filters)	990 37 04
Round filter set (consisting of 50 replacement filters)	990 37 05
Round filter set (consisting of 100 replacement filters)	990 37 06
O-ring	990 32 05

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14. Accessories - spare parts

Handle complete	990 31 00
hose connector "grey"	990 31 06
Microfibre-filter set for pump	990 41 04
Nut M 5	990 42 01
Rubber foot	990 42 03
Pump without filter, version EUR, 230 V	990 34 00
Pump with filter, version EUR, 230 V	990 44 00
Pump without filter, version UK, 230 V	990 34 10
Pump with filter, version UK, 230 V	990 44 10
Pump without filter, version EUR, 120 V	990 54 00
Pump with filter, version EUR, 120 V	990 64 00

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for unit ..... (Serial number)

Medium used: .....

Blood, bodily fluids, patholog. samples:

**YES**

**NO**

**Please define in case of Yes**

\_\_\_\_\_

Other hazardous biologic substances:

\_\_\_\_\_

Biologically-degradable materials which can pose a risk:

\_\_\_\_\_

Chemicals containing substances harmful to health:

\_\_\_\_\_

Radioactive materials:

\_\_\_\_\_

Other hazardous and toxic substances:

\_\_\_\_\_

### We hereby declare that

- the unit has been carefully cleaned and decontaminated prior to shipment
- no risks resulting from contamination exist
- we are aware that liability claims can be made against us in the event of damage being caused by contaminated units.

### Address of sender:

Name: \_\_\_\_\_

Date: \_\_\_\_\_

Signature: \_\_\_\_\_