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USER MANUAL

The Microsteril is a sterilizer of professional custom, booked to all the tools of work in metal. The sterilizing action, that develops to a temperature of 240°C, requires a time finite. The consumption of energy is extremely finite. His effectiveness has been butt and verified against: HIV - Vibrio Cholerae - Salmonella typhimurium - Escherichia coli - Bacillus subtilis - Bacillus stercatermophilus

ATTENTIONS

- ◆ We recommend to clean and dry carefully the material to sterilizer
- ◆ Before the start, please check that your load withstands without any damages at the set temperature (240°C +10% for safety).
- ◆ Don't touch the spheres of quartz during the operation of the sterilizer.
- ◆ Do not close the lid while the sterilizer is on.
- ◆ The apparatus could achieve elevated external temperatures.

C.B.M. is responsible for safety, reliability and performances provided that possible modifications or repairs are carried out by authorised people, the corresponding electric plant complies with IEC regulations and the equipment is used in compliance with these instruction for use.

INSTRUCTION FOR USE

- 1) Fill the pot with the quartz balls.
- 2) Insert the plug in a 230V socket and press the green switch. It will light itself and the orange lamp:
 - > Green lamp: power on
 - > Orange lamp: Heating elements working
- 3) After 45' minute the inner temperature gets 240°C (464°F)
The orange light will be off and on to keep steady the temperature.
- 4) Put inside the balls the cleaning instruments and leave them for the time they need to become sterile. (30" seconds)
- 5) For switch the sterilizer off, just press the green switch.
The sterilizer can remain on and ready for sterilizing till 10 - 12 hours.

IMPORTANT: in the sterilizer go inserted exclusively objects in metal, adequately washed.
Ascertaining that the sterilized instruments don't remain to the inland of the sterilizer beyond the necessary time

The C.B.M. S.r.l. guarantees his own products for one year against possible defects in workmanship, or against damages to the components not due to an improper use of the products.

CLEANING

ATTENTION: Before unscrewing any part of the oven be sure that this one is disconnected from the feeding network

We counsel of replace the quartz balls in the case they will be dirty

In order to do a good cleaning use neutral detergents only.

Do not use sponges or abrasive materials that can damage the surface.

To clean do not use detergents, disinfectants or other liquids containing chlorides as they are corrosive for stainless steel; in case of contact please rinse immediately.

WARRANTY

The C.B.M. guarantees his own products for one year against possible defects in workmanship, or against damages to the components not due to an improper use of the products.

C.B.M. S.r.l. is considered responsible concerning safety, reliability and performances on condition that possible modifications or repairs shall be effected by personnel authorised by C.B.M. , provided that the environment electric equipment is in compliance with IEC regulations and provided that the equipment is used according to these user's instructions.

DISPOSAL OF ELECTRIC AND ELECTRONIC REFUSE

Waste from Electrical and Electronic Equipment - Directive 2002/96/EC (WEEE)

The symbol (shown above) that appears on this product means that it should not be disposed of together with normal household refuse but should be treated separately.

Electric and electronic equipment may contain substances that are harmful to the environment and to the health of humans and should therefore be disposed of at special waste disposal sites.

There are harsh legal penalties in force against persons found disposing of such equipment in an irresponsible manner.

When the equipment is no longer of any use, it may be taken back to the outlet from which you intend to purchase a replacement.

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We: C.B.M. S.r.l., Via Castello 10, 26038 Torre de Picenardi (CR) Italy, in the person of its Sole Administrator and Legal Representative Mrs. Trusi Innocenta declare on our own responsibility, that the following electrical equipment:

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meet the provisions of following Directives which apply to them and its modifications:

- Directive CE 93/42 "Medical Devices"
- Directive CE 2006/95 "Low Voltage"
- Directive CE 2004/108 "Electro-Magnetic Compatibility"

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C.B.M. S.r.l.*L'Amministratore Unico – Sole Administrator