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## Instruction manual Vortex Mixer

### **Content:**

Introduction  
Service  
Warranty  
Safety instructions  
Proper use  
Inspection and scope of delivery  
Operation  
Maintenance and Cleaning  
Associated standards and regulations  
Technical datas  
Accessories

### **Introduction:**

Thank you very much for purchasing a Vortex Mixer from PHOENIX INSTRUMENT GmbH.

Please read the instruction manual carefully before you start to working with the instrument.

### **Service:**

When help is needed, you can always contact the service department of Phoenix Instrument.

Please provide the following information for Service:

- Serial number of the instrument (on the bottom panel)
- Description of the problem
- Methods and procedures adopted to resolve the problem.
- Your contact details

### **Warranty:**

The instrument is warranted to be free from defects in materials under normal use and service, for a period of 24 months from the date of invoice from Phoenix Instrument GmbH.

The warranty will not apply for the product or parts of it which have been damaged on account of improper installation and handling, misuse accident or abnormal conditions of operation.

For claims under warranty please contact your local distributor or Phoenix Instrument GmbH directly. Please always indicate article no., serial number and a description of the problem.

## **Safety instructions:**

- Read the operation instruction carefully before use.
- Ensure that only trained staff use the instrument.
- Make sure the socket is earthed before use.

When work, wear personal guard to avoid the risk from

- splashing of liquids
- mechanical vibrations can lead to glass breakage.

During operation

- hold the tube as vertical as possible
- mix the medium in the minimum pressure.
- use only original accessories from Phoenix Instrument.

Set the instrument on a stable, clean, dry, non-slip and fireproof surface. Do not operate in hazardous areas.

Put the "speed-control knob" to the lowest position. Gradually increase the speed (adjustable model). Reduce the speed if the medium splashes out of the vessel or the instrument is unstable on the stage.

Safe operation is only guaranteed with the original accessories from Phoenix Instrument. You will find the complete list in the appendix.

Accessories must be securely attached.

Check instrument and accessories before hand for damage each time you use it. Do not use damaged components.

Place a vessel at the centre of the mixing head. A sharp-edged vessel can wear out the mixing head.

Always disconnect the plug before fitting accessories.

The operation voltage mentioned on the label must match.

The instrument may only be opened by qualified and trained technician.

## **Proper use:**

The instrument is designed for mixing liquids in professional use in schools, laboratories, universities etc...

Do not use it in residential areas.

## **Inspection and scope of delivery:**

Unpack the equipment carefully and check for any damages which may have arisen during transport.

If it happens, please contact your local dealer or Phoenix Instrument GmbH promptly.

The packing includes:

- 1 Instrument with power supply
- 1 Instruction manual

Please retain all packing materials to guarantee a proper packing for warranty or service cases.

## **Operation:**

### **Permanent operation**

1. Place the mixer on the table, turn the speed control knob on the lowest position and plug-in the power supply.
2. Push the two-direction switch leftward to ON, power LED is lit. Then you can start to work.
3. Turn the speed control knob to set the rated speed (adjustable mode)
4. If you wish to finish, push the two-direction-knob rightwards to OFF, LED power is closed and instrument turned off.

### **Touch operation**

1. Place the mixer on the table and plug-in the power supply.
2. Push the two-direction knob rightward to the touch mode.
3. Press the test tube vertically on to the mixing head, the LED-lit is on and the instrument begins to work.. .
4. If you wish to finish, push the two-direction-knob rightwards to OFF, LED power is closed and instrument turned off.

## **Maintenance and cleaning:**

Proper maintenance and cleaning can keep instruments working in a good stage and will lengthen its lifetime.

Be careful not to spray the cleanser into the instrument. Unplug the power cord when cleaning. Please use only Isopropyl alcohol and water containing tenside. Before using other methods or cleaning agents, the user must ascertain with the manufacturer that this will not destroy the instrument.

Wear protective gloves during cleaning of the instrument.

Switch-off the instrument and put it in a dry, clean and stable place at room temperature in terms of long disuse.

## **Associated standards and regulations:**

The instruments are manufactured in accordance to the following standards and EMC-Richtlinien:

EN 61010-1

EN 61326-1

Associated EU-guidelines:

EMC-guideline: 89/336/EEG

Instrument guideline: 73/23/EEG

**Technical Datas:**

<b>Parameter</b>	<b>Datas</b>
Voltage (V)	220-230/110-120
Frequency (Hz)	50/60
Motor power (W)	60
Shaking movement	orbital
Orbital diameter	4mm
Motor type	Shaded pole motor
Motor rating input	58W
Motor rating output	10W
Speed range (rpm)	0-2500/ 2500(fix)
Dimension (mm)	127x130x160
Net weight	3,5 kg
Permitted ambient temperature	5-40°C
Relative humidity	80,00%
IP-protection according to DIN 60529	IP21

**Accessories (for both models):**

<b>Article number</b>	<b>Description</b>
RSV-E 10	Universal top plate for tubes and vessels up to 30mm Ø
RSV-E 44	Universal top plate for tubes and vessels up to 99mm Ø
RSV-E 12	Holding bar for RSV-E 20-40
RSV-E 15	100 mm plate with rubber mat
RSV-E 20	adapter for 48 tubes, 6mm Ø, to be used with RSV-E 15
RSV-E 25	adapter for 18 Tubes, 10mm Ø, to be used with RSV-E 15
RSV-E 30	adapter for 12 Tubes, 12mm Ø, to be used with RSV-E 15
RSV-E 35	adapter dor 8 Tubes, 16mm Ø, to be used with RSV-E 15
RSV-E 40	adapter for 8 Tubes, 20mm Ø, to be used with RSV-E 15
RSV-E 46	Rubber foot