
Cole-Parmer®

ST-200 Series

Portable Mini Stirrer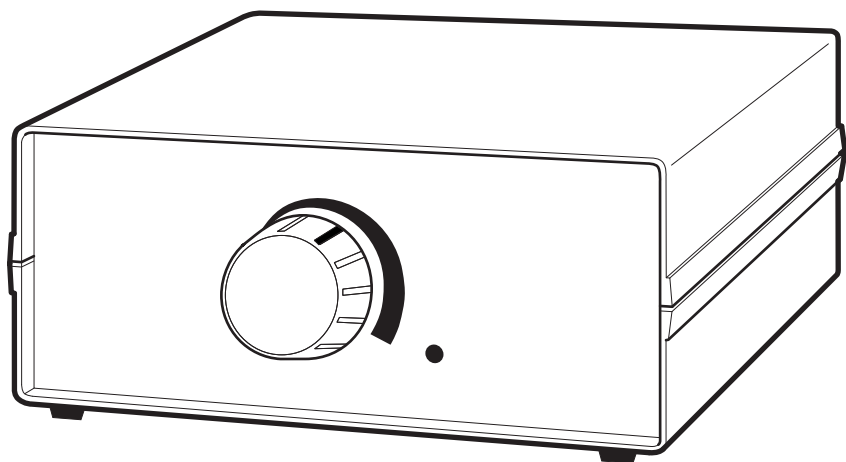

Instruction Manual

ISM1C5(S)-CPB Version 1.5

Cole-Parmer®
essentials

English

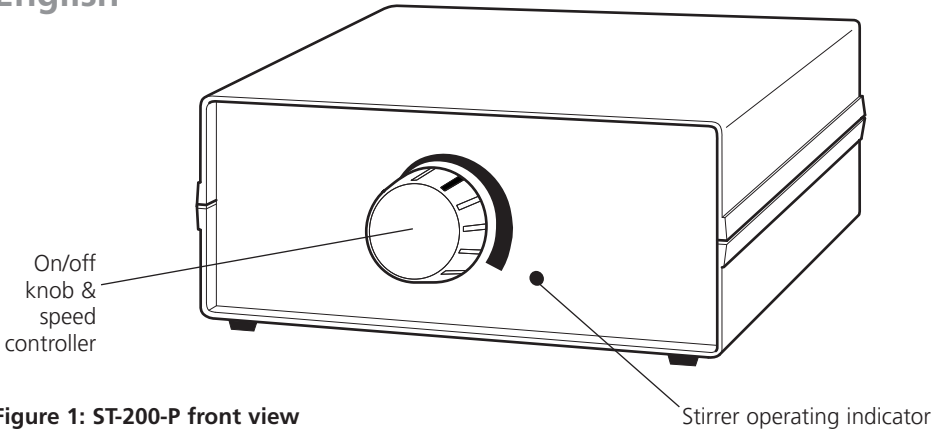
Figure 1: ST-200-P front view

ST-200-P PORTABLE MAGNETIC STIRRER

Thank you for purchasing this Cole-Parmer product. To get the best performance from the equipment, and for your own safety, please read these instructions carefully before use.

This equipment is designed to operate under the following conditions: -

- ❖ For indoor use only
- ❖ Use in a well ventilated area
- ❖ Ambient temperature +5°C to +40°C
- ❖ Altitude to 2000m
- ❖ Relative humidity not exceeding 80%
- ❖ Mains supply fluctuation not >10%
- ❖ Over voltage category II IEC60364-4-443
- ❖ Pollution degree 2 IEC664
- ❖ Use with a minimum distance all around of 200mm from walls or other items

If the equipment is not used in the manner described in this manual the protection provided by the equipment may be impaired.

Battery information

FOUR batteries, 1.5V, size D, IEC no. LR20

The batteries are fitted as shown in figure 2, with the positive terminals facing the rear of the instrument:

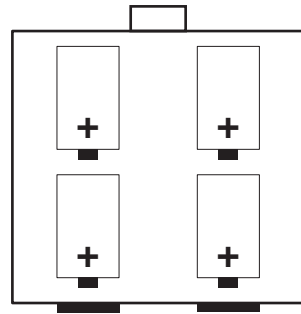
Figure 2: Battery orientation

Standard Alkaline Magnesium cells are recommended, but Zinc Chloride or rechargeable Nickel Cadmium batteries can also be used. Rechargeable batteries **MUST** be removed from the stirrer before recharging.

Main adapters

The ST-200-P stirrer can also be operated from an external 230Vac power supply PROVIDED THAT THE APPROPRIATE MAINS ADAPTER IS USED. Suitable adapters are detailed under 'Spares & Accessories'.

Adapters can be used with or without the batteries in place.

WARNING: The ST-200-P **MUST NOT** be connected directly to the mains electricity supply.

IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN

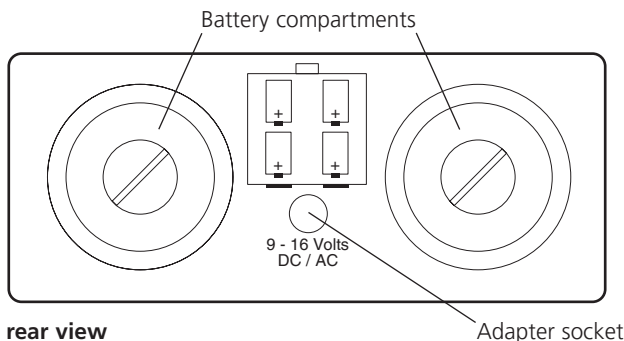
Figure 3: ST-200-P rear view

Safety Advice Before Use

- ❖ Never move or carry the unit with containers on the top plate.
- ❖ Do not leave discharged batteries in the unit as these may leak and cause corrosion.
- ❖ Remove rechargeable batteries before recharging
- ❖ Use only an appropriate adapter if utilising a power source other than batteries.
- ❖ Do not mix flammable liquids or use the equipment in hazardous atmospheres.
- ❖ In case of battery or mains interruption, a fault or mechanical failure, the unit will continue to operate on removal of fault.
- ❖ Mechanical energy can lead to breakage of glass vessels. Use with care.

ST-200-P General Description

The ST-200-P stirrer shown in figure 1 can be battery powered or used on mains supply. It is ideal for use inside incubators, glove boxes or in the field.

Operation

Stirring

Place the solution to be stirred together with a PTFE follower in a suitable vessel and place in the centre of the unit. Switch the stirrer 'ON' and adjust the speed of stirring using the speed control knob (figure 1). When the unit is

operating, the LED will be lit (figure 1).

Replacing the batteries

When the batteries are running low, the stirrer operating indicator (figure 1) will dim. To replace the batteries, open the battery compartments (figure 3) using a coin. Remove used batteries and dispose of safely. Replace new batteries with positive terminals facing the rear as shown in figures 2 & 3.

Using an adapter

Optional external power supply adapters are available for 230Vac and a car cigarette socket (see accessories overleaf).

Connect the stirrer to the adapter using the low voltage socket at the back of the unit (figure 3) before connecting the power supply.

Maintenance and Servicing

WARNING: Ensure the batteries are removed and /or the unit is disconnected from the mains adapter before attempting maintenance or servicing.

Cleaning

Periodically clean the instrument using a damp cloth and mild detergent solution. Do not use harsh abrasives or abrasive cleaning agents. Avoid strong acids and solvents on the ABS case.

Spares and Accessories

The following spares and accessories are available from Antylia Scientific or your laboratory dealer:

| Description | Catalogue Number |
|--|------------------|
| Mains adapter, 230V ac, UK plug | SM27/1 |
| Mains adapter, 230Vac, 2 pin European plug | SM27/2 |
| Adapter for car cigarette plug | SM27/3 |

For a comprehensive list of parts required by service engineers conducting internal repairs, please contact the sales department of Antylia Scientific quoting both the model and serial number.

Repairs

Only spare parts supplied or specified by Antylia Scientific or its agents should be used. Fitting of non-approved parts may affect the performance and safety features designed in to the instrument.

Any repairs or replacement of parts MUST be undertaken by suitably qualified personnel.

If in any doubt please contact the Technical Department at Antylia Scientific or the point of sale.

Technical Specification

| | |
|----------------------------------|---------------------------------------|
| Max Speed | 1300rpm |
| Capacity litres H ₂ O | 1.5 |
| Overall dimensions (w x d x h) | 150 x 160 x 70mm |
| Net weight | 1.3kg |
| Electricity supply | 4 x alkaline batteries or 230V, 50 Hz |
| Battery type | 1.5V, size D, ICE No. LR20 |
| Battery life | 6000 hours continuous use |
| Power | 1 Watt |

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Notes

Notes

This product meets the applicable CE Directives and UKCA Legislation for radio frequency interference and may be expected not to interfere with, or be affected by, other equipment with similar qualifications. We cannot be sure that other equipment used in its vicinity will meet these standards and so we cannot guarantee

that interference will not occur in practise. Where there is a possibility that injury, damage or loss might occur if equipment malfunctions due to radio frequency interference, or for general advise before use, contact the manufacturer.

Declaration of Conformity is available to view online at www.coleparmer.com

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Ordering Information

| Order No. | Series | Model | Legacy SKU |
|-----------|--------|----------|------------|
| 04801-05 | ST-200 | ST-200-P | SM27 |

Warranty Registration

UK

T: +44 (0) 1480 272279
E: uk.sales@antylia.com
W: coleparmer.co.uk

Germany

T: +49 (0) 9377 92030
E: de.sales@antylia.com
W: coleparmer.de

France

T: +33 (0) 1486 37800
E: fr.sales@antylia.com
W: coleparmer.fr

Italy

T: +39 (0) 284349215
E: it.sales@antylia.com
W: coleparmer.it

India

T: +9122 61394444
E: info@coleparmer.in
W: coleparmer.in

China

T: +1 847 549 7600
E: sales@antylia.com
W: coleparmer.com

USA

T: +1 847 549 7600
E: sales@antylia.com
W: coleparmer.com

Canada

T: +514 355 6100
E: info@antylia.ca
W: coleparmer.ca

Other

T: +1 847 549 7600

Cole-Parmer®
essentials

Antylia Scientific Ltd.
Beacon Road,
Stone,
Staffordshire,
ST15 0SA,
United Kingdom