KCL-Informationsbroschüre

Kat. III

Art. 866

D	Informationsbroschüre	946
GB	Information brochure	950
F	Brochure d'information	953
N	Informasjonsbrosjyre	954
DK	Informationsbrochure	955
FIN	Tiedotuslehtisen	966
H	Információs füzet	
S	Informationsbroschyr	
E	Folleto Informativo	
I	Opuscolo Informativo	
NL	Informatie brochure	
SK	Informačná brožúra	
HR	Informacije brošura	
EST	Infobrošüür	
CZ	Informační brožura	
P	Brochura de informaçaõ	
BG	Информационна брошура	
GR	Ευήμερωτικό φυλλάδιο	
LT	Informacinę brošiūrą	
PL	Broszura informacyjna	
LV	Direktivos	
RO	Broşuri de Informare	
SLO	Informacijska brošura	
TR	Direktif	
RUS	Информационная брошюра	

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The CE marking on this glove means that it meets the essential requirements for in the European Directive EEC 89/686 concerning Individual Protection Equipment (IPE): Harmlessness-Comfort-Dexterity-Sturdiness. Protective Gloves conform to the EN 420. Approved to EN 407 and EN 388 according to Article 10 of Directive 89/686/EEC: IFA, Alte Heerstraße 111, D-53757 St. Augustin, Identification number: 0121.

Surveillance after 11 B of Directive 89/686/EEC, chemical risks according to EN 374: same location as type test. The identification number for testing and certification centre 0121 for EC type testing and monitoring of quality assurance measures relates solely to the content of EN 407:2004-11 and the PPE Directive 89/686/EEC.

Article	Name	Size	EN 388	EN 407
866	866 K-MEX® L Spezial	9, 10	3441	X2423X
946	946 KarboTECT® LL	9, 10	3444	4344XX
950	KarboTECT® 950	7, 9, 10	1441	4241XX
953	953 K-TECT	9, 10	2540	4243XX
954	KarboTECT® L 954	7, 9, 10	1441	4243XX
955	Thermoplus® 955	8, 9, 10	244X	X1XXXX
966	966 Thermoplus® L	8, 9, 10	4343	X1XXXX

EN 388, mechanical Risks

1st digit Abrasion resistance (Min. 0; Max. 4)
3rd digit Tear strength resistance (Min. 0; Max. 4)

2nd digit Cut resistance (Min. 0; Max. 5)
4th digit Puncture resistance (Min. 0; Max. 4)

EN 407, thermal risks

1st digit Combustion characteristics (Min. 0; Max. 4)
3rd digit Convective heat (Min. 0; Max. 4)
5th digit Small spatters of molten metal (Min. 0; Max. 4)
(Keep protective gloves away from open flames.)

2nd digit Contact heat (Min. 0; Max. 4)
4th digit Radiated heat (Min. 0; Max. 4)
6th digit Large amounts of liquid metal (Min. 0; Max. 4)
X = not tested

Article 950, 953, 954: Gloves conform to Regulation (EC) No 1935/2004.
Protective gloves approved for short-term accidental contact with food.

! WARNING!

These protective gloves provide no protection against extreme cold (< -20 °C), electricity, chemical.

Do not use in the vicinity of moving machine parts. Risk of being drawn into the machine.

The specified protective index may be affected by factors such as temperature, abrasion, etc.

Unusable when the gloves are cracked, porous and stiff.

Article 866, 946, 955, 966: Safety glove not approved for food contact.

Article 950, 953, 954: Protective gloves only approved for a short-term accidental contact with food.

Allergy Notice: Contains in elastic ribbon natural rubber latex. Natural rubber latex can cause allergic reactions including anaphylactic reactions.

Article 866, 946, 954, 966: Gloves contains traces of chromate VI in the range permitted by law.

Use: The protective gloves you wear must be of the correct size. Note that using undergloves may result in some usage restrictions. Check the protective gloves for damage before you use them.

Cleaning: The specified gloves are not washable. Do not use any chemicals or sharp implements when cleaning (wire brushes, sandpaper, etc.). Dirt brush off.

Expiry date: With proper storage, no reduction in performance levels within 36 months.

Storage/Transport: Flat, dry, dark, with no additional weight load in its original packaging, at a temperature of 5 °C - 25 °C. Protect from sunlight and ozone.

Disposal: Gloves when used properly disposed of with household.

For further information please contact.

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KCL GmbH