# Poulsen & Graf

SUPERIOR LABORATORY PRODUCTS

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Operating manual  
Mode d'emploi  
Istruzioni per l'uso  
Instrucciones de  
manejo  
操作说明

FORTUNA<sup>®</sup>

<b>GER</b> -Bedienungsanleitung Zubehör	Seiten Seite	4 - 10 46 - 47
<b>ENG</b> -Operating Manual Accessories	Pages Page	11 - 17 46 - 47
<b>FRE</b> -Mode d'Emploi Accessoires	Pages Page	18 - 24 46 - 47
<b>ITA</b> -Istruzioni per l'uso Accessori	Pagine Pagina	25 - 31 46 - 47
<b>SPA</b> -Instrucciones de manejo Accesorios	Páginas Página	32 - 38 46 - 47
<b>CHI</b> -操作说明 配件	页 页	39 - 45 46 - 47

**Hersteller/Manufacturer/生产厂商:**

Poulsen & Graf GmbH, Karl-Carstens-Str. 10, 97877 Wertheim, Germany,  
Tel.: +49 - 9342/92290, Fax: +49 - 9342/922980, email: sales@poulsen-graf.com

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## Preface

Thank you for purchasing the POLYFIX® bottle top dispenser. You did a good choice. With 40 years of experience in making dispensers we guarantee a high standard of product quality and safety.

To guarantee quality and safety of the dispenser also for a long term, a correct operating and maintenance is necessary.

Therefore please read this operating manual carefully before using the dispenser and take notice of the warning instructions as well as the corresponding regulations.

Please pay attention to the colour code of our 3 POLYFIX bottle top dispensers:

■ 101.070, ■ 101.071, ■ 101.072

The delivery contents: 1 dispenser, 1 intake tube, 2 adapters made of PP (GL40+GL45), 1 volume setting system, 1 operating manual. Please check this before using the dispenser.

## Operating Manual Bottle Top Dispenser POLYFIX®

### I. Safety Instructions

1. Before using the dispenser please read the operating manual carefully and follow the safety instructions.
2. Observe general notes on hazards and safety regulations, e. g. wear protective clothing, gloves and glasses.
3. Observe any hazard symbols on reagent bottles.

4. Use only appropriate containers to dispense liquids. If using small bottles, please use a stand.
5. Always make sure that the discharge tube does not point at the user or any other person.
6. When transporting the dispenser please ensure that the piston is in zero-position and no liquid is in the cylinder.

**⚠ The stopper cap must close the discharge tube and the volume setting system must be set at zero-position!  
Never carry the dispenser at the finger guard, only at the PP valve block!**

7. Please check the correct assembling of the dispenser and the tight fit on the bottle before use.
8. To guarantee a trouble-free working please clean and use the dispenser according to the operating manual.
9. Repairs if necessary must be carried out by the manufacturer himself or authorized service stations only by using original parts.

## II. The World of POLYFIX®

**POLYFIX® 10 ml** Bottle top dispenser for aqueous liquids and acids

**Art. No. 101.070-38**

Piston made of precision ground glass. Integrated glass dispensing cylinder, plastic coated. Valve block made of PP. Valve system with Hastelloy metal spring in discharge channel. FIX adapter.

**Art. No. 101.071-38**

PTFE-coated piston. Integrated glass dispensing cylinder, plastic coated. Valve block made of PP. Valve system with Hastelloy metal spring in discharge channel. FIX adapter.

**Art. No. 101.072-48**

Piston made of glass. Integrated amber glass dispensing cylinder, plastic coated. Valve block made of PP. Valve system with Hastelloy metal spring in discharge channel. FIX adapter.

**POLYFIX® 50 ml** Bottle top dispenser for aqueous liquids and acids.

**Art. No. 101.070-48**

Piston made of precision ground glass. Integrated glass dispensing cylinder, plastic coated. Valve block made of PP. Valve system with Hastelloy metal spring in discharge channel. FIX adapter.

**Art. No. 101.071-48**

PTFE-coated piston. Integrated glass dispensing cylinder, plastic coated. Valve block made of PP. Valve system with Hastelloy metal spring in discharge channel. FIX adapter.

**Art. No. 101.072-38**

Piston made of glass. Integrated amber glass dispensing cylinder, plastic coated. Valve block made of PP. Valve system with Hastelloy metal spring in discharge channel. FIX adapter.

**⚠ Warning: Please see the technical information about polypropylene. This dispenser can only be used with chemicals against which Polypropylene is resistant.**

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#### Volume setting system

Easy-to-use volume setting system with fine adjustment for calibration. Suitable to be used by left- and right-handers.

#### Dispensing cylinder

Integrated glass cylinder, vacuum-calibrated and plastic coated, with stopper ring. Graduation black (101.070 + 101.071), white (101.072)

#### Valve block

Valve block made of PP. FIX Adapter.

#### Valves

Intake and discharge valve made of PP. Cristal valve balls. Hastelloy metal spring in discharge channel.

#### Dispensing piston

101.070: Dispensing piston made of precision ground glass, 101.071: PTFE-coated piston, 101.072: Dispensing piston made of precision ground glass.

#### Discharge tube

Integrated discharge tube made of PP. For safety reasons the discharge tube can not be taken apart. White stopper cap made of PP.

### IV. Handling

#### 1. Assembling / Filling

Before using the dispenser please check the dispenser for any damages or glass breakage.

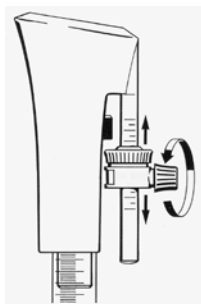
Push intake tube over the hose connection of the PTFE valve block.

Screw dispenser direct onto containers with thread GL 32 or use adapters to fix on any other threads or ground joint necks. Check the correct fitting. The dispenser can be turned on the flask and fixed in any direction by using the built-in FIX adapter.


For filling the dispenser the first time lift piston several times just a few until no air bubbles are seen in the cylinder.

## 2. Volume setting

Loosen the nut of the volume setting system, set the desired volume. The upper edge of the fine adjustment screw shows the volume set. Fix the screw.

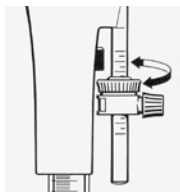For calibrating purposes turn the screw of the fine adjustment. Turn downwards to increase the volume, turn upwards to decrease the volume.

**⚠ Warning:**  
If the stopper cap is not removed when dispensing, there may be an over-pressure in the channel system and uncontrolled dispensing may be the consequence.

### 3. Dispensing procedure

- a) For safety reasons make sure that the discharge tube does not point at the user or any other person. By using the built-in FIX adapter it is possible to fit the dispenser accordingly. Remove stopper cap by holding the discharge tube tight with other hand.
- b) Set the desired volume with the volume setting system.
- c) Lift the dispensing piston up with the finger guard. Do not touch the piston itself to avoid contact with chemicals. The piston stops at the volume set. The graduation on the glass cylinder is an additional optical control of the volume set or air-bubbles.
- d) To dispense the liquid push the piston down.
- e) **Avoid uneven piston movements or hard pushes against the stopper ring.**
- f) Replace stopper cap after the dispensing procedure by holding the discharge tube tight with other hand.

#### **Warning:**

**The piston must be pushed down manually. Therefore chemicals may remain in the dispensing cylinder after dispensing. To avoid unintentional dispensing, always dispense the total volume set. The stopper cap must be placed on the discharge tube after every use. When dispensing observe safety precautions!**

### V. Operating Temperature

The operating temperature is from +15° C (+59° F) to +50° C (+122° F).

### VI. Sterilization

The dispenser can be autoclaved at +121° C (+248° F) / 1 bar. **Before sterilization loosen screw of the volume setting system.**

**POLYFIX 101.070 + 101.072:** the piston can remain in the cylinder when autoclaving  
**POLYFIX 101.071:** the dosing piston has to be sterilized separately.

 **Warning: Do not autoclave with higher temperatures. This will damage the dispenser. Wait before use until the dispenser has room temperature.**

**Sterilisation of POLYFIX® dispensers will put excessive stress on the dispensers both thermally and mechanically. As a consequence sterilization of POLYFIX dispensers will invalidate the warranty offered on the dispensers. Replacement for complaint after sterilization can be made only if sterilization protocol is available.**

### VII. Transport and Storage

1. **When transporting the dispenser never carry on the finger guard but always on the PP valve block to avoid unintentional dispensing of liquid! Replace the stopper cap before transporting and fix the volume setting system at the zero-position.**
2. Please transport and store the dispenser only in the original box to avoid damages. Storage temperature: -10° C (+14° F) to +50° C (+122° F).

### VIII. Maintenance and Cleaning

1. To guarantee a trouble-free function also for a long term it is necessary to clean the dispenser regularly.
2. Do not leave the dispenser empty without having it cleaned before to avoid blocking of the valve balls! If this happens all the same please put the dispenser for some hours in cold water or clean it by using ultrasound. Once the piston can be moved again, dismantle and remove any residual impurities by cleaning with a normal solution.
3. The dispenser can be left on the bottle all the time if the channel system is filled.
4. When changing the dispensing liquid the dispenser should be cleaned.

## IX. Repair Service

1. The POLYFIX® dispenser is mainly maintenance-free and reliable. If there is a problem or a defect and the user himself is not able to solve it, the faulty parts will be exchanged through our dealer or the manufacturer. In case a repair would be necessary, it must be done by the manufacturer or an authorized service station. This does not refer to parts which are available as spare parts, such as piston with finger guard, volume setting system, stopper cap, intake tube, adapters.

**2. Before returning the dispenser please do not forget to register it with the manufacturer in advance. Make sure that it is perfectly cleaned to avoid any contamination. Please also enclose a filled in safety declaration.**

3. Please operate the dispenser according to the operating manual to be entitled to the warranty. Repairs must be carried out by the manufacturer or authorized service stations.

## X. Disposal

When disposing the dispensers please pay attention to the national disposal regulations.

## XI. Technical Datas

### 1. Volume Range - Subdivisions - Accuracy

Volume range in ml	Graduation in ml	Accuracy A in %	Precision CV in %
2,0 - 10,0	0,2	$\leq \pm 0,6$	$\leq \pm 0,2$
10,0 - 50,0	1,0	$\leq \pm 0,6$	$\leq \pm 0,2$

The datas of accuracy and precision are established according to the German calibration ordinance (PTB) at a room temperature of +20° C (+68° F) by using distilled water. By using chemicals with a different viscosity or density a calibration on an analytical balance is necessary.

### 2. Packing

POLYFIX® dispensers 10 ml and 50 ml: single carton, 25,5 x 15,5 x 7 cm.

### 3. Weight, Dimensions

Volume range in ml	Weight in g appr.	Height in mm
2,0 - 10,0	132	160
10,0 - 50,0	237	220

## XII. Equipment

The dispenser comes in a break-proof single carton with 1 PTFE intake tube, 2 PP adapters (GL40+45), 1 volume setting system and 1 operating manual.

## XIII. Warranty

The POLYFIX® dispenser probably is the most reliable bottle top dispenser in the world. Therefore the manufacturer guarantees for manufacture and original parts for a 12 months period excluding all claims for damage, conversion and other claims. Under this warranty all parts that have been defective due to faulty material or production will be replaced. The manufacturer is not responsible for any damages or problems caused by transport or improper handling. All claims must be reported to the supplier. After a warranty repair the warranty period will not be extended. This warranty is not transferrable and void if repairs are carried out by unauthorized service stations.

Description of the malfunction	Possible reason	Solution
Air bubbles in the dispensing cylinder	<p>The dispensing piston has been pulled up too quickly.</p> <p>The system has not been vented.</p>	<p>Slowly and uniformly pulling out and pressing down the dispensing piston.</p> <p>Repeated filling and emptying of the dispensing cylinder</p>
The dispensed volume is too low.	<p>The dispensing piston has been pulled up too quickly.</p> <p>The system has not been vented.</p> <p>There are impurities in the intake tube.</p>	<p>Slowly and uniformly pulling out and pressing down the dispensing piston.</p> <p>Repeated filling and emptying of the dispensing cylinder</p> <p>Remove and clean the intake tube</p>
The dispensing piston is sluggish.	<p>There is contamination between the dispensing piston and the dispensing cylinder.</p> <p>The PTFE piston (<b>POLYFIX 101.071</b>) is swollen.</p>	<p>Disassembly and cleaning of dispensing piston and dispensing cylinder</p> <p>Disassembly of the dispenser, put the PTFE piston together with the finger protection cap in the freezer for a short time. Check the dosing liquid, liquids with Solvent-characteristics can only be dispensed with <b>101.070</b> or <b>101.072</b>.</p>
The dispensing cylinder can not be filled.	<p>The volume setting system sits at the top.</p> <p>The intake tube is missing.</p>	<p>Adjust the volume setting system to the desired volume</p> <p>Unscrew the dispenser from the bottle and insert the intake tube</p>

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All rights reserved, especially the rights to illustrations and original texts in their whole and in part, for any and all reproduction.
**POLYFIX Manual Ref 2018 01**, Copyright 2017 by Poulsen & Graf GmbH,