Bedienungsanleitung
Operating Instructions
Mode d'emploi

- » Auto-Exsikkatoren
- » Auto-Desiccators
- » Auto-Dessiccateurs

SICCO Auto-Star-Exsikkator, Artikelnummer: V 1877-07.

SICCO Auto-Star-Exsikkator Vitrum, Artikelnummer: V 1823-07.

SICCO Auto-Big-Star-Exsikkator, Artikelnummer: V 1894-07.

SICCO Auto-Maxi-Exsikkatoren, Artikelnummer: V 1914-01 und V 1916-01

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SICCO Auto-Maxi- Desiccator, Cat.-No.: V 1914-01 and V 1916-01

OPERATING INSTRUCTIONS

1. Supplied with	10
2. Security Advice	10
3. Taking into operation	10
4. Functionality	12
5. Service / Maintenance	13
6. Disturbances / Repairs	14
7. Waste Disposal	14
8. Specifications	14
9. Accessories	16

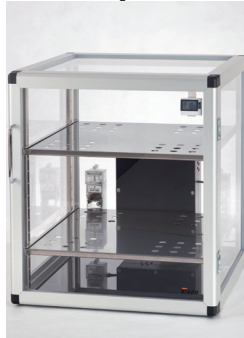
SICCO Auto-Star-Desiccators V 1877-07

SICCO Auto-Star-Desiccator Vitrum V 1823-07

SICCO Auto-Big-Star-Desiccators V 1894-07

SICCO Auto-Maxi-Desiccators V 1914-01 and V 1916-01

Sicco Auto-Desiccators

Please read the information listed below thoroughly. We urgently ask you to observe the instructions about security, use and maintenance.

SICCO Automatic Desiccators are made for long-term storage of reference samples, electronic components, photo equipment, documents, valuable assets and many more. The automatic drying unit reduces the relative humidity inside the desiccator to 20-30 %. The drying units work maintenance-free. The panels of SICCO Auto-Desiccators are made of high-quality and transparent plastic material respectively of Borosilicat glass 3.3 for model Vitrum with a strong aluminium frame and connectors made of polyamide.

1. Supplied with

- 1 Sicco Auto-Desiccator
- 4 Shelves made of acrylic glass incl. rails made of polyamide (Auto-Star-Desiccator)
- 4 Shelves made of stainless steel (Auto-Star Vitrum, Auto-Maxi 1- and Auto-Maxi 2-Desiccators)
- 2 Shelves made of stainless steel (Auto-Big-Star-Desiccator)
- 8 Rails for shelves (Auto-Star Vitrum, Auto-Maxi 1- and Auto-Maxi 2-Desiccators)
- 4 Rails for shelves (Auto-Big-Star-Exsikkator)
- 1 Hygrometer (fitted into the door)
- 1 Hygrometer per chamber (fitted into the door) for SICCO Auto-Maxi 2 Desiccator
- 1 Operating instructions

2. Security Advice

- » Do not take a damaged desiccator into operation.
- » SICCO Automatic-Desiccators are suitable only for 230 V AC.
- » Do not open the control unit.
- » Do not store any solvents or acids in the desiccator.
- » Do not put any explosive or inflammable products into the desiccator.
- » SICCO Desiccators are not suitable for overpressure or vacuum.
- » Provide a safe stand for the desiccator.
- » Place the desiccator in a reachable working height.
- » The ambient temperature range of -20 °C to $+70\text{ °C}$ respectively of -20 °C to $+120\text{ °C}$ shall not be exceeded.
- » SICCO Automatic-Desiccators are to be used in closed and dry rooms only.
- » Protect the Automatic-Desiccators from splash water.
- » Do not expose the desiccator to direct sunlight.
- » Remove only one shelf at once.

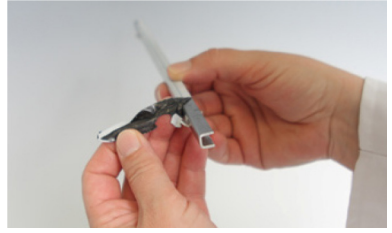
3. Taking into operation

- » Remove the protection cover.

Sicco Auto-Desiccators

» For versions with rails made of aluminium you need to remove the protection film from the rails.

» Remove the protection film from both sides of each shelf.

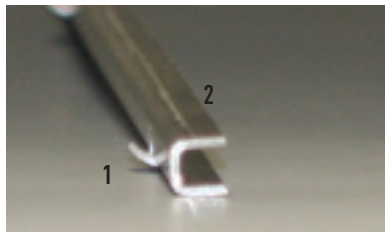
» Opening the door: Hold the desiccator housing with one hand while you open the door by pulling the door grip with the other hand.

» **Please note:** Due to the strong magnetic lock, it is harder to open the door the longer it was closed before. This is normal and can be compensated by pulling stronger than in the everyday use.

» Hook the rails made of aluminium (2) into the lateral mounted suspension arrangements (snap-in pin (1) up wards)

» Mount the rails (2) horizontally in pairs. Pay attention to mount them in the same height. A slight rotation of the rail helps to fix the snap-in pin safely in the suspension arrangement (3).

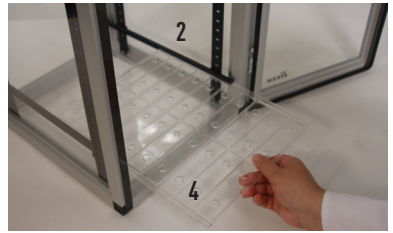
» For models with rail (2) and suspension arrangement (3) made of plastic: Mount the rails (2) horizontally in pairs on both sides of the suspension arrangement (3). The engagement of the rail (2) is clearly audible. For a better orientation, the levels are marked by numbers on the suspension arrangement (3).

Sicco Auto-Desiccators

- » The shelves (4) can then be slid in on the rails (2).
 - » Mount another pair of rails (2) and slide in another shelf (4).
- Please note:** Mount the rails in pairs and slide in immediately a shelf to avoid that other rails can be fall down accidentally.


- » Open the battery case on the back of the hygrometer by pushing the cover downwards.
- » Remove the battery safety strip and reinsert the battery. Close the battery case.

- » SICCO Auto-Star and Auto-Star Vitrum:
Connect the small plug (4) of the power supply (5) with the drying unit (6). Connect the power supply to a socket (230V AC). The green control lamp inside the drying unit starts to glow and you can hear a soft humming of the fan.

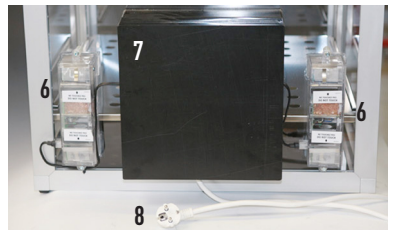
- » SICCO Auto-Maxi Desiccators/ SICCO Auto-Big-Star Desiccators:
The power supplies are connected with the switch box (7). Connect the grounding plug to a socket (230V AC). The green control lamps inside the drying units (6) start to glow and you can hear a soft humming of the fans.

4. Functionality


Drying is made by an automatic drying unit (6). This unit controls the alternation between drying of the air inside the cabinet and the regeneration of the integrated desiccant. The automatic dehumidification and desiccant regeneration cycle operates continuously every 20 minutes as follows:

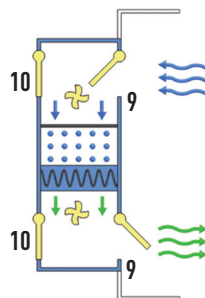- » During the drying cycle the circulating fan draws for approx. 14 minutes interior chamber air across the desiccant and thus dehumidifies the chamber. During this phase, the one-way flap valves are open to the interior of the desiccator (9) and closed to the exterior (10).

Sicco Auto-Desiccators

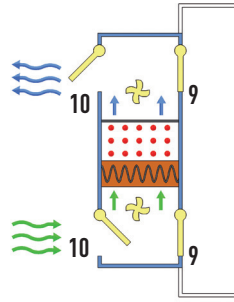» During the regeneration cycle the silica gel is heated electronically for about one minute. Afterwards the regeneration fan starts to operate and draws air in an opposite direction to the first fan. Due to the air flow, the one-way flap valves to the interior of the chamber (9) are closed and the valves to the exterior environment (10) are open. The desiccant is heated for another four minutes to release the captured moisture to the ambient air.

At the end of the cycle, the circulating fan operates for one minute with ambient air to cool down the desiccant to the ambient temperature. Then the cycle starts again.

Functions of the digital Thermometer/Hygrometer	
Measuring range:	<ul style="list-style-type: none"> » Temperature - 10 °C to + 60 °C » Air humidity 20 % to 90 % relative humidity
Max. error in measurement:	<ul style="list-style-type: none"> » +/- 1°C at -10 °C to + 50 °C » +/- 5 % relative humidity at 25 % to 75 % relative humidity and 0 °C to 50 °C » Attention: The mounted thermometer/hygrometer cannot display any values or data if the relative inside humidity of the desiccator is less than 25 % or higher than 75 %. To display exact values, we recommend using a thermometer/hygrometer of higher quality.
Battery:	<ul style="list-style-type: none"> » Please use a button cell LR44 or AG13.
Functionality of the Keys :	<p>Indication of maximum / minimum values:</p> <ul style="list-style-type: none"> » By pressing the MAX/MIN-key, the maximum value since the last reset will be displayed. » By pressing the MAX/MIN-key again, the minimum value since the last reset will be displayed. » The current values will be displayed after pressing MAX/MIN a third time. <p>Reset of the MAX/MIN-memory:</p> <ul style="list-style-type: none"> » Press CLEAR and the MAX/MIN-memory will be deleted. The current values will be stored instead. <p>Battery sign in the display:</p> <ul style="list-style-type: none"> » If the battery is low, a battery flag is shown on the left bottom of the display. Exchange the battery to ensure best possible operation of the device.

5. Service / Maintenance

- » Use only soft clothes for cleaning. Do not use any detergents which may contain solvents or which scrub.
- » Keep the magnetic door strip clean.

6. Disturbances / Repairs

The following information shall help you to recognise and, if possible, to eliminate possible reasons for disturbances by yourself. If you are not able to handle the disturbance, please give us a call. We will then try to help you this way. If then a solution to your problem cannot be found, we will ask you to return the cleaned desiccator into the original packing to us. Then we will check, whether a cost-effective repair is possible.

Objection	Reason and help
Humidity inside the desiccator does not fall.	<ul style="list-style-type: none"> » The desiccator is not connected to the mains. Connect the plug to a socket (230V AC). » The door is not closed tightly. Close the door. » The battery in the hygrometer is empty. Exchange the battery. » The built-in hygrometer might be faulty. Double-check the displayed values with an external hygrometer. Exchange the hygrometer.

7. Waste Disposal

The desiccators as well as their packing are made of recyclable materials. Please take further information on the materials used from the following specifications:

8. Specifications

SICCO Auto-Star-Desiccator

Cat.-No.:	V 1877-07
Drying Performance:	approx. 53 g moisture in 24 hours
Power Consumption:	max. 20 Watts
Mains Voltage:	230 Volt AC
Material panels:	PMMA (Polymethylmethacrylate), rear panel: Aluminium
Temperature resistance:	- 20° to + 70°C
Material frame:	Aluminium with connectors made of polyamide
Lock:	Magnetic tape
Overall dimensions (W x H x D):	310 x 525 x 375 mm
Excess length of drying unit:	60 mm (back side)
Usable interior space (W x H x D):	260 x 480 x 330 mm
Weight:	approx. 7,3 kg
Max. number of shelves:	26 pieces
Max. all-over load per shelf:	10 kg
Total all-over load:	30 kg

Sicco Auto-Desiccators

SICCO Auto-Star-Desiccator Vitrum

Cat.-No.:	V 1823-07
Drying Performance:	approx. 53 g moisture in 24 hours
Power Consumption:	Maximal 20 Watt
Mains Voltage:	230 Volt AC
Material panels:	Bororsilicate glass 3.3, rear panel: Aluminium
Temperature resistance:	- 20° to + 120°C
Material frame:	Aluminium with connectors made of polyamide
Lock:	Magnetic tape
Overall dimensions (W x H x D):	310 x 525 x 375 mm
Excess length of drying unit:	60 mm (back side)
Usable interior space (W x H x D):	260 x 480 x 330 mm
Weight:	approx. 13,5 kg
Max. number of shelves:	26 pieces
Max. all-over load per shelf:	10 kg
Total all-over load:	30 kg

SICCO Auto-Big-Star Desiccator

Cat.-No.:	V 1894-07
Drying Performance:	approx. 115g moisture in 24 hours per chamber
Power Consumption:	max. 80 Watts
Mains Voltage:	230 Volt AC
Material panels:	PMMA (Polymethylmethacrylate)
Temperature resistance:	- 20° to + 70°C
Material frame:	Aluminium with connectors made of polyamide
Lock:	Magnetic tape
Overall dimensions (W x H x D):	560 x 560 x 580 mm
Excess length of drying unit:	100 mm (back side)
Usable interior space (W x H x D):	495 x 500 x 540 mm
Weight:	approx. 13,5 kg
Max. number of shelves:	17 pieces per chamber
Max. all-over load per shelf:	30 kg
Total all-over load:	80 kg

Sicco Auto-Desiccators

SICCO Auto-Maxi 1 Desiccator

Cat.-No.:	V 1914-01
Drying Performance:	approx. 220g moisture in 24 hours
Power Consumption:	max. 80 Watts
Mains Voltage:	230 Volt AC
Material panels:	PMMA (Polymethylmethacrylate)
Temperature resistance:	- 20° bis + 70°C
Material frame:	Aluminium with connectors made of polyamide
Lock:	Magnetic tape
Overall dimensions (W x H x D):	560 x 1150 x 580 mm
Excess length of drying unit:	100 mm (back side)
Usable interior space (W x H x D):	495 x 1030 x 540 mm
Weight:	approx. 31 kg
Max. number of shelves:	34 pieces
Max. all-over load per shelf:	30 kg
Total all-over load:	160 kg

SICCO Auto-Maxi 2-Desiccator

Cat.-No.:	V 1916-01
Drying Performance:	approx. 115g moisture in 24 hours per chamber
Power Consumption:	max. 80 Watts
Mains Voltage:	230 Volt AC
Material panels:	PMMA (Polymethylmethacrylate)
Temperature resistance:	- 20° to + 70°C
Material frame:	Aluminium with connectors made of polyamide
Lock:	Magnetic tape
Overall dimensions (W x H x D):	560 x 1150 x 580 mm
Excess length of drying unit:	100 mm (back side)
Usable interior space (W x H x D):	495 x 500 x 540 mm per chamber
Weight:	approx. 35 kg
Max. number of shelves:	17 pieces per chamber
Max. all-over load per shelf:	80 kg per chamber
Total all-over load:	160 kg

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CE Declaration of Conformity

EC-Declaration of Conformity in compliance with Machinery Directive 2006/42/EC

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herewith declares that the following products:

Auto-Star-Desiccator, item number : V 1877-07

Auto-Star-Desiccator Vitrum, item number : V 1823-07

Auto-Big-Star Desiccator, item number : V 1894-07

Auto-Maxi 1-Desiccator, item number : V 1914-01

Auto-Maxi 2-Desiccator, item number : V 1916-01

Auto-Desiccator Wall, item number : V 1995-08

apply to the above mentioned directive including any amendments valid at the time this declaration was signed.

The low voltage directive 2006/95/EC has been fulfilled according to 2006/42/EC terms of protection objectives.

The following harmonized standards have been applied:

DIN EN ISO 12100

Safety of machinery

ISO 12100:2010

Risk assessment and risk reduction

DIN EN 61326-1: 1997

**Electrical equipment for measurement, control
and laboratory use, EMC requirements Part 1:
General Requirements**

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As a manufacturing company we can make products to your specifications. Send us a drawing or sketch together with answers to the questions listed below.

Return this form with contact details by fax or e-mail and we will be in touch to discuss details and give you a free, no obligation quote.

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BOHLENDER GmbH