Instruction Manual
Manuale di istruzioni
Manuel d'instructions
Manual de instrucciones
Bedienungsanleitung
操作指南

Rotilabo®-Heiz- und Magnetrührer MH 15

Roth-# AAN2.1

General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise / 基本信息

Before using the unit, please read the following instruction manual carefully.

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Attention, surface chaude!

Prudencia, superficie caliente!

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Do not dispose of this equipment as urban waste, in accordance with EEC directive 2012/19/UE.

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The product can be used with flammable liquids.

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This unit must be used for indoor laboratory applications only.

The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions. If the product is used in a way not specified by the manufacturer or with non specified accessories, product's safety may be compromised.

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This unit has been designed and manufactured in compliance with the following standards:

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Safety requirements for electrical equipment for measurement, control and for laboratory use

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IEC/EN 61010-2-051

IEC/EN 61010-2-010

Electrical equipment for laboratory use

UL 61010-1

General requirement - Canadian electrical code

CAN/CSA-C22.2 No.61010-1

VELP reserves the right to modify the characteristics of its products with the aim to constantly improve their quality.

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The plug disconnects the instrument. Therefore, place the instrument where it can be quickly disconnected.

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Use only the power cable provided with the instrument.

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The heated solution may release toxic, dangerous or poisonous gases. Adequate safety measures must be taken, in accordance with the safety regulations in force, including the presence of hood and personal protective equipment (masks, gloves, goggles, etc.).

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The vessel must be made of a suitable material to withstand the foreseen temperature.

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Magnetic field doesn't interfere with the function of cardiac pacemakers or data media if they are farther than 20 cm to the instrument.

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Position the instrument on a flat surface, with a distance from the wall of 30 cm (at least).

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It is responsibility of the user to appropriately decontaminate the instrument in case of dangerous substances fall on or in it accordingly to the safety datasheet of substances used and to the current laboratories safety standards. It is not possible to decontaminate the product under steam.

It is also responsibility of the user to use substances for cleaning or decontaminating which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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1. Introduction

EN

The MH 15 heating magnetic stirrer is used to heat and mix liquids inside a suitable container placed on the aluminium plate. The stirring is possible thanks to a magnetic drive stir bar placed into the vessel.

The instrument is used in the laboratory for general use and for all those applications that require precise regulation of the heating plate temperature and stirring speed.

The Aluminum alloy heating plate ensures:

- Optimum heat distribution and a high specific power thanks to the circular configuration
- Temperature homogeneity
- High resistance to thermal stress and thermal shock

The powerful brushless motor is able to stir from 100 up to 1500 rpm.

SpeedServo technology maintains constant speed as viscosity changes.

Note: it's important to choose the most suitable magnetic stirring bar related to the type and the quantity of the liquid to be stirred as well as to the type of the beaker. Size and shape of the magnetic stir bar determines the stirring efficiency at any given speed. The stirring bar which satisfies most applications is code A00000356 (8 x 40 mm).

2. Installation

- Unpacking
 - Check the integrity of the unit after unpacking
- The box includes
 - MH 15 heating magnetic stirrer
 - Instruction manual
 - Power cable
- Installation
 - Place the unit on a non-flammable surface
 - Make sure that the technical specification of the unit and of the power supply line are the same
 - Make sure that the electrical network is grounded
 - Make sure that the main switch is on position "0" (OFF)
 - Connect the instrument to an easily accessible socket (compliant with the current safety norms), using only the provided power cable

Note: keep the power cable far away from the hot plate.

- Place the container with the liquid and the stir bar on the stirring plate

3. Operation

Power-up

- Turn the unit on using the main switch
- The right led between the two knobs lights up

Stirring

- Rotate the right knob to set the stirring speed, the led over the speed knob lights up
- The real speed increases until setpoint achievement
- A microprocessor ensures constant speed even if the viscosity changes (SpeedServo)
- Turn the knob to the 0 positions to stop stirring

Heating

- Rotate the left knob to set the heating plate temperature
- If the set temperature is higher than the actual one, the led over the temperature knob lights up
- When the heating plate temperature exceeds 50 °C, the left led between the two knobs lights up until the temperature drops below 50 °C

Note: This warning is active even if the main switch is off but not if the instrument is not powered.

- Turn the knob to the 0 positions to stop heating

4. Error and warning messages

When the control panel shows one of the following error messages, the instrument functions automatically stop.

Display indicator	Malfunction:
Flashing Heating led and Hot plate led ON	Heating plate overtemperature ($T > 330^{\circ}\text{C}$)
Flashing Heating led and Hot plate led OFF	Heating plate safety probe overtemperature ($T > 370^{\circ}\text{C}$)
Flashing Stirring led	The stirring system doesn't run correctly

To remove the error message, restart the instrument from the main switch. For the overtemperature of the safety probe is necessary to disconnect the instrument from the power supply.

If the alarm persists on the control panel, please contact VELP Scientifica's technical assistance service.

5. Maintenance

No routine or extraordinary maintenance is required except periodic cleaning.

Cleaning

Disconnect the unit from the power supply and use a cloth dampened with a non-flammable non-aggressive detergent.

Repair

Repairs must be carried out by authorized Velp personnel only.

The transport of the instrument by freight forwarders, couriers or others must be carried out using the original shockproof packaging. Follow any instructions on the original packaging (e.g. palletizing).

It is the responsibility of the user to properly decontaminate the unit in case of hazardous substances remaining on the surface or interior of the device. If in doubt about the compatibility of a cleaning or decontamination product, contact the manufacturer or distributor.

6. Technical data

General features	Model	MH 15
	Voltage	230 V 50/60 Hz
	Power input	630 W
	Current consumption	2,8 A
	Dimensions (WxHxD)	160x85x270 mm (6,3x3,3x10,6 in)
	Weight	1,7 kg (3,7 lb)
	Construction material	Aluminum body –Technopolymer enclosure
	Working in continuous	Admitted
	Maximum load on the plate	25 kg
	Noisiness	<< 80 dBA
	Environmental temperature admitted	+5...+40 °C
	Storage temperature admitted	-10...+60 °C
	Level of electrical protection CEI EN60529	IP 42
	Max humidity	80%
	Overvoltage category	II
	Pollution degree CEI EN61010-1	2
	Max altitude	4000 m
Heating	Heating plate power output	600 W
	Heating plate dimensions	Ø 135 mm
	Temperature range	50 ÷ 310 °C
	Heat control type	Analogic
	Heating plate construction material	Aluminum
	Safety circuit	Separated with dedicated probe
	Hot plate alarm	Over 50°C
Stir	Overtemperature alarm	Over 330°C
	Stirring capacity	20 l H ₂ O
	Speed range	100 ÷ 1500 rpm
	Stir control type	Analogic
	Motor type	BLDC
	Motor rating output	10 W

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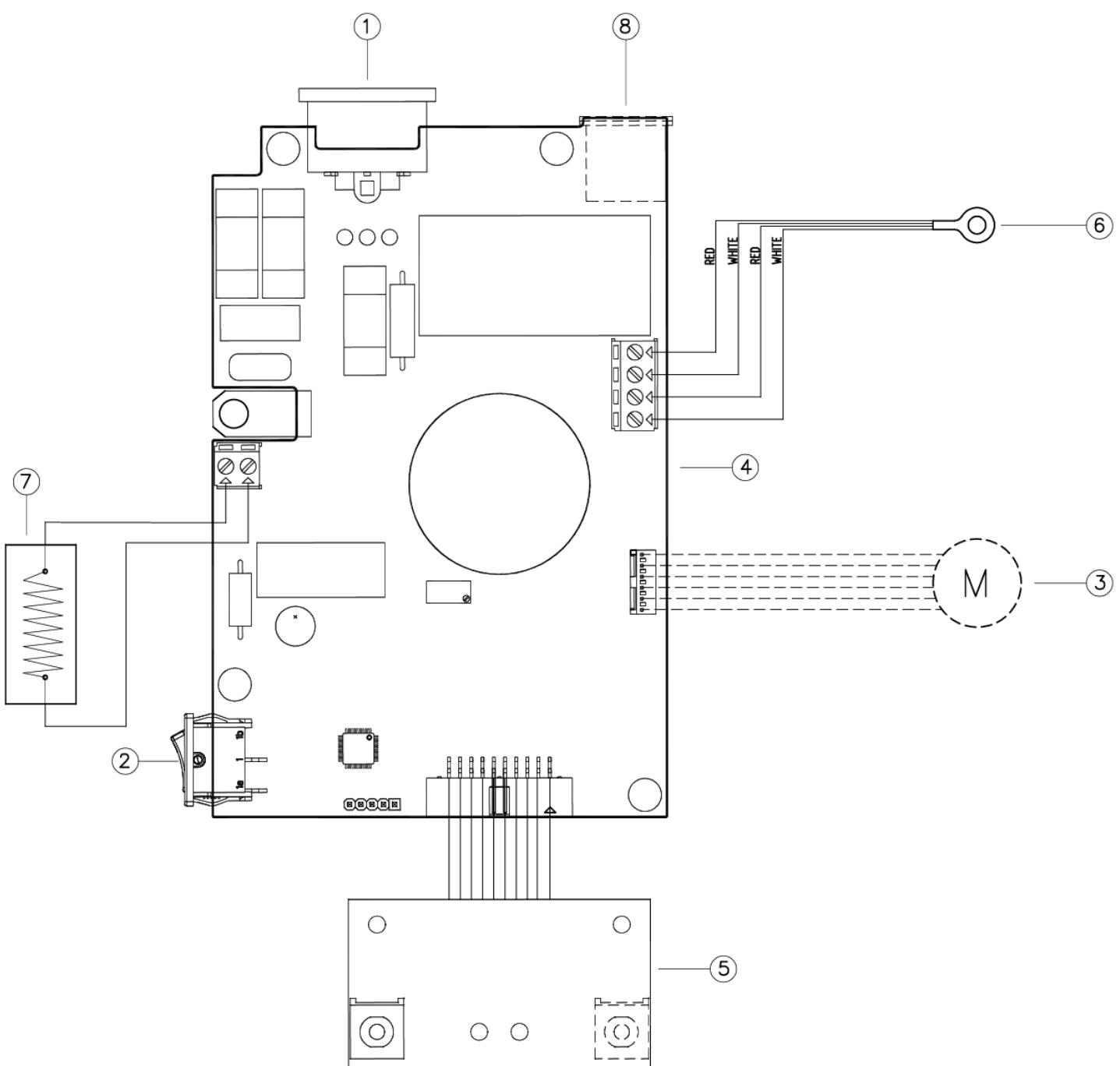
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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EN 61010-1	Safety requirements for electrical equipment for measurement, control, and laboratory use Part 1: General requirements
EN 61010-2-010	Particular requirements for laboratory equipment for the heating of material
EN 61010-2-051	Particular requirements for laboratory equipment for mixing and stirring
EN 61326-1	Electrical equipment for measurement, control and laboratory use - EMC requirements - Part 1: General requirements

and satisfies the essential requirements of the following directives:

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2006/42/EC	Machinery directive
2014/30/EU	Electromagnetic compatibility directive
2015/863/EU (RoHS III)	Restriction of the use of certain hazardous substances
2012/19/EU (WEEE)	Waste of electric and electronic equipment

plus modifications / più modifiche / plus modifications / más sucesivas modificaciones / in der jeweils gültigen Fassung /
加修改.

11. Declaration of conformity

We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following regulations:

S.I. 2016/1101 Electrical Equipment (Safety) Regulations 2016

S.I. 2016/1091 Electromagnetic Compatibility Regulations 2016

according to the relevant designated standards:

EN 61010-1 Safety requirements for electrical equipment for measurement, control, and laboratory use
Part 1: General requirements

EN 61010-2-010 Particular requirements for laboratory equipment for the heating of material

EN 61010-2-051 Particular requirements for laboratory equipment for mixing and stirring

EN 61326-1 Electrical equipment for measurement, control and laboratory use - EMC requirements -
Part 1: General requirements

and satisfies the essential requirements of regulations:

S.I. 2008/1597 Supply of Machinery (Safety) Regulations 2008

S.I. 2012/3032 Restriction of the Use of Certain Hazardous Substances in Electrical and Electronic
Equipment Regulations 2012

S.I. 2013/3113 Waste Electrical and Electronic Equipment Regulations 2013

plus modifications.

Thank you for having chosen VELP!

Established in 1983, VELP is today one of the world's leading manufacturer of analytical instruments and laboratory equipment that has made an impact on the world-wide market with Italian products renowned for innovation, design and premium connectivity. VELP works according to **ISO 9001**, **ISO 14001** and **OHSAS 18001** Quality System Certification.

Our instruments are manufactured in Italy according to the IEC 1010-1 and CE regulation.

Our product lines:

Analytical instruments

Elemental Analyzers
Digestion Units
Distillation Units
Solvent Extractors
Fiber Analyzers
Dietary Fiber Analyzers
Oxidation Stability Reactor
Consumables

Laboratory Equipment

Magnetic Stirrers
Heating Magnetic Stirrers
Heating Plates
Overhead stirrers
Vortex mixers
Homogenizers
COD Thermoreactors
BOD and Respirometers
Cooled Incubators
Flocculators
Overhead Shakers
Turbidimeter
Radiation Detector
Open Circulating Baths
Pumps

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Carl Roth GmbH + Co. KG