Intelligent Evaporation

Leading Safety Standards

Superior Ease of Use

Reduced Cost of Ownership

Operating Manual

2

Rotary Evaporator Hei-VAP Expert, Hei-VAP Expert Control, Hei-VAP Ultimate, Hei-VAP Ultimate Control

Operating manual must be read before initial start-up.

Please follow the safety instructions provided.

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About this Document	4
Version and variants	4
About this manual	4
Symbols and keywords	5
Safety instructions	6
General safety instructions.....	6
EU Declaration of Conformity.....	6
Intended use.....	6
Installation / electrical safety	7
Personnel training	7
Operator duties	7
Installation site	7
Modifications to the device	8
Employee safety	8
Safety during use	8
Disposal	9
Device description	10
Device overview	10
Overall view Hei-VAP Expert, Hei-VAP Expert Control	10
Overall view: Hei-VAP Ultimate, Hei-VAP Ultimate Control	11
Control box (for control variants only)	13
Control panel: Hei-VAP Expert, Hei-VAP Expert Control	13
Control panel: Hei-VAP Ultimate, Hei-VAP Ultimate Control	14
Start-up	15
Setting up the device	15
Connecting/disconnecting the power cord	15
Remove transportation lock.....	15
Switching the device on and off.....	16
Heating bath and flask lift.....	16
Basic Settings	19
Menu navigation.....	21
Overview of the menu tree.....	21
Operation of the menu	23
Changing the basic settings	26
Language and formats.....	26
Date and time	26
Settings	27
Applications	30
Symbols used	31
Operation	35
Process sequence	35
Selecting an application.....	35
Recording process times	37
Applications	38
Setting parameters manually ("Application Basic")	38
Using and managing favorites (Application "Extended")	42
Using and managing ramp profiles (Application "Extended")	46

Using automatic mode DAA (Application "Extended").....	51
Starting an application with a timer.....	53
Interfaces.....	54
Evaporating and receiving flasks.....	55
Assembling and removing the evaporating flask.....	55
Assembling and removing the receiving flask.....	56
Ventilation.....	56
Cleaning and Maintenance	57
Cleaning.....	57
Maintenance.....	58
Troubleshooting	59
Malfunctions and remedies.....	59
Resetting temperature limiter.....	66
Assembly	67
Electrical connections.....	67
Installation.....	68
Assembly parts.....	68
Transportation lock.....	70
Heating bath.....	70
Control panel.....	72
Vapor tube.....	73
Condenser.....	75
Evaporating flask: Angle and immersion depth.....	80
Control box.....	82
Vacuum manifold.....	82
Attaching the peripheral devices.....	83
Connecting vacuum.....	83
Cooling water connection.....	87
Optional accessories.....	88
Dismantling and Storage	92
Dismantling, transportation and storage.....	92
Dismantling.....	92
Transportation and storage.....	93
Spare parts and accessories	94
Scope of delivery.....	94
Accessories.....	94
Software copyright.....	95
Appendix	96
Technical data.....	96
Service	97
Contact / Technical Service.....	97
Warranty.....	98
Declaration of no objection.....	99

Copyright

This operating manual is protected by copyright and is exclusively intended for the internal use of the purchaser.

Handing over this operating manual to third parties, reproducing it in any type or form –in part or in full – and utilizing and/or communicating its contents is prohibited without prior written approval from the manufacturer.

Any violations will result in an obligation to provide compensation for damages. Other claims remain unaffected.

➤ Version and variants

Version

This manual describes the function, operation and maintenance of the Hei-VAP Expert and Hei-VAP Ultimate rotary evaporators.

Version	Alteration date
1.0	04/2019

Variants

The device is available in different variants. Certain features or functions are only available in specific product variants. These variants are described in this operating manual.

➤ About this manual

This operating manual is an integral part of the device described here.

- Carefully read through this operating manual and observe all of the safety instructions and warnings notices.
- Ensure that every user has carefully read the operating manual before using the device for the first time.
- Store this operating manual so that it is accessible to all users at all times.
- Pass the operating manual on to any subsequent owner.

The current version of this operating manual can also be found at:
[https://heidolph-instruments.com/en/service/downloads/operation-manuals#Rotary Evaporators \(Hei-VAP Expert, Hei-VAP Ultimate\)](https://heidolph-instruments.com/en/service/downloads/operation-manuals#Rotary-Evaporators-(Hei-VAP-Expert,-Hei-VAP-Ultimate))

In addition, also observe the local regulations at the installation site.

➤ Symbols and keywords

Standardized symbols and signal words are used in this operating manual to warn against any dangers and provide important instructions. These instructions must be strictly observed to avoid accidents and damage.

The following terms and basic symbols are used:

Symbol	Additional signal words / explanation
Warning signs 	<p>The yellow triangle indicates hazardous situations. It is used in combination with the following signal words:</p> <p>DANGER: Indicates a hazardous situation which, if not avoided, will result in serious injury or death.</p> <p>WARNING: Indicates a hazardous situation which, if not avoided, may result in serious injury or death.</p> <p>CAUTION: Indicates a hazardous situation which, if not avoided, may result in property damage and minor or moderate injury.</p>
Prohibitory signs 	<p>Prohibited:</p> <p>The red circle indicates a situation that should be avoided under all circumstances and which, if not avoided, may result in serious injury or death.</p>
Mandatory signs 	<p>Must be observed:</p> <p>The blue circle indicates important information. Please observe this information to avoid any property damage.</p>

Other symbols used:

Symbol	Description
→	Handling instruction, action required
✓	Result of action
▪	List of information
▪	
▪	
a.	List of variants
b.	
c.	

➤ General safety instructions

EU Declaration of Conformity

This device complies with the following EC-Directives*:

- 2006/42/EC Machinery Directive
- 2014/30/EU Electromagnetic Compatibility Directive

* See also the attached "EU Declaration of Conformity".

This device was tested according to the following Directives:

- UL 61010-1 :2012/R:2016-04
CAN/CSA-C22.2 No. 61010-1:2012/U2:2016-04
- UL 61010-2-010:2015
CAN/CSA-C22.2 NO. 61010-2-010:2015

The device has been manufactured according to state-of-the-art technology and in compliance with recognized safety regulations. However, risks may still arise during installation, operation and maintenance.

→ Please ensure that the operating manual is available at all times.

The device may only be used under the following conditions:

- Only operate the device if it is in a technically perfect condition.
- Only operate the device if you are properly aware of the risks and required safety measures.
- Only operate the device in accordance with the instructions given in this operating manual.
- If any information is ambiguous or missing, ask your superior or contact the manufacturer.
- Do not operate anything on the device without authorization.
- Only use the device in accordance with its intended use.

Intended use

The device may only be operated by authorized personnel.

The device is suitable for the following uses:

- Distillation to separate substances
- Concentration
- Crystallization
- Powder drying

The device may be operated in research laboratories, other laboratories and production facilities in the following sectors:

- Chemistry
- Pharmacy
- Biology
- Environmental analytics
- Basic research
- Similar research laboratories

Improper use / Foreseeable misuse

Any use which deviates from the device's intended use is considered to be improper. The manufacturer is not liable for damage that occurs as a result. The risk is borne by the operator alone.

In case of application in the medical and food sector, the operator is solely responsible for observing the regulations and standards pertaining in each case.

Installation / electrical safety

- The device may only be connected up if the available power supply voltage corresponds to the information stated on the rating plate for the device.
- The power supply connection must be easily accessible at all times.
- An RCD circuit breaker (residual current circuit breaker) must be fitted to the electrical system in the building so that the device is disconnected from the power circuit in the event of a fault.
- Repairs may only be performed by electricians authorized by Heidolph Instruments.
- Never operate the device with a damaged power cord.
- Always turn the device OFF and disconnect the power cord before carrying out any maintenance or repair work.

Personnel training

- The device may only be operated by trained personnel.
- The device may only be operated by persons who have been instructed and supervised in its proper use by trained specialist personnel.
- The device may only be operated by specialist personnel who are the legal minimum age or above.
- Other persons may only work on the device under the constant supervision of experienced and trained specialist personnel.
- This operating manual must be read and understood by all persons working with the device.
- Personnel must receive safety training that ensures responsible and safe work practices.

➤ Operator duties

Installation site

- The device must be installed in a suitable location.
- The device must be installed on a firm, stable and temperature-resistant surface.
- Ensure that the device and all of its components are easily accessible at all times.
- Maintain a sufficient safety clearance to, in particular, any moving and/or hot equipment components.
- It is not permitted to place or store any objects such as accessories, tools or chemicals within this distance area during operation.
- All screw connections must be securely tightened.
- It is not permitted to operate the device near to highly flammable or explosive substances.
- Operate the device in conjunction with a closed extractor hood if working with potentially harmful media (see DIN EN 14175 and DIN 12924).

Safety Instructions

- The device is designed for indoor use only under the following ambient conditions:

Ambient temperature	5 - 31 °C at 80% relative humidity 32 - 40 °C decreasing linearly to a maximum 50% relative humidity
Installation altitude	0 - 2,000 m above sea level
Contamination level	2
Overvoltage category	II
Permissible supply deviations	± 10 %
- If the device is operated in corrosive atmospheres, the service life of the device will decrease depending on the concentration, duration and frequency of the exposure to the corrosive atmosphere, e.g. concentrated hydrochloric acid (HCl)

Modifications to the device

- It is not permitted to make any unauthorized modifications or changes to the device.
- Do not attach or install any parts that have not been approved by the manufacturer.
- Unauthorized modifications or changes will void the EC Declaration of Conformity for the device and operation of the device will no longer be permitted.
- The manufacturer is not liable for any damage, dangers or injuries that result from unauthorized modifications and changes or due to the non-observance of the instructions in this manual.

Employee safety

- Ensure that the device is only operated by qualified specialist personnel and trained employees.
- Comply with the following instructions to avoid any personal injuries and property damage:
 - Laboratory regulations
 - Accident prevention regulations
 - Hazardous Substance Act
 - Other generally accepted occupational health and safety regulations
 - Local regulations

➤ Safety during use

- Wear the appropriate clothing when working on the device (protective clothing, safety glasses and, if necessary, safety gloves).
- Do not use the device in potentially explosive areas. The device is not protected against explosion. There is no explosion or ATEX protection available.
- Do not carry out work with exposed flames in the vicinity of the device (risk of explosion).
- Do not operate or assemble devices in the vicinity which are emission or radiation sources (electromagnetic waves) for the frequency range ($3 \cdot 10^{11}$ Hz to $3 \cdot 10^{15}$ Hz).
- Do not operate or assemble appliances in the vicinity of the device which constitute emission or radiation sources for ionizing radiation or in the ultrasonic range.
- Do not operate the device where adiabatic compression or shock waves might occur (shock wave combustion).

Safety Instructions

- Do not use materials which pose a potential risk of uncontrolled energy release which might result in pressure increase (exothermic reaction; spontaneous combustion of dusts).
- Do not work with hard, brittle materials (such as stones, soil samples, etc.) which could result in the destruction of the evaporating flask.
- Use only heating bath fluids which guarantee suitable heat transfer.
- Do not operate the device with excess pressure.
- Limit the pressure difference for glass components to a maximum of 2 bar.
- Do not increase the maximum excess pressure of the cooling fluid above 2 bar.
- Wear safety gloves and safety glasses when working with coolants in very low temperature ranges, such as with dry ice.
- Do not increase the flow speed of 1 m/s while siphoning liquids in with combustible materials (electrostatic charge; risk of ignition).
- Eliminate the production of gases classified in explosion group IIC during distillation by materials or due to chemical reactions, such as hydrogen.
- Do not use distillation material where the distillation residue might be explosive.
- Only connect clean ground joints.
- Attach ground joints securely in order to connect them leak-tight.
- Attach all tube connections securely to ensure safe operation.
- Connect all cables and tubes without kinks and locate them outside the operating and danger zone.
- Avoid putting pressure on the display when you are not operating the device.
- Do not spill liquids on the device or any parts of it.
- Wipe up any accidentally liquids spilled on the device immediately.
- Correct errors immediately.
- Do not use abrasive material to clean the glass surface. Wipe only with a damp cloth.
- Always switch the device OFF after use.

Disposal

- Check the device components for hazardous substances and solvents.
- Clean all components before disposal.
- Dispose of the device in accordance with the relevant national regulations.
- Dispose of the packaging material in accordance with the appropriate national regulations.

Caution: **Possible environmental damage!**

The device contains a CR2032 lithium button cell battery with the as per IEC-2. Batteries must be disposed of separately pursuant to the European Batteries Directive.

When disposing of the device, make sure that the waste management facility takes part in the joint return system for batteries or has specifically established and practices an approved return system according to the Battery Directive 2012/56/EU.

➤ Device overview

Overall view Hei-VAP Expert, Hei-VAP Expert Control

Overall view: Hei-VAP Ultimate, Hei-VAP Ultimate Control

Device Description

Back side of the device (control variants): valve-controlled pumps / house vacuum:


Vacuum manifold with components for vacuum control

Ventilation valve

Vacuum sensor

Vacuum valve

Control box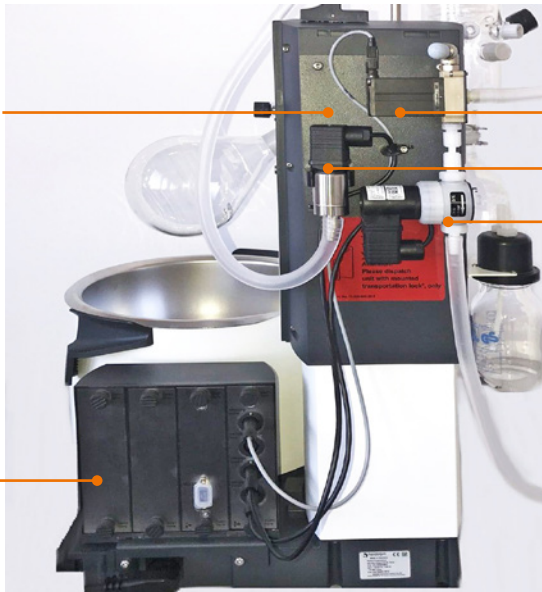


Back side of the device (control variants): speed-controlled pumps:

Vacuum manifold with components for vacuum control

Ventilation valve


Vacuum sensor

Control box

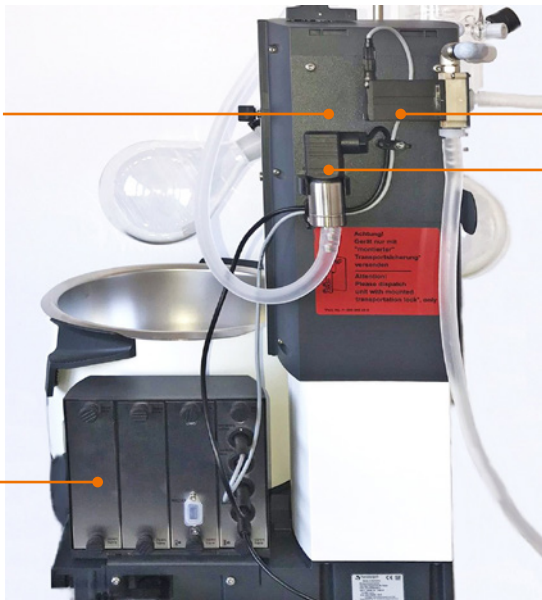Control box (for control variants only)

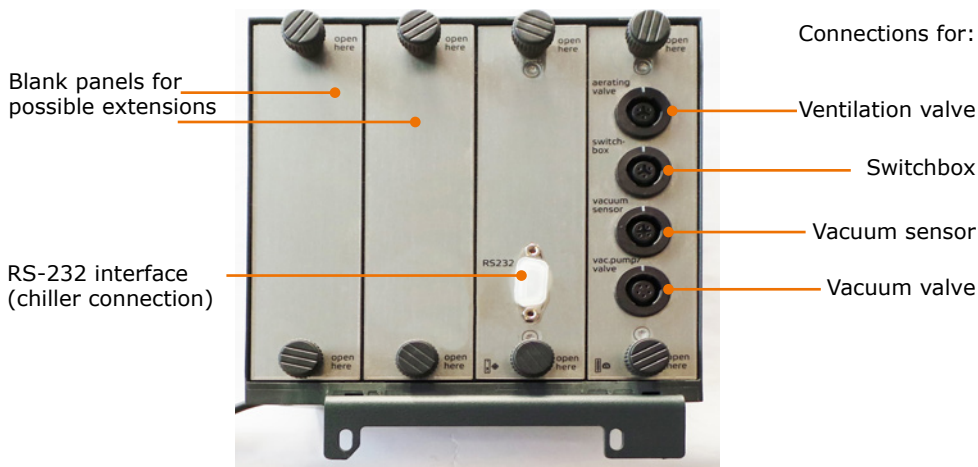Control panel: Hei-VAP Expert, Hei-VAP Expert Control

(Example display start page) The control panel is touch-sensitive

* Configuration of rotary knob assignments is possible.

Control panel: Hei-VAP Ultimate, Hei-VAP Ultimate Control

(Example display start page) The control panel is touch-sensitive

* Configuration of rotary knob assignments is possible.

➤ Setting up the device

Warning:
Risk of slipping for the device!

Due to an unsuitable surface or as a result of vibrations generated during the operation the device may tilt or start moving and slide off the table top.

The installation surface must be smooth, level, clean and temperature resistant. Avoid shaking and impact forces on the device as well as at the installation surface.

- Set up the device on a solid, smooth, level and temperature-resistant surface.
- Clean installation surface and feet of the device with a damp cloth and ethanol regularly.

➤ Connecting/disconnecting the power cord

Connecting the power cord

The power cord has a three-pole connector and a recess on its bottom side.


Power cord connector

Power cord

The device socket is located on the back side of the device.

- Insert the power cord connector into the device socket with the recess pointing downwards.
- ✓ The connector has been securely inserted when it has been pushed in up to the stop.
- Now insert the power cord plug into the power socket.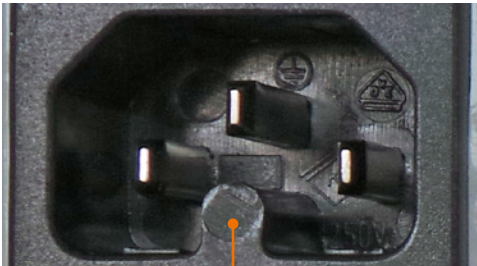

Device socket

Disconnecting the power cord

- First, remove the plug from the power socket.
- Then disconnect the power cord connector from the device.

➤ Remove transportation lock

Caution:
Possible damage to the device (motor lift variants only)

Switching the device on before the heating bath and the control panel are connected can cause malfunctioning of the electronics the moment the transportation lock is removed.

Prior to switching on the device for the first time, remove the transportation lock exactly in the order given in the instructions (see chapter "Assembly", section "Transportation lock").

➤ Switching the device on and off

The ON/OFF switch is located at the lower left side of the device.

Switching the device on

- ➔ Press the ON/OFF switch on the left side.
- ✓ The currently set values will appear on the display.
- ✓ The device is switched on.

ON/OFF switch: OFF

Switching the device off

- ➔ Press the ON/OFF switch on the left side.
- ✓ The display will go blank.
- ✓ The device is switched off.

➤ Heating bath and flask lift

The heating bath can be filled with different heat transfer fluids, e.g.,

20° C to 90° C	tap water
20° C to 170° C	polyethylene glycol (e.g. Budde heating bath fluid)*
150° C to 210° C	silicone oil with low viscosity (40 cP), flashpoint > 285 °C e. g. Wacker AP 200 silicone oil)*

* see chapter "Spare Parts and Accessories, Accessories".

Danger! **Risk of explosion!**

Heating bath fluid oil with a flash point < 285 °C could catch fire at high heating bath temperatures and lead to uncontrolled thermal reactions and possible explosion.

- ➔ Use only oil with a flash point > 285 °C.

The markings in the heating bath indicate the minimum and maximum liquid levels for the immersed evaporating flask.

The evaporating flask is moved up and down using a lift and is immersed in the heating bath in this manner.

When the lift moves downward, the evaporating flask displaces the already filled in water or oil.

To prevent the heating bath from overflowing:

- Prior to filling, move the lift with the mounted evaporating flask downward.
- Only then fill the water or oil up to at most the upper mark of the heating bath container.

1. Flask lift

Immersing the flask lift

Immerse the flask in the heating bath only so far that it does not touch the edge or the bottom.

For adjusting the immersion depth and flask angle, see Chapter "Assembly", Section "Evaporating Flask: Angle and immersion depth".

Immersing the flask: Manual lift

- Press the manual lift lever downward.
- While pressed, move the manual lift lever to the left for lowering or to the right for raising.
- Release the manual lift lever at the desired position.
- ✓ The manual lift locks in the corresponding position.


Manual lift lever: Lift position top

Immersing the flask: Motor lift

- ✓ The device is switched on.
- Press the up ▲ or down ▼ lift button.
- Release the lift button at the desired flask position.
- ✓ The motor lift will stop.

You can also set that the flask is automatically lowered at the start of the processes (see Section "Settings").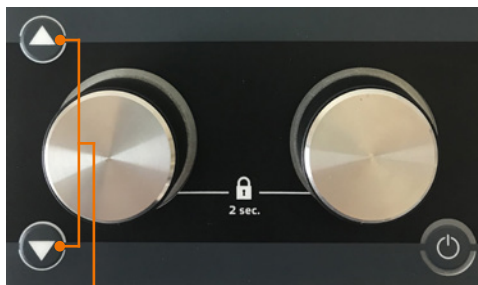

Lift buttons: up, down

2. Heating bath

Filling the heating bath

- ✓ The evaporating flask is immersed in the heating bath at the correct position.
- Fill the heating bath with the heating fluid at most up the upper mark.
- ✓ The liquid level is within the minimum and the maximum mark.

Connecting the heating bath

Warning: **Risk of crushing! Risk of burns!**

If the device is switched on, the heating or rotation can be started unintentionally. This may injure and burn your hands.

Connect the heating bath to the device base only when the device is switched off.

Start-up

The heating bath plug is located on the rear right of the heating bath.

- ✓ The device is switched off.
- ✓ The heating bath is locked in place on the base plate.

Heating bath connector socket

Heating bath plug

- Insert the 7-pin connector socket with the recess pointing upward into the heating bath plug.
- Tighten the union nut clockwise to the right.

Union nut

Heating bath plug

Recess

Feather

Emptying the heating bath

Warning: **Risk of scalding! Risk of slipping!**

When emptying the heating bath, you may burn yourself or slip on spilled fluid.

Before emptying heating bath:

- Wear protective clothing.
- Switch the device off.
- Wait until the heating bath fluid has cooled below 50 °C.
- Disconnect device base and heating bath.
- Remove evaporating flask.
- Lift the heating bath only at the insulated handles.

- ✓ The device is switched off.
- ✓ The evaporating flask has been removed.
- Turn the union nut on the heating bath connecting cable to the left in counter-clockwise direction until it is released.
- Disconnect the connector socket contact from the heating bath.
- Put your thumbs on the ribbed surfaces of the handles and then clasp them completely.
- Carefully lift the heating bath at both handles and empty it.

Protective lip

Heating bath handle


➤ Basic Settings

When the device is switched on for the first time, you will be asked to make several basic settings.

Setting the language

- ✓ A list of selectable languages is displayed.
- Swipe down on the screen to see more languages.
- Tap the desired language.
- ✓ The selected language is highlighted in orange.
- Confirm with
- ✓ The window for setting the formats is displayed.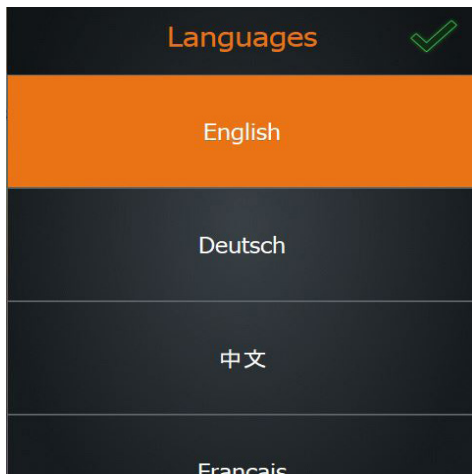

Determine formats and units


- Tap the desired format or the desired unit for:
 - Date
 - Clock
 - Decimal numbers
 - Heating bath fluid
 - Temperature
 - Pressure
- ✓ The corresponding selection is highlighted in orange.
- Confirm with
- ✓ The window for setting date and time is displayed.


Start-up

Setting date and time

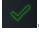- Tap the pair of digits for the date or the time which you want to change.
- Tap the number pad to change the corresponding setting highlighted in orange.
- Confirm with .

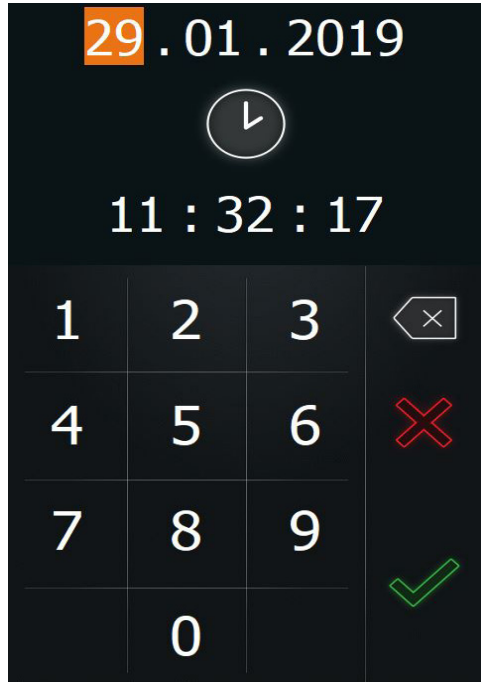





➤ Menu navigation

Overview of the menu tree

From the start page you can switch to the two menus "Basic Settings" and "Operating Modes")

Menu "Basic Settings"

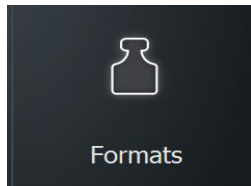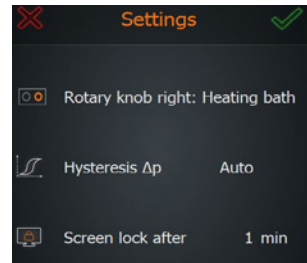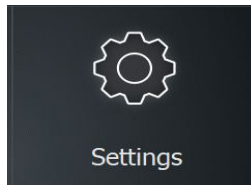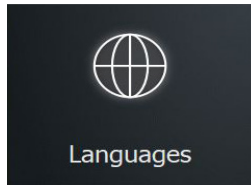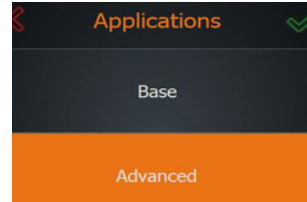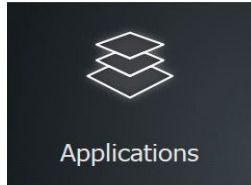
Start-up

Menu "Operating Modes" (only visible in "Applications extended"):

Favorites mode




3 menu pages

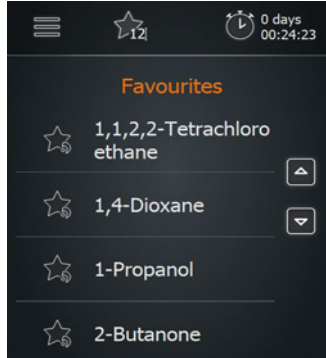Ramp mode

5 menu pages

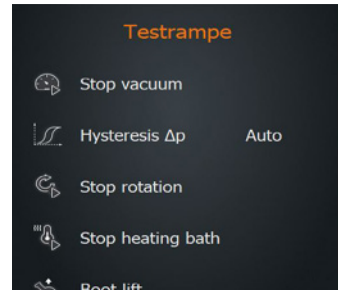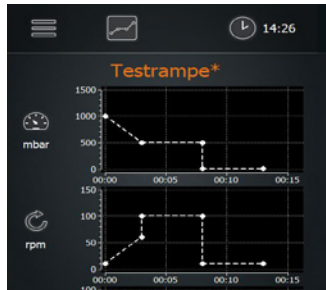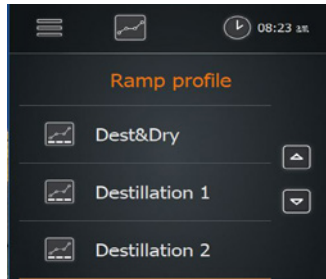AUTOaccurate mode

2 menu pages

Operation of the menu

All numbers and symbols are touch-sensitive with the exception of the temperature display for the condenser.


- The symbols of the menu bars direct you to more submenus or activate processes.
- The numbers on the right edge of the display show the selectable process parameters which can be activated and changed by tapping.
- Active elements are highlighted in color in the schematic drawing of the device. The colors are defined as follows:

White	indication
Orange	<ul style="list-style-type: none"> ▪ either: function activated ▪ or: setpoint has just been changed via the rotary knob
Grey	the indicated function is deactivated (cannot be used at the moment, e.g. because accessory is not attached)

Hei-VAP Expert and Hei-VAP Ultimate have the same menu navigation. In essence, only the arrangement of the control elements is different:

Hei-VAP Expert

The menu bars are located on the left and right edge of the control panel.

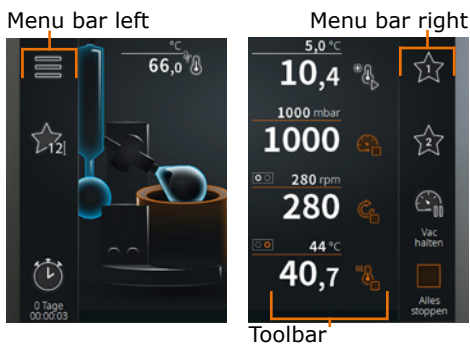Hei-VAP Ultimate

The menu bars are located on the top and bottom edge of the control panel.

Toolbar: Symbols

- In the toolbar, tap a symbol (such as "Rotation").
- ✓ The symbol changes its color to orange .
- ✓ The function starts.


The display varies depending on the chosen operating mode, the selected application and the attached peripheral equipment.

Start-up

Toolbar: Numbers


- Tap the number of a setpoint.
- The adjacent numbers menu is displayed.
- Tap the respective number to enter the desired parameter value.
- ✓ If you enter an invalid value, the valid range of values is indicated for a short time.
- Tap the green checkmark to confirm your entry.
- Tap the red cross to cancel your entry.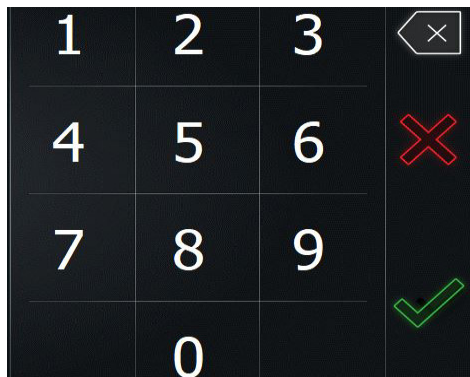

Open submenus

- Swipe the control panel from left to right (or the other way round) to switch to the next screen.

In all three operating modes, small squares appear at the bottom edge of the screen, the quantity of which indicates the page number of the submenus.

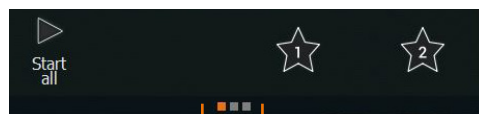Information on submenus (example)

Additionally active elements are highlighted in color in the schematic drawing of the device. For example, if the heating is switched on, the heating bath is lit orange.

If a vacuum pump or a chiller of series Hei-CHILL is attached, their setpoints and actual values are also displayed on the start page with the symbol for switching on and off. The setpoint can be changed by tapping on the corresponding page.


The values for rotation, temperature and vacuum can be adjusted both by means of the touch-sensitive control panel or by means of the rotary knob (see subsequent page).

Setpoints on the start page:

The setpoints are displayed in small print above the line. The actual values are displayed in large print below the line.

When adjusting a setpoint with the rotary knobs, this value is displayed in larger print and in orange for a short time.

Setpoint ——— Actual value

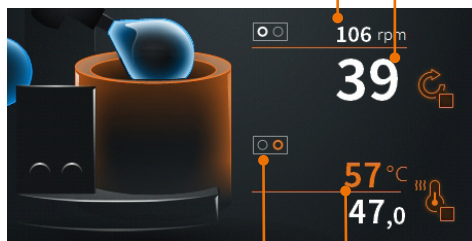Current assignment of the rotary knob

Setpoint during entry

Start-up

Adjustments via the rotary knobs:

- Turn the rotary knob to adjust the desired value.
- Press the rotary knob to activate the function.

Use the right rotary knob to adjust either the temperature or the vacuum.

* The assignment of the right rotary knob can be changed in the settings (see chapter "Start-up, Basic settings", section "Setting the process sequence parameters").

- ✓ When rotation is switched on, the left rotary knob lights up with a white illuminated ring.
- ✓ If the right rotary knob is assigned to "temperature", the rotary knob lights up with an orange illuminated ring when the heating is on.
- ✓ If the right rotary knob is assigned to "vacuum", the rotary knob lights up with a blue illuminated ring when the vacuum is on.

- ✓ Press the rotary knobs again to switch off the heating resp. the vacuum or the rotation.

You can protect the adjusted setpoints (rotation speed and temperature/vacuum) against unintentional adjustment by keeping the appropriate rotary knob pressed for approx. 2 sec.

You determine the increment by means of the rotation speed. You can adjust the desired value precisely or quickly by turning the rotary knob slowly (increment of 1), medium-fast (increments of 25) and quickly (increments of 50).


➤ Changing the basic settings

Language and formats

You can change the language and the formats in the main menu "Basic settings" at any time.


- Tap the symbol for basic settings  on the start page.
- In the menu "Basic settings", tap "Languages".
- ✓ An overview of the selectable languages is shown.
- In the menu "Basic settings", tap "Formats".
- ✓ An overview of the selectable formats is shown.

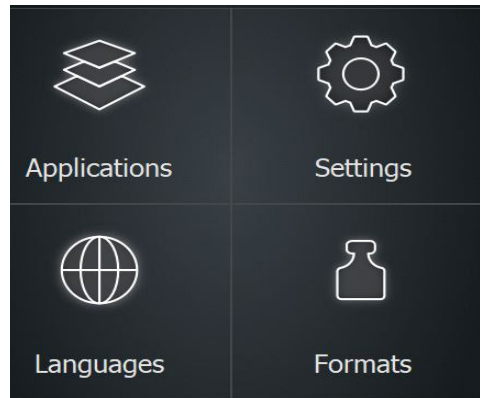Date and time

Three different time symbols can be displayed in the menu bar of the start page.


- Tap the displayed time symbol.
- ✓ Three different time symbols are shown.
- Tap to select the clock symbol.
- Keep the clock symbol pressed for approx. 3 sec.
- ✓ The window for setting date and time is displayed.

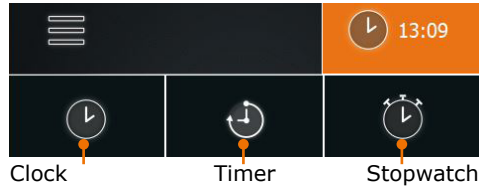Settings



- Tap the symbol "Basic settings"  on the start page.
- In the menu "Basic settings", tap "Settings".
- ✓ A list of configurable settings is displayed.
- Swipe down to see more settings.
- ✓ The following settings can be made:
 - Rotary knob assignment
 - Pressure hysteresis
 - Period to screen lock
 - Brightness of display
 - Maximum bath temperature
 - Excess pressure
 - Leakage pressure
 - Max. pump capacity
 - Pressure sensor calibration
 - Lift activation
 - Factory settings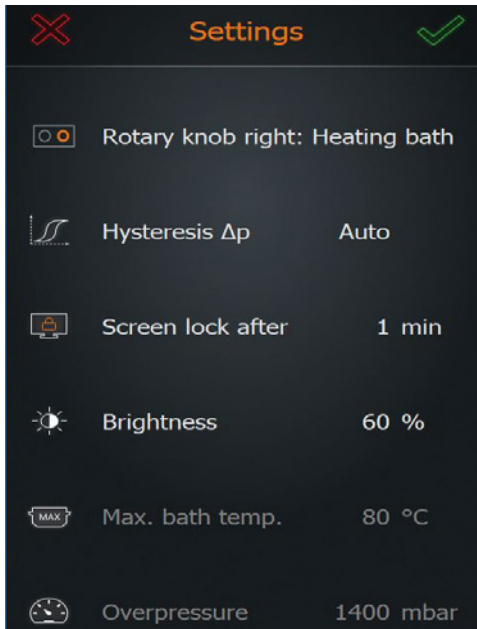

1. Assignment of the right rotary knob

It is possible to assign two different functions to the right rotary knob. You can choose if you want to use it to regulate either the temperature or the pressure (vacuum).

- Tap the menu item to change the rotary knob assignment.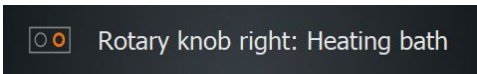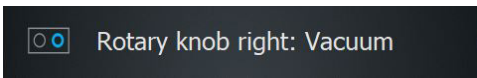


2. Pressure hysteresis

You can set by how many millibar the pressure is allowed to rise before the vacuum valve opens again.

- Tap "Hysteresis Δp ".
- Choose between automatic and manual hysteresis.

"Automatic hysteresis" contains predefined hysteresis values for several pressure ranges.

In "Manual hysteresis", you can enter a value between 1 and 50 mbar which applies uniformly over the whole vacuum pressure range.

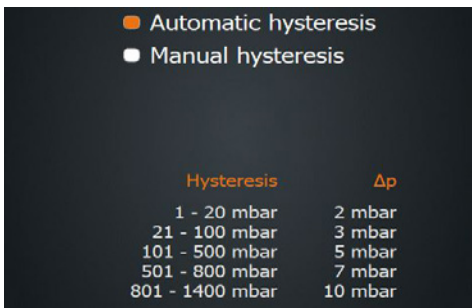
Start-up

3. Screen lock

The screen lock can be set to last between 1 to 60 min.


- Tap "Lock screen" and set the desired time.
- To unlock, tap the control panel.
- Swipe the orange ball from left to right to unlock.
- ✓ The screen is unlocked.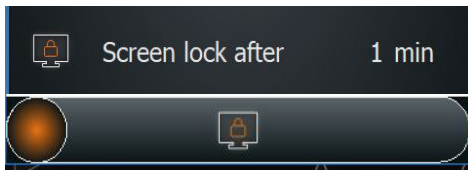

4. Brightness of display

The brightness of the screen can be adjusted from 10% to 100%.


- Tap "Brightness" and adjust the desired percentage value in increments of 10.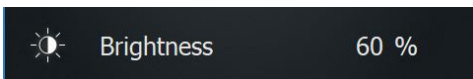

5. Maximum bath temperature

In order to protect, for example, sensitive samples, you can set a maximum temperature here. It will then not be possible to manually enter values above this temperature.

- Tap "Max. bath temp." and set the desired maximum temperature (max. 90°C for water and max. 210°C for oil).

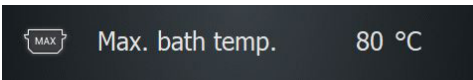When loading favorites and ramp profiles, the system checks if the set maximum bath temperature could influence the process and will display a message, if needed.


6. Excess pressure

Here you specify the vacuum pressure at which the device will automatically cancel a running process for safety reasons.

Default factory settings:	1400 mbar
Range of maximum values:	900 mbar - 1400 mbar

- Tap "Excess pressure" and set the desired maximum pressure for vacuum.

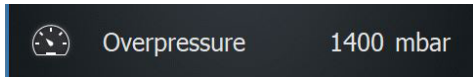- ✓ Once the set pressure is reached, vacuum, rotation and heating are stopped, the flask lift is moved up, and a corresponding warning message appears.

When loading favorites and ramp profiles, the system checks if the set maximum pressure will influence the process, and displays a message, if needed.

Start-up

7. Leakage pressure

Here, you define at which leakage pressure the device will automatically cancel a running process for safety reasons.

Default factory settings:	1399 mbar
Range of maximum values:	1 mbar - 1399 mbar

→ Tap "Leakage pressure" and set the desired maximum leakage pressure.

Leakage pressure

1399 mbar

- ✓ Once the pressure during pumping out falls below the leakage pressure, the safety function is activated.
- ✓ If the value is exceeded during pumping out - for example because of glass breakage or leakages - vacuum, rotation and heating are stopped, the flask lift is moved up, and a corresponding warning message appears.

8. Max. pump capacity

A reduction in the suction capacity leads to a higher recovery rate of the solvent.

When using a speed-controlled pump (Rotavac Vario control and Rotavac Vario tec), here you can reduce the max. pump capacity down to 20% by increments of 10% in order to prevent foaming or delay in boiling.

→ Tap "Max. pump capacity" and set the desired capacity in percent.

Max. pump performance!

100 %

If a valve-controlled pump is connected, the inscription is grayed out and it is not possible to adjust the pump capacity.

9. Calibrating the pressure sensor

For calibration, we recommend to switch on the rotation in order to ensure seal integrity of the system. If other processes are running apart from the rotation, all processes will be stopped for calibration.

When using the vacuum sensor, you can calibrate it. For this purpose, you need an calibrated external pressure gauge.

The basic calibration of the sensor is done by Heidolph Instruments. Normally, a recalibration is not required.

- ✓ The vacuum sensor is attached and connected.
 - ✓ The system is completely ventilated (no vacuum!).
 - ✓ During calibration, a calibrated external pressure gauge is attached to the evaporator system.
- Measure the atmospheric pressure on the calibrated external pressure gauge.
- Tap "Calibrate pressure sensor".
- Following the instructions of the information texts on the display:

Calibrate pressure sensor

Start-up

- Enter the atmospheric pressure determined with the external pressure gauge.
- Close all openings on the evaporator system.
- ✓ When the atmospheric pressure has been entered and confirmed, the vacuum control will start automatically. When the system has been evacuated to a low value and keeps it stable, the second calibration value can be entered.
- Wait until the attached vacuum pump cannot reduce the pressure any further and consequently the lowest possible vacuum pressure has been reached. Enter this value.
- Confirm this value.

The calibration has now been corrected on the basis of the entered values.

If calibration is cancelled, the entered values are discarded. If you restore the factory settings, the entered calibration values are also discarded and replaced by the default values.

10. Lift activation

For motor lift devices, you can set that the flask is automatically lowered into the heating bath upon starting the processes. The lift will move completely downwards. If you use a foaming inhibitor or a different individual arrangement with a bigger glass flask, you can also suppress the automatic downward movement of the lift at the start to prevent potential glass breakage.

- Tap "Start all: Lift deactivated" / "Start all: Lift activated".
- ✓ If you have selected "Start all: Lift deactivated", you have to immerse the flask by means of the lift buttons.

Start all: Lift inactive

11. Factory settings

You can restore the factory settings at any time.

When you tap "Factory settings" and confirm the security prompt, your favorites and ramp profiles are deleted!

- Tap "Factory settings".
- ✓ The factory settings or the settings of the last update are restored.


Factory settings

Applications

Changing the operating mode

- In the menu bar, tap the symbol for the current operating mode (example).
- The adjacent options menu is displayed.
- Tap the symbol to switch to the appropriate operating mode.

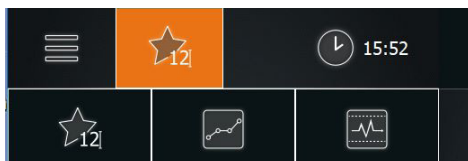








› Symbols used

Symbol	Designation	Function or display
	Menu	Open the "Basic Settings" menu
	Favorites mode	Activate favorites mode
	Ramp profile mode	Activate ramp profile mode
	Automatic mode DAA	Activate automatic mode DAA
	Clock/timer/stopwatch	Display/measure/stop time
	Vacuum off/on	Start/stop vacuum control
	Maintain/start vacuum	Stop/restart vacuum control
	Heating bath on/off	Switch the heating on/off
	Rotation on/off	Start / stop rotation
	Vapor temperature	Monitor vapor temperature
	Rotary knob right: Heating bath	Heating bath switched on
	Rotary knob left: Vacuum	Rotation switched on
	Rotary knob right: Vacuum	Vacuum switched on
	Rotary knob locked	Rotary lock
	Start all	Start all processes

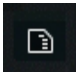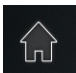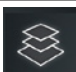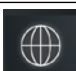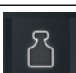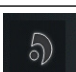Start-up

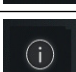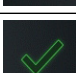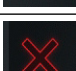Symbol	Designation	Function or display
	Stop all	Stop all processes
	USB port	USB stick is available
	SD port	SD card is available
	Start page	Return to start page
	Applications	Select basic or extended
	Settings	Open settings
	Language selection	Open language selection
	Format selection	Open format selection
	System check	Display list of all connected components
	System update	Update system
	Error	Open list of errors
	Information	Open software information
	OK	Confirm entry
	Cancel	Cancel process
	Hysteresis	Set hysteresis

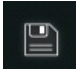
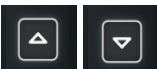
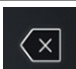




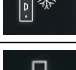
Start-up

Symbol	Designation	Function or display
	Lock screen	Activate screen lock
	Display brightness	Adjust display brightness
	Max. heating bath temperature	Indicate max. heating bath temperature
	Calibrated pressure	Calibrate the pressure sensor
	Factory settings	Restore factory settings
	Date format German/ English	Select date format
	Heating bath H ₂ O, oil	Select heating bath
	Temperature units	Select temperature units
	Error/Warning/Information	List of errors: Error/Warning/ Information
	Favorite Heidolph/Favorite 1 Heidolph/Favorite 2 Heidolph	Default favorites, Favorite 1 and Favorite 2 are on quick-selection buttons
	Favorite/Favorite 1/ Favorite 2	Own favorites, Favorite 1 and Favorite 2 are on quick-selection buttons
	Favorite 1 activated	Loaded favorites
	Ramp profile/Ramp profile 1/Ramp profile 2	Own ramp profiles, profile 1 and profile 2 are on quick-selection button
	Ramp profile 1 activated	Loaded ramp profile

Start-up
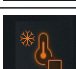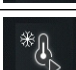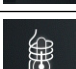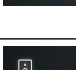
Symbol	Designation	Function or display
	Delete	Delete selection
	Save	Save selection
	Add/remove line	Add/remove ramp profile line
	Line or page up/down	Move or scroll ramp profile line up/down
	Backspace	Delete character
	Stepwise ramp	Ramp profile: Value adjustment in steps
	Linear ramp	Ramp profile: Continuous value adjustment
	Move lift down	Ramp profile settings: Stop lift
	Raise lift	Ramp profile settings: Move lift up
	Chiller on	Ramp profile setting; maintain cooling
	Chiller off	Ramp profile settings; stop chiller
	Temperature display	Current value indication of the external temperature sensor (vapor temperature sensor or AUTOaccurate sensor)
	Vacuum module	Plug-in card present in control box
	Chiller module	Plug-in card present in control box
	List on USB	Export list of favorites or ramp profiles to USB stick

➤ Process sequence


Performing a system check

After switching on the device, you can check in the overview list if all components are connected.

- Open the menu "Basic settings" with the symbol .
- In the menu "Basic settings", tap the symbol "System check" at bottom left.
- ✓ The list of attached components is displayed:
 - first, components which are not attached (tagged with a yellow triangle)
 - followed by the attached components (tagged with a green checkmark).

If an attached component is not indicated:

- check the connections and restart the device.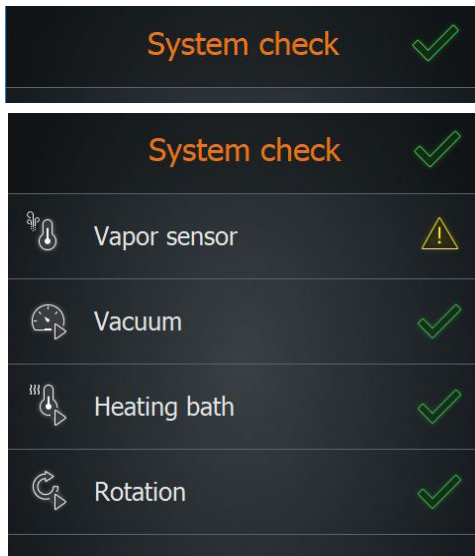


If peripheral devices, such as chillers and vacuum pump, are not attached, the corresponding indications are grayed out in all menu windows. Consequently, it is not possible to adjust their parameters.

Selecting an application

Depending on the desired process sequence, you can choose between two different applications:

- **Basic:** Here, you manually adjust the basic parameters (vacuum, rotation and temperature) before starting the evaporation.
 - **Extended:** Here, you can create and use individual favorites or ramp profiles or select the automatic process control.
- Tap "Applications" in the menu "Basic settings".
 - Select the desired application and confirm with the green checkmark.

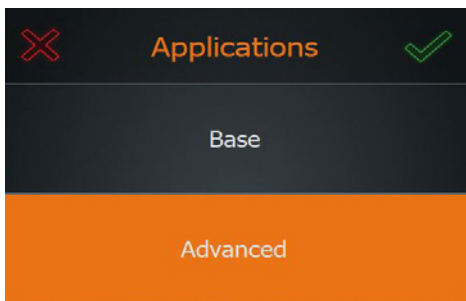
When you have selected the application "Extended", you can choose one of the following modes:

Operation

- Favorites mode (selected as a standard)
- Ramp profile mode
- Automatic mode DAA
- In the menu bar, tap the symbol of the operating mode used last.
- ✓ The symbols of the three operating modes are displayed.
- Tap the symbol of the desired operating mode.

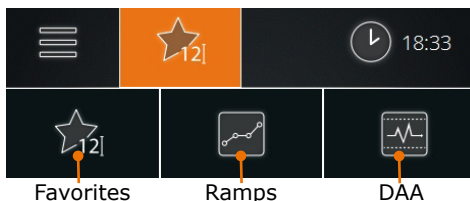To switch applications, the current processes have to be stopped.

Detailed descriptions are found in the following chapters:

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
Any chiller attached can be operated in all applications and modes.


Starting / stopping all processes

It is not only possible to start individual functions but also to start all functions at once.

- ✓ The desired settings have been made, parameters, favorites or ramp profiles have been selected.
- ✓ Ensure that no glass breakage will occur.
- In the menu bar of the current application, tap "Start all" (or "Start profile", "Start DAA").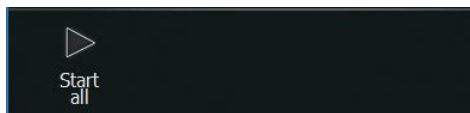

- ✓ The functions rotation, heating bath, pump (vacuum) and, if needed, stopwatch are started.
- ✓ If a chiller is attached, it will not be started.
- ✓ The lift will move down all the way if it is activated in the settings.
- ✓ With "Start DAA", the AUTOaccurate temperature is determined prior to starting the processes.
- In the menu bar of the current application, tap "Stop all" (or "Stop profile", "Stop DAA").


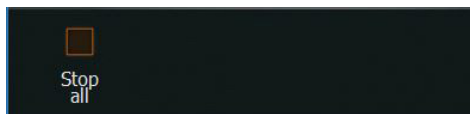
- ✓ All started processes are stopped.
- ✓ The possibly attached chiller is not stopped.
- ✓ The stopwatch stops and the value is displayed until the stopwatch is restarted with "Start all".

Recording process times

The stopwatch is automatically measuring the time of all processes which lets you keep track of the times of any process and enables you to record them if needed.

- Start any desired process.
- ✓ The time display switches to the stopwatch, and the stopwatch starts running.
- ✓ Thus, you know at any time how long the process has taken up until then.
- Stop the process.
- ✓ The time display switches to the clock.
- Switch the time display to "Stopwatch".
- ✓ On the display, the elapsed time is still shown next to the stopwatch symbol.

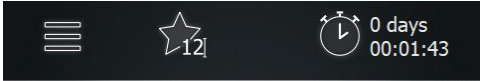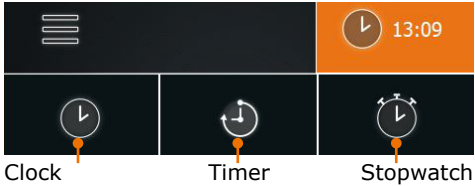The stopwatch shows the duration of the last process until a new process begins or until the device is switched off.

➤ Applications

Warning: **Possible unintentional parameter changes**

As long as the screen and the rotary knobs are activated, there is the danger that parameters are changed unintentionally.

If the device is to operate unattended, do not go away before the rotary knobs have been locked and the screen lock has been activated. (See chapter "Operation, Setting parameters manually", section "5. Protecting setpoints" as well as chapter "Start-up, Settings", section "3. Locking the screen".)

Setting parameters manually ("Application Basic")

1. Setting the rotation speed

The distillation speed can be increased by increasing the rotation speed. The shortened distillation time contributes to the thermal protection of the distillate.

However, depending on the size of the evaporating flask and the selected heating fluid, different rotation speeds are appropriate.

Rotation speed range and adjustment

The rotation speed can be set between 10 rpm and 280 rpm.

Warning: **Risk of scalding! Risk of slipping!**

At high rotation speeds, the heating bath fluid can splash and/or spill over, resulting in slippery floors and severe scalding.

If you operate the device without a safety hood:

- Always test the fluid used before a process and document the spraying behavior!
- Make this documentation available to all employees.
- Wear protective clothing near the device.
- Make sure the floor is clean.

- ➔ Test the fluid selected by you:
 - at slowly increasing temperatures
 - with your different evaporating flask sizes
 - with increasing rotation speeds
- ➔ Note the results and make them available to all employees.

The viscosity and volume of the oil change depending on the oil selected, the age of the oil and the setting of the bath temperature.

For this reason, document both the exact name/manufacturer specification and the period of use for oil and repeat the documentation at specified time intervals.

- ✓ The base device is switched on.
- ✓ The evaporating flask is immersed in the heating bath.

Operation

→ Adjust the desired rotation speed via the rotary knob.

Alternatively, you can also enter the rotation speed setpoint on the control panel:

- On the display, tap the indication of the rotation speed and enter the desired value.
- To start the rotation press the rotary knob or tap the symbol "rotation" on the display.

- ✓ The rotary knob and the symbol of the evaporating flask on the start page are lit white as long as rotation is switched on.
- Switch off rotation by pressing the rotary knob again or by tapping the symbol on the display again.

2. Setting the bath temperature

Temperature range available:

The temperature range can be set between 20 °C and 210 °C.

- Permissible temperature range when using water: 20 °C - 90 °C
- Permissible temperature range when using oil: 20 °C - 210 °C

Starting at a desired bath temperature of 75 °C, we recommend to use polyethylene glycol or silicone oil as heating fluid (see also chapter "Spare Parts and Accessories, Accessories").

Warning: **Risk of death! Risk of explosion!**

If you use an unsuitable oil as heating bath fluid, the oil can ignite and possibly cause an explosion.

→ Use only oil with a flash point > 285 °C.

Warning: **Risk of scalding!**

If you operate your device without a safety hood, severe scalding may occur at high bath temperatures.

Be sure to wear protective clothing (safety glasses, protective clothing, safety gloves) when you are near the device while a process is running.

Caution: **Possible device damage**

The heater responds very quickly. If the heating bath is empty and heating is switched on, the heating bath may develop dark discolorations.

Only switch heating on when the heating bath is sufficiently filled.

Operation

Adjusting the heating bath temperature on the display:

- ✓ The base device is switched on.
- ✓ The heating bath is filled with oil or water.
- ✓ The evaporating flask is immersed in the heating bath.
- On the display, tap the indication of the bath temperature and enter the desired value.

Adjusting the heating bath temperature with the rotary knob:

- ✓ The right rotary knob is assigned to temperature adjustment.
- Set the desired temperature with the right rotary knob.
- To start the heating, press the right rotary knob or tap the "Heating bath" symbol on the display.
- ✓ The right rotary knob and the heating bath symbol in the control panel are lit orange as long as heating is switched on.

- To switch off the heating, press the right rotary knob again or tap the "Heating bath" symbol on the display.
- ✓ The orange illuminated ring flashes if there is still residual heat.

The frequency of flashing depends on the level of the current residual heat. The flashing stops once the temperature drops below 50 °C.

If the vacuum setting is assigned to the right rotary knob, a message in the control panel will indicate if the temperature due to residual heat is above 50°C.

The display of the residual heat is only available when the device is switched on and in standby mode. The residual heat is not displayed if the device has been switched off via the main switch (ON/ OFF switch).

3. Setting the vacuum

Available pressure range:

The vacuum can be adjusted between 1 and 1400 [mbar], independent of the set safety pressure. When switching on the vacuum control, a notice about the safety pressure is displayed if needed.

- ✓ The base device is switched on.

Adjusting the vacuum on the display


- On the display, tap the pressure indication and enter the desired value.
- On the display, tap the "Start vacuum" symbol
- ✓ With valve-controlled pumps, the vacuum valve is switched on or off; with speed-controlled pumps, the rotation speed is set to zero.
- ✓ The schematic drawing of the glass flask and the condenser are lit blue.
- ✓ If the vacuum setting is assigned to the right rotary knob, the ring will be lit blue.
- Tap the symbol on the display again to switch off the vacuum.
- ✓ The blue light will be extinguished.

Operation



Setting the vacuum with the rotary knob

- ✓ The right rotary knob is assigned to the vacuum setting.
- Set the desired pressure with the right rotary knob.
- To start the vacuum, press the right rotary knob or tap the symbol "Start vacuum" on the display.
- ✓ The right rotary knob and the symbols of the glass flask/condenser are lit blue.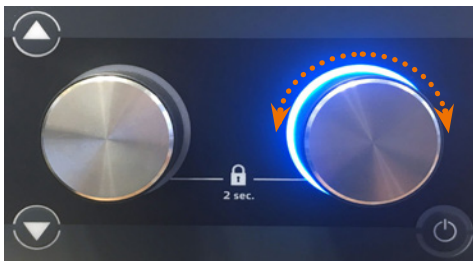

Maintaining the vacuum

You can maintain the current vacuum at any time, for example, if the setpoint has not been reached yet but you would like to stop the process for a short time.

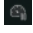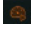- On the display, tap the "Maintain vacuum" symbol  (only available in favorites mode and in automatic mode DAA).
- ✓ The current actual value of the pressure is temporarily kept as setpoint.
- ✓ The symbol changes to the symbol for "Start vacuum".
- Tap the "Start vacuum" symbol  to stop maintaining the vacuum at the paused level.
- ✓ Now, the system will continue to the configured setpoint.

4. Display of vapor temperature / AUTOaccurate temperature

(optional accessory, see chapter "Spare Parts and Accessories").

The vapor temperature is displayed in the control panel when the vapor temperature sensor is attached.

The AUTOaccurate temperature is displayed, when the AUTOaccurate sensor is attached.


- ✓ The base device is switched on.
- ✓ You are in favorites mode or in ramp profile mode.
- ✓ When the vapor temperature sensor is attached, the vapor temperature is displayed at the top left of the control panel, next to the condenser.
- ✓ You are in automatic mode DAA.
- ✓ When the AUTOaccurate sensor is attached, the AUTOaccurate temperature is displayed at the top left of the control panel, next to the condenser.

5. Protecting setpoints

You can protect the setpoints for rotation speed and heating bath (or vacuum, depending on the button assignment) from unintentional maladjustment.

- Press the respective rotary knob for approx. 2 sec. to lock the setpoint.
- ✓ The locked value is signified by a lock.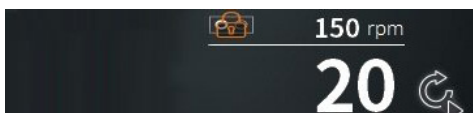

- Press the rotary knob again for approx. 2 sec. to unlock the setpoint.

Using and managing favorites (Application "Extended")

All parameters for the distillation of a substance are stored in a favorite with a fixed value. In this way, saving favorites allows quick access to different processes for your various samples.

A favorite is suited for distilling solvents and for concentrating or purifying substances.

Since the parameters are stored with a single value, the favorite has to be switched off manually at the end of the process or be stopped using the timer.


You access the favorites mode in the application "Extended". In the favorites mode, you can:

- load favorites from a list
- generate new favorites
- open a favorite directly by means of two quick-selection buttons

1. Loading favorites from a list

The factory settings already include various favorites for solvents. They are marked with a little Heidolph logo at the bottom right of the favorites symbol.

- ✓ You are in the favorites mode in the application "Extended".
- Swipe the screen from right to left to switch to the list of favorites.
- Tap the desired solvent.
- ✓ The solvent is highlighted in orange.
- Tap the "Load" symbol.
- The start page with the selected favorites is displayed, and the process can be started with the default parameters.

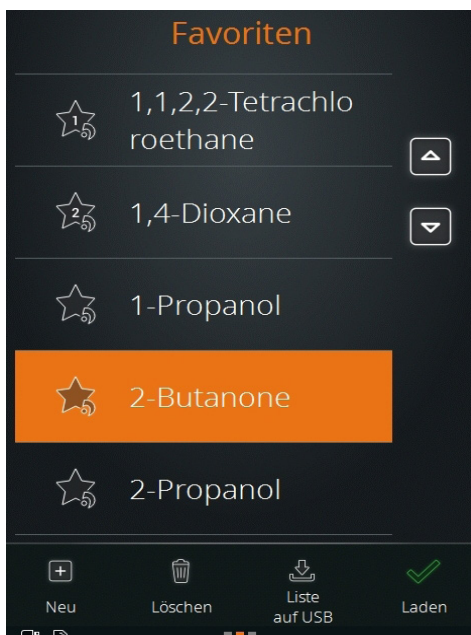It is also possible to delete and modify the Heidolph favorites. Once you have changed and saved their parameters, the favorite is stored without a Heidolph logo on the symbol.

2. Generating new favorites

You can add as many of your own favorites as you want to the existing favorites list.

✓ You are in the favorites mode in the application "Extended".

→ Swipe the screen from right to left to switch to the list of favorites.

Creating a favorite based on an existing favorite

→ Tap a solvent.

→ Swipe the screen from right to left.

✓ The parameters for pressure, hysteresis, rotation and temperature are displayed.

→ Tap the fields of the parameters you want to adapt and enter the values on the number pad.

→ Tap the "Save" symbol.

→ Enter a new name and confirm with the green checkmark.

✓ The new favorite is saved and will be inserted at the top of the favorites list.

✓ The Heidolph logo is removed from the favorites symbol.

✓ The original favorite still exists in the Heidolph list.

Creating new favorites

→ Tap the "New" symbol.

✓ Parameter suggestions are displayed for pressure, hysteresis, rotation and temperature.

→ Tap the fields of the parameters which you want to adapt and enter the values on the number pad.

→ Tap the "Save" symbol.

→ Enter a new name and confirm with the green checkmark.

✓ The new favorite is saved and added to the favorites list.

If no chiller of the Hei-CHILL series (Hei-CHILL 250-1200) or a vacuum valve or a speed-controlled vacuum pump by Heidolph Instruments is attached, the appropriate parameter is grayed out. You can still enter and save a value for them.

The screenshots illustrate the steps to create a new favorite:

- Top Screenshot:** Shows a list of favorites: Benzene, Cyclohexane, and Dichloromethane. A star icon is next to each name.
- Second Screenshot:** Shows the 'Cyclohexane' favorite selected. Parameters are displayed: Vacuum (152 mbar), Hysteresis Δp (Auto), and a 'Favourites' header.
- Third Screenshot:** Shows the 'Cyclohexane01' favorite created and highlighted in orange. The original 'Cyclohexane' is still present below it.
- Bottom Screenshot:** Shows the 'Name of favourite*' screen with a list of parameters: Vacuum (1000 mbar), Hysteresis Δp (Auto), Rotation (10 rpm), Heating bath (20 °C), and Chiller (5,0 °C). A clock icon shows '0 days 00:28:05'.


3. Assigning the quick-selection buttons


You can assign two favorites to the quick-selection buttons (asterisk symbol) on the start screen. This enables you to load these favorites directly.

Re-assigning the quick-selection buttons

- ✓ You are in the favorites mode in the application "Extended".
- ✓ The quick-selection buttons are not yet assigned to a favorite.
- Tap the symbol for quick-selection 1 or 2.
- ✓ The favorites list is displayed.
- Select a favorite and confirm with "Load".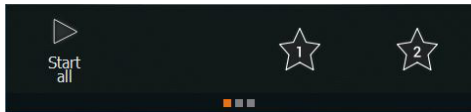

- Keep the desired quick-selection button pressed for approx. 3 sec.
- ✓ The selected favorite is stored on the quick-selection button.

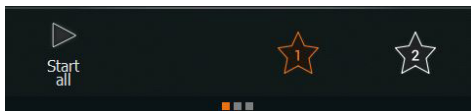
- ✓ The favorite in the list of favorites is tagged with the button assignment.


Changing the assignment of the quick-selection button

- ✓ One of the quick-selection buttons is assigned to a favorite.
- Swipe the screen from right to left to switch to the list of favorites.
- Load the desired favorite.
- Keep the quick-selection button pressed for approx. 3 sec. to store the new favorite.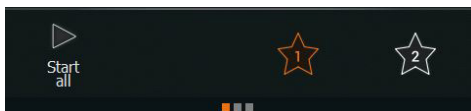

Loading a favorite with the quick-selection button

- ✓ You are in the favorites mode in the application "Extended".
- ✓ You have assigned favorites 1 and 2.
- Tap the 1 or 2 symbol.
- The symbol turns orange, the selected favorite is loaded, and you can start the process immediately.

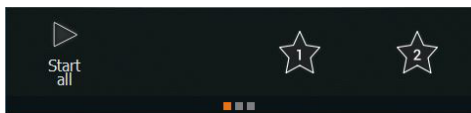
- If no favorite is currently loaded, you can jump to page 3 by keeping the symbol pressed for some time in order to save the current parameters. In doing so, the favorite is not stored on the quick-selection button.
- If a favorite is already loaded, you can save the current parameters by keeping the selected quick-selection button pressed for some time.

4. Deleting a favorite

- ✓ You are in the favorites mode in the application "Extended".
- Swipe from right to left to switch to the list of favorites.
- Mark the favorite you want to delete.
- Tap the "Delete" symbol and confirm the following prompt with the green checkmark.
- ✓ The favorite is removed from the list.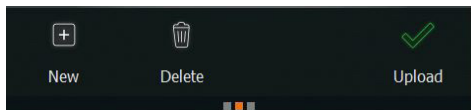

5. Exporting the favorites list

It is possible to save all favorites on a USB stick.


- ✓ You are in the favorites mode in the application "Extended".
- ✓ The USB stick is inserted in the USB port, and the USB stick symbol  appears at the bottom edge of the screen.
- Swipe the screen from right to left to switch to the list of favorites.
- Tap the "List to USB" symbol and confirm the prompt with the green checkmark.
- ✓ The list is saved.

6. Importing the favorites list

If you have saved a favorites list on a USB stick, you can import the list.

- ✓ You are in the favorites mode in the application "Extended".
- Insert the USB stick with the favorites list into the USB port.
- ✓ A message pops up asking if you want to import the favorites list.
- Confirm the prompt.
- ✓ The favorites are imported.

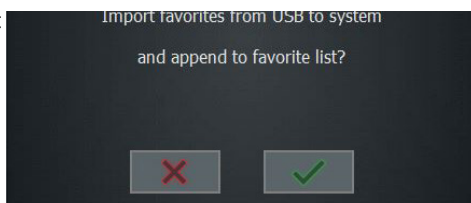Using and managing ramp profiles (Application "Extended")

A ramp profile consists of a sequence of several individual process steps.

Just like saving favorites, the saving of ramp profile allows fast access to suitable process sequences for your various samples.

A ramp profile is suited for the distillation of mixtures. At the end of a process, it is possible to determine for each parameter individually if it is to be maintained or stopped. You access the ramp profile mode in the application "Extended". In the ramp profile mode you can:

- generate new ramp profiles
- open an existing ramp profile directly by means of two quick-selection buttons


Warning: Risk of injury near the device!

When operating with ramp profiles, the rotation speed, the heating bath temperature and, if needed, the lift position are automatically changed.


Maintain a safe distance from all parts of the device when it is switched on and especially when a ramp profile is in operation.

- In order to start the ramp profile mode, tap the symbol in the menu bar depicting the operating mode that was used last.
- ✓ The symbols of the three modes are displayed.
- Tap the "ramp profile mode" symbol.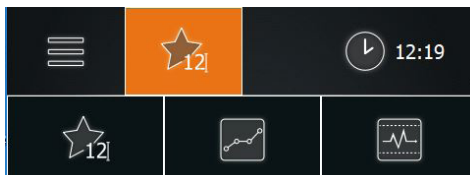

1. Generating ramp profiles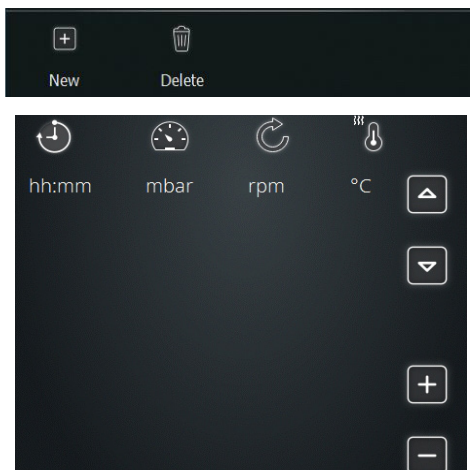

- ✓ You are in the ramp profile mode of the application "Extended".
- Swipe the screen from right to left to switch to the list of ramp profiles.
- Tap the "New" symbol.
- ✓ You switch to the ramp profile view.




Operation

Generating and adding lines

- Tap the + symbol.
 - ✓ The first two lines of your ramp profile are created.
 - ✓ The default parameters of the first line are grayed out.
 - ✓ The parameters of the second line are highlighted in orange.
 - Tap the second line.
 - ✓ The selected line is highlighted in orange.
 - ✓ Tap the time indication 00:01.
-
- ✓ The number pad is displayed.
 - Enter the desired duration in the format hh:mm.
 - Enter the desired values for all parameters this way.
 - Tap to select the step symbol  if the value has to be reached as quickly as possible.
 - Tap to select the linear symbol  if the value has to be achieved by a continuous increase.

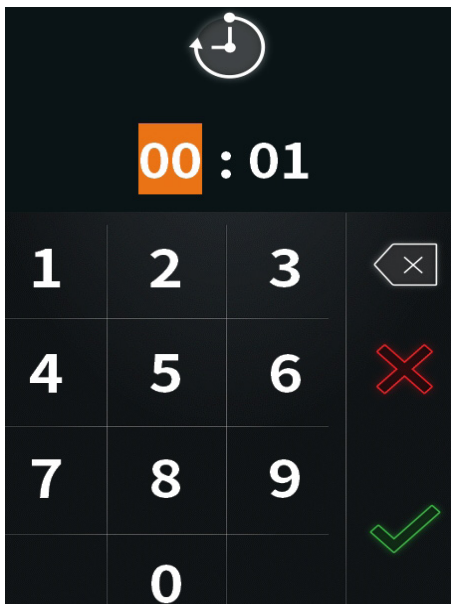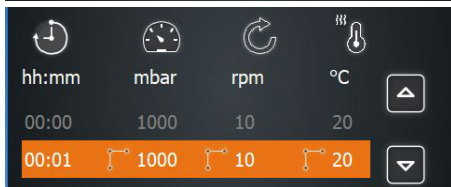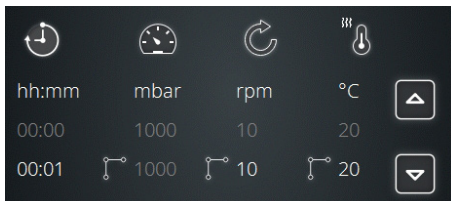Operation

- Tap the symbol + to add more lines to the profile.
- The next line of your ramp profile is created with the values of the previous step.

Changing the order of the lines


- To change the order of the lines, tap a line and use the arrows to alter its position in the profile.

Curve display

- Swipe the screen from right to left to switch to the curve display of the ramp profiles.
- ✓ The curves of pressure, rotation and temperature are plotted versus time.
- Tap a diagram to enlarge it and tap it again to shrink it.

Once the processes have started, the current values are also shown as an orange curve.

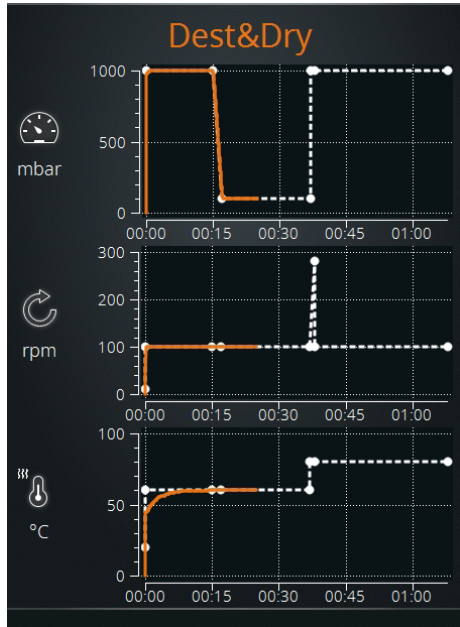
Operation


Parameters after the end of the ramp profile

- Swipe the screen from right to left to switch to the list of parameters which apply after the ramp profiles have been completed.
- Determine:
 - Maintain or stop vacuum
 - Automatic or manual hysteresis
 - Continue or stop rotation
 - Continue or stop heating bath
 - Hold or raise the lift
 - Continue or stop the chiller
- ✓ When parameters have been changed, the name of the profile gets an "*".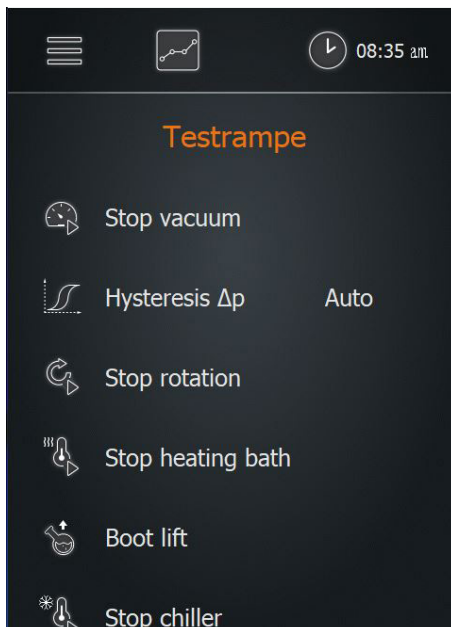

Saving the ramp profile


- Return to the ramp profile view.
- Tap the "Save" symbol.
- Enter a new name and confirm with the green checkmark.
- ✓ The new ramp profile is saved and can be loaded.

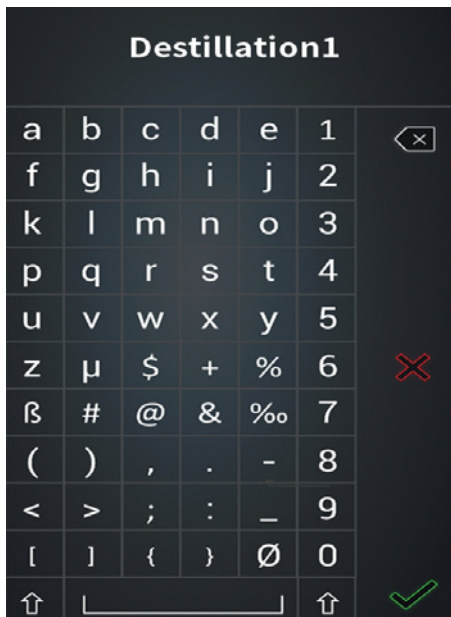
2. Loading ramp profiles from the list

- ✓ You are in the ramp profile mode of the application "Extended".
- Swipe the screen from right to left to switch to the list of ramp profiles.
- Tap the desired ramp profile.
- ✓ The selected ramp profile is highlighted in orange.
- Tap the "Load" symbol.
- ✓ The start page is displayed with the selected ramp profile, and the process can be started with the pre-set parameters.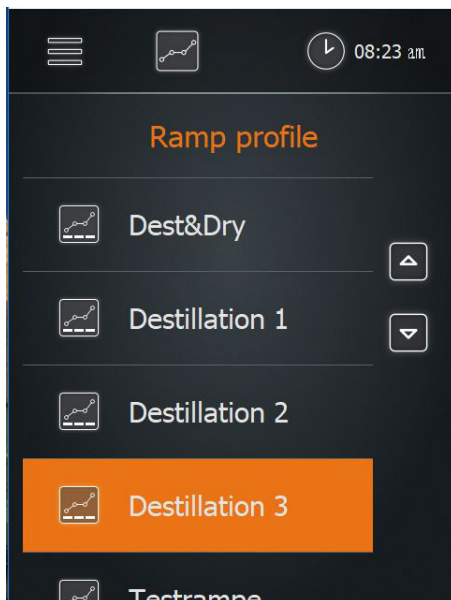


3. Loading existing ramp profiles 1 and 2

- ✓ You are in the ramp profile mode of the application "Extended".
- ✓ You assigned ramp profiles 1 and 2.
- Tap the 1 or 2 symbol (quick-selection buttons).
- The symbol turns orange, the selected ramp profile is loaded, and you can start the process immediately.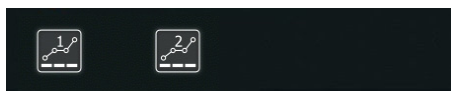

4. Assigning ramp profiles to a quick-selection button

- ✓ You are in the ramp profile mode of the application "Extended".
- ✓ You have selected and loaded a ramp profile from the list.
- Tap the 1 or 2 symbol for a couple of seconds.
- ✓ The message "Ramp 1 saved" or "Ramp 2 saved" appears, and 1 or 2 symbol is highlighted in orange.
- ✓ The favorites 1 and 2 are tagged in the list of ramp profiles.

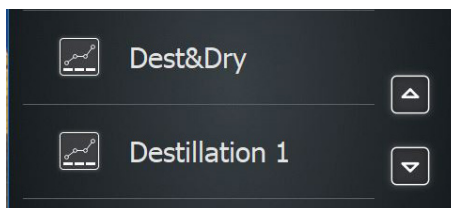You can assign two profiles to the quick-selection buttons (ramp symbol) on the start screen. This allows you to load these profiles directly.

- If the quick-selection button have not yet been assigned to a ramp profile, you can jump to page 2 by tapping on the symbol in order to load a favorite from the list.
- If a ramp profile has already been loaded, you can save the current parameters by keeping the selected quick-selection button pressed.

5. Exporting the ramp profiles to a USB stick

It is possible to save the complete list of ramp profiles on a USB stick.

- ✓ You are in the ramp profile mode of the application "Extended".
- ✓ The USB stick is inserted in the USB port.
- Swipe the screen from right to left to switch to the list of ramp profiles.
- Tap the "List to USB" symbol and confirm the prompt with OK.
- ✓ The list is saved.

6. Importing the ramp profile list

If you have saved a ramp profile list on a USB stick, you can import the list.

- ✓ You are in the ramp profile mode of the application "Extended".
- Insert the USB stick with the ramp profile list into the USB port.
- ✓ A message pops up asking if you want to import the ramp profile list.
- Confirm the prompt.
- ✓ The ramp profiles are imported.

Using automatic mode DAA (Application "Extended")

In automatic mode DAA, the vacuum is automatically controlled by the AUTOaccurate sensor. Therefore, this mode can only be used with the optional AUTOaccurate sensor.

This mode offers the following advantages:

- Only low quantities of solvent are sucked in by the vacuum pump.
- The boiling point is detected automatically and followed up for mixtures.
- Mixture distillation is possible without manual adjustments.
- Knowledge of the thermodynamic substance data is not required.
- It is not necessary to preheat the heating bath: cold starting is possible.

The system controls the vacuum in such a way that the produced vapor quantity regulates the temperature increase on the AUTOaccurate sensor. The temperature increase refers to the starting temperature during the deactivated process. It is thus necessary to have a constant temperature in the chiller prior to the start of the process. The evaporator checks autonomously for constant temperature and informs you if necessary.

Operation

Since the AUTOaccurate sensor is inserted into the condenser from above, the automatic mode DAA can only be operated with glassware sets G3, G3 XL and G6.


- ✓ You are in the automatic mode DAA of the application "Extended".
- ✓ The AUTOaccurate sensor is installed.
- Tap "Heat" if you want to heat up the bath.

→ Swipe the screen from right to left and select either "single solvent distillation" or "mixture distillation".

→ Define the "intensity", i.e. the temperature difference which will close the vacuum valve.

→ For mixture distillation, define the "final pressure".

The value of the switch point T_s is displayed after process start. T_s is the sum of the measured basic temperature and the indicated intensity.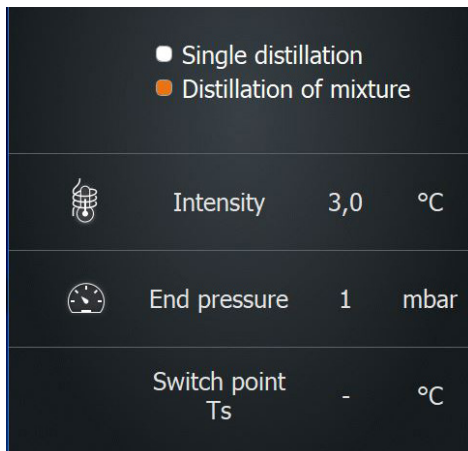


→ On the start page, tap "Start DAA".

✓ The AUTOaccurate sensor measures the temperature in the condenser for at least one minute.

✓ The message with the progress indication appears, and the stopwatch is started.

If you consider the temperature to be sufficiently stable, you can end this measuring process at any time by tapping on "Cancel". The current actual value is adopted.

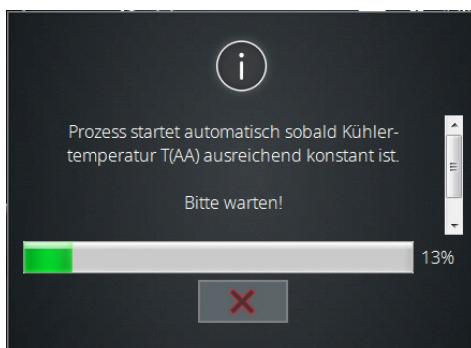- ✓ If the basic temperature remains stable ($\pm 0.5 \text{ °C}$) for 1 minute or if you have tapped on "Cancel", the DAA control will start.
- ✓ The pressure is lowered until the temperature at the AUTOaccurate sensor has risen by the specified intensity and has reached the switch point T_s . The vacuum valve will be closed, vapor condenses in the condenser.
- ✓ The pressure regulation stops and will restart when the temperature falls below the value of the switch point T_s by 0.5 °C .

You can also regulate the pressure at any time by tapping on the symbol "Maintain vacuum".

Operation

→ To stop the process, tap "Stop DAA".



If you do not tap "Stop DAA", the evaporation process of the single solvent distillation will stop automatically after the pressure regulation has been carried out three times.

The evaporation process of the mixture distillation stops upon reaching the specified final pressure.

Starting an application with a timer

Use the timer to specify the duration of a process. As soon as the process has started, the timer will show the elapsing time.

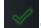- Tap the timer symbol for about two seconds and set the desired duration.
- Confirm with .

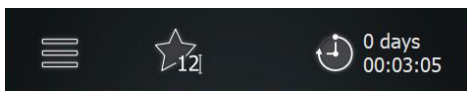The timer can be adjusted for all processes with the exception of the ramp profile mode.

Interfaces

RS-232 interface

The RS-232 interface is located on the back side of the control box.


The external chiller uses this interface to communicate with the rotary evaporator. (see chapter "Assembly", "Chiller connection")

USB port

You can use a USB stick to update the system.

- Insert the USB stick with the new software into the USB port on the left side of the control panel.
 - Tap the "Basic settings" symbol  on the start page.
 - Tap "System update" at the bottom in the "Basic settings" menu.
 - ✓ The current version is shown.
 - ✓ Tap the green checkmark to update the software version.
 - ✓ The installation is started, and the progress of the installation is displayed.
- After finishing the update, remove the USB stick from the device and cover the ports with the protective cap.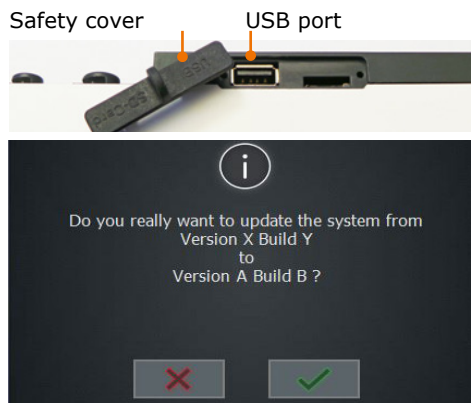

In accordance with the USB 2.0 specification, use the USB port only up to a maximum of 500 mA.

The SD card port is meant for our service technicians only.

Caution: Possible damage to the device

If the ports on the device are not covered, liquid may enter and touch parts carrying current. Penetrating liquids can lead to malfunctions of the interfaces and result in short circuits.

The interface ports on the back side of the device must be covered at all times.

Accessing system information

- In the "Basic settings" menu, tap "Information" at the bottom edge of the screen.
- Tap "Firmware" to see the version in use.
- Tap "Licenses" to open the license details.

➤ Evaporating and receiving flasks

Warning: **Risk of injury!**

If the device is switched on during assembly of the evaporating flask, it is possible for the device to start rotation due to unintentionally touching of the control panel; this could result in injuries.

Assemble and remove the evaporating flask only when the device is switched off.

Warning: **Risk of glass breakage!**


If one of the flasks is damaged, glass may break, and you may be seriously injured.

Only work with glass which has no defects. Inspect the glass regularly for damage.

Assembling and removing the evaporating flask

Assembling the evaporating flask

- ➔ Loosen the coupling of the Easy Clip and open the swiveling locking part of the Easy Clip.
- ➔ Now guide the evaporating flask over the ground joint of the vapor tube.
- ➔ Close the Easy Clip by sliding the swiveling part of the Easy Clip over the flange at the evaporating flask so that the Easy Clip engages.

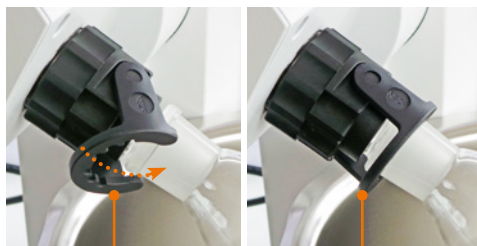Easy Clip open

Easy Clip closed

- ✓ The evaporating flask now hangs securely in the Easy Clip lock.
- ➔ Tighten the Easy Clip on the vapor tube.

Removing the evaporating flask

To remove the evaporating flask:

- Heating must be switched off
- Rotation must be switched off
- The system must be ventilated, and ambient pressure must be established
- The lift with the evaporating flask must be in highest position (see chapter "Start-Up", section "Heating bath and flask lift")
- The device must be switched off

Warning: **Risk of burns!**

Contact with the evaporating flask can cause severe burns if it has not cooled down sufficiently.

- Always allow the evaporating flask to cool down before removing it.
- The maximum temperature must not exceed 50 °C / 122 °F.
- Wear protective clothing.
- Wear safety gloves to ensure a secure grip.

- Loosen the coupling of the Easy Clip flask clamp from the vapor tube using 1-2 counterclockwise turns.
- Grip the evaporating flask with one hand and open the Easy Clip.
- Carefully remove the evaporating flask.

Assembling and removing the receiving flask

Assembling the receiving flask

- Hold the receiving flask with the opening facing up against the flange of the condenser.
- Slide the ground joint clamp over the flange on the condenser and the receiving flask.
- Carefully tighten the knurled screw of the ground joint clamp.
- ✓ The receiving flask is attached.

Removing the receiving flask

- Grip the receiving flask with one hand.
- Loosen the knurled screw of the ground joint clamp.
- Remove the ground joint clamp.

Knurled screw

Ground joint clamp

▶ **Ventilation**

The system can be ventilated manually using the ventilation cap:

- For ventilation, turn the ventilation cap half a turn or a maximum of 1 turn counterclockwise to the left.
- ✓ The ventilation speed increases with the rotation.
- Slowly turn the ventilation cap clockwise to the right.
- ✓ The air feed will close slowly.

Ventilation cap

> Cleaning

The housing and surfaces of the device can be cleaned using a damp cloth and, if required, a mild soap solution.

Warning:
Risk of electric shock!

Liquids penetrating into the device could cause the device to become electrically live if it is still connected to the power supply.

Before cleaning the device:

- Switch the device off.
- Disconnect the power cord plug.

Caution:
Damage to the device

Liquids penetrating into the device could damage the electronics.
Wipe the surfaces only with a damp cloth and not a wet one.

Caution:
Damage to surfaces

The surfaces could be damaged by improper cleaning.

Do not use the following under any circumstances:

- Chlorine bleach or cleaning agents containing chlorine
- Substances containing solvents (e.g. acetone)
- Ammonia
- Abrasive cleaning agents such as cleaning wool, scouring agents or cleaning agents with metallic components

Cleaning the heating bath

Rust spots on the bottom of the heating bath after shorter or longer use are caused by the deposition of foreign substances (iron particles or rust particles).

Rust spots can be removed by cleaning with a commercial available stainless steel polish.

Lime deposits can be removed easily by cleaning with a commercially available lime solvent, for example based on citric acid or acetic acid.

Please follow the instructions of the cleaning agent manufacturer.

Cleaning of coated glassware sets

If the cleaning temperature is too high, the coating of the glassware sets can be damaged.

Rinse coated glass parts only up to a maximum temperature of 50 °C.

➤ Maintenance

Caution:
Damage to the device

When operating with a damaged PTFE seal, the desired vacuum cannot be achieved, your sample can be contaminated, and your device can be damaged.

Never work with a damaged or worn PTFE seal. Examine the PTFE seal regularly and replace it as needed.

The device requires no maintenance. Any repairs that may be required must be performed by a specialist authorized by Heidolph Instruments.

In this case, contact your authorized Heidolph dealer or a representative of Heidolph Instruments.

➤ Malfunctions and remedies

Errors on the device	Possible cause	Remedy
<ul style="list-style-type: none"> The device cannot be switched on Display remains dark after switching on 	No power supply	Check that the plug is connected properly to the power socket and the device
	Defective fuse	Replace device fuses (see chapter "Assembly, Electrical Connection")
	ON/OFF switch off	Switch ON/OFF switch on
	ON/OFF switch defective	Contact local dealer
Device does not heat	Connecting cable to heating bath is not attached	Attach connecting cable
	Bath heating defective	Contact local dealer
	Overheating protection activated	<ul style="list-style-type: none"> Let the heating bath cool down Emptying the heating bath Reset overheating protection (see the following section "Resetting the temperature limiter")
Drive does not turn	Rotation is switched off.	Press rotation knob
	Drive defective	Contact local dealer
Motor lift does not run	Lift is at end stop	Move lift in the opposite direction
	Lift is at height stop	Change height stop
	Mechanics / Motor defective	Contact local dealer
No vacuum	ON/OFF switch at the vacuum pump off	Switch ON/OFF switch on
	Vacuum valve defective	Replace vacuum valve
Vacuum insufficient	Leak in system	<ul style="list-style-type: none"> Check seals and connections Check ground glass joints, grease if necessary
	Vacuum pump defective	Follow information from the vacuum pump manufacturer
Device switches suddenly off	Fuses defective	Replace fuse (see chapter "Assembly, Electrical Connection")
	When pressure is detected, the device is emergency stopped, see the settings for excess pressure and leakage.	<ul style="list-style-type: none"> Check system and switch device on again Calibrate the vacuum sensor

Troubleshooting

Vapor temperature is not displayed	Vapor temperature sensor not attached	<ul style="list-style-type: none"> ▪ Attach vapor temperature sensor or check connection ▪ If needed, switch the device off and on again
	Vapor temperature sensor defective	<ul style="list-style-type: none"> ▪ Inspect vapor temperature sensor for defects ▪ Contact service, if needed, replace sensor
	Temperature is outside of measuring range	Let sensor warm up or cool down
Automatic mode (DAA) cannot be selected	AUTOaccurate sensor not attached	<ul style="list-style-type: none"> ▪ Attach AUTOaccurate sensor or check connection ▪ if needed, switch the device off and on again
	AUTOaccurate sensor defective	<ul style="list-style-type: none"> ▪ Inspect AUTOaccurate sensor for defects ▪ Contact service, if needed, replace sensor
	Temperature outside of measuring range	Let sensor warm up or cool down
Control box not connected or defective	Control box not connected or defective	<ul style="list-style-type: none"> ▪ Check connection between control box and evaporator ▪ Contact service
Ventilation valve defective	Short circuit in the ventilation valve	Contact service
Vacuum valve not attached or defective	Vacuum valve is not attached or the connection between control box and vacuum valve is interrupted	<ul style="list-style-type: none"> ▪ Connect vacuum valve ▪ Check connection between control box and vacuum valve ▪ Contact service
	Short circuit in the vacuum valve	Contact service
Vacuum pump not attached or defective	Vacuum pump is not attached or the connection between control box and vacuum pump is interrupted	<ul style="list-style-type: none"> ▪ Attached vacuum pump ▪ Check connection between control box and vacuum pump

Troubleshooting

After the rotary evaporator has been switched on, it checks the attached components. The scope of functions depends on the devices attached.

Missing component	Description / Response of the device	Remedy
Control box	Control box is not attached: <ul style="list-style-type: none"> ▪ Vacuum value is not shown in display ▪ Vacuum control deactivated 	<ul style="list-style-type: none"> ▪ Attach control box and components ▪ Switch the device off and then on again
Vacuum valve or vacuum pump	Vacuum valve or vacuum pump is not attached: <ul style="list-style-type: none"> ▪ Vacuum control is deactivated 	<ul style="list-style-type: none"> ▪ Attach components ▪ Switch the device off and then on again
Vapor temperature sensor:	The sensor for determining the vapor temperature is not attached: <ul style="list-style-type: none"> ▪ Steam temperature is not displayed 	<ul style="list-style-type: none"> ▪ Attach components ▪ Switch the device off and then on again
AUTOaccurate sensor	The sensor for controlling the automatic mode DAA is not attached: <ul style="list-style-type: none"> ▪ "Automatic (DAA)" mode is deactivated 	<ul style="list-style-type: none"> ▪ Attach components ▪ Switch the device off and then on again

Troubleshooting

Analyzing the list of errors

You can view a compilation of the following warnings and errors, which appear on the display during operation, at a later time in order to analyze your process.

All in all, a maximum of 32 errors is stored. After that, older messages are overwritten by newer ones.

→ In the menu bar of the “Basic settings” menu, tap “Error list” to view information, errors and warnings.

✓ An error history is displayed showing the corresponding date and time.

→ To view details, tap the individual entries.

The following messages/errors are displayed.

Error messages on display	Cause	Remedy
Warning: Transportation lock activated! Please move lift upward and switch the device off and then on again.	The transportation lock is still activated.	<ul style="list-style-type: none"> Remove transportation lock if needed Move lift upward Switch the device off and then on again
Warning: The vapor temperature sensor is defective or not attached!	Sensor is not attached or defective	Check vapor temperature sensor and replace it if needed
Warning: The AUTOaccurate sensor is defective or not attached!	Sensor not attached or defective	Check AUTOaccurate sensor and replace it if needed
Warning from chiller: Temperature has fallen below the lower limit!	Chiller temperature too low	Increase chiller temperature
Warning from chiller: Temperature has exceeded the upper limit!	Chiller temperature too high	Reduce chiller temperature
Warning from chiller: Filling level of thermal fluid too low!	Not enough cooling fluid	Add cooling fluid
Warning: Heating bath temperature Careful! Hot!	Heating bath temperature > 50 °C	Let the heating bath cool down
Error The AUTOaccurate sensor is defective or the temperature is below the measuring range.	Defective sensor or temperature below -40 °C	<ul style="list-style-type: none"> Check AUTOaccurate sensor and replace it if needed or increase temperature
Error The AUTOaccurate sensor is defective or the temperature is above the measuring range.	Defective sensor or temperature above 173 °C	<ul style="list-style-type: none"> Check AUTOaccurate sensor and replace it if needed or reduce temperature
Error The vapor temperature sensor is defective or the temperature is below the measuring range.	Defective sensor or temperature below -40 °C	<ul style="list-style-type: none"> Check vapor temperature sensor and replace it if needed or increase temperature

Troubleshooting

Error messages on display	Cause	Remedy
<p>Error A data bus error has occurred. Switch the device off and on again</p>	<p>Error in data bus, e.g. a bus node was disconnected during operation.</p>	<ul style="list-style-type: none"> Switch the device off and on again!
<p>Error A temperature above the safety temperature was measured in the heating bath. The temperature measurement unit is defective or the medium filled in was too hot. Switch the device off and on again!</p>	<p>Heating bath temperature is too high or the temperature measurement unit is defective.</p>	<ul style="list-style-type: none"> Let the heating bath cool down Switch the device off and on again Contact service
<p>Error An implausible temperature difference was measured between the heating bath sensors. The temperature measurement unit is defective. Switch the device off and then on again</p>	<p>Error in temperature measurement of heating bath</p>	<p>Switch the device off and then on again</p>
<p>Error The temperature sensor for the heating bath control is defective (hardware monitoring unit). Switch the device off and on again</p>	<p>Heating bath control is defective, perhaps due to short circuit</p>	<ul style="list-style-type: none"> Switch the device off and on again Contact service
<p>Error The temperature sensor for the heating bath control is broken or the filled in bath fluid is too hot. Switch the device off and on again</p>	<p>Heating bath control is defective or temperature in the heating bath >217°C</p>	<ul style="list-style-type: none"> Let the heating bath cool down Switch the device off and on again Contact service
<p>Error The temperature sensor for the heating bath control has shorted or the bath fluid was added too cold. Switch the device off and on again</p>	<p>Heating bath control is defective, perhaps due to short circuit or temperature <0 °C</p>	<ul style="list-style-type: none"> Heat up heating bath Switch the device off and then on again Contact service
<p>Error The lift motor is defective or has jammed! Switch the device off and on again</p>	<p>Defective or jammed lift motor</p>	<ul style="list-style-type: none"> Switch the device off and on again Contact service
<p>Error The pressure exceeds the maximum permissible excess pressure!</p>	<p>Pressure in the vacuum system is too high</p>	<ul style="list-style-type: none"> Reduce pressure in the vacuum system
<p>Error The set safety pressure was exceeded.</p>	<p>Pressure in the vacuum system is too high</p>	<ul style="list-style-type: none"> Reduce pressure in the vacuum system If required, adjust the set maximum pressure

Troubleshooting

Error messages on display	Cause	Remedy
Error The pressure sensor is not attached or defective.	Pressure sensor not attached or defective or control box is not connected correctly.	<ul style="list-style-type: none"> ▪ Check and re-attach vacuum sensor ▪ Check connection of control box
Error Defective communication with the pressure sensor!	Hardware fault on the vacuum plug-in board	<ul style="list-style-type: none"> ▪ Contact service ▪ Replace vacuum plug-in board
Error The ventilation valve is not connected or is defective.	Ventilation valve not connected or defective or control box is not connected correctly.	<ul style="list-style-type: none"> ▪ Check and re-connect the ventilation valve ▪ Check connection of control box
Error The ventilation valve is defective.	Short circuit in the ventilation valve	<ul style="list-style-type: none"> ▪ Contact service
Error The vacuum valve is not attached or defective.	Vacuum valve not attached or defective or control box is not connected correctly.	<ul style="list-style-type: none"> ▪ Check and re-attach vacuum valve ▪ Check connection of control box
Error The vacuum valve is defective.	Short circuit in the vacuum valve	<ul style="list-style-type: none"> ▪ Contact service
Error The vacuum pump is not attached or defective.	Vacuum pump not attached or defective or control box is not connected correctly.	<ul style="list-style-type: none"> ▪ Check and re-attach vacuum pump ▪ Check connection of control box
Error Vacuum module: Connection error between base plate and vacuum plug-in board. Please check the cable connection!	Bus error of the component, e.g. due to plugging/unplugging during operation.	<ul style="list-style-type: none"> ▪ Check cable connection ▪ Switch the device off and on again!
Error Chiller module: Connection error between base plate and chiller plug-in board. Please check the cable connection!	Bus error of the component, e.g. due to plugging/unplugging during operation.	<ul style="list-style-type: none"> ▪ Check cable connection ▪ Switch the device off and on again!
Error A communication error between base plate and control panel has occurred!	Bus error of the component, e.g. due to plugging/unplugging during operation.	<ul style="list-style-type: none"> ▪ Check cable connection ▪ Switch the device off and on again!
Error The rotary drive is overloaded or defective! Switch the device off and on again!	Drive motor defective or overloaded	<ul style="list-style-type: none"> ▪ Switch the device off and on again! ▪ Contact service
Error Alarm from the chiller: Circuit board is too hot	Chiller overloaded	Switch chiller off and let it cool down

Troubleshooting

Error messages on display	Cause	Remedy
Error Alarm from the chiller: Pump is overheated	Chiller overloaded	Switch chiller off and let it cool down
Error Alarm from the chiller: Hardware fault!	Fault in the chiller hardware	<ul style="list-style-type: none"> ▪ Switch the device off and on again! ▪ Contact service
Error Update unsuccessful! Use the main power switch to switch the device off and on again!	Software update unsuccessful	Switch the device off and on again!
Error Update unsuccessful!	Software update unsuccessful	Switch the device off and on again!
Information Reset system to factory settings? The device must be restarted afterward.	Reset to factory setting opened	Switch the device off and on again
Information The process will start automatically as soon as the chiller temperature T(AA) is sufficiently constant. Please wait!	Automatic mode DAA was started	<ul style="list-style-type: none"> ▪ Wait ▪ or cancel and start without constant temperature
Information The bath temperature exceeds the set max. bath temperature. The max. bath temperature must be adjusted!	The heating bath temperature exceeds the set max. heating bath temperature.	Adjust the max. heating bath temperature
Information The pressure exceeds the max. permissible excess pressure. Please adjust the pressure!	The pressure detected exceeds the maximum excess pressure value	Reduce pressure below the excess pressure value.
Information A calibrated external pressure gauge is required. Please enter the atmospheric pressure indicated on the external pressure gauge and confirm the value with OK!	Pressure sensor calibration has started	Enter/correct the actual atmospheric pressure value
Information Calibration will stop the running processes. Do you really want to calibrate the p-sensor?	Pressure sensor calibration was started	<ul style="list-style-type: none"> ▪ Confirmation will stop all running processes. ▪ Exception: if only rotation has started, it will continue running

Troubleshooting

Error messages on display	Cause	Remedy
Information Please close all openings on the evaporator system and evacuate it. When the pressure value has become stable, enter the value indicated on the external pressure gauge and confirm it with OK!	The actual atmospheric pressure value has been entered	Enter/correct the actual value of the current minimum vacuum pressure
Information Vacuum sensor and/or AUTOaccurate sensor not available! Automatic mode (DAA) cannot be started.	Vacuum sensor and/or AUTOaccurate sensor are not attached, automatic mode (DAA) cannot be started.	Attach vacuum sensor and/or AUTOaccurate sensor
Information No ramp loaded Ramp mode cannot be started!	No ramp profile loaded	<ul style="list-style-type: none"> ▪ Load ramp profile ▪ Start ramp profile mode

➤ Resetting temperature limiter

- ➔ Switch the device off Remove the cooled evaporating flask
- ➔ Disconnect the heating bath plug and remove the heating bath.
- ➔ If necessary, empty the heating bath and turn it over.
- ➔ Use a suitable object to briefly press the red button on the bottom until it clicks audibly into place.
- ✓ Reset the temperature limiter

Temperature limiter

If you experience a malfunction that cannot be resolved using the information listed above, please inform your authorized Heidolph dealer immediately.

Electrical connections

- Fuses must **only** be replaced by a **professional electrician**.
- **Repairs** to the device must **only be carried out by a qualified electrician approved by Heidolph Instruments**. Improper repairs could result in serious dangers. Please contact your specialist dealer if the device needs to be repaired.

Warning: Risk of electric shock

The device must be connected to an earthed power socket.



The rated voltage for the device and the supply voltage must be the same. The rated voltage for the device can be found on the rating plate on the rear of the device.

The device is supplied with a grounded power cord. In case the original plug is replaced, the new plug must have a protective conductor!

If the device is still connected to power while changing the fuse, you might get in contact with live parts.

Prior to replacing the fuse ensure that the unit is switched off and pull the plug from the electric socket. Only use original spare fuses from Heidolph Instruments.

After changing a fuse check the device for safe operating condition according to IEC 61010-1.

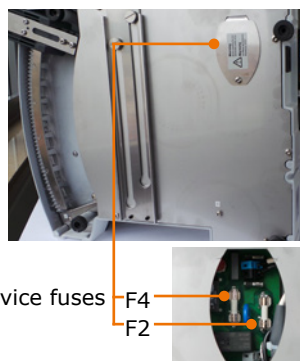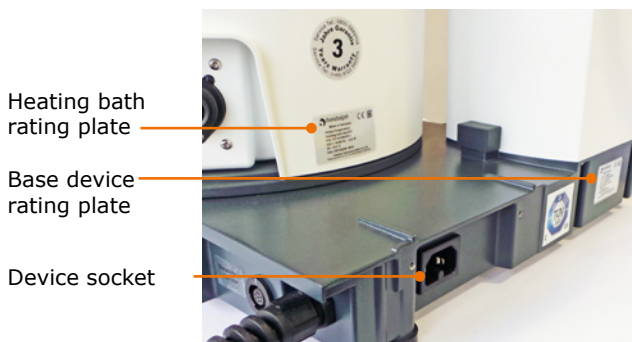Color code for electrical connections:

Europe

GREEN/YELLOW	PE: Protective conductor (Earth)
BLUE	N: Neutral conductor
BROWN	P: Phase

North America

GREEN	PE: Protective conductor (Earth)
WHITE	N: Neutral conductor
BLACK	P: Phase

The device is connected to the mains with the power cord supplied. The power connection is located on the rear of the housing.

For countries where a different plug than the standard plug supplied is used:

- The plug supplied may only be changed by a professional electrician.
- If you wish to use an adapter, ensure that it is grounded and approved in accordance with the local regulations.
- Wiring, adapter and plug must at least meet the electrical power stated on the rating plate.


Installation

Assembly parts


Base device

Control panel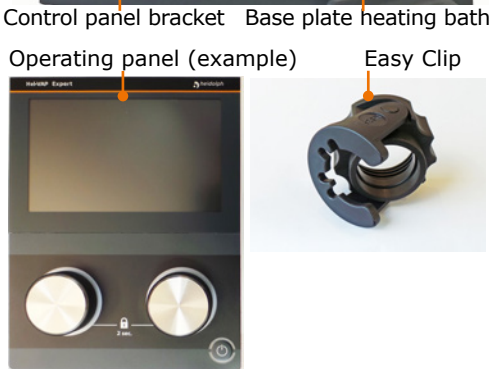

Heating bath

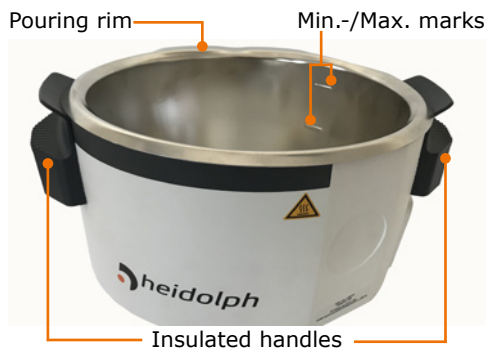
Assembly

Control box

Vacuum manifold

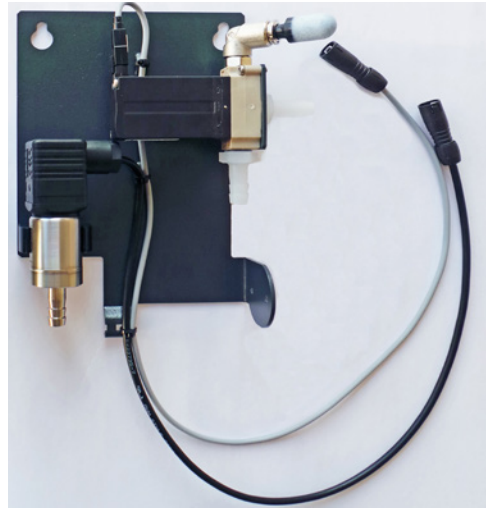Small parts for assembly

- 1x tube nipple with screw cap (for vacuum connection)
- 2x metric screws
- 2x threaded screws
- 1x torx wrench

Screws: Plastic - metric

Torx wrench

Tube nipple with screw cap

Warning:
Risk of crushing! Risk of burns!

If the device is switched on, the heating or rotation can be started unintentionally. This may injure and burn your hands.

Install the device only when the device is switched off.

Transportation lock

(Motor lift only)

The transportation lock must be removed before start-up of the device. Retain the transportation lock for a possible return.

The transportation lock consists of three M5x8 screws and a connecting plate with holes for fixing the position.

Remove transportation lock prior to start-up

- ✓ The base device is positioned on a stable, level surface.
- Connect the heating bath
- Connect the control panel.
- Switch the device on.
- Use the provided Allen key to loosen all 3 screws and remove the metal sheet of the transportation lock.
- Move the flask lift to the highest position.
- Switch the device off.
- ✓ You can now switch the device back on and put it back into operation.
- Retain all parts of the transportation lock for later use.

Transportation lock


To deactivate the additional electronic transportation lock, the device must be switched off once when the lift is in the highest position.

Heating bath

Insert the heating bath

- ✓ A sliding base plate with a raised guide rail is located on the right side of the base device.

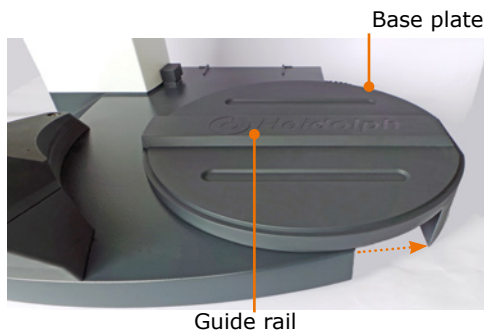
Guide rail

Assembly

- Place the heating bath with the recess to the left on the guide rail.
- Move the heating bath slightly to the right and left, so that it locks securely.

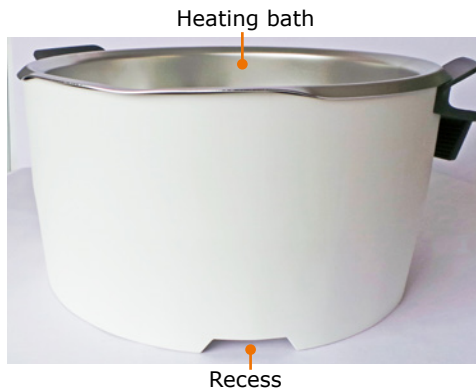Moving the heating bath

When using larger evaporating flasks or adapters between evaporating flask and vapor tube, it may be necessary to increase the distance between the heating bath and the drive.

- ✓ The base device is positioned on a stable, level surface.
- ✓ There is about 20 cm space for movement on the same base to the right of the heating bath.
- Hold the heating bath by both handles and pull it to the right to the desired position.

Caution:
Risk of crushing!

When moving the heating bath towards the device, your fingers may get caught between heating bath and device.


When moving the heating bath, always hold it at the insulated handle.

Control panel

The control panel support is located in front of the flask drive.

Control panel bracket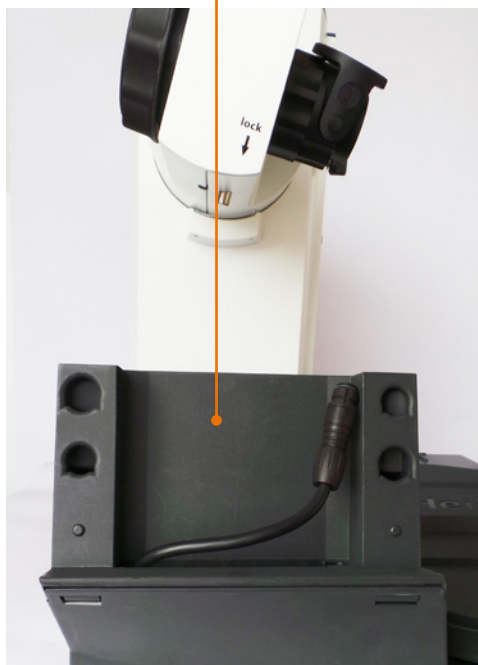

Warning: **Risk of poisoning!**


If you operate the device inside a fume hood and have to reach the control panel during a running process, you may be endangering your health when you open the fume hood.

Remove the control panel and position it outside the fume hood. (see also chapter "Scope of supply and accessory")

Mounting the control panel on the bracket

- Connect the cable from the back side of the control panel to the connection cable in the support.
- Place the control panel on the support so that the buttons on the back of the panel are seated in the recesses of the support.
- Then slide the control panel down to the stop.
- ✓ The control panel is now ready for use.

Connection cable

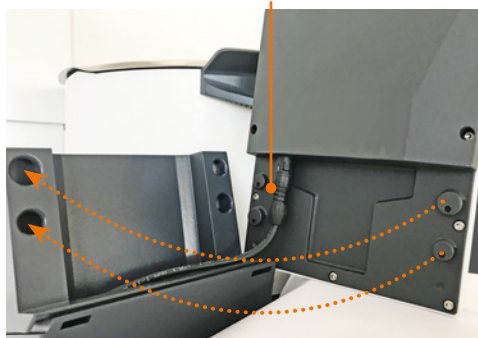Vapor tube



Assembly parts

- Vapor tube
- Easy Clip flask clamp
- Coupling sleeve
- Clamping sleeve


The Easy Clip flask clamp is pre-mounted in the NS29 size. The Easy Clip NS24 is included for smaller flasks.

When the device is delivered, the coupling for the vapor tube and the condenser are pre-assembled.

- Press and hold the locking knob below "lock" for a quarter turn.
- Unscrew the coupling of the vapor tube together with the Easy Clip to the right of the drive.
- Unscrew the coupling for the condenser to the left of the drive.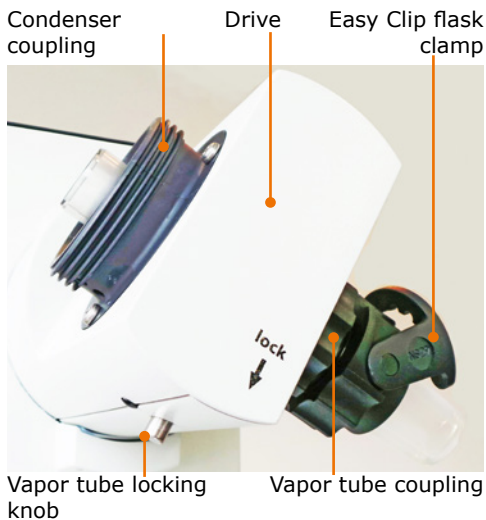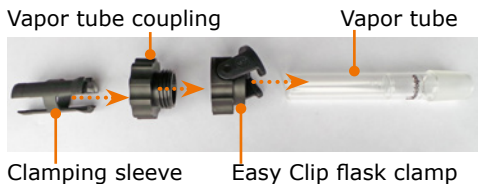

- Remove the tension spring and the PTFE seal on the condenser side.

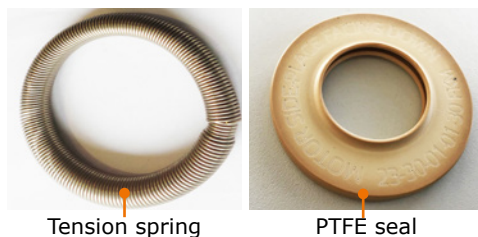Caution:
Contamination of your sample, device damage

A damaged seal can prevent reaching the desired vacuum, can contaminate your sample or damage your device.

Never work with a damaged or worn PTFE seal. Inspect the PTFE seal regularly and replace it if necessary.

Assembly

- ✓ Coupling and Easy Clip are screwed together
- Push the coupling of the vapor tube together with the Easy Clip onto the vapor tube. (Direction and sequence as shown under "Assembly parts")
- Push the clamping sleeve with the bead toward the glass onto the vapor tube until it clicks into place.
- Insert the vapor tube with the clamping sleeve first from below into the drive and screw it tight.
- Press the locking knob during the last quarter turn.

Vapor tube, Easy Clip and coupling

Condenser

Warning: **Risk of glass breakage!**

If a glass component is damaged, glass may break, and you may be seriously injured.

Only work with glass which has no defects. Inspect the glass regularly for damage.

The coated glassware sets from Heidolph Instruments reduce the risk of injury in the event of glass breakage. However, they are only suitable for a temperature range up to max. 80 °C.

PTFE seal

- Slide the PTFE seal with the mark "Motor Side" first forward onto the vapor tube.
- ✓ The PTFE seal has been pushed to the stop.

PTFE seal

Condenser: **Glassware set G1, G3 and G3 XL**

- Slide the coupling for the condenser over the chiller flange.
- Slide the tension spring over the chiller flange.
- Hand-tighten the chiller with the coupling on the drive.

Tension spring Coupling

- ✓ The chiller is mounted on the drive.

Assembly

Condenser:

Glassware set G5 dry ice cooler

- ✓ The chiller is mounted on the drive.
- Screw the ventilation cap onto the opening opposite the vapor tube.
- If you are working with inert gas, screw the inlet hose with inert gas valve onto the opening above the ventilation cap.

Inert gas valve

Ventilation cap

- Place the red seal ring in the groove at the top of the chiller flange.
- Place the white PTFE ring above it with the collar facing upwards in the chiller flange.
- Now insert the cold trap into the condenser.
- ✓ The flange is drawn in by the vacuum and the system is closed vacuum-tight by the installed seal.
- Place the black lid on top to protect your dry ice.


Assembly

Condenser: Glassware set G6 (chiller for reflux distillation)

Additional assembly parts at the vapor tube:

- PTFE seal ring
- Spacer sleeve
- Additional coupling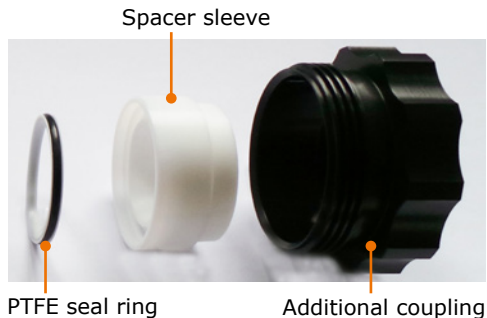

- ✓ The vapor tube is installed together with the PTFE seal.
- Install the PTFE seal ring in the spacer sleeve
- Install the spacer sleeve together with the PTFE seal ring with the narrow side first in the vapor tube on the condenser side
- Screw the additional coupling over it with the male thread pointing left.
- Tighten the distributor with the coupling including the tension spring onto the male thread.
- ✓ The distributor is installed.

Distributor


- Insert the chiller into the upper ground glass joint on the distributor.
- Screw the ventilation cap onto the opening opposite the vapor tube.

Assembly

- Turn reflux valve clockwise until the seal ring is clearly visible and white plug blocks the output to the receiving flask
- ✓ The valve is closed, and the sample is guided back to the evaporating flask.
- Turn reflux valve counterclockwise until the white plug does not block the outlet to the receiving flask.
- ✓ The valve is open, and the sample is guided into the receiving flask.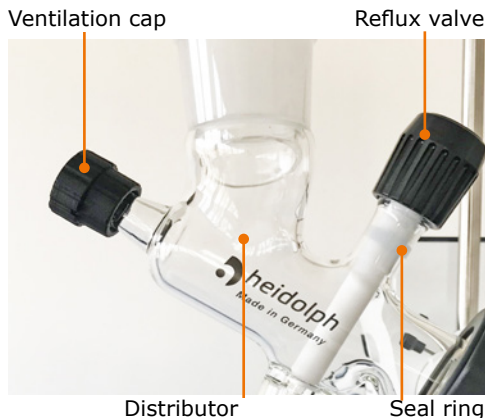

The glassware set is no longer vacuum-tight if the reflux valve is opened too far. Open the valve only so far that the white seal ring underneath the knurled screw remains visible about 0.5 cm wide.

Support system for vertical condenser: Glassware sets G1, G3, G3 XL and G6

Assembly parts:




- Support rod
- Cross clamp with condenser clamp

A hole is located on the left side of the device for receiving the support rod.

- ✓ First mount the condenser and then slide the condenser clamp over the condenser.

- Insert the support rod with the flattened side downward into the hole.
- Secure the support rod with the knurled screw.

- Slide the cross clamp over the support rod.
- Use the knurled screw to secure the cross clamp on the support rod.
- Secure the condenser clamp with the knurled screw around the condenser.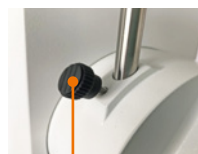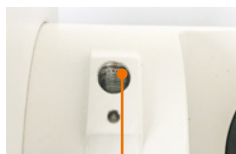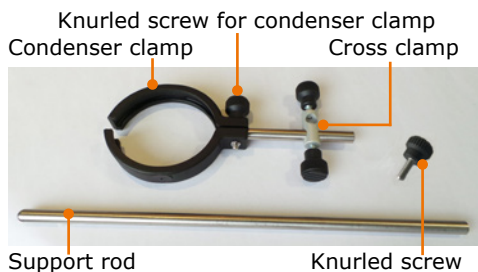


Support system for vertical condenser: Glassware set G5

Assembly parts:

- Support rod
- Cross clamp with tension strap

- Mount the support system like the support for G3, G3 XL and G6.
- Use the tension strap to secure the condenser.

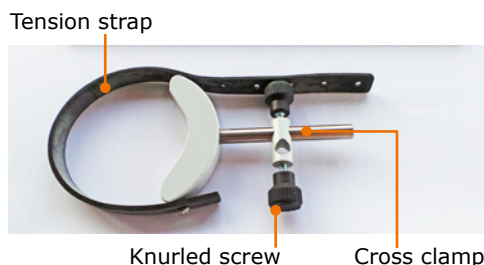
Evaporating flask: Angle and immersion depth

Flask angle

The evaporating flask angle can be adjusted by turning the drive.

- ✓ The rotation is switched off.
- ✓ The evaporating flask is mounted (see chapter "Operation", section "Mounting the evaporating flask").
- Slide the locking mechanism on lift column to the right and hold it against the spring force.
- Turn the condenser carefully until the connected evaporating flask has the correct angle. Ensure that the flask can rotate freely and does not touch anything.
- Release the locking mechanism.
- If necessary, turn the condenser slightly until the drive unit locks in place.

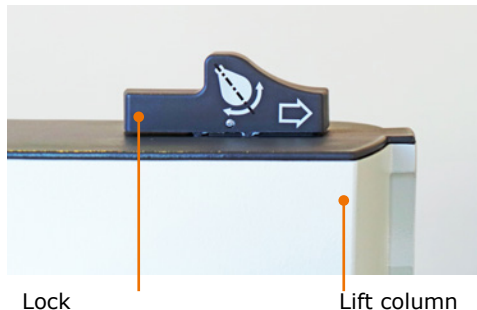Flask immersion depth (height stop)

The stop band of the height stop is located on the right side of the flask column that can be used to adjust a constant immersion depth for the evaporating flask.

- ✓ The rotation is switched off.
- ✓ The evaporating flask is mounted (see chapter "Operation", section "Mounting the evaporating flask").
- ✓ The height stop is at the top position.

Adjusting the lift stop without the safety hood:

- Lower the lift until the desired immersion depth is reached (see chapter "Start-Up", section "Heating bath and flask lift").
- Turn the locking screw on the stop band of the height stop to the right of the lift column by about 180° counterclockwise.
- Slide the stop band down as far as it will go using the locking screw.
- Tighten the locking screw clockwise.
- Return the lift to the top.
- ✓ The locking screw holds the stop in position.
- ✓ The lift moves down no further than the stop and ensures a constant immersion depth of the evaporating flask

Adjusting the lift stop with the safety hood:

- Hold the safety hood with one hand.
- Guide the safety hood by hand when moving the safety hood (see chapter "Assembly, Safety hood").

Control box

The control box is mounted to the base plate behind the heating bath.

- Attach the control box to the base plate with the two included metric screws.

Fixing bolt

The control box plug is on the rear of the heating bath to the right, next to the heating bath plug.

- Plug the cord of the control box into the connection on the base plate.

Control box connection


Vacuum manifold

Assembly parts

- 1x vacuum manifold (holding plate with vacuum sensor, ventilation valve and dirt filter)
- 2x threaded screws
- 1x torx wrench

The sheet of the vacuum manifold is mounted onto the back side of the flask lift.

- Use the torx wrench to screw the two threaded screws into the back side of the flask lift by a few turns.
- Hang the sheet of the vacuum manifold onto the two threaded screws.
- Tighten the two screws.
- ✓ The vacuum manifold is installed.

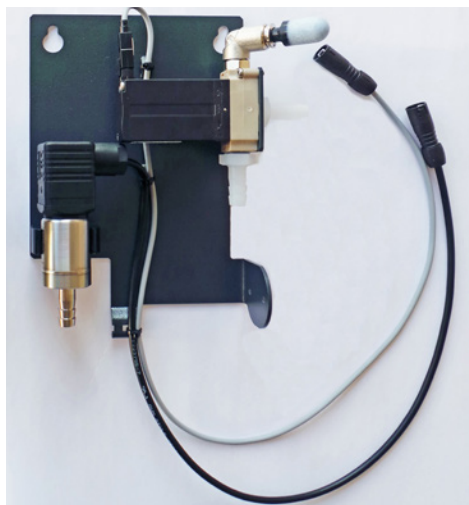Threaded screws

➤ Attaching the peripheral devices

Connecting vacuum

Warning:
Risk of implosion, risk of poisoning

Glass flasks that show even the slightest damage can burst and, depending on the sample, can poison the surrounding air.

Do not increase the vacuum pressure unless necessary.

Never exceed the permissible pressure of your distillate.

To connect the vacuum, you need:

- a vacuum source
- a vacuum tube (Ø 7-8 mm inside)

Glassware sets G1, G3, G3 XL and G6

The opening for the vacuum connection is located on the back and top of the condenser.

- Unscrew a screw cap with the tube nipple.
- Push the vacuum tube onto the tube nipple.
- Tighten the screw cap with the tube nipple and the tube.
- Attach the other end of the tube to the ventilation valve on the back side of the evaporator.

Vacuum connections G1, G3, G3 XL and G6

Vacuum connection screw caps

Glassware set G5

The opening for the vacuum connection is located on the back and top of the condenser.

Vacuum connection G5

Vacuum connection opening


Woulfe's bottle

(optional accessories: see chapter "Spare Parts and Accessories")

The Woulfe's bottle is attached between ventilation valve and condenser.

- Unscrew the glass bottle from the lid.
- Use the two included knurled screws to fix the retainer bracket to the side of the flask lift.
- Hand-tighten the glass bottle on the lid again.
- Lay a short tube from the ventilation valve to the outlet of the Woulfe's bottle.
- Lay a slightly longer tube from the inlet of the Woulfe's bottle to the opening for the vacuum connection on the back side of the condenser at the top.

Retainer bracket Knurled screws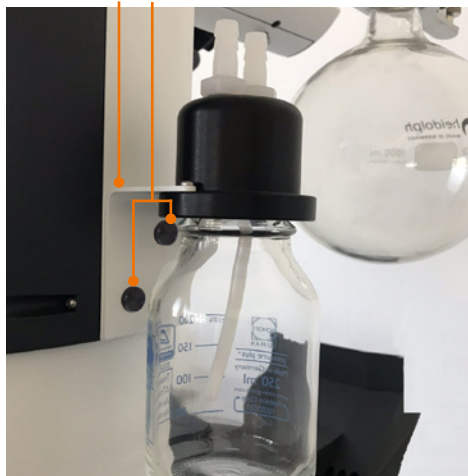


Connecting inert gas

The inert gas connection is situated at the ventilation valve on the vacuum manifold. Prior to connecting the inert gas, the dirt filter must be removed from the ventilation valve.

- Pull back the release ring around the opening of the ventilation valve and pull off the dirt filter.
- Now connect the inert gas.
- When removing the inert gas connection, put the dirt filter back onto the ventilation valve.

Dirt filter Release ring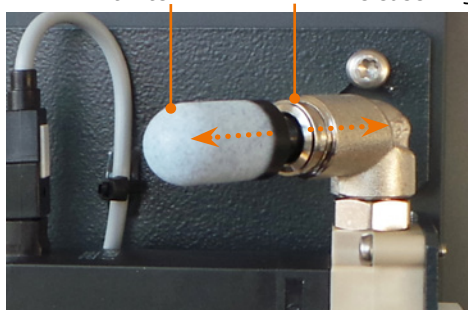

When the device has been in operation for some time, the ventilation valve will get very hot. Before connecting the inert gas, switch off all processes and let the ventilation valve cool down for a bit.


Assembly

Attaching the vacuum control to the control box


The components for the vacuum control are situated on the vacuum manifold at the back side of the flask lift. The connections for these components are situated on the control box.

- Attach the cable of the ventilation valve to the connection "aerating valve". The ventilation valve is situated on the top right of the vacuum manifold.
 - Attach the cable of the vacuum sensor to the "vacuum sensor" connection.
- Only for valve-controlled pumps:
- Attach the cable of the vacuum valve to the connection "vacuum valve".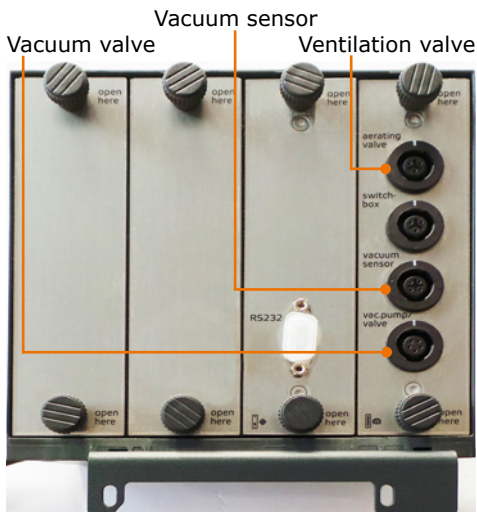

If you operate with a valve-controlled vacuum pump, you require the vacuum valve. This is available as optional accessories (see chapter "Spare Parts and Accessories")

Connection diagram for vacuum: Example valve-controlled pump/house vacuum (View: evaporator back side)

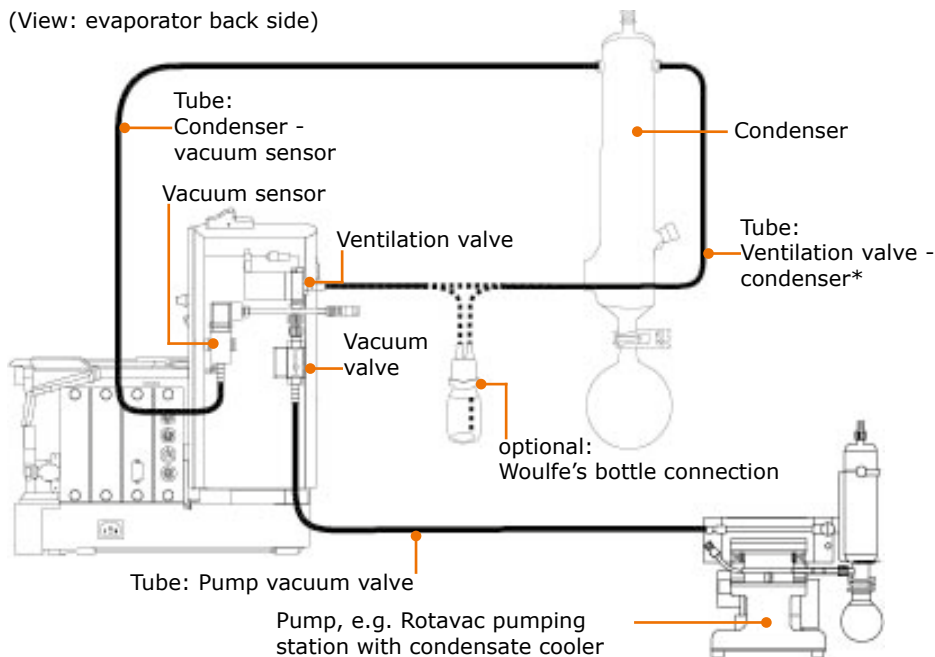
* The optionally available Woulfe's bottle is attached between condenser and ventilation valve, if required.

Assembly

Connection diagram for vacuum: Example speed-controlled pump


(View: evaporator back side)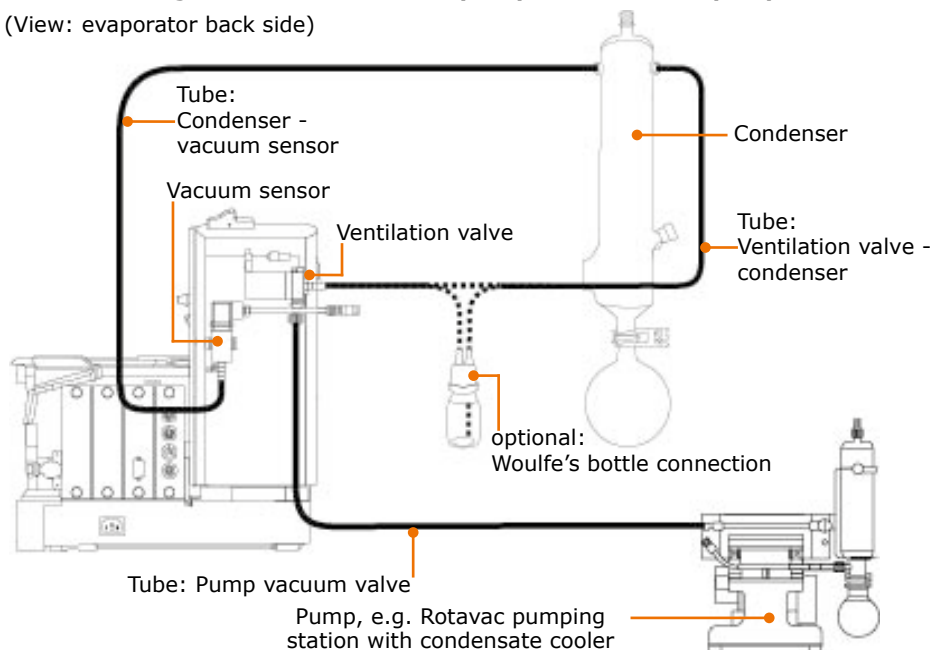

* The optionally available Woulfe's bottle is attached between condenser and ventilation valve, if required.

Connection diagram for vacuum: Example pumping station with integrated vacuum control

(View: evaporator front)

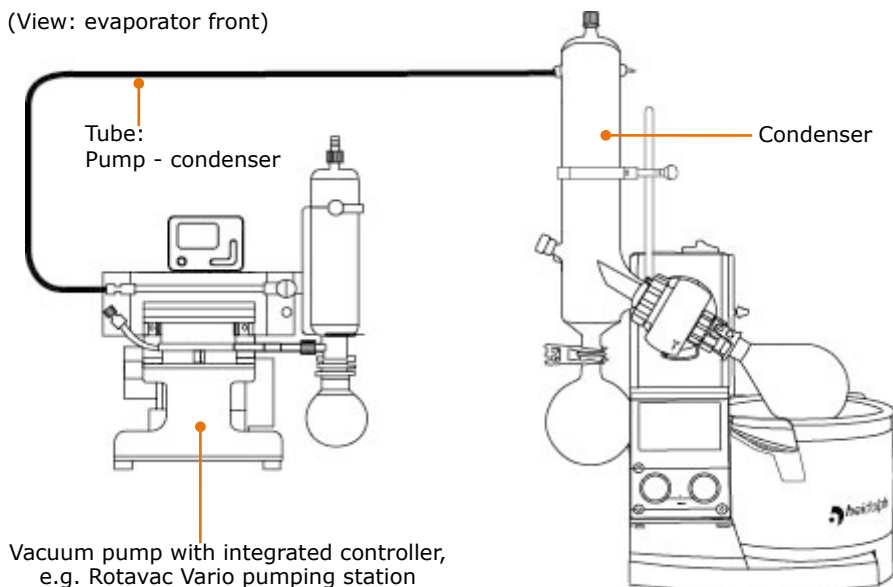Cooling water connection

The external chiller Hei-CHILL is attached to the RS-232 interface on the back side of the control box. The cooling temperature is then displayed in the control panel and can be adjusted there.


Warning: Risk of explosion

A closed valve in the return line will build up high pressure in the glass coils of condensers and may result in glass bursting.

The cooling coils of the condenser may be subjected to a maximum cooling water pressure of 2 bar. To prevent a higher pressure from ever occurring, no valve or throttle must be installed in the cooling water return line.

Assembly parts:


- 2x tube nipple (included in scope of delivery)
- 2x tube clamps (included in scope of delivery)
- 2x tube clamps
- 1x tube, inside Ø 7-8 mm (included in the optionally available tube set)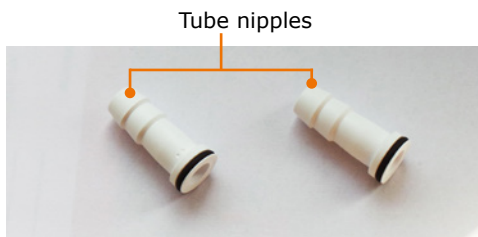

Cooling fluid inlet and outlet are located on the back and bottom of the condenser.

- Remove the screw caps from the cooling water inlet and outlet.
- Push the screw caps over the tube nipples so that the black seal is located inside the cap.
- Push the respective feed and outlet hoses for the cooling water through the screw cap over a tube nipple.
- Tighten the screw caps with the seal first on one of the respective openings of the condenser.
- Secure the tubes additionally with the tube clamps included in the delivery.
- ✓ Cooling water feed and return line are connected.

Cooling water inlet and outlet

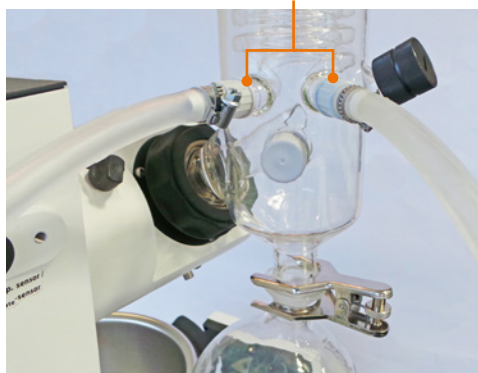Optional accessories

Safety hood

Caution:
Possible damage to the safety hood


If temperatures are too high and/or if the wrong heating bath fluids are used, the material will become soft and may be subject to deformations and possibly tearing.

The safety hood is only suitable up to heating bath temperatures of <130°C.


(For suitable heating bath fluids see chapter "Start-up, Heating bath and flask lift" as well as chapter "Spare Parts and Accessories, Accessories")

Before you can use the safety hood, you need to screw the support of the safety hood to the height stop.

- If needed, slide the safety hood off the carrier.
- Unscrew the locking screw from the height stop band.
- Tighten the safety hood carrier on the height stop band using the two supplied Allen screws.
- ✓ The knurled screw on the safety hood carrier takes over the function of the locking screw for the height stop band.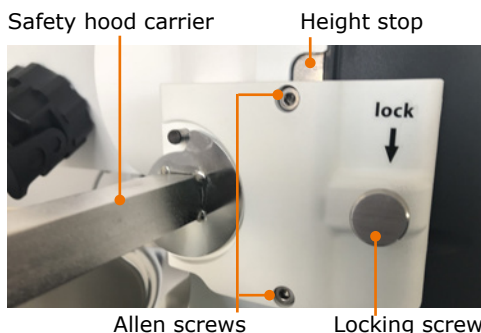


- Loosen the clamping screw and slide the safety hood horizontally onto the carrier until the hood sits above the heating bath.
- Tighten the clamping screw.
- ✓ The safety hood is installed. The safety hood can be lifted open at the handle until it clicks into place.

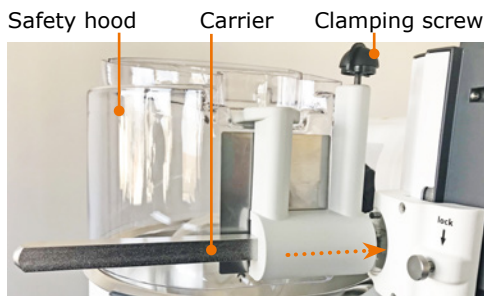Safety shield

- Move the flask lift upward
- Hang the safety shield to the edge of the heating bath and position it so that the (potentially present) safety hood can be lifted open easily.
- Turn the eccentric screw 180° and use to secure the safety shield on the heating bath.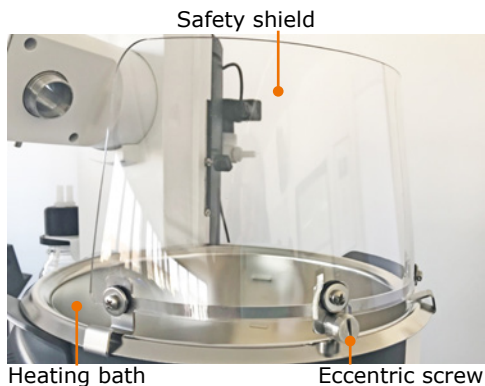

Inlet valve


(Optional accessories) Assembly parts:

- 2x inlet hose
- 1x inlet valve

During a running process, the sample flow for the distillation is metered via the inlet valve.

- ✓ The glassware set is under vacuum.
- Turn the inlet valve counterclockwise.
- ✓ The in-feed is opened, and the distillate is drawn into the evaporating flask.
- ✓ The more you open the inlet valve, the faster your sample will be supplied.
- Turn the inlet valve clockwise.
- ✓ The in-feed is reduced, or the inflow of the sample is stopped.

Inlet valve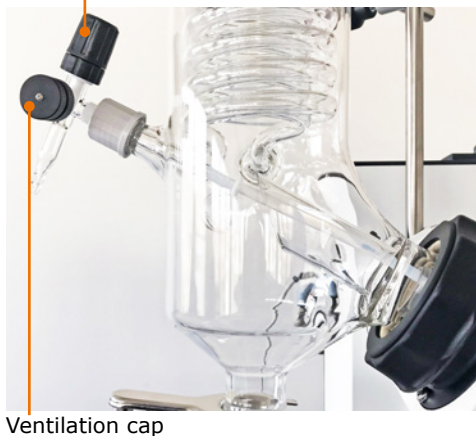

Mounting the inlet valve with the inlet hoses


All glassware sets (with the exception of glassware set G1) have a screw opening on the lower left side of the condenser to accommodate the inlet hose with the inlet valve.

Assembly

Inlet hose 1


- Shorten the supplied PTFE tube so that the end of the tube ends in the evaporating flask.
- Slide the drain disk onto the PTFE tube so that it is positioned before the vapor tube of the evaporating flask.
- ✓ This prevents condensed fluid from flowing back into the evaporating flask via the inlet hose.
- Unscrew the screw cap from the condenser.
- Slide the seal ring and screw cap about 3 cm onto the PTFE tube.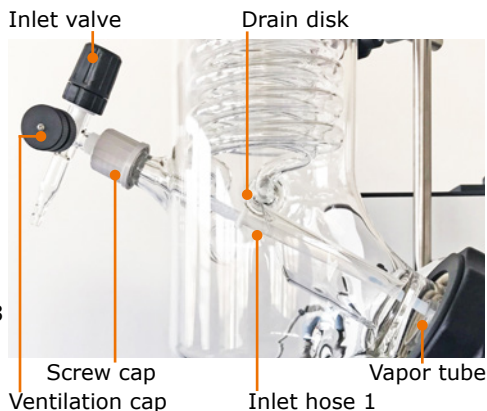

- Insert the inlet valve with the inlet hose into the inlet opening of the condenser and tighten the screw cap.

Inlet hose 2

- Slide the PTFE tube onto the feed opening.
- ✓ The inlet hoses are installed.

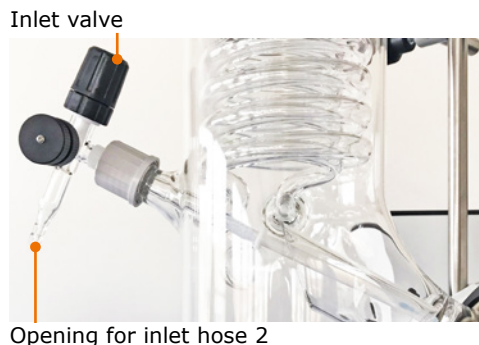Glassware set G1 inlet hoses

The opening for the inlet hoses is located at the top of the condenser.

Vapor temperature sensor, AUTOaccurate sensor

(optional accessory, see chapter "Spare Parts and Accessories").

The opening for the vapor temperature sensor is located at the bottom on the side of the condenser.

The opening for the AUTOaccurate sensor is located at the top center of the condenser.

- Remove the sealing cap from the respective opening of the condenser.
- Insert the appropriate temperature sensor.
The vapor temperature sensor is inserted in the condenser from the side whereas the AUTOaccurate sensor is inserted from the top.
- Tighten the temperature sensor with the screw connection cap.
- Plug in the connector plug of the temperature sensor with the white mark pointing upwards into the plug socket on the left side of the device.

Opening for AUTOaccurate sensor

Opening for vapor temperature sensor
Temperature sensor connection

The AUTOaccurate sensor can only be used with glassware sets G3, G3 XL and G6.

➤ Dismantling, transportation and storage

Dismantling

Warning

Risk of scalding!

Hot glass surfaces and/or heating bath fluids may result in scalding.

Allow all components to cool down before dismantling the device.

Risk of poisoning!

Leakage of hazardous substances and solvents at leak points can cause hazardous substances to be released into the environment.

Risk of electric shock!

If the device is still connected to the power supply, you may come in contact with electrically live parts when you dismantle the device.

Before dismantling the device, switch it off and disconnect the mains plug from the socket. Do not reconnect the device to power under any circumstances.

- ✓ The device is switched off.
- ✓ The heating bath is emptied and cleaned.
- ➔ Remove all tube and cable connections to cooling water, vacuum and other devices.
- ➔ Remove all glassware and remove the heating bath.

Motor lift version: Install the transportation lock

- ✓ The device is switched off.
- ➔ Press and hold the lift stop on the flask lift.
- ➔ Switch the device on and keep the lift stop pressed until the device is fully raised.
- ➔ Release the lift stop.
- ➔ Move the flask lift to the bottom position.

Lift stop

Transportation lock

- ✓ "Transportation lock activated" will appear in the control panel.
- ➔ Use the provided Allen key to attach the metal sheet of the transportation lock with the 3 screws.
- ➔ Switch the device off.
- ➔ Disconnect the power cord from the socket and the device.

Transportation and storage

- Only store and transport the device and its components after emptying and cleaning.
- Store and transport the device and its components in the original packaging or in another suitable container to prevent any damage.
- Carefully seal the packaging against any unauthorized or unintentional opening.
- Store the device in a dry and frost-free location.

Improper handling during storage and transportation may result in damage to the device and the mechanical components of the device.

Avoid jolts and vibrations when transporting the device.

➤ Scope of delivery

Component	Quantity	Product number
Hei-VAP Expert: Motor lift device with glassware set G3*	1	574-01000-00
Hei-VAP Ultimate: Motor lift device with glassware set G3*	1	576-01000-00
Control box	1	569-00200-00
PTFE/FKM vacuum seal	1	23-30-01-01-30
Vapor tube	1	514-00000-01
Clamping sleeve	1	23-30-01-05-31
Motor lift transportation lock	1	11-300-006-28
Coupling	1	23-09-03-01-03
Tension spring	1	22-03-02-01-05
Easy Clip flask clamp, NS 29	1	23-30-01-05-29
Easy Clip flask clamp, NS 24	1	23-30-01-05-57
Ground joint clamp	1	515-42000-00
Operating manual German / English	1	01-005-005-86
Warranty registration / declaration of no objection	1	01-006-002-78
EU Declaration of Conformity	1	01-001-025-03
Power cord	1	country-dependent

* The indicated product numbers are valid for 230 V devices in the European Community. Please contact our sales department or www.heidolph-instruments.com for product numbers of country-specific varieties.

➤ Accessories

Component	Quantity	Product number
Device fuse, 230 V	2	14-002-028-22
Device fuse 115 V	2	14-002-028-29
Vapor temperature sensor	1	569-00030-00
AUTOaccurate sensor	1	569-00040-00
Safety hood	1	569-00010-00
Safety shield	1	569-00020-00
5 L Budde heating bath fluid (20° C to 170°C)	1	515-31000-00
5 L Wacker AP 200 silicone oil (150° C to 210° C)	1	569-00600-00
Tube (suitable for vacuum and water)	1	591-35000-00
PTFE/FKM vacuum seal	1	23-30-01-06-70
Extension cord for control panel 1.5 m	1	14-300-020-12
Vacuum valve	1	14-300-005-28
Woulfe's bottle	1	569-00071-00
Glassware sets		see general catalog
Vacuum pumps and chillers type and size vary		see general catalog

Additional accessories can be found in our general catalog or on our website at www.heidolph-instruments.com

Software copyright

© Copyright 2018 - Copyright notice

The software contained in this product is protected by copyright.

The copyright is held by Heidolph Instruments GmbH & Co. KG, Walpersdorfer Strasse 12, 91126 Schwabach. Any Open Source software components which may be included in the software are exempt. For details about the Open Source software included in the product please see the supplementary sheet "Details about Open Source software included in products" which is enclosed to the product as well as the respective information on the also enclosed data medium.

Copyright violations (e.g. the unauthorized use or alterations of the software) can prompt claims according to civil law (such as omission, compensation of damages, German Copyright Act [UrhG] 97 et seq.) and/or indictable consequences (UrhG 106 et seq.).

➤ Technical data

Hei-VAP Core

Standard supply voltage	230 V 50/60 Hz or 115 V 50/60 Hz
Power consumption (W)	1400
Protection class (IEC 60529)	<ul style="list-style-type: none"> ▪ Device: IP 20 ▪ Heating bath cable: IP 67 ▪ Control box: IP 42 ▪ Control panel: IP 42
Noise pressure level (dB(A)) (based on IEC 61010)	< 85 dB(A)
Flask lift vertical travel (mm)	155
Motor	EC motor
Rotation speed range (rpm)	10-280
Speed control	electronic
Heating bath container	<ul style="list-style-type: none"> ▪ Inner Ø: 253 mm, Ø outer: 291 mm ▪ Material: stainless steel (1.4404) X2CrNiMo 17-12-2 ▪ Volume 4.5 l
Heating capacity (W)	1,300
Heating bath temperature range (°C)	<ul style="list-style-type: none"> ▪ 20-100 H₂O ▪ 20-210 oil
Heating bath temperature control	electronic
Heating bath control accuracy (°C)	±1
Excess temperature protection	cut off at 5 °C difference to set temperature
Overheating protection (°C)	250
Condensing area (cm ²)	1,400
Evaporation rates (L/h) ΔT* = 40°C (continuous operation)	<ul style="list-style-type: none"> ▪ Toluene 8.5 ▪ Acetone 5.8 ▪ Ethanol 3.5 ▪ Water 1.2
Digital displays	<ul style="list-style-type: none"> ▪ Bath temperature display (°C) ▪ Vapor temperature display (°C) (only with optional vapor temperature sensor) ▪ Rotation speed display ▪ Vacuum display

Lift operation

	Manual lift	Motor lift
Weight (without glassware set and control box) (kg)	14.7	15
Lifting speed (mm/s)	-	25
Dimensions with glassware set G3 (w x d x h) (mm)	739 x 532 x 887	739 x H 477 x 887

* ΔT = Difference between bath temperature and vapor temperature

➤ **Contact / Technical Service**

Questions and repair work

If any aspect of installation, operation or maintenance is not covered in this manual, please get in touch with the following address.

For repairs, please call Heidolph Instruments directly or your local authorized Heidolph Instruments distributor beforehand.

Warning: Risk of poisoning!

Contaminated devices can lead to severe injury or death of our employees!

When shipping devices for repair that may have been contaminated by substances harmful to health, please:

- advise exact substance
- take adequate protective action towards our parts receiving and service personnel
- mark the pack in accordance with Ordinance on Hazardous Substances

A declaration of no objection can be found at the end of this operating manual.

→ Prior to shipping the device for repair, complete a copy of this declaration of no objection form and submit it in advance.

Our contact details for Germany, Austria, Switzerland

Please only send devices to this address after prior consultation with us:

Heidolph Instruments GmbH & Co. KG

Technical Service
Walpersdorfer Str. 12
91126 Schwabach/Germany

Phone: +49 9122 9920-74

Fax: +49 9122 9920-84

E-mail: service@heidolph.de

Local distributors

You will find the contact details of your local Heidolph distributor at www.heidolph.com

➤ Warranty

Heidolph Instruments provides a 3 year warranty for the products described here (excluding glass and wearing parts) if you register using the warranty card enclosed or online (www.heidolph.com). The warranty period begins with the date of purchase. The device serial number is also valid without registration.

This warranty covers material and manufacturing defects.

In the event of a material or manufacturing defect, the device shall be repaired or replaced free of charge under the terms of the warranty.

Heidolph Instruments shall not assume liability for any damage incurred as a result of improper handling or transport.

→ Please notify Heidolph Instruments if you wish to make a warranty claim.

➤ Declaration of no objection

→ In the case of repair, copy and complete the declaration of no objection form and send it to Heidolph Instruments

1. Details about the device

Product number _____

Serial number _____

Reason for repair _____

2. Has the device been cleaned, decontaminated/sterilized?

Yes _____ No _____

3. Is the device in a condition which does not represent any health threats for our service department staff?

Yes _____ No _____

If not, which substances has the device come into contact with?

4. Legally binding declaration

The customer is aware of being legally liable to Heidolph Instruments for any damages arising from incomplete and incorrect information.

Date

Signature

Company stamp

Please note

The shipper is responsible for the return of the goods in well packed condition, suitable for the mode of transport.

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The screenshot shows the application's interface for creating a new favorite. At the top, there is a list of existing favorites: Benzene, Cyclohexane, and Dichlormethane. Below this, the 'Cyclohexane' favorite is selected, and its parameters are displayed: Vakuumpumpe (152 mbar), Hysterese Δp (Auto), and Rotation (10 rpm). Below this, the 'Favoriten' screen is shown, where a new favorite 'Cyclohexane01' has been created. The parameters for this new favorite are: Vakuumpumpe (1000 mbar), Hysterese Δp (Auto), Rotation (10 rpm), Heizbad (20 °C), and Umlaufkühler (5,0 °C). At the bottom, there are buttons for 'Abbruch' (Cancel) and 'Speichern' (Save).

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© Copyright 2018 - Urheberrechtshinweis
This is a translation of the original operating manual. We reserve the right to make any technical modifications without prior notification. This documentation is not subject to revision; the latest version of this documentation can be found on our homepage. The copyright for all texts and images is held by Heidolph Instruments GmbH & Co. KG.
© Heidolph Instruments GmbH & Co. KG