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## ROTI®ImmunoBlock

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# Instructions for use

## ROTI®ImmunoBlock

Blocking reagent for immunofluorescent applications

### A. Introduction

ROTI®ImmunoBlock is an innovative, ready to use blocking reagent made on a polymer basis without detergent and following advantages:

- Your detection system is not contaminated by any foreign proteins. The signal intensity remains constant.
- Specific binding sites are preserved whilst nonspecific reactions are suppressed. Signal intensity increased.
- The ready to use solution saves you time and labour by eliminating the need for dissolving.

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### B. Application

#### ROTI®Block working solution

Dilute ROTI®ImmunoBlock (10 x concentrate) before use 1:10 with H<sub>2</sub>O (distilled). The diluted solution can be used directly.

#### ROTI®ImmunoBlock and samples

When blocking please cover your entire sample with ROTI®ImmunoBlock working solution.

#### Blocking time and temperature

Blocking should occur under gentle shaking at room temperature. A blocking period from 15 mins to 1 h has proved successful.

#### ROTI®ImmunoBlock and antibodies

You can also incubate your sample with primary and secondary antibody diluted in ROTI®ImmunoBlock working solution. Alternatively incubation can occur in ROTI®ImmunoBlock working solution diluted 1:10 in PBS or TBS, respectively, or PBST or TBST, respectively.<sup>14</sup>

TBST: ROTI®Stock 10 x TBST (Art. No. 1061.1)  
(25 mM Tris, pH 7.4, 150 mM NaCl, 3 mM KCl, 0.05% Tween 20)

PBST: ROTI®Stock 10 x PBST (Art. No. 1059.1)  
(137 mM NaCl, 2.7 mM KCl, 10 mM Na<sub>2</sub>HPO<sub>4</sub>, 2 mM KH<sub>2</sub>PO<sub>4</sub>, pH 7.4, 0.05 % Tween 20)

TBS: ROTI®Stock 10 x TBS (Art. No. 1060.1)  
(25 mM Tris, pH 7.4, 150 mM NaCl, 3 mM KCl)

PBS: ROTI®Stock 10 x PBS (Art. No. 1058.1)  
(137 mM NaCl, 2.7 mM KCl, 10 mM Na<sub>2</sub>HPO<sub>4</sub>, 2 mM KH<sub>2</sub>PO<sub>4</sub>, pH 7.4)

<b>ROTI®ImmunoBlock</b>	T144.1	250 ml
-------------------------	--------	--------

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#### **Carl Roth GmbH+Co. KG**