Temperature

Controlled Shaking

Leading Safety Standards

Superior Ease of Use

Reduced Cost of Ownership

Operating Manual

Page 2

Test Tube Shakers Hei-MIX Reax top/Hei-MIX Reax control
Operating manual must be read before initial start-up.
Please follow the safety instructions provided.
Please keep for future reference.

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About this Document	4
Version and variants	4
About this manual	4
Symbols and signal words.....	5
Safety Instructions	6
General safety instructions.....	6
EU Declaration of Conformity.....	6
Intended use.....	6
Installation / electrical safety	7
Qualifications of personnel	7
Obligations of the operator	7
Installation site	7
Modifications to the device.....	8
Safety of personnel.....	8
Safety during use	9
Disposal	9
Device Description	10
Device overview	10
Hei-MIX Reax top	10
Hei-MIX Reax control	10
Start-Up	11
Setting up device	11
Connect / disconnect power cord	11
Switch unit on and off	12
Operation	13
Shaking mode.....	13
Automatic mode	13
Continuous mode.....	13
Speed / Shaking intensity	14
Set Speed.....	14

Cleaning and Maintenance	15
Cleaning.....	15
Maintenance	15
Troubleshooting	16
Possible errors and how to resolve	16
Assembly	17
Electrical connection	17
Mounting test tube tray	18
Disassembly and Storage	19
Dismantling, transportation and storage	19
Dismantling	19
Transportation and storage	19
Accessories and Spare Parts	20
Scope of delivery.....	20
Accessories.....	20
Attachments	21
Technical data	21
Service	22
Contact / Technical Service	22
Warranty	23
Confirmation of condition.....	24

About this Document

Copyright

This operating manual is protected by copyright and is exclusively intended for the internal use of the purchaser.

Handing over this operating manual to third parties, reproducing it in any type or form –in part or in full – and utilizing and/or communicating its contents is prohibited without prior written approval from the manufacturer.

Any violations will result in an obligation to provide compensation for damages. Other claims remain unaffected.

➤ Version and variants

Version

This manual describes the function, operation and maintenance of the test tube shakers Hei-MIX Reax top and Hei-MIX Reax control.

Version	Alteration date
1.3	05/2018

Variants

The device is available in different variants. Certain features or functions are only available in specific product variants. These variants are described in this operating manual.

➤ About this manual

This operating manual is an integral part of the device described here.

- ➔ Carefully read through this operating manual and observe all of the safety instructions and warnings notices.
- ➔ Ensure that every user has carefully read the operating manual before using the device for the first time.
- ➔ Store this operating manual so that it is accessible to all users at all times.
- ➔ Pass the operating manual on to any subsequent owner.

The current version of this operating manual can also be found at:
[https://heidolph-instruments.com/en/service/downloads/operation-manuals#Shakers & Mixers \(Reax top and Reax control\)](https://heidolph-instruments.com/en/service/downloads/operation-manuals#Shakers & Mixers (Reax top and Reax control))

In addition, also observe the local regulations at the installation site.

➤ Symbols and signal words

Standardized symbols and signal words are used in this operating manual to warn against any dangers and provide important instructions. These instructions must be strictly observed to avoid accidents and damage.

The following terms and basic symbols are used:

Symbol	Additional signal words / explanation
Warning signs 	The yellow triangle indicates hazardous situations. It is used in combination with the following signal words: DANGER: Indicates a hazardous situation which, if not avoided, will result in serious injury or death. WARNING: Indicates a hazardous situation which, if not avoided, may result in serious injury or death. CAUTION: Indicates a hazardous situation which, if not avoided, may result in property damage and minor or moderate injury.
Prohibitory signs 	Prohibited: The red circle indicates a situation that should be avoided under all circumstances and which, if not avoided, may result in serious injury or death.
Mandatory signs 	Must be observed: The blue circle indicates important information. Please observe this information to avoid any property damage.

Other symbols used:

Symbol	Description
→	Handling instruction, action required
✓	Result of action
▪	List of information
▪	
▪	
a.	List of variants
b.	
c.	

➤ General safety instructions

EU Declaration of Conformity

This device complies with the following EC-Directives*:

- 2006/42/EC Machinery Directive
- 2014/30/EU Electromagnetic Compatibility Directive

*Also see the attachment "EU Declaration of Conformity".

The device has been manufactured according to state-of-the-art technology and in compliance with recognized safety regulations. However, risks may still arise during installation, operation and maintenance.

→ Please ensure that the operating manual is available at all times.

The device may only be used under the following conditions:

- Only operate the device if it is in a technically perfect condition.
- Only operate the device if you are properly aware of the risks and required safety measures.
- Only operate the device in accordance with the instructions given in this operating manual.
- If any information is ambiguous or missing, ask your superior or contact the manufacturer.
- Do not operate anything on the device without authorization.
- Only use the device in accordance with its intended use.

Intended use

The device may only be operated by authorized personnel.

The device is suitable for the following uses:

- Shaking
- Mixing
- Emulsifying
- Suspending
- Dissolving

The device may be operated in research laboratories, other laboratories and production facilities in the following sectors:

- Chemistry
- Pharmacy
- Biology
- Environmental analytics
- Basic research
- Similar research laboratories

Reasonable foreseeable misuse

Any use which deviates from the device's intended use is considered to be improper. The manufacturer is not liable for damage that occurs as a result. The risk is borne by the operator alone.

Installation / electrical safety

- The device may only be connected up if the available power supply voltage corresponds to the information stated on the rating plate for the device.
- The power supply connection must be easily accessible at all times.
- An RCD circuit breaker (residual current circuit breaker) must be fitted to the electrical system in the building so that the device is disconnected from the power circuit in the event of a fault.
- Repairs may only be performed by electricians authorised by Heidolph Instruments.
- Never operate the device with a damaged power cord.
- Always turn the device OFF and disconnect the power cord before carrying out any maintenance or repair work.

Qualifications of personnel

- The device may only be operated by trained personnel.
- The device may only be operated by persons who have been instructed and supervised in its proper use by trained specialist personnel.
- The device may only be operated by specialist personnel who are above the legal minimum age.
- Other persons may only work on the device under the constant supervision of experienced and trained specialist personnel.
- This operating manual must be read and understood by all persons working with the device.
- Personnel must receive safety training that ensures responsible and safe working practices.

➤ Obligations of the operator

Installation site

- The device must be installed in a suitable location.
- The device must be installed on a firm and stable surface.
- Ensure that the device and all of its components are easily accessible at all times.
- Maintain a sufficient safety clearance to, in particular, any moving and/or hot equipment components.
- It is not permitted to place or store any objects such as accessories, tools or chemicals within this clearance area during operation.
- All screw connections must be securely tightened.
- It is not permitted to operate the device near to highly flammable or explosive substances.
- The device may only be operated inside buildings that have been equipped with appropriate laboratory equipment (e.g. a fume cupboard).
- Only operate the device (depending on the substances being used) in conjunction with an extractor hood (see DIN EN 14175 and DIN 12924).

Safety Instructions

- It is only permitted to operate the device in dry and enclosed rooms and under the following ambient conditions:

Ambient temperature	5 - 31 °C at 80 % relative humidity
	32 - 40 °C decreasing linearly to a maximum 50 % relative humidity
Installation altitude	0 - 2,000 m above sea level
Contamination level	2
Overvoltage category	II
Permissible supply deviations	± 10 %

- If the device is operated in corrosive atmospheres, the service life of the device will decrease based on the concentration, duration and frequency of the exposure to the corrosive atmosphere e.g. concentrated hydrochloric acid (HCl).

Modifications to the device

- It is not permitted to make any unauthorised modifications or changes to the device.
- Do not attach or install any parts that have not been approved by the manufacturer.
- Unauthorized modifications or changes will void the EC Declaration of Conformity for the device and operation of the device will no longer be permitted.
- The manufacturer is not liable for any damage, dangers or injuries that result from unauthorized modifications and changes or due to the non-observance of the instructions in this manual.

Safety of personnel

- Ensure that the device is only operated by qualified specialist personnel and trained employees.
- Observe the following instructions to avoid any personal injuries and property damage:
 - Laboratory regulations
 - Accident prevention regulations
 - Hazardous Substance Act
 - Other generally accepted occupational health and safety regulations
 - Local regulations

➤ Safety during use

- Wear the appropriate clothing when working on the device (protective clothing, protective glasses and, if necessary, safety gloves).
- Do not use the device in potentially explosive areas. The device is not protected against explosion. There is no explosion or ATEX protection available.
- Do not carry out work with naked flames in the vicinity of the device (risk of explosion).
- Do not operate or assemble devices in the vicinity which are emission or radiation sources (electromagnetic waves) for the frequency range ($3 \cdot 10^{11}$ Hz to $3 \cdot 10^{15}$ Hz).
- Do not operate or assemble appliances in the vicinity of the device which constitute emission or radiation sources for ionizing radiation or in the ultrasonic range.
- Do not operate the device where adiabatic compression or shock waves might occur (shock wave combustion).
- Do not use substances, where the energy input due to mixing might pose a hazard.
- Do not spill liquids over the device or any parts of it.
- Remove any accidentally spilled liquids immediately.
- Eliminate errors immediately.
- Always switch the device OFF after use.
- Avoid putting pressure on the display when you are not operating the device.
- Do not use abrasive material to clean the surface. Only wipe with damp cloth.
- Observe the Directive 2002-44 EC regarding the exposure of workers to the risks arising from physical agents (vibration).

➤ Disposal

- Check the device components for hazardous substances and solvents.
- Clean all components before disposal.
- Dispose of the device in accordance with the relevant national regulations.
- Dispose of the packaging material in accordance with the appropriate national regulations.

➤ Device overview

Hei-MIX Reax top

Hei-MIX Reax control

➤ Setting up device

Warning: **Risk of slipping device!**

As a result of vibrations generated during operation the device might slide off the table top.

The surface must be smooth and clean and should be properly leveled.

- Locate the shaker on a stable, horizontal surface.
- Clean surface and feet with a damp cloth and ethanol regularly.

➤ Connect / disconnect power cord

Connect power cord

The power cord comes with a three-wire plug and a recess on the bottom side.

Unit plug

Power cord

The socket is located on the back side of the unit.

- ✓ The device is switched OFF.
- Connect plug to the socket with the recess facing down.
- Push plug up against socket until securely attached.
- Connect plug to power socket.

Unit socket

Disconnect power cord

- Disconnect plug from power socket first.
- Then disconnect plug from socket on unit.

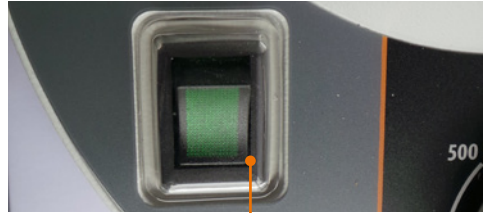
➤ Switch unit on and off

The ON/OFF switch is situated on the front panel.

Switch unit on


- Press the On/Off switch on the lower side.
- ✓ The green control light switches on.
- ✓ The unit is switched on.

Switch On: device switched on

Switch unit off

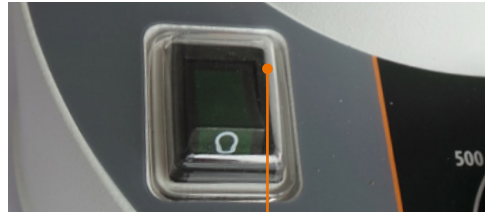- Press the On/Off switch on the upper side.
- ✓ The green control light switches off.
- ✓ The unit is switched off.

Switch Off: device switched off

➤ Shaking mode

You can choose from two shaking options:

- Automatic mode: The unit starts shaking as soon as you put pressure on the test tube tray.
- Continuous mode: The unit shakes at an ongoing vibration mode until you switch off the continuous mode or the complete unit.

Warning: **Risk of injury!**

Returning power after a power cut off will restart shaking mode. There is a risk of injury in the vicinity of moving parts of device.

Before working with the device or in the vicinity ensure all controls are switched to "0" or "OFF" mode, when device is not moving.

Warning: **Risk of injury!**

High speeds in combination with high loads and/or a tall assembly may lead to an unstable device. As result the device may start moving and slide off the table top.

When working with with high loads and/or a tall assembly, increase the speed with care and watch for the devices stability. Ensure the device is installed in such a way as it can not collide with the surrounding environment or slip off its work surface.

Automatic mode

- ➔ Turn the slide switch to the right as far as it will go.
- ✓ A delta icon ▼ appears on the left side of the slide switch.
- ➔ Place your tube in the test tube tray and push the tube with soft pressure downwards.
- ✓ The shaking operating starts as soon as you position a tube on the test tube tray.
- ➔ Remove the tube.
- ✓ The shaking operating stops.

Symbol automatic mode

Slide switch

Continuous mode

- ➔ Turn the slide switch to the left side as far as it will go.
- ✓ A circle icon ● appears on the right side of the slide switch.
- ✓ The shaking operating starts.
- ➔ Turn the slide switch to automatic mode or switch off the unit.
- ✓ The shaking operating stops.

Slide switch

Symbol continuous mode

➤ Speed / Shaking intensity

Set Speed

Warning: **Risk of splashing!**

Open vessels and too high vibration frequencies may result in samples splashing.

Wear safety glasses and appropriate protective clothing.

Select low speed when starting the unit to avoid splashing from tubes.

If necessary, increase the speed carefully.

Caution: **Possible risk of exertion**

There is a possible risk of exertion due to physical force required to holding test tube in place.

Do not use the device any longer than recommended in the directive 2002-44 EC on vibration strain.

REAX top

- ➔ Turn the speed rotation knob clockwise to the right.
- ✓ The shaking intensity increases.
- ➔ Turn the speed rotation knob counter-clockwise to the left.
- ✓ The shaking intensity decreases.

The Reax top is equipped with a guideline to set the speed.

REAX control

The Reax control is equipped with an analog dial speed control with accurate rpm numbers setting.

> Cleaning

All surfaces of the unit allow for cleaning with a damp cloth and if required a mild soap lotion.

Warning:
Risk of electric shock!

If the device is still connected to power during cleaning, you might get in contact with live parts due to accidentally entering liquids.

Prior to cleaning:

- ensure that the device is switched OFF
- ensure that the plug is pulled from the electric socket.

Caution:
Damage to the device

Accidentally entering liquids will cause the device to fail .
Only wipe with a damp cloth. Never use a soaked cloth.

Attention:
Risk of damage to the surfaces

All of the surfaces may be damaged by improper cleaning.

Never use:

- Chlorine bleach or any cleaning agent containing chlorine
- Solvent-bearing agents (e.g. acetone)
- Ammonia
- Abrasive cleaning agents such as scouring agents, cleaning wool or agents with metallic particles

As a result of vibrations generated during operation the device might slide off the table top.

To ensure a stable position clean surface and feet with a damp cloth and ethanol regularly.

> Maintenance

The unit is maintenance-free. Any necessary repair must be performed by an authorized Heidolph distributor.

Please contact Heidolph Instruments or your local Heidolph Instruments distributor.

➤ Possible errors and how to resolve

Malfunctions	Possible reason	Troubleshooting
On-/Off switch does not enlighten	No power	<ul style="list-style-type: none"> ▪ Check size and shape of plug and compatibility to your electrical socket ▪ Check circuit breakers
	Light diode defective	Contact your local Heidolph Instruments distributor
No shaking movement in tilt automatic mode	Slide switch not in correct position	Push slide switch to the far right position
	Not enough pressure on the tray	Increase pressure by hand
	Pressure on tray not applied vertical	<ul style="list-style-type: none"> ▪ Apply vertical pressure ▪ Switch over to continuous operation mode
Shaking movement stops in continuous operation mode	<ul style="list-style-type: none"> ▪ Motor defective ▪ Switch defective ▪ Control electronics defective 	Contact your local Heidolph Instruments distributor

If you experience a malfunction which can not be resolved, please contact your authorized Heidolph Instruments distributor immediately.

Electrical connection

Repairs to the device must **only be carried out by a qualified electrician approved by Heidolph Instruments**. Improper repairs could result in serious dangers. Please contact your specialist dealer if the device needs to be repaired.

Warning: Risk of electric shock

The device must be connected to an earthed power socket.

The rated voltage for the device and the supply voltage must be the same. The rated voltage for the device can be found on the rating plate on the rear of the device.

The device is earthed via the supplied power cord when delivered. If the original plug is replaced, it is essential that the new plug is connected to a protective earth!

Rating plate

Unit socket

Colour coding for the electrical connection:

Europe

**GREEN/
YELLOW**

PE: Protective earth
(earth)

BLUE

N: Neutral conductor

BROWN

P: Phase

North America

GREEN

PE: Protective earth
(earth)

WHITE

N: Neutral conductor

BLACK

P: Phase

The device is connected to the power supply with the supplied power cord. The device socket is located on the rear of the device.

If you want to use the device in a country with a different type of plug:

- The supplied plug may only be replaced by a qualified electrician.
- If you want to use an adapter, it must be earthed and approved in accordance with local regulations.
- The cable, adapter and plug must correspond at least to the power stated on the rating plate.

➤ Mounting test tube tray

Take off test tube tray

The test tube tray can easily be attached and pulled off the unit manually without using a tool.

- Grab the test tube tray and lift it upwards to pull it off the unit.

Attach test tube tray

A groove is located on the bottom of the test tube tray.

Groove

- Place the test tube tray with groove in line with the tongue of the unit.
- Push the test tube tray with soft pressure onto the unit.

Tongue

➤ Dismantling, transportation and storage

Dismantling

Warning: Risk of injury!

While the device is plugged in you might be at harm if the device is accidentally switched to ON mode.

Prior to dismantling device ensure that the device is switched off and the plug is pulled from the electric socket.

- ✓ The device is switched off.
- Disconnect plug from power socket.
- Disconnect plug from socket on the device.
- Remove all vessels.

Transportation and storage

- Store and transport the device and its components only if they were emptied, cleaned and dismantled.
- To carry the device wear safety gloves and lift the unit from below.
- Store and transport the device and parts of device in the original packing material or alternatively in an adequate container to prevent damage.
- Seal the package carefully against unauthorized or accidental opening.
- Store the device in a dry and frost-free place.

Improper storage and transportation may result in damages to the system and the mechanical components.

Avoid any kind of shocks during transportation.

➤ Scope of delivery

Component	Variant	Quantity	Product number
Hei-MIX	Reax top	1	541-10000-00
	Reax control	1	541-11000-00
Tray, small		1	23-07-06-05-01
Operating Manual	English / German	1	01-005-002-74
Warranty registration / Confirmation of condition		1	01-006-002-78
EC Declaration of Conformity		1	01-001-025-06
Power cord		1	depending on country

➤ Accessories

Component	Quantity	Product number
Tray, large	1	549-19000-00

Additional accessories are shown in the general catalog or on our website at

- www.heidolph-instruments.com.

➤ Technical data

Hei-MIX Reax top, Hei-MIX Reax control

Standard supply voltage	230 V/50 Hz or 115 V/60 Hz
Supply power (W)	51
Protective class (IEC 61140)	1 ⊕
Protection class (IEC 60529)	IP22
Sound pressure level (dB(A)) (based on IEC 61010)	< 45
Drive	shaded pole motor
Overheat protection	self-resetting
Speed control	electronic
Speed range (U/min)	0 - 2,500
Motion	circular vibration
Orbit (mm)	5
Operating mode	automatic or continuous
Weight (kg)	2.8
Dimensions (w x h x d) (mm)	134 x 105 x 172

Model Hei-MIX	Reax top	Reax control
Speed setting	analog, +/- scale	analog, numbered scale
Control accuracy	unregulated	± 2%

➤ Contact / Technical Service

Questions / Repair work

If any aspect of installation, operation or maintenance remains unanswered in the present manual, please get in touch with the following address.

For repairs please call Heidolph Instruments or your local authorized Heidolph Instruments distributor.

Warning: Danger of poisoning!

Contaminated devices can lead to severe injury or death of our employees!

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise exact substance
 - take adequate protective action towards our parts receiving and service personnel
 - mark the pack in accordance with Ordinance on Hazardous Substances
-

A "Confirmation of Condition" form can be found at the end of this operating manual.

→ Prior to shipping the device for repair, complete a copy of this form and submit it in advance.

Contact details:

Heidolph Instruments Germany

Heidolph Instruments GmbH & Co. KG
Technical Service
Walpersdorfer Str. 12
D-91126 Schwabach / Germany

Tel.: +49 - 9122 - 9920-74
Fax: +49 - 9122 - 9920-84

E-Mail: service@heidolph.de
www.heidolph.de

Heidolph Instruments North America

Phone: 1-866-650-9604
E-mail: service@heidolph.com
www.heidolphNA.com

Heidolph Instruments United Kingdom

Phone: 01799 - 5133-20
E-mail: service@radleys.co.uk
www.heidolph-instruments.co.uk

All other countries

You will find contact details of your local Heidolph distributor at www.heidolph.com

➤ Warranty

Heidolph Instruments provides a 3 year warranty for the products described here (excluding glass and wearing parts) if you register using the warranty card enclosed or online (www.heidolph.com). The warranty period begins with the date of registration. The serial number is also valid without registration.

This warranty covers material and manufacturing defects.

In the event of a material or manufacturing defect, the device shall be repaired or replaced free of charge under the terms of the warranty.

Heidolph Instruments shall not assume liability for any damage incurred as a result of improper handling or transport.

Warranty claim?

- ➔ Please notify Heidolph Instruments or your local Heidolph distributor if you wish to make a warranty claim.

➤ Confirmation of condition

→ In the case of repair, copy and complete the Confirmation of condition form and send it to Heidolph Instruments.

1. Details about the device

Product number _____

Serial number _____

Reason for repair _____

2. Has the device been cleaned, decontaminated/sterilized?

Yes _____ No _____

3. Is the device in a condition which does not represent any health threats for the staff of our service department?

Yes _____ No _____

If not, which substances has the device come into contact with?

4. Legally binding declaration

The customer is aware of being legally liable to Heidolph Instruments for any damages arising from incomplete and incorrect information.

Date

Signature

Company stamp

Please note

The shipper is responsible for the return of the goods in well packed condition, suitable for the mode of transport.

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This is a translation of the original operating manual. We reserve the right to make any technical modifications without prior notification. This documentation is not subject to revision; the latest version of this documentation can be found on our homepage. The copyright for all texts and images is held by Heidolph Instruments GmbH & Co. KG.
© Heidolph Instruments GmbH & Co. KG