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Instructions for All-layer-samplers

Art. No. 5316-2055, 5316-2085, 5316-2150 (AISI 316 L); 5317-2055, 5317-2085, 5317-2150 (aluminium);
5319-1085, 5319-1150, 5319-1250 (Jumbo/aluminium)

Read carefully before use and keep for later reference

Safety Regulations

Every user must know these instructions and have them at hand at all times. The sampler may only be used for easy-flowing granulates. Do not use the sampler for explosive powders/granulates. This device can only be used within the limits set by the resistance of the materials it is made out of. Sampler parts coming into contact with the sample material are

Sampler No. 5316- stainless steel 1.4404/AISI 316 L

Sampler No. 5317- anodised aluminium/stainless steel 1.4404/AISI 316 L

Sampler No. 5319- anodised aluminium/stainless steel 1.4404/AISI 316 L

Always use the sampler in such a way as not to endanger yourself or others. When using the sampler, never use force.

Hazardous materials

If hazardous materials are used during operation or maintenance or while the device is being cleaned, observe the current safety data sheets. Wear personal protective equipment when working with hazardous materials. The operator must determine requirements for specific cases and ensure the requirements are implemented (safety data sheet, hazard analysis). Devices that are contaminated with hazardous materials must not be returned to the device manufacturer for repairs.

Packaging/Contents

Please check the contents to see if they are complete or if they have been damaged during transport. Damage to the packaging which can be seen on the outside should be immediately shown to the deliverer/forwarding agent who in turn must confirm it.

Function, Precautionary Measures, Use Restrictions

The device has been designed to take samples of easy-flowing powders or granulates from sacks, big-bags, canisters, barrels. The operator is responsible for ensuring that only persons who have read and understood the operating instructions are permitted to use the device.

The sampler cannot be used for chemicals that attack 1.4404 stainless steel or the anodised aluminium face.

Sampler Instructions

Set the sampler in the "closed" position by rotating the handle clockwise to the right. Insert the sampler to the desired depth in the powder/granulate. Rotate the handle anti-clockwise to the left (now the powder/granulate will flow into the sampler). Now turn the handle clockwise to the right, thus sealing the sample chambers. Remove the sampler carefully, turn it around and empty the sample through the open inner tube into a suitable container.

Warning!

Please note that when turning the handle, no parts of the body (finger, palm of the hand) are in the opening. The opening has sharp corners that can cause injury.

Maintenance and Care

The samplers need no maintenance, but after sampling, they should be cleaned as follows:

1. Pull out inner tube completely. Unscrew tip.
2. Wipe both tubes inside and out. With the cleaning brush (accessory), the sampling tube can be cleaned thoroughly.

The samplers can be sterilised with steam in the autoclave at temperatures of up to 121 °C.

The samplers should not be lubricated, because this would contaminate the samples.

Accessories/spare parts

Cleaning brush

for all-layer-samplers
for Jumbo samplers

5304-0100
5304-5100

Protective clothing

Safety goggles
Protective gloves
Plastic apron

2502-1001
2507-0001
2507-1001

Self-adhesive seals for closing punctured sacks

close-it control seal, printed, black
close-it control seal, printed, red
close-it control seal, printed, yellow
close-it control seal, printed, green
close-it control seal, unprinted, white

5303-0011
5303-0012
5303-0014
5303-0016
5303-0002

Sample container, sterilisable, made of PP

Screw-cap tubs 100 ml
Screw-cap tubs 60 ml
Screw-cap tubs 180 ml

7103-0100
6203-0060
6203-0180

More accessories at www.buerkle.de