# Temperature

## Controlled Shaking

Leading Safety Standards

Superior Ease of Use

Reduced Cost of Ownership

### **Operating Manual**

**Page 2**

Overhead Shaker Hei-MIX Reax 20  
Operating manual must be read before initial start-up.  
Please follow the safety instructions provided.  
Please keep for future reference.

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<b>About this Document</b>	<b>4</b>
Version and variants .....	4
About this manual .....	4
Symbols and signal words.....	5
<b>Safety Instructions</b>	<b>6</b>
General safety instructions.....	6
EU Declaration of Conformity.....	6
Intended use.....	6
Installation / electrical safety .....	7
Qualifications of personnel .....	7
Obligations of the operator .....	7
Installation site .....	7
Modifications to the device.....	8
Safety of personnel.....	8
Safety during use .....	9
Disposal .....	9
<b>Device Description</b>	<b>10</b>
Device overview .....	10
Overall view.....	10
Control panel .....	10
<b>Start-Up</b>	<b>11</b>
Setting up device .....	11
Emergency button/stop .....	11
Switch unit on and off .....	11
Connect/disconnect power cord .....	12
<b>Operation</b>	<b>13</b>
Installing bottles .....	13
Starting rotation.....	14

<b>Cleaning and Maintenance</b>	<b>16</b>
Cleaning.....	16
Maintenance .....	16
<b>Troubleshooting</b>	<b>17</b>
Possible errors and how to resolve .....	17
<b>Assembly</b>	<b>18</b>
Electrical connections .....	18
Attachment for smaller bottles .....	19
<b>Disassembly and Storage</b>	<b>20</b>
Dismantling, transportation and storage .....	20
Dismantling .....	20
Transportation and storage .....	20
<b>Accessories and Spare Parts</b>	<b>21</b>
Scope of delivery.....	21
Accessories.....	21
<b>Attachments</b>	<b>22</b>
Technical data .....	22
<b>Service</b>	<b>23</b>
Contact / Technical Service .....	23
Warranty .....	24
Confirmation of condition.....	25

# About this Document

## Copyright

This operating manual is protected by copyright and is exclusively intended for the internal use of the purchaser.

Handing over this operating manual to third parties, reproducing it in any type or form –in part or in full – and utilizing and/or communicating its contents is prohibited without prior written approval from the manufacturer.

Any violations will result in an obligation to provide compensation for damages. Other claims remain unaffected.

## ➤ Version and variants

### Version

This manual describes the function, operation and maintenance of the Hei-MIX Reax 20 overhead shaker.

Version	Alteration date
1.2	05/2018

### Variants

The device is available in different variants. Certain features or functions are only available in specific product variants. These variants are described in this operating manual.

## ➤ About this manual

This operating manual is an integral part of the device described here.

- ➔ Carefully read through this operating manual and observe all of the safety instructions and warnings notices.
- ➔ Ensure that every user has carefully read the operating manual before using the device for the first time.
- ➔ Store this operating manual so that it is accessible to all users at all times.
- ➔ Pass the operating manual on to any subsequent owner.

The current version of this operating manual can also be found at:  
[https://heidolph-instruments.com/en/service/downloads/operation-manuals#Shakers & Mixers \(Reax 20\)](https://heidolph-instruments.com/en/service/downloads/operation-manuals#Shakers & Mixers (Reax 20))

In addition, also observe the local regulations at the installation site.

## ➤ Symbols and signal words

Standardized symbols and signal words are used in this operating manual to warn against any dangers and provide important instructions. These instructions must be strictly observed to avoid accidents and damage.

The following terms and basic symbols are used:

Symbol	Additional signal words / explanation
<b>Warning signs</b> 	The yellow triangle indicates hazardous situations. It is used in combination with the following signal words: <b>DANGER:</b> Indicates a hazardous situation which, if not avoided, will result in serious injury or death. <b>WARNING:</b> Indicates a hazardous situation which, if not avoided, may result in serious injury or death. <b>CAUTION:</b> Indicates a hazardous situation which, if not avoided, may result in property damage and minor or moderate injury.
<b>Prohibitory signs</b> 	Prohibited: The red circle indicates a situation that should be avoided under all circumstances and which, if not avoided, may result in serious injury or death.
<b>Mandatory signs</b> 	Must be observed: The blue circle indicates important information. Please observe this information to avoid any property damage.

### Other symbols used:

Symbol	Description
→	Handling instruction, action required
✓	Result of action
▪	List of information
▪	
▪	
a.	List of variants
b.	
c.	

## ➤ General safety instructions

### EU Declaration of Conformity

This device complies with the following EC-Directives\*:

- 2006/42/EC Machinery Directive
- 2014/30/EU Electromagnetic Compatibility Directive

\*Also see the attachment "EU Declaration of Conformity".

The device has been manufactured according to state-of-the-art technology and in compliance with recognized safety regulations. However, risks may still arise during installation, operation and maintenance.

→ Please ensure that the operating manual is available at all times.

#### **The device may only be used under the following conditions:**

- Only operate the device if it is in a technically perfect condition.
- Only operate the device if you are properly aware of the risks and required safety measures.
- Only operate the device in accordance with the instructions given in this operating manual.
- If any information is ambiguous or missing, ask your superior or contact the manufacturer.
- Do not operate anything on the device without authorization.
- Only use the device in accordance with its intended use.

### Intended use

The device may only be operated by authorized personnel.

The device is suitable for the following uses:

- Shaking
- Mixing
- Emulsifying
- Suspending
- Separating

The device may be operated in research laboratories, other laboratories and production facilities in the following sectors:

- Chemistry
- Pharmacy
- Biology
- Environmental analytics
- Basic research
- Similar research laboratories

#### **Reasonable foreseeable misuse**

Any use which deviates from the device's intended use is considered to be improper. The manufacturer is not liable for damage that occurs as a result. The risk is borne by the operator alone.

### Installation / electrical safety

- The device may only be connected up if the available power supply voltage corresponds to the information stated on the rating plate for the device.
- The power supply connection must be easily accessible at all times.
- An RCD circuit breaker (residual current circuit breaker) must be fitted to the electrical system in the building so that the device is disconnected from the power circuit in the event of a fault.
- Repairs may only be performed by electricians authorised by Heidolph Instruments.
- Never operate the device with a damaged power cord.
- Always turn the device OFF and disconnect the power cord before carrying out any maintenance or repair work.

### Qualifications of personnel

- The device may only be operated by trained personnel.
- The device may only be operated by persons who have been instructed and supervised in its proper use by trained specialist personnel.
- The device may only be operated by specialist personnel who are above the legal minimum age.
- Other persons may only work on the device under the constant supervision of experienced and trained specialist personnel.
- This operating manual must be read and understood by all persons working with the device.
- Personnel must receive safety training that ensures responsible and safe working practices.

### ➤ Obligations of the operator

#### Installation site

- The device must be installed in a suitable location.
- The device must be installed on a firm, stable surface.
- Ensure that the device and all of its components are easily accessible at all times.
- Maintain a sufficient safety clearance to, in particular, any moving and/or hot equipment components.
- It is not permitted to place or store any objects such as accessories, tools or chemicals within this clearance area during operation.
- All screw connections must be securely tightened.
- It is not permitted to operate the device near to highly flammable or explosive substances.
- The device may only be operated inside buildings that have been equipped with appropriate laboratory equipment (e.g. a fume cupboard).
- Only operate the device (depending on the substances being used) in conjunction with an extractor hood (see DIN EN 14175 and DIN 12924).

## Safety Instructions

- It is only permitted to operate the device in dry and enclosed rooms and under the following ambient conditions:

Ambient temperature	5 - 31 °C at 80 % relative humidity
	32 - 40 °C decreasing linearly to a maximum 50 % relative humidity
Installation altitude	0 - 2,000 m above sea level
Contamination level	2
Overvoltage category	II
Permissible supply deviations	± 10 %

- If the device is operated in corrosive atmospheres, the service life of the device will decrease based on the concentration, duration and frequency of the exposure to the corrosive atmosphere e.g. concentrated hydrochloric acid (HCl).

## Modifications to the device

- It is not permitted to make any unauthorised modifications or changes to the device.
- Do not attach or install any parts that have not been approved by the manufacturer.
- Unauthorized modifications or changes will void the EC Declaration of Conformity for the device and operation of the device will no longer be permitted.
- The manufacturer is not liable for any damage, dangers or injuries that result from unauthorized modifications and changes or due to the non-observance of the instructions in this manual.

## Safety of personnel

- Ensure that the device is only operated by qualified specialist personnel and trained employees.
- Observe the following instructions to avoid any personal injuries and property damage:
  - Laboratory regulations
  - Accident prevention regulations
  - Hazardous Substance Act
  - Other generally accepted occupational health and safety regulations
  - Local regulations

### ➤ Safety during use

- Wear the appropriate clothing when working on the device (protective clothing, protective glasses and, if necessary, safety gloves).
- Do not use the device in potentially explosive areas. The device is not protected against explosion. There is no explosion or ATEX protection available.
- Do not carry out work with naked flames in the vicinity of the device (risk of explosion).
- Do not operate or assemble devices in the vicinity which are emission or radiation sources (electromagnetic waves) for the frequency range ( $3 \cdot 10^{11}$  Hz to  $3 \cdot 10^{15}$  Hz).
- Do not operate or assemble appliances in the vicinity of the device which constitute emission or radiation sources for ionizing radiation or in the ultrasonic range.
- Do not operate the device where adiabatic compression or shock waves might occur (shock wave combustion).
- Do not use substances, where the energy input due to mixing might pose a hazard.
- Do not spill liquids over the device or any parts of it.
- Remove any accidentally spilled liquids immediately.
- Eliminate errors immediately.
- Always switch the device OFF after use.
- Do not use abrasive material to clean the surface. Only wipe with damp cloth.

### ➤ Disposal

- Check the device components for hazardous substances and solvents.
- Clean all components before disposal.
- Dispose of the device in accordance with the relevant national regulations.
- Dispose of the packaging material in accordance with the appropriate national regulations.

## ➤ Device overview

### Overall view

### Control panel

### ➤ Setting up device

**Warning:**  
**Risk of slipping device!**

As a result of vibrations generated during operation the device might slide off the table top.

The surface must be smooth and clean, should be properly leveled and has to feature an anti-skid surface coat.

- Locate the shaker on a stable, horizontal surface.
- Clean surface and feet with a damp cloth and ethanol regularly.

### ➤ Emergency button/stop

The emergency stop is located on the front side of the unit above the ON/OFF switch.

- ✓ The unit is switched on.
- Push the emergency stop.
- ✓ The ON/OFF switch jumps to OFF and is blocked.
- ✓ The rotation stops.
- If necessary resolve the failure.
- Turn the red cap of the emergency stop 30° clockwise to the right.
- ✓ The emergency stop returns to the initial position.
- ✓ The ON/OFF switch is enabled.
- ✓ Start the unit via the ON/OFF switch.

Emergency button

### ➤ Switch unit on and off

The ON/OFF switch is situated on the front panel.

**Switch unit on**

- Press the ON/OFF switch on the bottom side.
- ✓ The unit is switched on.

**Switch unit off**

- Press the ON/OFF switch on the upper side.
- ✓ The unit is switched off.

ON/OFF switch: device switched OFF

### ➤ Connect/disconnect power cord

The three-wire socket is firmly connected to the unit.

The connection is at the rear side of the unit

- ✓ The ON/OFF switch is in the "Off" position.
- Connect the power cord with a grounded socket.
- Disconnect the plug if the unit is not in use.

## ➤ Installing bottles

The standard version of the unit is designed for 2000 ml bottles. There are 2 attachments available for smaller bottles:

- Attachment for 4 x 1.0 l bottles
- Attachment for 4 x 0.5 l bottles

Set the rotor in the proper position in order to install the bottles.

- Set rotation speed to 1 rpm.
- Switch unit on and rotate until proper position is reached to install bottles.

Analog knob to set speed

- Release clamping lever by turning it from back to front.

Clamping lever released

- Pull knurled knob upwards to raise tension plate.

Knurled knob pulled upwards

- Insert the bottle and take care that the tension plate centers exactly above the screw cap of the bottle.
- Push knurled knob down in the direction of the bottle.
- Tighten clamping lever by turning it from front to back.
- Repeat the procedure according to the number of bottles.
- Check the stability of the bottles.
- Re-tighten them, if necessary.

A single flask should be arranged in the middle of the shaker plate. Several flasks should be distributed equally on the plate in front as well as in the back.

When using shakers inside conditioning cabinets, make reference to ambient conditions as stipulated in the "Technical data" section.

## Operation

Rotation angle of the clamping lever is limited to 180°. In case this is not enough to tighten the bottle holder, it is possible to reposition the lever using a ratchet mechanism.

Pull the lever in the direction shown in diagram (inwards, towards the center) and hold, pull the lever in the anti-clockwise direction, release the inward pressure on the lever, and then tighten the lever in the clockwise direction. Always load the REAX 20 in a symmetrical way so as to distribute the weight equally.

- Pull clamping lever in direction of arrow and hold.
- Turn clamping lever anti-clockwise.
- Release clamping lever.
- The clamping lever locks into the new position.
- Tighten clamping lever by turning it from front to back.

## ▶ Starting rotation

### **Warning:** **Risk of injury!**

Returning power after a power cut off will restart shaking mode. There is a risk of injury in the vicinity of moving parts of device.

Before working with the device or in the vicinity ensure all controls are switched to "0" or "OFF" mode, when device is not moving.

### **Caution:** **Risk of breakage and spillage**

If the vessels are not attached securely they may fall out and spill the sample.

Prior to switching the unit on ensure that all vessels are securely fastened.

### **Warning:** **Risk of poisoning**

Open vessels will result in samples splashing.

Close all vessels tightly especially if sample is dangerous or toxic.

Adjust rotation speed step by step until you have reached the required rpm settings.

Wear safety glasses and especially with dangerous and toxic samples appropriate protective clothing.

A power cut will stop the unit.

With power return the device is switched on automatically. Shaking must be started manually.

**Warning:  
Risk of injury!**

High speeds in combination with high loads and/or a tall assembly may lead to an unstable device. As result the device may start moving and slide off the table top.

When working with with high loads and/or a tall assembly, increase the speed with care and watch for the devices stability. Ensure the device is installed in such a way as it can not collide with the surrounding environment or slip off its work surface.

A single flask should be arranged in the middle of the shaker plate. Several flasks should be distributed equally on the plate.

Rotation of the unit is steplessly adjustable in the range from 1 - 16 rpm. An electronic overspeed monitor provides for constant rotation.

Rotation starts as soon as unit is switched on.

- Turn unit on with the ON/OFF switch.
- Set rotation speed with analog knob.
- To remove bottles wait for proper rotation position.
- Switch unit OFF and remove bottles from front side.
- Switch unit ON and let rotate into position for second side.
- Switch unit OFF and remove bottles from opposite side.

### ➤ Cleaning

All surfaces of the unit allow for cleaning with a damp cloth and if required a mild soap lotion.

**Warning:**  
**Risk of electric shock!**

If the device is still connected to power during cleaning, you might get in contact with live parts due to accidentally entering liquids.

Prior to cleaning:

- ensure that the device is switched OFF
- ensure that the plug is pulled from the electric socket.

**Caution:**  
**Damage to the device**

Accidentally entering liquids will cause the device to fail .  
Only wipe with a damp cloth. Never use a soaked cloth.

**Attention:**  
**Risk of damage to the surfaces**

All of the surfaces may be damaged by improper cleaning.

Never use:

- Chlorine bleach or any cleaning agent containing chlorine
- Solvent-bearing agents (e.g. acetone)
- Ammonia
- Abrasive cleaning agents such as scouring agents, cleaning wool or agents with metallic particles

As a result of vibrations generated during operation the device might slide off the table top.

To ensure a stable position clean surface and feet with a damp cloth and ethanol regularly.

### ➤ Maintenance

The unit is maintenance-free. Any necessary repair must be performed by an authorized Heidolph distributor.

Please contact Heidolph Instruments or your local Heidolph Instruments distributor.

## ➤ Possible errors and how to resolve

<b>Malfunctions</b>	<b>Possible reason</b>	<b>Troubleshooting</b>
ON/OFF switch does not enlighten	No power	<ul style="list-style-type: none"> <li>▪ Check size and shape of plug and compatibility to the mains</li> <li>▪ Check circuit breakers</li> <li>▪ Check device fuse (see below)</li> </ul>
	Light diode defective	Contact your local Heidolph Instruments distributor
Unit stops rotating	Thermal motor circuit breaker triggered by motor overload	Wait 20 minutes, decrease overall load
	<ul style="list-style-type: none"> <li>▪ Mechanical parts broken</li> <li>▪ Electronic failure</li> </ul>	Contact your local Heidolph Instruments distributor

If you experience a malfunction which can not be resolved, please contact your authorized Heidolph Instruments distributor immediately.

## Electrical connections

- Fuses must **only** be replaced by a **professional electrician**.
- **Repairs** to the device must **only be carried out by a qualified electrician approved by Heidolph Instruments**. Improper repairs could result in serious dangers. Please contact your specialist dealer if the device needs to be repaired.

### Warning: Risk of electric shock

The device must be connected to an earthed power socket.

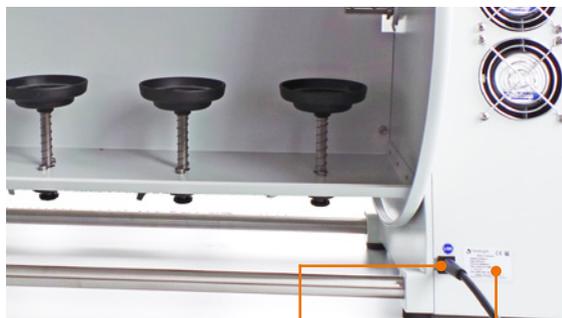
The rated voltage for the device and the supply voltage must be the same. The rated voltage for the device can be found on the rating plate on the rear of the device.

The device is earthed via the supplied power cord when delivered. If the original plug is replaced, it is essential that the new plug is connected to a protective earth!

If the device is still connected to power while changing the fuse, you might get in contact with live parts.

Prior to replacing the fuse ensure that the unit is switched off and pull the plug from the electric socket. Only use original spare fuses from Heidolph Instruments.

After changing a fuse check the device for safe operating condition according to IEC 61010-1.

Power cord

Rating plate

Device fuse

### Color code for electrical connections:

#### Europe

**GREEN/YEL-  
LOW**

PE: Protective conductor (**Earth**)

**BLUE**

N: Neutral conductor

**BROWN**

P: Phase

#### North America

**GREEN**

PE: Protective conductor (**Earth**)

**WHITE**

N: Neutral conductor

**BLACK**

P: Phase

The device is connected to the mains with the power cord supplied. This plugs into the socket on the rear of the housing.

For countries where a different plug than the standard plug supplied is used:

- The plug supplied may only be changed by a professional electrician.
- If you wish to use an adapter, ensure that it is grounded and approved in accordance with the local regulations.
- Wiring, adapter and plug must at least meet the electrical power stated on the rating plate.

## ➤ Attachment for smaller bottles

Components attachment for smaller bottles: Example: attachment 4 x 1 l bottles

- 4 extension pieces
- 1 allen key

4 x extension pieces

1 x allen key

- ➔ Tighten clamping lever by turning backwards.
- ➔ Detach tension plate by removing allen screw with allen key supplied.
- ➔ Remove tension plate.

Tension plate

Allen key

- ➔ Tighten extension piece with thread upwards.  
Attention: Do not clamp spring.

Thread

Extension piece

- ➔ Mount tension plate by tightening allen screw with allen key supplied.

Allen screw

- ➔ Secure smaller bottle.

# ➤ Dismantling, transportation and storage

## Dismantling

**Warning:**  
**Risk of injury!**

While the device is plugged in you might be at harm if the device is accidentally switched to ON mode.

Prior to dismantling device ensure that the device is switched off and the plug is pulled from the electric socket.

- ✓ The device is switched off.
- Disconnect plug from power socket.
- Disconnect plug from socket on the device.
- Remove all vessels.
- Disassemble attachments.

## Transportation and storage

- Store and transport the device and its components only if they were emptied, cleaned and dismantled.
- To carry the device wear safety gloves and lift the unit from below.
- Store and transport the device and parts of device in the original packing material or alternatively in an adequate container to prevent damage.
- Seal the package carefully against unauthorized or accidental opening.
- Store the device in a dry and frost-free place.

Improper storage and transportation may result in damages to the system and the mechanical components.

Avoid any kind of shocks during transportation.

## ➤ Scope of delivery

Component	Variants	Quantity	Product number
Hei-MIX	REAX 20/4	1	541-20004-00
	REAX 20/8	1	541-20008-00
	REAX 20/12	1	541-20012-00
Operating Manual English / German		1	01-005-002-18
Warranty registration / Confirmation of condition		1	01-006-002-78
EC Declaration of Conformity		1	01-001-025-06
Power cord		1	depending on country

## ➤ Accessories

Component	Quantity	Product number
Device fuses 230 V	1	14-002-015-46
Device fuses 115 V	1	14-002-015-49
Attachment for 4 x 1,0-l-bottles	1	549-26000-00
Attachment for 4 x 0,5-l-bottles	1	549-27000-00
Tension plate for caps Ø 94 mm	1	11-001-001-81

Additional accessories are shown in the general catalog or on our website at [www.heidolph-instruments.com](http://www.heidolph-instruments.com)

## ➤ Technical data

### Hei-MIX Reax 20

Standard supply voltage	230 V 50/60 Hz or 115 V 60 Hz		
Supply power (W)	280		
Device fuses	T 3,15 A (230 V) or T 6,25 A H (115 V)		
Protective class (IEC 61140)	1 ⊕		
Protection class (IEC 60529)	IP 21		
Sound pressure level (dB(A)) (based on IEC 61010)	< 45		
Drive	asynchronous motor		
Overload	overheat protection, self reset		
Speed control	electronic		
Motion	overhead rotation		
Speed range (rpm)	1 - 16		
Rotation speed setting	analog		
Load capacity (kg)	30		
Bottles max. Ø (mm)	136		
Bottles max. h (mm)	270		
<b>Model</b>	<b>Reax 20/4</b>	<b>Reax 20/8</b>	<b>Reax 20/12</b>
Dimensions (mm) (w x h x d)	490 x 520 x 465	770 x 520 x 465	1050 x 520 x 465
Weight (kg)	23	28	33

## ➤ Contact / Technical Service

### Questions / Repair work

If any aspect of installation, operation or maintenance remains unanswered in the present manual, please get in touch with the following address.

For repairs please call Heidolph Instruments or your local authorized Heidolph Instruments distributor.

---

#### **Warning: Danger of poisoning!**

Contaminated devices can lead to severe injury or death of our employees!

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise exact substance
  - take adequate protective action towards our parts receiving and service personnel
  - mark the pack in accordance with Ordinance on Hazardous Substances
- 

A "Confirmation of Condition" form can be found at the end of this operating manual.

→ Prior to shipping the device for repair, complete a copy of this form and submit it in advance.

### Contact details:

---

#### **Heidolph Instruments Germany**

Heidolph Instruments GmbH & Co. KG  
Technical Service  
Walpersdorfer Str. 12  
D-91126 Schwabach / Germany

Tel.: +49 - 9122 - 9920-74  
Fax: +49 - 9122 - 9920-84

E-Mail: [service@heidolph.de](mailto:service@heidolph.de)  
[www.heidolph.de](http://www.heidolph.de)

#### **Heidolph Instruments North America**

Phone: 1-866-650-9604  
E-mail: [service@heidolph.com](mailto:service@heidolph.com)  
[www.heidolphNA.com](http://www.heidolphNA.com)

#### **Heidolph Instruments United Kingdom**

Phone: 01799 - 5133-20  
E-mail: [service@radleys.co.uk](mailto:service@radleys.co.uk)  
[www.heidolph-instruments.co.uk](http://www.heidolph-instruments.co.uk)

#### **All other countries**

You will find contact details of your local Heidolph distributor at [www.heidolph.com](http://www.heidolph.com)

---

## ➤ Warranty

---

Heidolph Instruments provides a 3 year warranty for the products described here (excluding glass and wearing parts) if you register using the warranty card enclosed or online ([www.heidolph.com](http://www.heidolph.com)). The warranty period begins with the date of registration. The serial number is also valid without registration.

This warranty covers material and manufacturing defects.

---

In the event of a material or manufacturing defect, the device shall be repaired or replaced free of charge under the terms of the warranty.

Heidolph Instruments shall not assume liability for any damage incurred as a result of improper handling or transport.

Warranty claim?

- ➔ Please notify Heidolph Instruments or your local Heidolph distributor if you wish to make a warranty claim.

## ➤ Confirmation of condition

→ In the case of repair, copy and complete the Confirmation of condition form and send it to Heidolph Instruments.

### 1. Details about the device

Product number \_\_\_\_\_

Serial number \_\_\_\_\_

Reason for repair \_\_\_\_\_

\_\_\_\_\_

### 2. Has the device been cleaned, decontaminated/sterilized?

Yes \_\_\_\_\_ No \_\_\_\_\_

### 3. Is the device in a condition which does not represent any health threats for the staff of our service department?

Yes \_\_\_\_\_ No \_\_\_\_\_

If not, which substances has the device come into contact with?

\_\_\_\_\_

\_\_\_\_\_

### 4. Legally binding declaration

The customer is aware of being legally liable to Heidolph Instruments for any damages arising from incomplete and incorrect information.

Date \_\_\_\_\_

Signature \_\_\_\_\_

Company stamp \_\_\_\_\_

Please note

The shipper is responsible for the return of the goods in well packed condition, suitable for the mode of transport.

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This is a translation of the original operating manual. We reserve the right to make any technical modifications without prior notification. This documentation is not subject to revision; the latest version of this documentation can be found on our homepage. The copyright for all texts and images is held by Heidolph Instruments GmbH & Co. KG.
© Heidolph Instruments GmbH & Co. KG