KCL-Informationsbroschüre

Kat. II

Art. 470+

D	Informationsbroschüre	471+
GB	Information brochure	472+
F	Brochure d'information	623+
N	Informasjonsbrosjyre	624+
DK	Informationsbrochure	625+
FIN	Tiedotuslehtisen	
H	Információs füzet	
S	Informationsbroschyr	
E	Folleto Informativo	
I	Opuscolo Informativo	
NL	Informatie brochure	
SK	Informačná brožúra	
HR	Informacije brošura	
EST	Infobrošüür	
CZ	Informační brožura	
P	Brochura de informaçaõ	
BG	Информационна брошура	
GR	Ευήμερωτικό φυλλάδιο	
LT	Informacinę brošiūra	
PL	Broszura informacyjna	
LV	Direktivos	
RO	Broşuri de Informare	
SLO	Informacijska brošura	
TR	Direktif	
RUS	Информационная брошюра	

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**Information brochure in section 1.4 of Annex II of Directive 89/686/EEC
Personal Protective gloves Cat. II, Article 470+, 471+, 472+, 623+, 624+, 625+**

The CE marking on this glove means that it meets the essential requirements for in the European Directive EEC 89/686 concerning Individual Protection Equipment (IPE): Harmlessness-Comfort-Dexterity-Sturdiness.
Protective Gloves conform to the EN 420 . Approved to EN 388 according to

Article 10 of Directive 89/686/EEC:
TÜV Rheinland LGA Products GmbH, Tillystrasse 2, 90431 Nürnberg,
Identification number: 0197

Article	Name	Size	EN 388
470+ / 471+	CovaSpec®	6, 7, 8, 9, 10	113X
472+	CovaSpec®	6, 7, 8, 9, 10, 11	3131
623+	Camapur® Comfort Antistatik 623+	7, 8, 9, 10	X13X
624+	Camapur® Comfort Antistatik 624+	5, 6, 7, 8, 9, 10	X13X
625+	Camapur® Comfort Antistatik 625+	6, 7, 8, 9, 10	3131

These gloves are suitable for activities where protection from mechanical injury is required (eg assembly, dirty work).

EN 388, mechanic risks

1st digit Abrasion resistance (min. 0; max. 4)

3rd digit Tear strength resistance (min. 0; max. 4)

X = Cannot be tested

2nd digit Cut resistance

4th digit Puncture resistance

(min. 0; max. 5)

(min. 0; max. 4)

Antistatic properties according to prEN 16350 tested by KCL laboratory

Anti-static properties of the above items according to prEN 16350 tested by the laboratory KCL. Products meet the strict requirements of the normative type A: Vertical resistance RV von < 1,0 x 10⁸ Ω.

Testing atmosphere: air temperature (23 ± 2) °C, relative humidity (25 ± 4) %

More data about the electrostatic properties can be requested at KCL. There is no existing European Standard that currently can be used for testing of electrostatic Properties (also ESD) of gloves.

There is no existing European Standard that currently can be used for testing of electrostatic Properties (also ESD) of gloves.

Warning!

These protective gloves provide no protection against extreme cold (<-5 °C), heat (>50 °C), electricity, Chemical.

Do not use in the vicinity of moving machine parts. Risk of being drawn into the machine.

The specified protective index may be affected by factors such as temperature, abrasion, etc.

472+, 625+:- The specified level of protection is only present on the coated surfaces.

Unusable when the gloves are cracked, porous and stiff.

Safety glove not approved for food.

The person wearing the electrostatic dissipative protective gloves shall be properly earthed. The resistance between the person on the earth shall be less than 10⁸ Ω e.g. by wearing adequate footwear.

Electrostatic dissipative protective gloves shall not be adjusted or removed whilst in flammable or explosive atmospheres or while handling flammable or explosive substances.

The electrostatic properties of the protective gloves might be adversely affected by ageing, wear, contamination and damage.

The gloves lose by washing their anti static properties.

Allergy Notice: Article contain natural rubber latex which may cause allergic reactions including anaphylactic reactions. Protective gloves contain traces of Dimethyl formamide in legally permitted ppm concentrations.

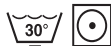
Use: The protective gloves you wear must be of the correct size.

Note that using undergloves may result in some usage restrictions. Check the protective gloves for damage before you use them. Damaged protective gloves must not be used under any circumstances.

Expiry date: With proper storage, no reduction in performance levels within 36 months.

Cleaning: Delicate wash cycle up to 30 °C with greatly reduced mechanical action. Can be dried on care programme. The certified index values are not guaranteed after washing/drying. Do not use any chemicals or sharp implements when cleaning (wire brushes, sandpaper, etc.).

Storage/Transport: Flat, dry, dark, with no additional weight load in its original packaging, at a temperature of 5 °C - 25 °C. Protect from sunlight and ozone.

Disposal: Gloves when used properly disposed of with household.

For further information please contact.

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KCL GmbH