Bürkle GmbH

Rheinauen 5 | D - 79415 Bad Bellingen
Telefon +49 (0) 76 35 8 27 95-0
Telefax +49 (0) 76 35 8 27 95-31
info@buerkle.de | www.buerkle.de

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Do not use any force when operating the instrument.

WARNING!

Damage to the respiratory tract and other organs due to escaping or outgassing media.

- ▶ When sampling hazardous substances, follow the safety data sheet of the medium to be sampled.
- ▶ Wear the appropriate protective clothing.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

Safety instructions during and after use

- ▶ The device can be used in the temperature range of 0 - 60 °C.
- ▶ Pay attention to have sufficient space above during the application (minimum height: height of the maximum extended piston + 20 cm).
- ▶ Do not perform any lever movements during the application. The instrument may be deformed.

Layout and function

- ▶ The instrument is made of PP (item no. 5331-1xx0) or PTFE/FEP (item no. 5331-2xx0).
- ▶ The instrument is suitable to take samples of viscous substances such as pastes, mud, creams and oils.
- ▶ The instrument is only suitable to take samples and not for continuous operation.
- ▶ The instrument is not suitable for the storage of samples.
- ▶ The instrument is maintenance-free and suitable for steam sterilisation (up to 121°C).

Parts in contact with medium

- ▶ PP (item no. 5331-1xx0)
- ▶ PTFE/FEP (item no. 5331-2xx0)

Point out any damage in the packaging to the freight forwarder/supplier immediately.

Safe sampling

WARNING!

Bodily injury due to uncontrolled escaping liquid under pressure.

- ▶ Wear the appropriate protective clothing.

Point sampling/soil sample

1. Slide the wiping ring ① to the desired sampling depth.
2. Slide the piston rod ② with the aid of the handle ③ into the instrument up to the stop.
3. Submerge the instrument into the medium to be sampled up to the desired depth / to the ground.
4. Pull the handle. Thereby, the medium is sucked in by the piston.
5. Remove the instrument from the medium to be sampled.
6. Empty the sample into a suitable collecting container by sliding in the piston rod.

All-layer sample

1. Slide the wiping ring ① to the desired sampling depth.
2. Slide the piston rod ② with the aid of the handle ③ into the instrument up to the stop.
3. Pull the handle while submerging the instrument into the medium to be sampled up to the desired depth / to the ground. Thereby, the medium is sucked in by the piston.

4. Remove the instrument from the medium to be sampled.
5. Empty the sample into a suitable collecting container by sliding in the piston rod.

After taking the sample

1. Slide the wiping ring with the aid of the supplied pusher down to wipe any remaining liquids.
2. Screw off the discharge funnel ④.
3. Screw off the guide nut ⑤.
4. Pull the handle and pull out the piston rod completely.
5. For cleaning purposes, the hexagon screw of the suction piston can be opened by using the supplied key.
6. Clean all parts. Use the cleaning brushes (see accessories) for the inside of tubes.

Troubleshooting

Problem	To eliminate the error
The instrument doesn't fit through the container opening.	Screw off the discharge funnel and take the sample without it. Screw the discharge funnel on again before emptying the sample.
Poor suction power or suction piston is leaky.	Increase the contact pressure of the suction piston. Screw off the discharge funnel and adjust the o-rings clockwise with a hexagon key (the o-rings must seal tightly and should have a low friction).

Accessories

Type	Specification	Number	Item No.
Cleaning brush	PVC/stainless steel, 100 cm	1	5304-0100
Drain safety stand	Total length 117 cm, Ø 35 mm	1	5618-1000
Sample bottle HDPE	250 ml	1	0342-0250
Sample bottle HDPE	750 ml	1	0342-0750

More accessories can be found at www.buerkle.de

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