***Instruction Manual***  
***Manuale di istruzioni***  
***Manuel d'instructions***  
***Manual de instrucciones***  
***Bedienungsanleitung***

## **MULTISTIRRER 6 & 15 DIGITAL Magnetic Stirrers**

F203A0179, F203A0180

### **General Information / Informazioni Generali / Informations Générales / Información General / Allgemeine Hinweise**

Before using the unit, please read the following instruction manual carefully.

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Do not dispose of this equipment as urban waste, in accordance with EEC directive 2002/96/CE.

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#### **This unit must be used for laboratory applications only.**

The manufacturer declines all responsibility for any use of the unit that does not comply with these instructions. If the product is used in a not specified way by the manufacturer or with non specified accessories, product's safety may be compromised.

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**This unit has been designed and manufactured in compliance with the following standards:**

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**IEC/EN 61010-1  
IEC/EN 61010-2-051**

Electrical equipment for laboratory use

**UL 61010-1**

General requirement - Canadian electrical code

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Do not use with explosive and dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres.

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It is responsibility of the user appropriately decontaminate the instrument in case of dangerous substances fall on or in it.

It is also responsibility of the user to use safety substances for cleaning or decontaminating, which do not react with internal parts of the instrument or with the material contained in it. In case of doubts on the compatibility of a cleaning solution, contact the manufacturer or local distributor.

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In case of black-out, once the power is back the device will restart automatically.

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Do not use with explosive and dangerous materials for which the equipment is not designed. The stirrer must not be used in explosive atmospheres, in bain-marie and to stir combustible liquids that have a low combustion temperature. The minimum fire point of flammable solution is 200 °C. Only small amounts (< 50 ml) of flammable liquid can be used with the device.

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## 1. Introduction

MULTISTIRRER DIGITAL is able to stir up to 6 beakers having a maximum diameter of 85 mm (MULTISTIRRER DIGITAL 6) or up to 15 beakers having a maximum diameter of 64 mm (MULTISTIRRER DIGITAL 15) at the same time. The stirrer stays cold even after several days of continuous operation. This feature makes it particularly appreciated in microbiology and biochemistry. Epoxy painted metal structure, with stainless steel cover, are studied and tested in order to give to the instrument an excellent resistance to the attack of chemical agents and to corrosion in general. The stirring speed can be set within a range of 80 - 1500 rpm using the knob on the front panel. The Stirrer display shows when the speed set.

**NOTE:** it's important to choose the most suitable magnetic stirring bar related to the quantity and the liquid to be stirred as well as to the type of the beaker. For MULTISTIRRER DIGITAL 6 is suggested the stirring bar A00001056 (6x35 mm); for MULTISTIRRER DIGITAL 15 the A00001057 (6x20 mm).

## 2. Assembly and installation

Upon receipt and after having removed the packaging, please check the integrity of the instrument. The box includes:

- Magnetic stirrer MULTISTIRRER DIGITAL
- Switching VELP 100-240V/12V without plug
- Plug for power supply
- Instruction manual

### 2.1 Electrical connection

After having unpacked the instrument, place the unit on the laboratory bench.

Before connecting the instrument to the power supply, make sure that the values on the rating plate correspond to those of the power supply. Connect the unit to the power supply using the transformer supplied. Ensure that the socket and the relative cut-off device conform to current safety norms and easy to reach. When the instrument is connected, the display shows the software version.

### 2.2 Start-up

Place the flasks containing the sample and a suitable magnetic stirring bar on the stirring plate. Then, press the knob and set the speed by turning it. The display shows the set value of speed.

## 3. Operating controls

Turn the stirring speed knob on the front panel in order to start the stirring function. If the display is off, press the speed knob. It is possible to select speeds between 80 and 1500 rpm.

With the instrument on, pressing the knob, it is possible select the Autoreverse time (sec.). With this option it is possible invert the rotation direction automatically every set seconds (from 5 till 900 sec.). Select - - - in order to remove it.

Press again the knob, in order to select the Timer mode (min). With this option it is possible turn off automatically the instrument after the minutes set. Select - - - in order to remove it.

Keep the knob pressed in order to turn off the unit.

## 4. Maintenance

No routine or extraordinary maintenance is necessary apart from periodically cleaning the unit as described in this manual. In compliance with the product guarantee law, repairs to our units must be carried out in our factory, unless previously agreed otherwise with local distributors. The instrument must be transported in its original packaging and any indications present on the original packaging must be followed (e.g. palletized).

### 4.1 Cleaning

Disconnect the unit from the power supply and use a cloth dampened with an non-inflammable non-aggressive detergent.

## 5. Technical data

Power	MULTISTIRRER DIGITAL 6: 4,1 W    MULTISTIRRER DIGITAL 15: 9 W
Admitted power supply	100-240 V / 50-60 Hz
Dimensions (WxHxD)	230x51.5x370 mm
Weight (with plug)	MULTISTIRRER DIGITAL 6: 1,75 Kg MULTISTIRRER DIGITAL 15: 2,1 Kg
Speed range	80 ÷ 1500 rpm
Stirring capacity	MULTISTIRRER DIGITAL 6: 6x400 ml MULTISTIRRER DIGITAL 15: 15x250 ml
Operating mode	Continuous
Construction material	Epoxy painted metal structure and stainless steel
Environmental ambient temperature	5 - 40 °C (41 - 104°F)
Storage temperature range	-10... + 60 °C (14 - 140°F)
Pollution degree	2
Overvoltage category	II
Sound level	35 dBa
Max humidity	80%
Altitude	Up to 2000 m.

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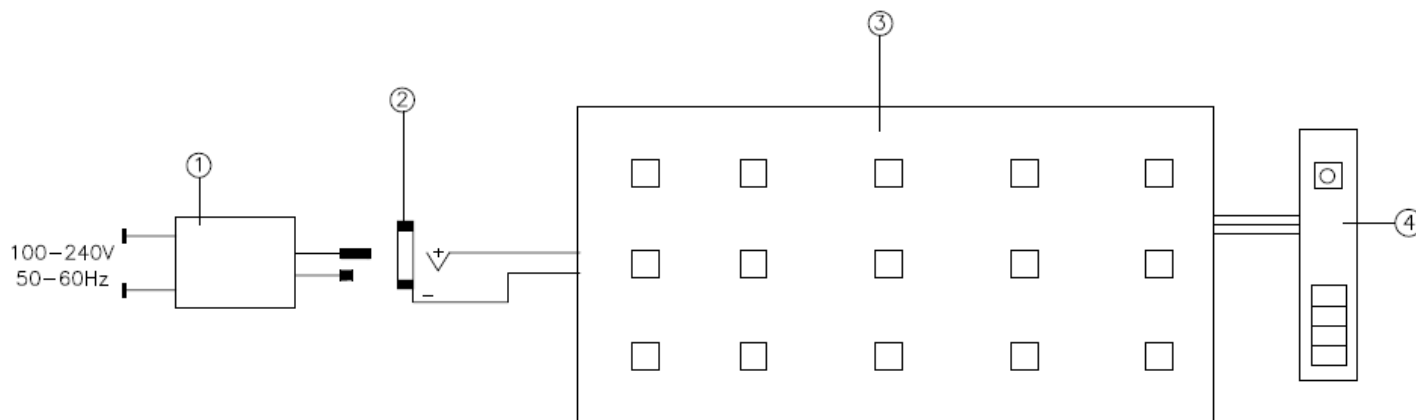
We, the manufacturer VELP Scientifica, under our responsibility declare that the product is manufactured in conformity with the following standards:

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## Thank you for having chosen VELP!

Since 1983 VELP has offered to professionals in the sector a range of sophisticated and reliable equipment.

VELP works according to **ISO 9001**, **ISO14001** and **OHSAS 18001** Quality System Certification.

Instruments are built according to the International norms IEC 1010-1 and to the rules of the CE mark.

Our product lines:

### Food&Feed Line

DK and DKL Digesters  
JP Recirculating water pump for fumes aspiration  
SMS Scrubber  
UDK Distillation units  
SER Solvent extractors  
FIWE Raw fiber extractors  
Dietary fiber extractors  
Oxitest Oxidation test reactor  
NDA 701 Elemental analyzer  
Consumables

### Environment Line

ECO Thermoreactors  
BOD. Determination systems  
Refrigerated thermostats and incubators  
Flocculators  
Overhead mixer  
Mineralization unit for trace heavy metals determination  
Turbidimeter  
Radiation detector

### Stirring Line

Heating magnetic stirrers  
Vertex digital thermoregulator  
Ultraflat magnetic stirrer  
Magnetic stirrers  
Overhead stirrers  
Heating plates  
Vortex mixers  
Homogenizer

### Other Lab Solution

Open circulating bath  
Recirculating water vacuum pump  
Peristaltic pump

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