# Kaiser slimlite plano


LED-Leuchtplatte

LED Light Box

Table lumineuse LED





Mesa de luz LED

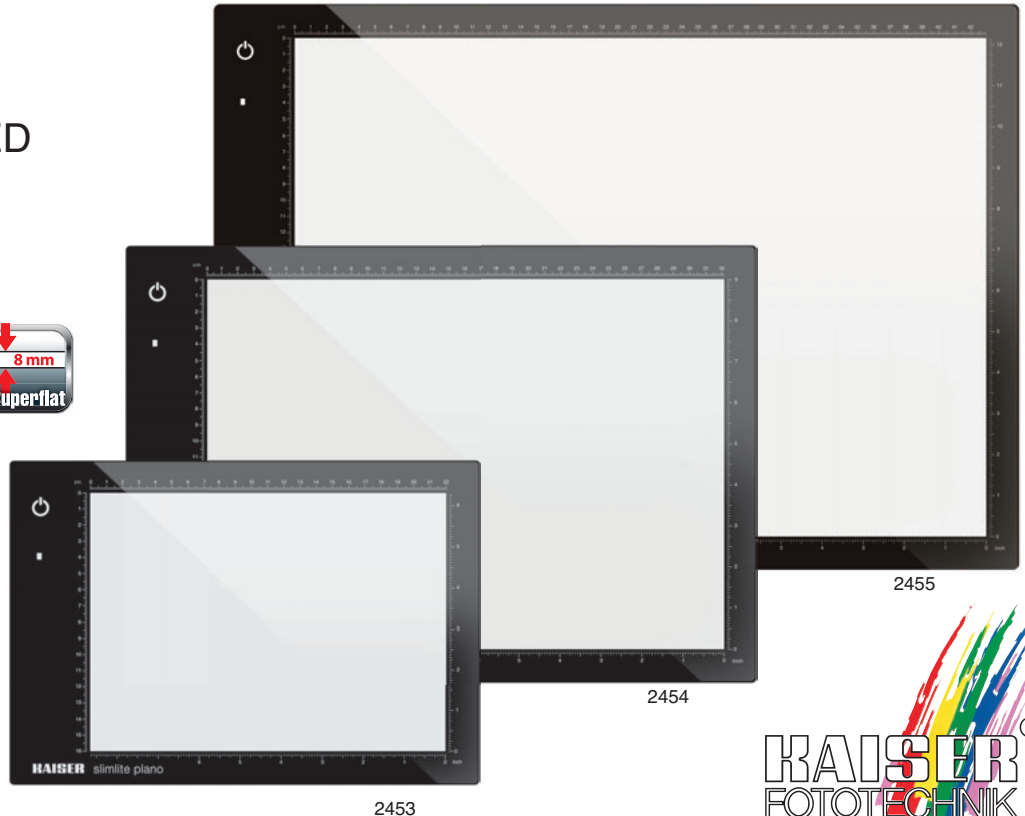Visore LED

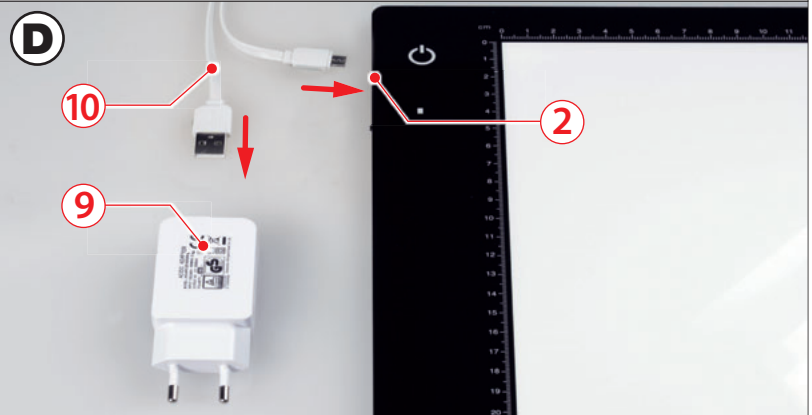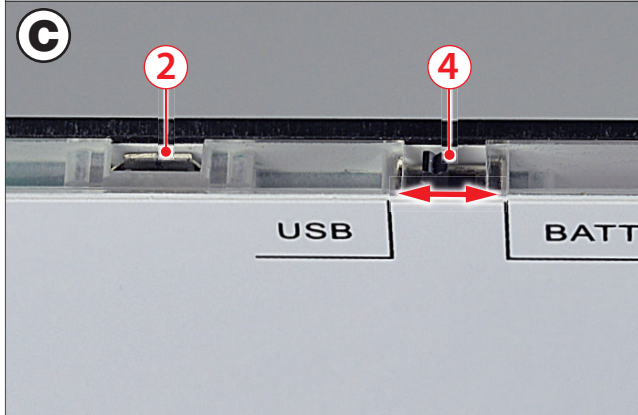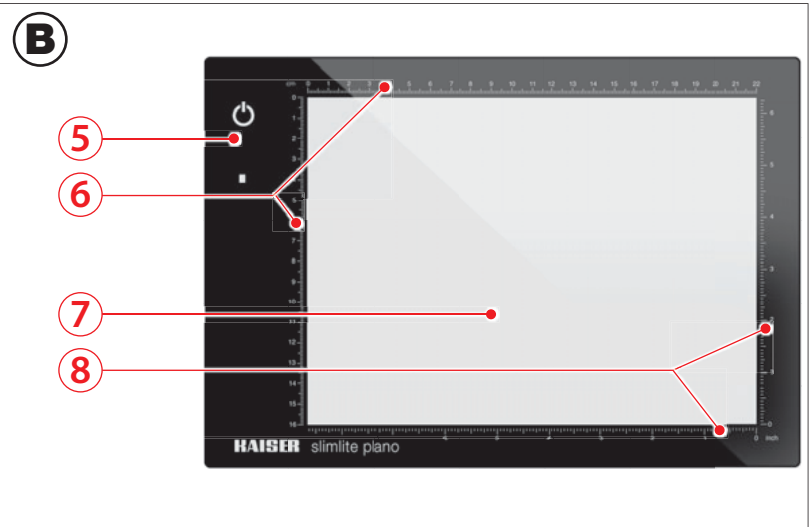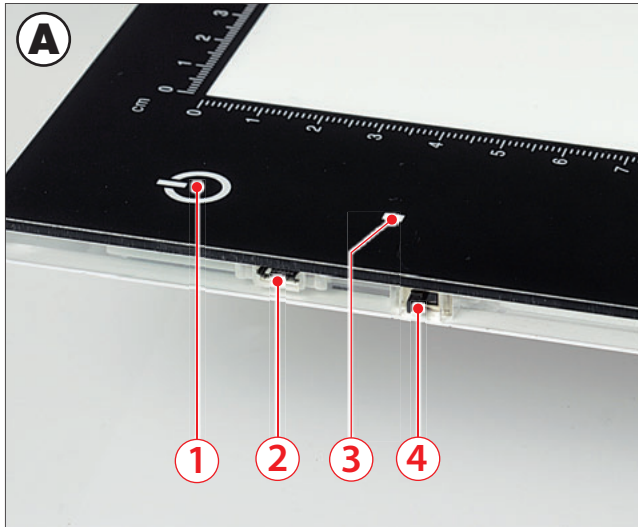Bedienungsanleitung  
Operating instructions  
Mode d'emploi  
Instrucciones  
Istruzioni

# DEUTSCH

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# ENGLISH

- ① ON/OFF and brightness control touch switch
- ② Micro-USB socket
- ③ Standby (red) and charging indicator (orange)
- ④ Mode switch USB/battery operation
- ⑤ Light box
- ⑥ cm scales
- ⑦ Illuminated area
- ⑧ inch scales
- ⑨ Mains/charging adapter
- ⑩ USB cable

Prior to using your light box, please read these instructions carefully. Keep them safe to have them at hand at any time. If you pass the light box on to somebody else, always include these instructions.

## 1. Intended use

The "slimlite plano" LED light box is intended to be used primarily for transillumination of transparent originals (like slides or foils for example). It is not intended for household lighting.

Use this light box only as described in these instructions. Any other use has to be considered as not as intended and can result in material damage or physical injury. This light box is not a toy.

Kaiser takes no responsibility for damages caused by not intended or incorrect use.

## 2. Safety instructions

Please observe the following safety instructions:

## WARNING!

Disconnect the devices (light box, mains/charging adapter, USB cable) from mains before cleaning and maintenance.

Keep the products away from children.

Do not put the devices into operation, if housing, plugs or cables are damaged.

Do not disassemble or modify the products. Repairs must only be carried out by qualified persons.

## CAUTION!

Only use the light box with parts designed for it.

Use the devices only in dry rooms.

Keep the devices away from water and other liquids and from combustible materials.

To avoid deep discharge of the battery, the light box must not be stored in an uncharged condition. If you do not use the light box for a longer period of time, the battery has to be recharged every 6 months.

Do not use or store the devices at places with extremely high air humidity or at extremely hot places. Keep the devices away from direct solar radiation.

Prior to boxing the light box, switch it off completely (see section 4 below) and let all devices cool down.

Do not place metal parts or subjects with sharp edges on the surface of the light box to avoid scratches.

## 3. Check devices and scope of supply

When unpacking the devices for the first time, please check if the scope of supply is complete. It comprises

- Light box ⑤
- Mains/charging adapter ⑨
- USB cable ⑩

Check also, if the devices are damaged by extraneous causes. If this is the case, do not take the devices into operation.

#### 4. Initial operation

The light box ⑤ can be operated with the combined mains/charger adapter ⑨ or with the built-in Li-ion battery.

Push the mode switch ④ to the "USB" position for mains operation or to "BATTERY" for battery operation (fig. C). In the latter case the standby indicator ③ might light up red.

Place the light box ⑤ on a flat and dry surface.

Now connect the mains/charging adapter ⑨ to the light box, no matter if you have selected mains or battery operation, because the battery is to be charged prior to the initial operation. Please proceed as follows:

Check that the mains voltage corresponds to the operating voltage given on the mains/charger adapter ⑨.

Connect the mains/charging adapter ⑨ to the light box ⑤ using the USB cable ⑩ (fig. D). Pay attention to the orientation of the respective plug, before you insert it into the USB and micro-USB socket. Misalignment may lead to damage of the plug or the socket. Use the USB cable only when uncoiled.

Now connect the mains/charging adapter ⑨ to the mains.

If mains operation ("USB") is selected, the standby indicator ③ will light up red.

If battery operation ("BATTERY") is selected, the standby indicator ③ will light up orange and the battery is charged. When charging is completed, the standby indicator ③ will light up red and the USB cable ⑩ can be disconnected from the light box ⑤. Depending on the state of charge of the battery, charging may last up to 6 hours. Nevertheless you may use the light box during battery charging. However it is recommended not to operate the light box during charging for longer battery life.

*Note: The light box ⑤ can also be operated or charged via the included USB cable ⑩ from any other live USB terminal with 5 V output voltage (e.g. a computer, notebook or a power bank).*

*The light can be switched on and changed in brightness by using the touch switch ①.*

Switching on: Keep your finger on the touch switch ① until the desired or maximum brightness is reached.

Switching off: A short touch of the switch ① will cause the brightness to go down until the light is switched off.

Adjusting brightness: If the light is switched on, keep your finger on the touch switch ① until the desired brightness is reached. The next time you touch the switch ① the brightness will change into the opposite direction.

If you switch off the light later by a short touch of the switch ①, another short touch will switch on the light with the previously selected brightness.

When operated from the mains ("USB"), the light box is suitable for unlimited continuous operation.

If there is a considerable decline of brightness with battery operation ("BATTERY"), the battery has to be recharged (see above).

Complete switch-off: If you do not use the light box for a certain time (temporarily), move the mode switch ④ to the "USB" position and disconnect the light box from the mains/charging adapter ⑨ if applicable. The standby indicator ③ goes off and a previously saved brightness level will be deleted. Now the light box is completely switched off.

#### 5. Maintenance and care

To avoid deep discharge of the battery, the light box must not be stored in an uncharged condition. If you do not use the light box for

a longer period of time, the battery has to be recharged every 6 months.

Make sure that the illuminated area ⑦ is free of stains and dirt. If necessary remove these before operation. Clean the devices with a soft dry cloth. On no account use petrol or other solvents.

## 6. Specifications

slimlite plano	2453	2454	2455
Illuminated area	22 x 16 cm (8.7 x 6.3 in.)	32 x 22.8 cm (12.6 x 9 in.)	42,9 x 30.9 cm (16.9 x 12.2 in.)
Light source	23 LEDs	33 LEDs	90 LEDs
Color temperature	5000 K	5000 K	5000 K
Color rendition index	CRI = 95	CRI = 95	CRI = 95
Light density	750 cd/m <sup>2</sup>	600 cd/m <sup>2</sup>	650 cd/m <sup>2</sup>
Power consumption	3.6 W	4.1 W	6.6 W
Operating voltage	5 V	5 V	5 V
Li-ion battery	1050 mAh	1050 mAh	1050 mAh
Operating time*	5 hours	3.5 hours	1.5 hours
Charging time	5-6 hours	5-6 hours	5-6 hours
Operating temperature	-5 to +40 °C	-5 to +40 °C	-5 to +40 °C
Outside dimensions (cm/in.)	29 x 20 x 0.8 (11.4 x 7.9 x 0.3)	39 x 27 x 0.8 (15.4 x 10.6 x 0.3)	49.9 x 35 x 0.8 (19.6 x 13.8 x 0.3)
Weight	570g / 20.1oz.	1010g / 35.6oz.	1690g / 59.5oz.
Length of USB cable	2 m (6.6 ft.)	2 m (6.6 ft.)	2 m (6.6 ft.)

\*at full power, longer when dimmed

Mains/charging adapter: AC 100-240 V, 50/60 Hz, DC 5 V / 2 A

Right reserved to make technical modifications.  
All specifications are approximate.

## Disposal Instructions

Dispose unserviceable products in accordance of the relevant statutory regulations.

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