# pH 3310

EN English .....	2	KO 한국어 .....	24
DE Deutsch .....	4	NL Nederlands .....	26
CS Český .....	6	NO Norsk .....	28
DA Dansk .....	8	PL Polski .....	30
EL Ελληνικά .....	10	PT Português .....	32
ES Español .....	12	RO Română .....	34
FI Suomi .....	14	RU Русский .....	36
FR Français .....	16	SK Slovenský .....	38
HU Magyar .....	18	SV Svenska .....	40
IT Italiano .....	20	TR Türkçe .....	42
JA 日本語 .....	22	ZH 中文 .....	44

## pH 3310

### Overview


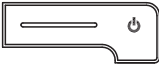
The pH 3310 meter enables you to perform measurements (pH) quickly and reliably.

The pH 3310 meter provides a maximum degree of operating comfort, reliability and measuring certainty.

### Keypad

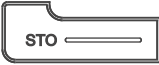Keys are indicated by brackets <..>.

The key symbol (e.g. <ENTER>) generally indicates a short keystroke (under 2 sec). A long keystroke (approx. 2 sec) is indicated by the underscore behind the key symbol (e.g. <ENTER\_\_>).


	<F1>: <F1__>:	Situation dependent function (is shown on the display) e.g.: <F1>/[Menu]: Opens the menu for measurement settings <F1__>/[Menu]: Opens the menu for system settings
	<F2>: <F2__>:	
	<On/Off>:	Switches the meter on or off
	<M>:	Selects the measured parameter
	<CAL>: <CAL__>:	Calls up the calibration procedure Displays the calibration data
	<STO>: <STO__> :	Saves a measured value manually Opens the menu for the automatic save function
	<RCL>: <RCL__>:	Displays the manually stored measured values Displays the automatically stored measured values
	<▲><▼>: :	Menu control, navigation
		
	<ENTER>: <ENTER__>:	Opens the menu for measurement settings / confirms entries Opens the menu for system settings
	<AR>:	Freezes the measured value (HOLD function) Switches the AutoRead measurement on or off

## Display

The diagram shows the display screen of the pH 3310 meter. It features a large digital readout of '6.093' with a 'pH' label to its left. Above the readout is a scale from 0 to 14. Below the readout, the temperature is shown as '25.0 °C'. At the bottom left, there are softkeys labeled 'HOLD' and 'AR'. Below these are 'Menu', a date/time display '15.11.2008 08:37', and a 'USB output' indicator. A vertical bar on the right side of the display is labeled with numbers 1 through 8, corresponding to the legend table.

1	Status information
2	Measured value (with unit)
3	Measured parameter
4	Continuous measurement control (CMC function)
5	Sensor symbol (calibration evaluation)
6	Measured temperature (with unit)
7	Further status information
8	Softkeys and date/time
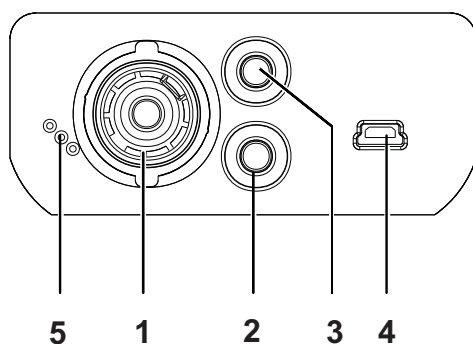
## Jack field /connections

1	pH electrode
2	Reference electrode
3	Temperature sensor
4	USB B interface (device)
5	Service interface

## Initial commissioning

- 1 Insert the supplied batteries. Make sure that the poles are in the correct position.
- 2 Switch on the meter with the <On/Off> key.

### Caution

Follow the safety instructions of the used sensors. The operating manuals of the sensors are available on the supplied CD and on the Internet under [www.WTW.com](http://www.WTW.com).

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
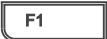
## Visuel

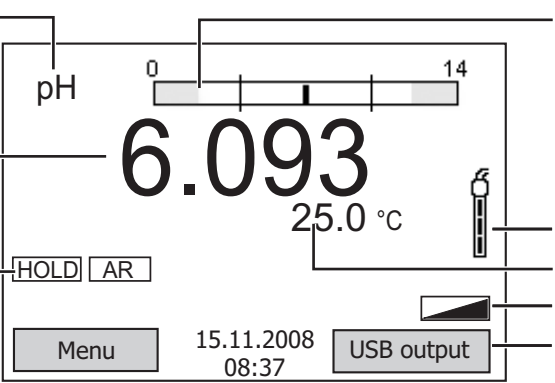The diagram shows the display screen of the pH 3310 meter. It features a large digital readout of '6.093' with a pH symbol to its left. Above the readout is a scale from 0 to 14. Below the readout, the temperature is shown as '25.0 °C'. At the bottom left, there are 'HOLD' and 'AR' softkeys. At the bottom center, the date and time '15.11.2008 08:37' are displayed. At the bottom right, there is a 'Menu' softkey and a 'USB output' indicator. A vertical bar on the right side of the screen indicates the calibration status. Numbered callouts 1 through 8 point to these various elements.

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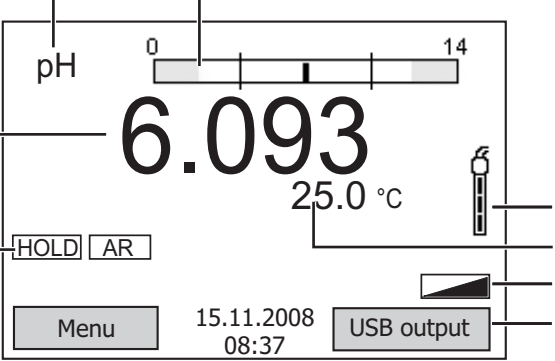
## Visor

The diagram shows the display screen of the pH 3310 meter. It features a large digital readout (DRO) showing '6.093' with a 'pH' label to its left. Above the DRO is a horizontal scale from 0 to 14. Below the DRO, the temperature is displayed as '25.0 °C' with a thermometer icon to its right. At the bottom left, there are 'HOLD' and 'AR' softkeys. At the bottom center, the date and time are shown as '15.11.2008 08:37'. At the bottom right, there is a 'Menu' softkey and a 'USB output' indicator. A legend on the right side of the diagram maps numbers 1 through 8 to specific elements on the screen.

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## **Wissenschaftlich-Technische Werkstätten GmbH**