pH 3310

EN English	2	KO 한국어	24
DE Deutsch	4	NL Nederlands	26
CS Český	6	NO Norsk	28
DA Dansk	8	PL Polski	30
EL Ελληνικά	10	PT Português	32
ES Español	12	RO Română	34
FI Suomi	14	RU Русский	36
FR Français	16	SK Slovenský	38
HU Magyar	18	SV Svenska	40
IT Italiano	20	TR Türkçe	42
JA 日本語	22	ZH 中文	44

pH 3310

Overview

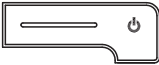
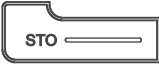
The pH 3310 meter enables you to perform measurements (pH) quickly and reliably.

The pH 3310 meter provides a maximum degree of operating comfort, reliability and measuring certainty.

Keypad

Keys are indicated by brackets <..>.

The key symbol (e.g. <ENTER>) generally indicates a short keystroke (under 2 sec). A long keystroke (approx. 2 sec) is indicated by the underscore behind the key symbol (e.g. <ENTER__>).

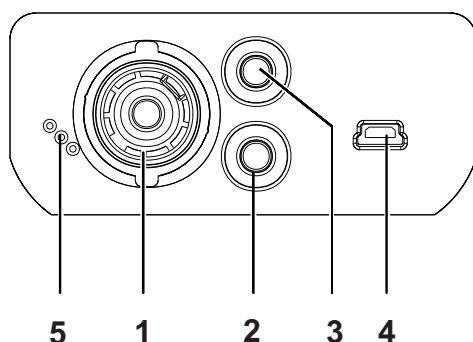
	<F1>: <F1__>:	Situation dependent function (is shown on the display) e.g.: <F1>/[Menu]: Opens the menu for measurement settings <F1__>/[Menu]: Opens the menu for system settings
	<F2>: <F2__>:	
	<On/Off>:	Switches the meter on or off
	<M>:	Selects the measured parameter
	<CAL>: <CAL__>:	Calls up the calibration procedure Displays the calibration data
	<STO>: <STO__> :	Saves a measured value manually Opens the menu for the automatic save function
	<RCL>: <RCL__>:	Displays the manually stored measured values Displays the automatically stored measured values
	<▲><▼>: :	Menu control, navigation
		
	<ENTER>: <ENTER__>:	Opens the menu for measurement settings / confirms entries Opens the menu for system settings
	<AR>:	Freezes the measured value (HOLD function) Switches the AutoRead measurement on or off

Display

The diagram shows the display screen of the pH 3310 meter. It features a large digital readout of '6.093' with a pH scale bar above it ranging from 0 to 14. Below the main reading, the temperature is shown as '25.0 °C'. At the bottom, there is a date and time '15.11.2008 08:37' and a 'USB output' indicator. On the left side, there are 'HOLD' and 'AR' softkeys. A 'Menu' button is located at the bottom left. A vertical bar on the right side of the display indicates the sensor's status.

1	Status information
2	Measured value (with unit)
3	Measured parameter
4	Continuous measurement control (CMC function)
5	Sensor symbol (calibration evaluation)
6	Measured temperature (with unit)
7	Further status information
8	Softkeys and date/time

Jack field /connections

1	pH electrode
2	Reference electrode
3	Temperature sensor
4	USB B interface (device)
5	Service interface

Initial commissioning

- 1 Insert the supplied batteries. Make sure that the poles are in the correct position.
- 2 Switch on the meter with the <On/Off> key.

Caution

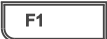
Follow the safety instructions of the used sensors. The operating manuals of the sensors are available on the supplied CD and on the Internet under www.WTW.com.

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Wissenschaftlich-Technische Werkstätten GmbH